Eutech EcoTestr pH 1 pocket tester

Instruction Manual

EcoTestr pH1
Revision Date: 13-11-2015

Language Index	
English	Page 2
Deutsch	Page 3
Français	Page 4
Nederlands	Page 5
Portuguese	Page 6
Español	Page 7

EcoTestr pH 1

Note: DO NOT BE ALARMED if white crystals form around the cap, this is normal.

CONDITIONING before first use. Fill the cap with water, close the cap and soak the sensor for 1 hour. This will activate the sensor and dissolve the crystals.

Note: Keep the tester upright during soaking, as the cap is not leak proof.

CALIBRATION: Select pH 7 buffer for general testing or pH 4 or 10 buffer when measuring acidic/basic samples • Press on/off button to power on. Immerse sensor in chosen buffer about 20 mm deep and stir gently. Wait for displayed value to stabilize at or near the pH buffer chosen • Next, press **cal** button to enter calibration sequence. The display will flash continuously • To cancel calibration at this point, press **cal** button to escape • To complete calibration, press **hold/ent** button to confirm. You have successfully calibrated the instrument.

Note: If no button is pressed for 15 seconds, tester will automatically confirm calibration.

pH TESTING: Remove cap and press on/off button to turn on the tester • Dip the sensor about 20 mm into the test solution • Stir once and let the display stabilize. Note the pH value • Press **hold/ent** button if you wish to hold the reading. Press the same button again to release the held reading • Press the on/off button to shut off • Tester automatically shuts off after 8.5 minutes of non-use to conserve batteries!

MAINTENANCE IS EASY: Rinse the sensor in tap water after use. In harsh samples, take readings quickly, rinse sensor immediately afterwards with de-ionized water to remove residues and prevent electrode contamination. To maximize sensor life, fill the cap with 20 mm clean tap water (NOT DE-IONIZED WATER), and replace cap. This will keep the sensor conditioned and be ready for use.


CHANGING BATTERIES: Lift up front battery cover and two sides of the pocket clip to remove battery cover. Remove old batteries and replace with fresh ones noting polarity as shown in battery compartment. No need to re-calibrate after battery change since tester's non-volatile memory function retains calibration.

USEFUL NOTE: To avoid cross contamination, rinse between samples and buffer with de-ionized water. Calibration should be done regularly or daily to ensure good tester accuracy.

ERROR MESSAGES: **Lb** - Weak batteries. Need replacement. **Er** - Wrong or bad buffer value (out of range) or the sensor is failing. **Or / Ur** – Over and Under range signal or sensor is contaminated / damaged.

RESET OPTION: Reset option allows you to restore the calibration back to factory default settings • Turn off the tester • Press and hold **cal** button and then switch on using the on/off button. The display shows a blinking 'rS' (reset) • Press **hold/ent** button to confirm reset or press the **cal** button to cancel reset option.

WARRANTY: This tester is supplied with a warranty against manufacturing defects for a period of ONE year from date of purchase.

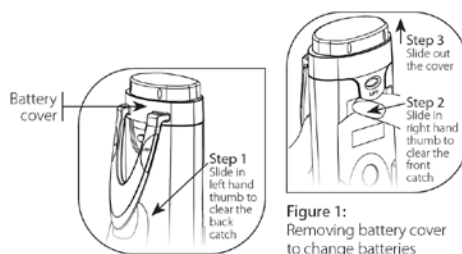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