# VWR® Timer

## INSTRUCTION MANUAL

**North American Catalog Numbers: 61161-340**

**European Catalogue Numbers: 609-0090**

Version: 1  
Issued: 2<sup>nd</sup> May 2017

## Legal Address of Manufacturer

### United States

VWR International  
Radnor Corporate Center  
Building One, Suite 200  
100 Matsonford Road, P.O. Box 6660  
Radnor, PA 19087  
800-932-5000  
<http://www.vwr.com>

### Europe

VWR International bvba  
Researchpark Haasrode 2020  
Geldenaaksebaan 464  
B-3001 Leuven  
+ 32 16 385011  
<http://be.vwr.com>

Country of origin - China

### Intended use

This item is intended for general laboratory use to monitor time duration.

## Product Specifications

|                              |                                                                           |
|------------------------------|---------------------------------------------------------------------------|
| <b>Display:</b>              | LCD, 4-Digit 1"                                                           |
| <b>Resolution:</b>           | 1 second                                                                  |
| <b>Accuracy:</b>             | ± 0.01%                                                                   |
| <b>Power:</b>                | AAA battery                                                               |
| <b>Size:</b>                 | 3¾ x 5½ X ½ inches                                                        |
| <b>Accessories Supplied:</b> | Wall-mount bracket, stand, battery, Traceable® Certificate, instructions. |

## Getting Started

Before using this timer, open the battery compartment and remove the battery tab.

## Operation

### Setting the Clock:

1. Insert battery in battery compartment located on the back side of unit
2. To set clock, slide the CLOCK/TIMER switch to CLOCK.
3. Select 12 hour or 24-hour format
4. Slide SET/LOCK switch located on the back of unit to SET. Press HR, MIN, and SEC keys until desired time is reached. Slide SET/LOCK switch to LOCK position.

### Counting up

1. Place the unit in the clock mode by sliding the CLOCK/TIMER switch to CLOCK.
2. Press the START/STOP key once to begin counting up. Press the START/STOP key a second time to pause the count (time-out). Press the START/STOP key again to resume counting. Any number of time-outs may be taken. Press CLEAR while the count is paused to reset the unit to 00:00:00.

3. The count-up is shown on the lower portion of the display. When 19:59:59 is reached, the timer will return to 00:00:00 and begin counting up again.

### **Counting Down**

1. Place the unit in the timer mode by sliding the CLOCK/TIMER switch to TIMER. Slide the switch to T1 to set channel 1 and to T2 to set channel 2. Slide SET/LOCK switch located on the back of unit to SET.
2. Press HR, MIN, and SEC keys until desired time is reached. Press MEMORY key once. "MEMORY" will appear on the display. Slide SET/LOCK switch to LOCK. Recall stored time period, press MEMORY key.
3. Press the START/STOP key once to start timing. Press the START/STOP key a second time to pause the timing (time-out). Press the START/STOP key again to resume timing. Any number of time-outs may be taken. Press CLEAR while the timing is paused to reset the unit to 00:00:00.
4. The upper portion of the display will show channel one. The lower portion of the display will show channel two. To display both channels simultaneously, slide the CLOCK/TIMER switch to the TIMER 1 & 2 position.
5. When 00:00:00 is reached, the timer will alarm and begin counting up. "TIME'S UP" will flash in the upper left portion of the display for the appropriate channel. To silence the alarm and stop the count-up, press the START/STOP key.

### **Low Battery**

Erratic readings, a faint display, or no display are all indicators that the battery is low and needs to be replaced. To replace the battery, remove the battery cover located on the back of the unit and insert a new AAA battery. Incorrectly installed batteries may damage electronics. Replace the Battery Cover.

### **All Operational Difficulties**

If this timer does not function properly for any reason, please properly replace the battery with a new 1.5V AAA alkaline battery (see Low Battery section). Low battery power can occasionally cause any number of "apparent" operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

### **Equipment disposal**

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste. Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment. For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

### **Compliance with local laws and regulations**

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

### **Technical service**

#### **Web Resources**

Visit the VWR website at [vwr.com](http://vwr.com) for:

- Complete technical service contact information
- Access to the VWR Online Catalogue, and information about accessories and related products
- Additional product information and special offers

**Contact us** For information or technical assistance contact your local VWR representative or visit [vwr.com](http://vwr.com).

### **Warranty**

**VWR** warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

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VWR International AB
VWR International AG
Teknolojileri Ltd.Şti.
VWR International Ltd