Instruction manual

VWR® Standard and Advanced Orbital Shakers

EU cat. no

3500 Standard Orbital Shakers: 444-2903

3500 Advanced Orbital Shakers: 444-2906

3750 Advanced Reciprocating Shakers: 444-2909

5000 Standard Orbital Shakers: 444-2912

5000 Advanced Orbital Shakers: 444-2915

10000-1 Advanced Orbital Shakers: 444-2918

10000-2 Advanced Orbital Shakers: 444-2921

15000-1 Advanced Orbital Shakers: 444-2924

15000-2 Advanced Orbital Shakers: 444-2927

EN	- English		1
FR	- Français		18
ES	- Español		36
DE	- Deutsch		54
IT	- Italiano		72
PT	- Português		90
BG	- Български		108
CZ	- Čeština		112
DK	- Dansk		116
EE	- Eesti		120
FI	- Suomi		124
GR	- Ελληνικά		128
HR	- Hrvatska		132
HU	- Magyar		136
IS	- Íslenska		140
LT	- Lietuvių		144
LV	- Latviešu		147
NL	- Nederlands		152
NO	- Norsk		156
PL	- Polski		160
RO	- Română		164
SI	- Slovenščina		168
SK	- Slovenčina		172
SV	- Svenska		176

Legal address of Manufacturer

Europe

VWR International bv,
Researchpark Haasrode 2020
Geldenaaksebaan 464 B-3001
Leuven
+3216385011
be.vwr.com

UK Importer:

VWR International Ltd
Hunter Boulevard, Magna Park
Lutterworth, Leicestershire, LE17 4XN
uk.vwr.com

Country of Origin

United States

Table of Contents

Safety Instructions	4
Package Contents	4
Installation	4
Intended Use	4
Maintenance & Service	5
Environmental Conditions	5
Product Specifications	
<i>Model 3500</i>	6
<i>Model 3750</i>	6
<i>Model 5000</i>	6
<i>Model 10000</i>	7
<i>Model 15000</i>	7
Operating Instructions	
<i>Standard Orbital Shakers</i>	7
<i>Advanced Orbital Shakers</i>	9
Control Panel - Advanced	11
Troubleshooting	11
Accessories	
<i>Platform Usage Chart</i>	12
<i>Test Tube Rack Platform Capacity</i>	13
<i>Flask Clamp Platform Capacity</i>	14
<i>Microplate Clamp</i>	15
Technical Service	15
Warranty	15
Disposal	16
Distributor Contact Information	17

Safety Instructions

Please read the entire instruction manual before operating the Orbital Shaker.

WARNING! DO NOT use the Orbital Shaker in a hazardous atmosphere or with hazardous materials for which the unit was not designed. Also, the user should be aware that the protection provided by the equipment may be impaired if used with accessories not provided or recommended by the manufacturer, or used in a manner not specified by the manufacturer.

Always operate unit on a level surface for best performance and maximum safety.

DO NOT lift the unit by the tray.

CAUTION! To avoid electrical shock, completely cut off power to the unit by disconnecting the power cord from the unit or unplug from the wall outlet. Disconnect unit from the power supply prior to maintenance and servicing.

Spills should be removed promptly. **DO NOT** immerse the unit for cleaning.

DO NOT operate the unit if it shows signs of electrical or mechanical damage.

Earth Ground - Protective Conductor Terminal

Alternating Current

Pinch Point - Keep fingers clear during operation

Package Contents

Orbital or Reciprocating Shaker

Description	Quantity
Non-skid rubber mat	1
234cm detachable power cords (EU and UK) CH adapter	3

Installation

Upon receiving the VWR Orbital or Reciprocating Shaker, check to ensure that no damage has occurred during shipment. It is important that any damage that occurred in transport is detected at the time of unpacking. If you do find such damage the carrier must be notified immediately. After unpacking, place the shaker on a level bench or table, away from explosive vapors. Ensure that the surface on which the unit is placed will withstand typical heat produced by the unit and place the unit a minimum of 15cm from vertical surfaces. Always place the unit on a sturdy work surface. The Orbital or Reciprocating Shaker is supplied with a power cord that is inserted into the IEC connector on the back of the unit first, then it can be plugged into a properly grounded outlet. The 230V unit plugs into a 230 volt, 50/60 Hz source. To prevent the rubber mat from sliding, remove backing from adhesive on the underside of mat, reposition in tray, and firmly press.

Intended use

These Shakers are intended for general laboratory use.

Maintenance & Servicing

The Orbital or Reciprocating Shaker is built for long, trouble-free, dependable service. No lubrication or other technical user maintenance is required. It needs no user maintenance beyond keeping the surfaces clean. However at least every three (3) months you should:

- Unplug the unit.
- Remove any accumulated dirt from the base and tray.
- Check all accessible items to make sure they are properly tightened.

The unit should be given the care normally required for any electrical appliance. Avoid wetting or unnecessary exposure to fumes. **DO NOT** use a cleaning agent or solvent on the front panel which is abrasive or harmful to plastics, nor one which is flammable. Always ensure the power is disconnected from the unit prior to any cleaning. If the unit ever requires service, contact your VWR representative.

Environmental Conditions

Standard Models

Operating Conditions: Indoor use only.

* For use in dry CO₂ environments, incubators, or cold rooms. Not for use in CO₂-incubators with high humidity.

Temperature: 0 to 40°C

Humidity: maximum 80% relative humidity, non-condensing

Altitude: 0 to 2000 M above sea level

Non-Operating Storage:

Temperature: -20 to 65°C

Humidity: maximum 80% relative humidity, non-condensing

Installation Category II and Pollution Degree 2 in accordance with IEC 664.

Advanced & Reciprocating Models

Operating Conditions: Indoor use only.

* For use in dry CO₂ environments, incubators, or cold rooms. Not for use in CO₂-incubators with high humidity.

Temperature: -10 to 60°C

Humidity: maximum 80% relative humidity, non-condensing

Altitude: 0 to 2000 M above sea level

Non-Operating Storage:

Temperature: -20 to 65°C

Humidity: maximum 80% relative humidity, non-condensing

Installation Category II and Pollution Degree 2 in accordance with IEC 664.

* **Avoid cold starts:** Unit is not designed to start after being in a cold room environment. Bring unit into cold room from a room temperature environment, operate and remove unit from cold room as soon as operation is complete.

Product Specifications

Model 3500

Overall dimensions (LxWxH): 41,3 x 35,6 x 14,6cm

Tray dimensions (LxW): 33 x 28cm

Electrical: 230 volts, 2,5 amps, 75 watts

Fuses: 5mm x 20mm, 5 amp quick acting

Speed Range:

(Advanced) 15 to 500 rpm

(Standard) 25 to 500 rpm

Speed Accuracy: (Advanced)

Above 100rpm $\pm 1\%$ of set speed

Below 100rpm ± 1 rpm

Stroke: 19mm

Capacity: ~16kg @ 75 rpm

~2,3kg @ 500 rpm

Timer:

(Advanced) 1 second to 160 hours

(Standard) 1 to 120 minutes

Controls:

(Advanced) See page 8

(Standard) See page 7

Ship Weight: 22,2kg

Model 3750

Overall dimensions (LxWxH): 41.3 x 35.6 x 14.6cm

Tray dimensions (LxW): 33 x 28 cm

Electrical: 230 volts, 2,5 amps, 40 watts

Fuses: 5mm x 20mm, 5 amp quick acting, 250V

Speed Range: 20 to 300 rpm

Speed Accuracy:

Above 100rpm $\pm 2\%$ of set speed

Below 100rpm ± 2 rpm

Stroke: 19mm

Capacity: ~6kg @ 300 rpm

Timer: 1 second to 160 hours

Controls: See page 8

Ship Weight: 22,2kg

Model 5000

Overall dimensions (LxWxH): 59,9 x 67,8 x 15,2cm

Tray dimensions (LxW): 45,7 x 61cm

Electrical: 230 volts, 2,5 amps, 75 watts

Fuses: 5mm x 20mm, 5 amp quick acting

Speed Range:

(Advanced) 20 to 500 rpm

(Standard) 25 to 500 rpm

Speed Accuracy: (Advanced)

Above 100rpm $\pm 1\%$ of set speed

Below 100rpm ± 1 rpm

Stroke: 25.4mm

Capacity: ~22,3kg @ 125 rpm

~4,5kg @ 500 rpm

Timer:

(Advanced) 1 second to 160 hours

(Standard) 1 to 120 minutes

Controls:

(Advanced) See page 8

(Standard) See page 7

Ship Weight: 49,5kg

Model 10000

Overall dimensions (LxWxH): 71,8 x 67,8 x 17,8cm

Tray dimensions (LxW): 61 x 61cm

Electrical: 230 volts, 2,5 amps, 80 watts

Fuses: 5mm x 20mm, 5 amp quick acting

Speed Range:

10000-1: 15 to 500 rpm

10000-2: 15 to 300 rpm

Speed Accuracy: (Advanced)

Above 100rpm $\pm 1\%$ of set speed

Below 100rpm ± 1 rpm

Stroke :

10000-1: 25mm

10000-2: 50mm

Capacity: ~45,3 @ 100 rpm

~18,1 @ 300 rpm

Timer: 1 second to 160 hours

Controls: See page 8

Ship Weight: 90,7kg

Model 15000

Overall dimensions (LxWxH): 71,8 x 67,8 x 17,8cm

Tray dimensions (LxW): 61 x 91,4cm

Electrical: 230 volts, 2,5 amps, 80 watts

Fuses: 5mm x 20mm, 5 amp quick acting

Speed Range:

15000-1: 15 to 500 rpm

15000-2: 25 to 500 rpm

Speed Accuracy: (Advanced)

Above 100rpm $\pm 1\%$ of set speed

Below 100rpm ± 1 rpm

Stroke :

15000-1: 25mm

15000-2: 50mm

Capacity: ~68,0kg @ 100 rpm

~18,1kg @ 300 rpm


Timer: 1 second to 160 hours

Controls: See page 8

Ship Weight 104,3kg

Operating Instructions

*Standard
Orbital Shaker
with Optional
platform and
flask clamps*

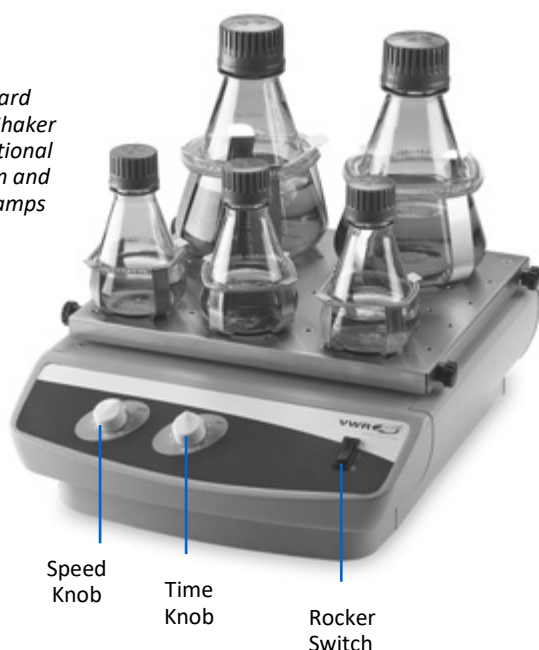Standard Orbital Shaker

The Standard Orbital Shaker is used for general laboratory shaking needs.

Models 3500 and 5000

1. Getting ready:

- The speed and time knobs have a built in on-off switch at their extreme counter-clockwise position. Turn both knobs to the off position.
- Make sure the rocker switch is in the standby position.
- Plug the cord into a properly grounded outlet.

2. Setting speed:

- To run in continuous mode, push rocker to the run position. Set speed knob to desired setting and adjust if necessary. Unit will run until you move the rocker switch to the standby position. The microprocessor speed control slowly ramps to set speed to avoid splashing.

3. Setting time:

- To run in timed mode, set timer and the speed knob to the desired settings. Press the rocker switch to the time position and release. The shaker will now run for the set time.
- To exit the timed mode turn the time knob to the extreme counter-clockwise position.

4. Turning unit off:

- To stop shaking function, turn the speed knob to the extreme counterclockwise position and push the rocker switch to the standby position. The shaker should be kept in the standby position when not in use. To completely cut off power to the unit, disconnect the power cord from the unit or unplug from the wall outlet.

OPERATING TIPS

Centering your sample and even weight distribution on the tray helps with balance and stability. As a safety feature, a built-in program will shut power off to the motor if the tray is prevented from rotating, or the unit is overloaded beyond its recommended weight capacity. To reset the unit, press the rocker switch to standby and then press the rocker switch to the on position. The shaker will automatically restart after a power interruption. Built-in memory maintains the last used speed and time settings during a power interruption. To prevent the rubber mat from sliding, remove backing from adhesive on the underside of mat, reposition in tray, and firmly press.

Advanced Orbital Shaker

1. Getting ready:

- Plug the cord into a properly grounded outlet. The standby indicator light will illuminate, verifying power to the shaker.
- Press the standby button to move the unit from standby mode. The standby indicator light will turn off and the speed and time displays will illuminate, displaying the previously used settings.

2. Setting speed:

- Press the up/down arrow below the speed display until you reach the desired speed. When you release the button, the display will blink off and then on indicating the new set speed has been accepted.
- Press the on/off button to start the shaking function. The indicator light below the speed display will illuminate and blink until the setpoint is reached. Once the set-point is reached the light will stop blinking and remain lit until shaking has ceased. The microprocessor controlled ramping feature slowly increases speed until the set-point is reached which helps to avoid splashing, and provides excellent low end control.
- Speed adjustments can be made without interrupting shaking by using the up/down arrows below the speed display. After the change has been made and you release the button, the display will blink off and then on indicating the new set speed has been accepted.
- To stop the shaking function press the on/off button below the speed display. The speed indicator light will turn off.

*Advanced Orbital Shaker
with optional platform*

3. Setting time to zero (0:00) and continuous mode: Accumulated time.

- Press and hold the on/off button below the time display. After three (3) seconds, the display will indicate the previous set time.
- Simultaneously press both the up and the down arrows, the display will indicate zero (0:00). The unit time is now set to zero (0:00) minutes. Alternately, you can use the up/down arrows to get to zero (0:00).
- Press the on/off button below the time display. The display will indicate accumulated time. The up/down arrows will become inactive. To stop timer, press the on/off button again. **IMPORTANT:** This will **NOT** interrupt the shaking function. Press the on/off button below the speed display to interrupt the shaking function.
- To reset, press and hold the on/off button below the time display. After three (3) seconds the display will indicate the previous set time, which was zero (0:00).

4. Setting timed mode: Programmed time.

- Press the up/down arrows below the time display until you reach the desired time.
- Start this function by pressing the on/off button below the time display. The unit will run for the selected time, the up/down arrows will become inactive while the timer is running. The unit will stop shaking when time display reaches zero (0:00). Four (4) audible beeps will indicate the count down function is complete. The time display will default back to the set time. To repeat for the same time, simply depress the on/off button again.
- To interrupt an automatic timing cycle before it is completed, press the on/off button below the time display. The display will flash off and on to indicate the time function is on "hold". **IMPORTANT:** This will **NOT** interrupt the shaking function. Press the on/off button below the speed display to interrupt the shaking function. Restart the timer by pressing the on/off button below the time display. Unit will continue counting down to zero (0:00). When the display reaches zero (0:00), you will hear the four (4) audible beeps that indicate the countdown function is complete and the shaking function will cease.

5. Turning unit off:

- To turn the unit off, press the standby button. The speed and time displays will be blank, the standby indicator light will illuminate. The Orbital Shaker should be kept in standby mode when not in use. To completely cut off power to the unit, disconnect the power cord from the unit or unplug from the wall outlet.

Load Sensing Function

The Advanced Orbital Shaker Models 3750 and up is equipped with a load sensing function that can be activated by the user. This function provides protection against improper positioning of load and maximum load being exceeded. When activated, the unit will automatically sense improper load conditions and slow to a safe running speed, then display that speed followed by E04 on the speed display. The unit will also beep three (3) times every 60 seconds until the error is reset by pressing the speed on/off button.

To activate the load sensing function use the following steps:

1. Place the unit in standby mode.
2. Press and hold the speed on/off button and press the standby button. The unit will beep two (2) times, confirming the function is enabled.
3. To restore normal operation, remove AC power to the unit for ten (10) seconds and then restore. If the E04 error occurs be sure the load is within the maximum specification and properly balanced (centered on tray) and/or reduce sample size/speed before restarting the unit. If the E04 occurs due to acceptable sample vibration or another vibration source, the vibration sensing function can be disabled as described above.

Additional Load Sensing Feature

The Advanced Orbital Shaker Models 3750 and up, are equipped with an additional overload protection feature that lets the user know when the unit's set speed is higher than the achievable speed of the unit. The unit speed display will show an E7. The unit will also beep three (3) times every 60 seconds until the error is reset by pressing the speed on/off button.

To disable this feature, perform the following steps:

1. Place the unit in standby mode by pressing the standby button.
2. Press and hold the speed up button and press on/off button. The unit will beep two (2) times, confirming the function is disabled.
3. To restore unit to normal operation, remove AC power to unit (unplug power cord from wall outlet) for ten (10) seconds and restart.

Beeper Preference

To silence beeper operation (except for error codes), with the unit in standby mode, press and hold the time on/off button and press the standby button. To restore normal beeper operation, remove AC power to unit for ten (10) seconds and then restore.

Calibration Procedure (Advanced Orbital Shaker Models 3500 and up)

This procedure is used to self-calibrate the Advanced Orbital Shaker. The tray should be free of any samples, vessels, and accessories prior to calibrating.


1. Turn unit on. Speed and time displays will be illuminated.
2. Press and hold the standby button and momentarily press the speed on/off button. The speed display should read "CAL".
3. The unit will run for approximately one (1) minute and automatically calibrate.

RS-232 Serial Port (Advanced Orbital Shaker Models 3500 and up)

RS-232 serial port provides two-way communications for data logging and unit control. If you need additional details, please contact your local VWR After Sales Service Department.

Operating Tips

Centering your sample and even weight distribution on the tray helps with balance and stability. As a safety feature, a built-in program will shut power off to the motor if the tray is prevented from rotating, or the unit is overloaded beyond its recommended weight capacity. The shaker will automatically restart after a power interruption. Built-in memory maintains the last used speed and time settings during a power interruption.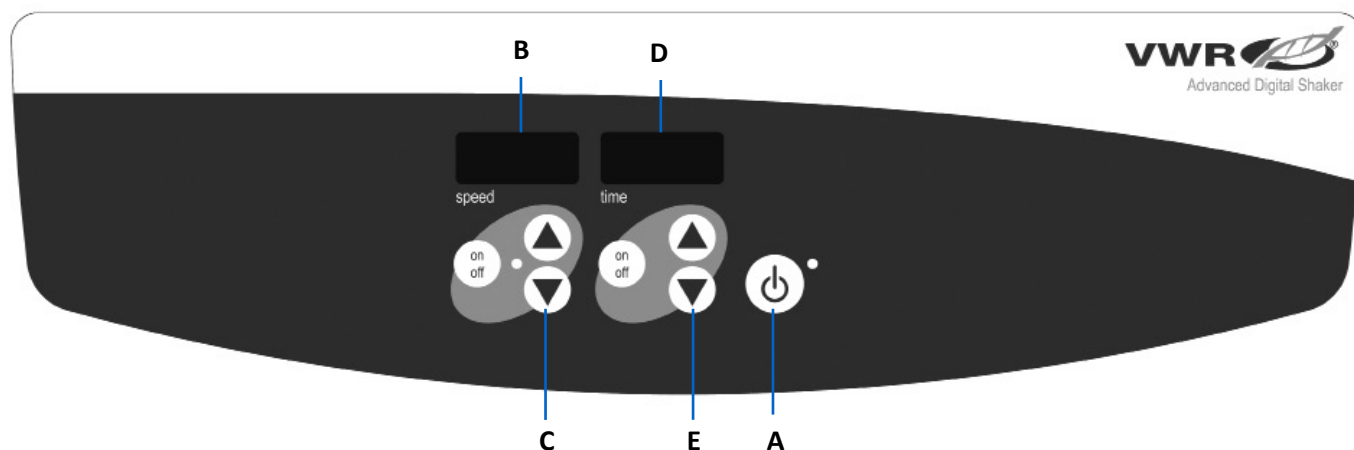

Control Panel - Advanced Orbital Shaker

The front panel of the Advanced Orbital Shaker contains all the controls and displays needed to operate the unit.

- A. Standby button/standby indicator light:** The standby indicator light will illuminate when the unit is plugged in. The unit will be in standby mode. Press the standby button to start the speed and time functions. The standby indicator light will shut off. Press the standby button again and the unit will once again be in standby mode.
- B. Speed display:** Displays the speed of the shaker.
- C. Up/down arrows:** For set-point control. On/off button starts/stop shaking function.
- D. Time display:** Displays accumulated time (continuous mode) or how much time is remaining (timed mode). The display range is from 0 to 9999 minutes in one (1) second increments. The display will indicate minutes and seconds until the timer reaches 99 minutes and 59 seconds (99:59), then the display will automatically display minutes up to 9999.
- E. Up/down arrows for setpoint control.** On/off button starts/stops the timer function.


Troubleshooting

Review the information in the table below to troubleshoot operating problems.

Problem	Cause	Solution
Unit will not run	Mechanical obstruction Motor obstruction	Add or replace fuse as necessary. If problem persists, please contact your VWR representative for repair.
Unit is excessively noisy	Sensor fan misaligned Motor misaligned	Ensure that tray is secured tightly. If problem persists, please contact your VWR representative for repair.
Unit not shaking at proper speed	-	Perform speed calibration test on page 10 If problem persists, please contact your VWR representative for repair.
E3	Mechanical obstruction Drive system failure Ceased bearing Drive belt broken	Remove mechanical obstruction. If problem persists, the reason may be the drive system and should not be addressed by the end user. Contact your VWR representative for repair.
E4	Improper positioning of load Maximum load exceeded	Ensure the load is evenly distributed and does not exceed the maximum load capacity for the unit. See "Load Sensing Function" on page 10. If problem persists, please contact your VWR representative for repair.
E7	Unable to reach set speed	Reduce the speed setting or weight. This function can be disabled by following the instructions listed under "Additional Load Sensing Feature" on page 10. (NOTE: This function is only available on models 3750 and up and is not available on Model 3500 shakers.)
E8	Electronics error	This error cannot be fixed by the end user. Please contact your VWR representative for repair.

Accessories

Platform Usage Chart

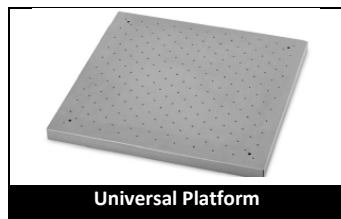Universal Platform

Rubber Mat

Dedicated Platform

Tier Braces with Universal Platform

Universal Platform Size	Description	Cat. No.	Compatible Shaker Models
28 x 33cm	Universal platform with holes for mounting accessories	444-7087	Model 3500, Model 3750
33 x 33cm	Universal platform with holes for mounting accessories	444-2930	Model 3500, Model 3750
46 x 46cm*	Universal platform with holes for mounting accessories	444-2931	Model 3500, Model 3750**
46 x 61cm*	Universal platform with holes for mounting accessories	444-2932	Model 3500, Model 3750**, Model 5000**
46 x 76cm*	Universal platform with holes for mounting accessories	444-2934	Model 5000
61 x 61cm	Universal platform with holes for mounting accessories	444-2933	Model 10000
61 x 91cm	Universal platform with holes for mounting accessories	444-2935	Model 15000
28 x 33cm	Rubber Replacement Mat	444-2966	Model 3500, Model 3750
46 x 61cm	Rubber Replacement Mat	444-2967	Model 5000
61 x 61cm	Rubber Replacement Mat	444-2968	Model 10000
61 x 91cm	Rubber Replacement Mat	444-2969	Model 15000

Dedicated Platform Size	Dedicated Platforms with PVC Flask Clamp Capacities	Cat. No.	Compatible Shaker Models
33 x 33cm (125mL)	16 x 125mL	444-2945	Model 3500, Model 3750
33 x 33cm (250mL)	12 x 250mL	444-2946	Model 3500, Model 3750
33 x 33cm (500mL)	8 x 500mL	444-2947	Model 3500, Model 3750
33 x 33cm (1L)	4 x 1L	444-2948	Model 3500, Model 3750
46 x 46cm (125mL)*	27 x 125mL	444-2949	Model 3500, Model 3750**
46 x 46cm (250mL)*	20 x 250mL	444-2950	Model 3500, Model 3750**
46 x 46cm (500mL)*	13 x 500mL	444-2951	Model 3500, Model 3750**
46 x 46cm (1L)*	9 x 1L	444-2952	Model 3500, Model 3750**
Tier Braces	Provides 25.4cm of clearance between stacking platforms	444-2965	

Culture Platform

Adjustable Platform

Large Vessel Carrier

Separatory Funnel Platform

Platform Size	Platform Type	Cat. No.	Compatible Shaker Models
46 x 46cm*	Culture Platform with anti-slip mat for petri dishes	444-2936	Model 3500, Model 3750**
46 x 61cm*	Culture Platform with anti-slip mat for petri dishes	444-2937	Model 3500, Model 3750**, Model 5000**
46 x 46cm	Adjustable Platform 4 Roller Bars	444-2938	Model 3500
46 x 61cm	Adjustable Platform 4 Roller Bars	444-2939	Model 5000
46 x 76cm	Large Vessel Carrier with 4 Roller Bars	444-2940	Model 5000
61 x 61cm	Large Vessel Carrier with 4 Roller Bars	444-2941	Model 10000
61 x 91cm	Large Vessel Carrier with 5 Roller Bars	444-2942	Model 15000
46 x 46cm	Separatory Funnel Platform accommodates 3 funnels 500mL to 2L	444-2953	Model 3500, Model 3750
Tier Braces	Provides 25.4cm of clearance between stacking platforms	444-2965	
46cm	Replacement Rollerbar	444-2943	444-2938, 444-2939, 444-2940
61cm	Replacement Rollerbar	444-2944	444-2941, 444-2942

* Two Tier Ready

**Platform Stacking is NOT Available on Model 3750 and Model 5000

Test Tube Rack Platform Capacity

Full Size Stationary

Rack Size: 9.5 x 41.9 x 10.2 cm

Platform Size	Cat. No.	10 to 14mm Test Tube Rack capacity = 48 444-7037	16 to 20 mm Test Tube Rack capacity = 33 444-7038	21 to 25mm Test Tube Rack capacity = 21 444-7039
28 x 33cm	444-7087	N/A	N/A	N/A
33 x 33cm	444-2930	N/A	N/A	N/A
46 x 46cm	444-2931	3	3	3
46 x 61cm	444-2932	5	5	5
46 x 76cm	444-2934	6	6	6
61 x 61cm	444-2933	7	7	7
61 x 91cm	444-2935	5	5	5

Full Size Pivoting

Internal Rack Size: 12.9 x 27.4 x 9.9cm

External Rack Size: 12.7 x 27.6 x 12.7cm

Platform Size	Cat. No.	13mm Test Tube Rack Capacity = 90 444-2955	16mm Test Tube Rack Capacity = 60 444-2956	20mm Test Tube Rack Capacity = 40 444-2957	25mm Test Tube Rack Capacity = 24 444-2958	30mm Test Tube Rack Capacity = 21 444-2959
28 x 33cm	444-7087	1	1	1	1	1
33 x 33cm	444-2930	2	2	2	2	2
46 x 46cm	444-2931	2	2	2	2	2
46 x 61cm	444-2932	3	3	3	3	3
46 x 76cm	444-2934	4	4	4	4	4
61 x 61cm	444-2933	4	4	4	4	4
61 x 91cm	444-2935	6	6	6	6	6

Half Size Stationary

Rack Size: 12.7 x 17.8 x 10.2cm

Microtube Rack: 4.3cm


Platform Size	Cat. No.	1.2 to 2 mL Microtube Rack Capacity = 70 444-0276	10 to 13mm Test Tube Rack Capacity = 63 444-0278	14 to 16mm Test Tube Rack Capacity = 48 444-0279	18 to 20mm Test Tube Rack Capacity = 35 444-0280	22 to 25mm Test Tube Rack Capacity = 24 444-0281	15mL Centrifuge Test Tube Rack Capacity = 35 444-0283	50mL Centrifuge Test Tube Rack Capacity = 12 444-0284
28 x 33 cm	444-7087	2	2	2	2	2	2	2
33 x 33 cm	444-2930	2	2	2	2	2	2	2
46 x 46 cm	444-2931	4	4	4	4	4	4	4
46 x 61 cm	444-2932	6	6	6	6	6	6	6
46 x 76 cm	444-2934	8	8	8	8	8	8	8
61 x 61 cm	444-2933	8	8	8	8	8	8	8
61 x 91 cm	444-2935	7	7	7	7	7	7	7

Full Size Stationary

Full Size Pivoting

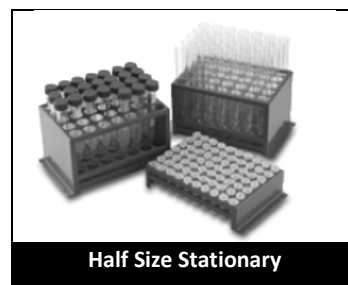Half Size Stationary

PVC Erlenmeyer Flask Clamps

Stainless Steel Erlenmeyer Flask Clamps

Stainless Steel Media Bottle Clamps

Flask Clamp Platform Capacity

PVC Erlenmeyer Flask Clamps

Platform Size	Cat. No.	125 mL 444-2960	250mL 444-2961	500 mL 444-2963	1L 444-2963	2L 444-2964
28 x 33 cm	444-7087	10	8	5	2	N/A
33 x 33 cm	444-2930	12	10	6	4	3
46 x 46 cm	444-2931	20	18	12	8	4
46 x 61 cm	444-2932	28	25	16	10	6
46 x 76 cm	444-2934	36	33	20	14	8
61 x 61 cm	444-2933	41	35	24	13	9
61 x 91 cm	444-2935	61	35	38	22	13

Stainless Steel Media Bottle Clamps

Platform Size	Cat. No.	500mL 444-7036	1L 444-7036
28 x 33 cm	444-7087	5	2
33 x 33 cm	444-2930	6	5
46 x 46 cm	444-2931	16	10
46 x 61 cm	444-2932	20	13
46 x 76 cm	444-2934	28	18
61 x 61 cm	444-2933	25	18
61 x 91 cm	444-2935	40	30

Stainless Steel Erlenmeyer Flask Clamps

Platform Size	Cat. No.	10mL 444-7040	25mL 444-7041	50mL 444-7042	125mL 444-7043	250mL 444-7044	500mL 444-7045
28 x 33 cm	444-7087	60	25	13	10	9	7
33 x 33 cm	444-2930	60	30	15	12	12	8
46 x 46 cm	444-2931	113	64	32	20	20	13
46 x 61 cm	444-2932	158	88	44	28	28	20
46 x 76 cm	444-2934	203	112	56	36	36	26
61 x 61 cm	444-2933	221	121	61	41	41	25
61 x 91 cm	444-2935	336	160	94	64	64	40

Stainless Steel Erlenmeyer Flask Clamps

Platform Size	Cat. No.	1L 444-7030	2L 444-7031	2.8L 444-7032	4L 444-7033	6L 444-7034
28 x 33 cm	444-7087	4	N/A	N/A	N/A	N/A
33 x 33 cm	444-2930	4	3	1	1	1
46 x 46 cm	444-2931	8	5	2	4	2
46 x 61 cm	444-2932	12	6	3	4	3
46 x 76 cm	444-2934	15	8	3	6	4
61 x 61 cm	444-2933	16	9	5	5	5
61 x 91 cm	444-2935	24	14	7	9	7

Microplate Clamp

Microplate Clamp		
Platform Size	Cat. No.	Microplate Clamp 444-2954
28 x 33 cm	444-7087	4
33 x 33 cm	444-2930	6
46 x 46 cm	444-2931	12
46 x 61 cm	444-2932	18
46 x 76 cm	444-2934	21
61 x 61 cm	444-2933	24
61 x 91 cm	444-2935	36

Microplate Clamp

Technical service

Web Resources

Visit the VWR website at www.vwr.com for:

- Complete technical service contact information
- Access to the VWR Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit. www.vwr.com.

Warranty


VWR warrants that this product will be free from defects in material and workmanship for a period of five (5) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorisations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorisation, unless any refusal is due to a defect of the product.

Equipment disposal

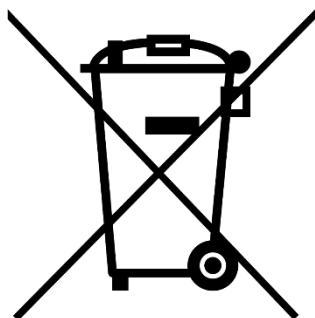This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

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India

VWR Lab Products Private Limited
No.139. BDA Industrial Suburb,
6th Main, Tumkur Road, Peenya Post,
Bangalore, India – 560058
Tel.: +91 80 28078400
Email: vwr_india@vwr.com

Ireland

VWR International Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 01 88 22 222
Email sales.ie@vwr.com

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