

Bone Clones®

800-914-0091



www.boneclones.com

OSTEOLOGICAL REPRODUCTIONS

Assembly Instructions:

SC-027-A Articulated Komodo Dragon Skeleton



Bone Clones, Inc. 9200 Eton Ave. Chatsworth, CA 91311

Phone: 818-709-7991 Fax: 818-709-7993

email: info@boneclones.com website: www.boneclones.com

Assembly Instructions
SC-027-A Articulated Komodo Dragon Skeleton
All photos © Bone Clones, Inc.



Parts Included:




- Skull
- Front Legs
- Back Legs
- Pelvis
- Shoulders
- Ribcage
- Acrylic Posts (3)
- Pins (2)
- Tail (4 pieces)



Assembly Instructions
SC-027-A Articulated Komodo Dragon Skeleton
All photos © Bone Clones, Inc.



- Step 1

Attach the shoulder assembly to the ribcage. Attach the wire hook on each side to the rib.

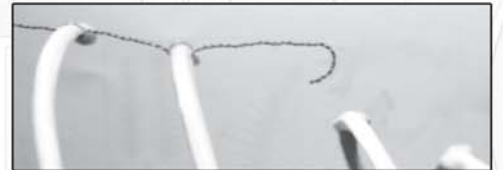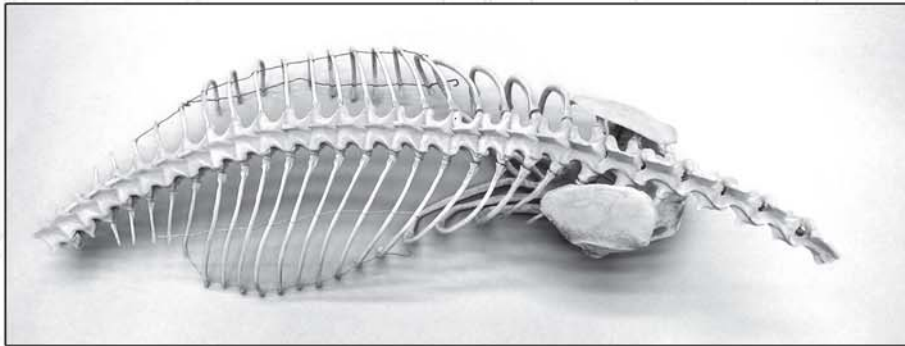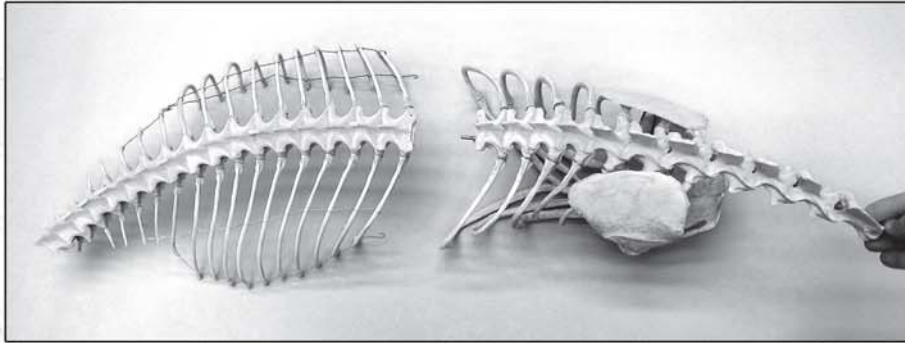


- Step 2

Slide the torso onto one of the acrylic posts.

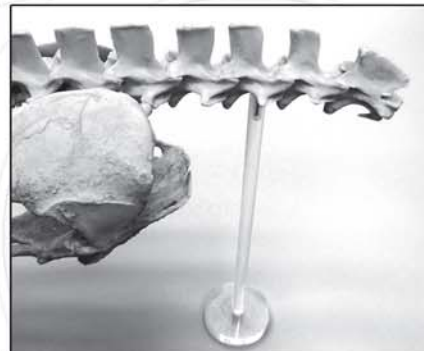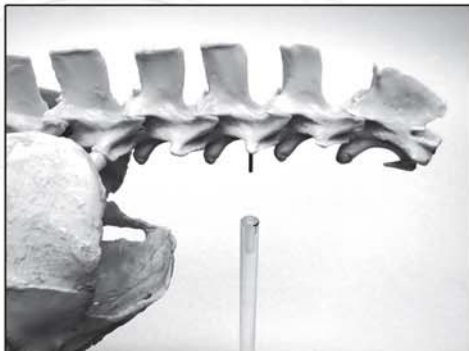


- Step 3

Repeat for the remaining 2 posts.


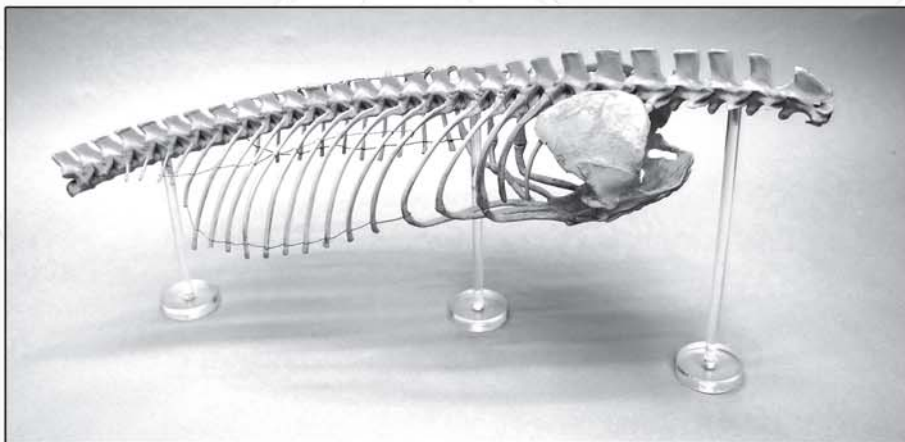


Assembly Instructions
SC-027-A Articulated Komodo Dragon Skeleton
All photos © Bone Clones, Inc.

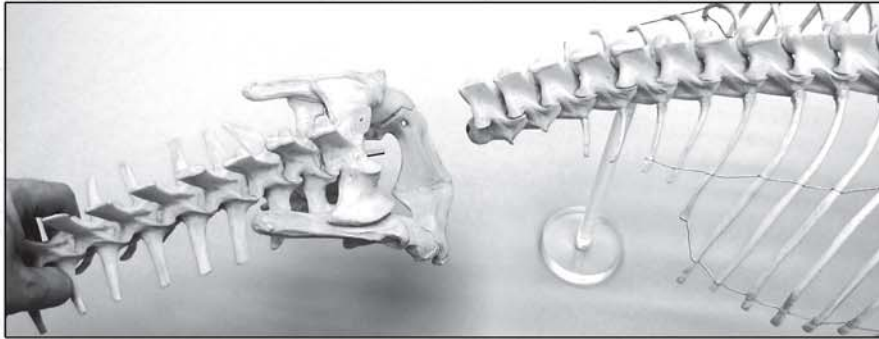

- Step 4

Attach the pelvis assembly to the torso.

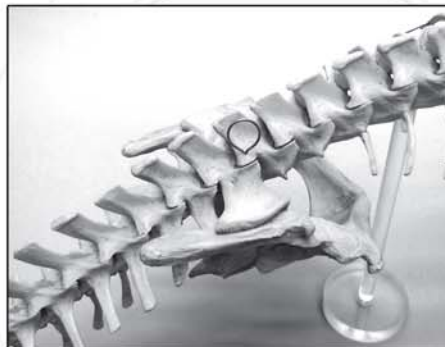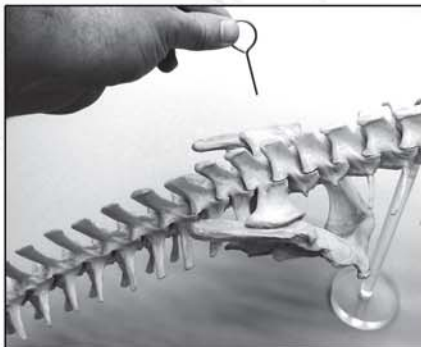

- Step 5


Slide a pin through the pelvis to secure it.



- Step 6


Repeat with second pin on the other side.

Assembly Instructions
SC-027-A Articulated Komodo Dragon Skeleton
All photos © Bone Clones, Inc.

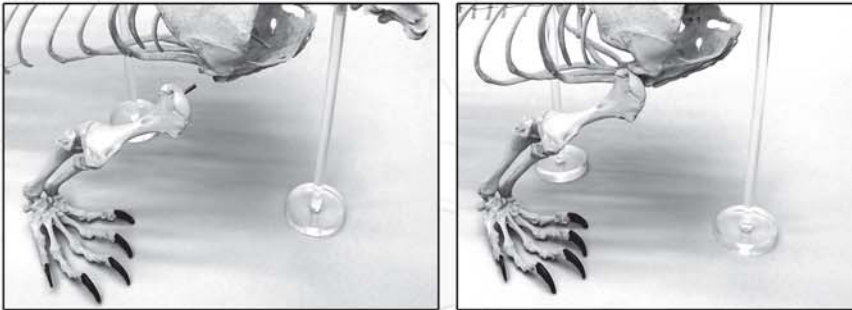


- Step 7

Attach the front-right leg to the torso.

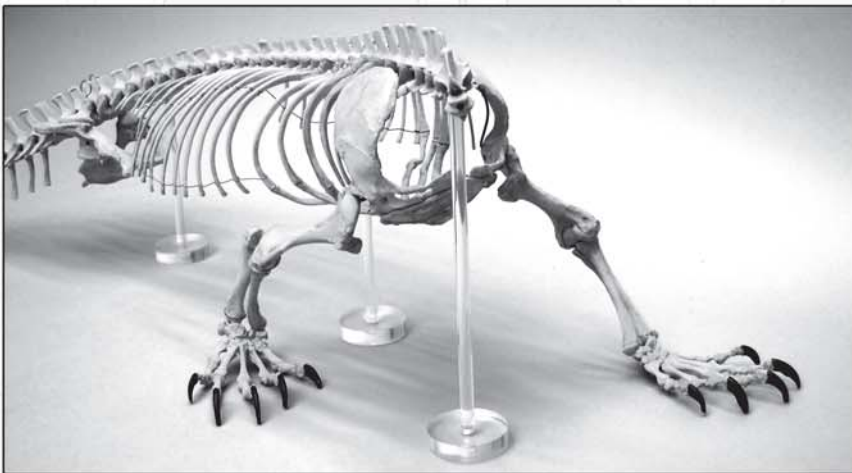


- Step 8

Repeat for the other side.

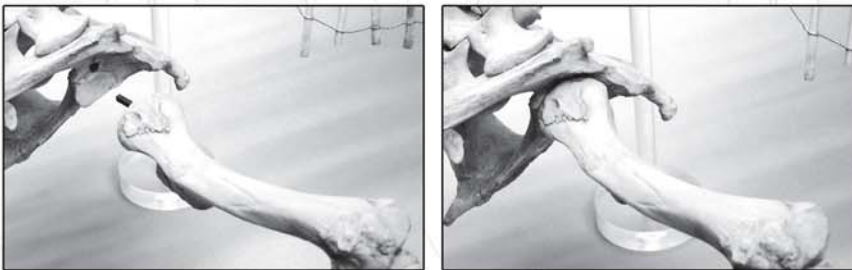

- Step 9

Attach the back-right leg to the torso.


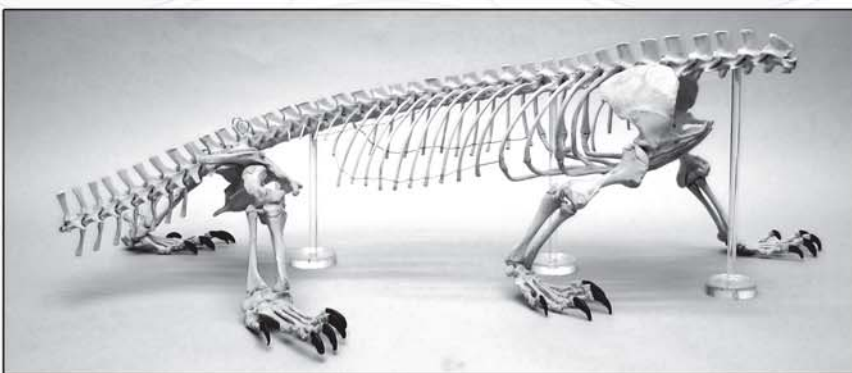


- Step 10

Repeat for the other side.



Assembly Instructions
SC-027-A Articulated Komodo Dragon Skeleton
All photos © Bone Clones, Inc.

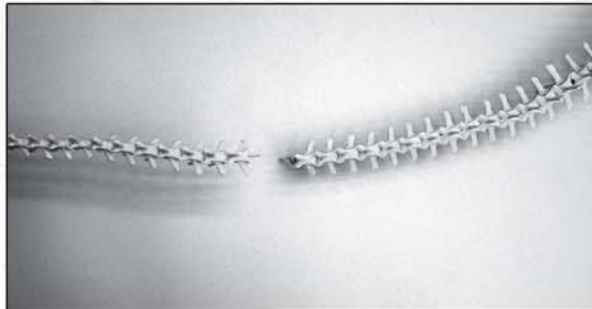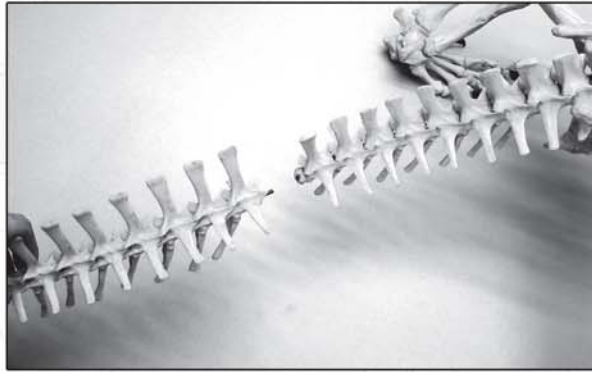

- Step 11

Attach the tail to the torso (4 pieces).



Assembly Instructions
SC-027-A Articulated Komodo Dragon Skeleton
All photos © Bone Clones, Inc.



- Step 12

Attach the head to the torso.

