# Instruction manual

## VWR® INCU-Line Tower

EU cat. no 390-0963 INCU-Line Tower

CE

UK  
CA

## **Legal address of Manufacturer**

### **Europe**

VWR International bv,  
Researchpark Haasrode 2020  
Geldenaaksebaan 464  
B-3001 Leuven  
+3216385011  
be.vwr.com

UK Importer:  
VWR International Ltd  
Hunter Boulevard, Magna Park  
Lutterworth, Leicestershire, LE17 4XN  
uk.vwr.com

## **Country of Origin**

SLOVENIA

# Table of Contents

|                                                   |    |
|---------------------------------------------------|----|
| Warning .....                                     | 4  |
| Safety information .....                          | 4  |
| Package Contents .....                            | 4  |
| Unpacking .....                                   | 4  |
| Installation.....                                 | 5  |
| Intended Use .....                                | 5  |
| Symbols and conventions .....                     | 5  |
| Product specifications .....                      | 5  |
| Overview .....                                    | 6  |
| Description of front panel .....                  | 6  |
| Instructions for use .....                        | 7  |
| Getting started .....                             | 7  |
| Operation .....                                   | 7  |
| Temperature re-adjustment.....                    | 7  |
| Troubleshooting .....                             | 8  |
| Cleaning and maintenance .....                    | 8  |
| Accessories .....                                 | 8  |
| Technical Service .....                           | 9  |
| Warranty .....                                    | 9  |
| Compliance with local laws and regulations.....   | 9  |
| Equipment disposal .....                          | 9  |
| Local VWR offices in Europe and Asia Pacific..... | 10 |

## Warning

Configuring and operating this equipment requires **precise knowledge** and **observance** of the instructions in this **user manual**. The equipment is designed only for its intended application.

**Important safety symbols** used in this manual:

**Warning!**

This is a warning of risk situations and dangers. Failure to observe this warning could be life-threatening. These warnings must be heeded.

This equipment is designed for 230V (50/60Hz).

Make sure it is connected to the correct power supply.

Disconnect the power cord before carrying out maintenance or cleaning work.

## Safety Information

The responsibility for operating the equipment lies exclusively with the owner or user if said equipment is incorrectly serviced, maintained or altered by persons not employed by an authorized dealer, or if the equipment is used in a manner contrary to its specified purpose.

The equipment must be maintained and operated in accordance with this user manual.

Ensure that this user manual is accessible at all times and has been read and understood.

To avoid electric shock do not spill any liquids on the equipment.

## Package Contents

| Description                      | Quantity |
|----------------------------------|----------|
| Cooled Incubator INCU-Line Tower | 1        |
| Stainless steel shelf            | 4        |
| Power cord                       | 1        |
| Test certificate                 | 1        |

## Unpacking

Carefully inspect the unit and accessories thoroughly upon receipt for any damage that may have occurred during shipping. Severe damage to the packaging container may indicate damage to the unit. If you suspect damage to the unit, immediately file a claim with the carrier in accordance with their instructions before contacting VWR. Also verify that all accessories are included and that the unit is in good working order before discarding the shipping packaging.

## Installation

Your satisfaction and safety require a complete understanding of this unit. Read the instructions thoroughly and be sure all operators are given adequate training before attempting to put the unit in use.

**NOTE:** This equipment must be used only for its intended application; any alterations or modifications will void your warranty and may cause injury.

## Intended use

VWR INCU-Line Tower is a general purpose convection incubator with two separate chambers, capable of uniformly cooling and heating samples from 10 °C below ambient up to a maximum of 50 °C. The unit has two transparent doors, so samples can be observed without opening the doors, thus preventing temperature changes in the chambers.

For research use only. Not intended for any animal or human therapeutic or diagnostic use.

## Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

|  |                                                                                                  |
|--|--------------------------------------------------------------------------------------------------|
|  | <b>CAUTION</b><br>This symbol indicates a potential risk and alerts you to proceed with caution. |
|--|--------------------------------------------------------------------------------------------------|

## Product Specifications

| Specifications                           | INCU-Line Tower                   |
|------------------------------------------|-----------------------------------|
| Electrical                               | 230 V, 50/60 Hz, 0.7 A            |
| Maximum total power consumption          | 160 W                             |
| Power consumption per chamber at 15 °C   | 65 W                              |
| Power consumption per chamber at 37 °C   | 35 W                              |
| Power consumption per chamber at 50 °C   | 60 W                              |
| Fuses                                    | 2 x 2 AT 250 V                    |
| Controls                                 | PID control                       |
| Temperature range                        | Ambient -10 °C to +50 °C          |
| Temperature resolution (display)         | 0.1 °C                            |
| Temperature accuracy at 37 °C            | ± 0.5 °C                          |
| Temperature fluctuation (time) at 37 °C  | ± 0.2 °C                          |
| Temperature variation (spatial) at 37 °C | ± 1.0 °C                          |
| Environment temperature                  | +15 °C to +40 °C                  |
| Interior volume                          | 2 x 23 l                          |
| Maximum load per shelf                   | 3 kg                              |
| Maximum load per chamber                 | 9 kg                              |
| Exterior dimensions W x D x H            | 380 x 470 x 790 mm                |
| Interior dimensions W x D x H            | 294 x 300 x 270 mm (each chamber) |
| Weight                                   | 34 kg                             |

## Overview

The INCU-Line Tower cooled incubator has two chambers - upper and lower. Temperature is individually controlled.


Access port (Ø 7 mm) for external temperature probe

## Description of front panel

The front panel is divided in two parts by a wave line. Display and keys on the left side (above the wave line) are used to control the upper chamber. Display and keys on the right side (below the wave line) are used to control the lower chamber.

Power switch turns the power ON and OFF for both chambers. If you use only one chamber, you can turn the other chamber OFF by simultaneously pressing the corresponding UP and DOWN keys.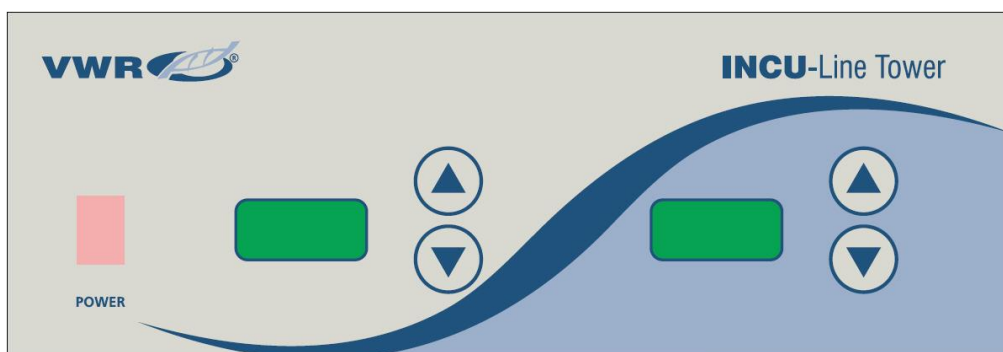

**POWER** - power switch: For turning the incubator ON and OFF. It lights when the incubator is switched ON.

**▲** - UP key: For increasing the set-temperature.

**▼** - DOWN key: For decreasing the set-temperature.

**Display**: For displaying the actual temperature (dot blinks when you change the set temperature).

## Instructions for use

### Getting Started

Check the data plate for voltage, cycle, wattage and ampere requirements. If matched to your power source, plug the power cord into a grounded outlet. Voltage should not vary more than 10 % from the data plate rating.

In selecting a location, consider all conditions which might affect performance, such as heat from radiators, autoclaves, etc. Avoid direct sun, fast-moving air currents, heating/cooling ducts and high traffic areas. Allow a minimum of 50 mm between the unit and walls or partitions which might obstruct free airflow.

### Operation

**Do not use this incubator in explosive or flammable environments. Do not heat or incubate flammable, explosive or highly reactive materials in this incubator as serious injury may result.**

Place the unit on a level surface and plug the unit into a properly grounded outlet of appropriate voltage.

Install the shelves into both chambers, at the desired height, with the shelf edges up. If you intend to place the samples on the bottom of the chamber, install the shelf so that it sits with the edges down on the chamber bottom, thus creating an air gap between the chamber bottom and your samples.

Place the samples in one or both chambers and close the doors. Make sure that both doors are securely closed.

Turn the power switch to the ON position. The switch should illuminate. On both displays all "888" are displayed at first, then software version "VX.X", and then the current temperatures for both chambers. Incubator starts to cool or to heat individual chambers independently, regarding to the set temperature for each chamber.

Display and keys on the left side (above the wave line) are used to control the upper chamber. Display and keys on the right side (below the wave line) are used to control the lower chamber.

Set desired temperature for each chamber with the corresponding UP or DOWN key. You must press and hold the UP or DOWN key for more than 2 seconds, until the DOT on the DISPLAY starts to blink. When the DOT starts to blink (it blinks for 3 seconds), set the temperature to the desired value, between 10 °C below ambient and up to a maximum of 50 °C. When the desired temperature is set, the display will automatically show the current temperature after 3 seconds (DOT stops to blink).

You can always display the SET temperature by pressing and holding the relevant UP or DOWN key for more than 2 seconds.

If you want to use only one chamber, you can turn the other chamber OFF by simultaneously pressing (short press) the corresponding UP and DOWN keys. Message "OFF" appears on the display to indicate that the chamber is turned OFF. When you want to use the chamber again, turn it ON by simultaneously pressing (short press) the corresponding UP and DOWN keys. Current temperature appears on display and cooling or heating of the chamber starts.

### Temperature readjustment

Perform the following procedure individually for each chamber.

The incubator has software for temperature readjustment. You must first measure the temperature in the middle of the chamber or in the sample for a minimum of 2 hours. You must measure the temperature with a certificated reference temperature measuring instrument with 0.1 °C precision or better. After 2 hours or more read the temperature on a thermometer and compare it with the temperature on the LED display. Difference between thermometer and display is the value that you will enter in the software.

You can insert the thermometer into the upper chamber through the opening on the top of the incubator, and into the lower chamber through the opening on the left side of the incubator.

Example 1: Reference temperature is 37.9 °C, temperature on display is 37 °C. Difference is  $37.9 - 37 = 0.9$   
Value is 0.9 (value which you enter in the software).

Example 2: Reference temperature is 36.2 °C, temperature on display is 37 °C. Difference is  $36.2 - 37 = -0.8$   
Value is - 0.8 (value which you enter in the software).

#### Procedure for temperature readjustment:

Press the UP and DOWN keys simultaneously for more than 5 seconds (while current temperature is displayed on display) - message "CO" (correction) appears on display and the DOT starts to blink.

With the relevant UP or DOWN key set the value (while DOT is blinking). Pay attention to the sign of the value (plus or minus). After 3 seconds, when you have set the temperature, display will automatically show the current temperature.

## Troubleshooting

Review the information in the table below to troubleshoot operating problems.

| Display | Cause                                                                               | Solution                                                                                             |
|---------|-------------------------------------------------------------------------------------|------------------------------------------------------------------------------------------------------|
| Er1     | Temperature probe fault, electronic board fault or incorrect setting of parameters. | Contact your distributor.                                                                            |
| Er4     | Overheating of the electronics.                                                     | Turn the main switch off and wait for the device to cool down. Then you can use the incubator again. |

## Cleaning and maintenance

**Prior to any cleaning or maintenance of the unit, disconnect the power cord from the wall outlet.**

When you are using the unit for cooling, some drops of water could condense on the inside of the chambers. Use a soft cloth to wipe the inside of the chambers, if necessary.

Clean the unit internally and externally with a damp cloth and mild soap. Do not immerse the unit or spill liquids into or on the unit. Use of alcohol or ammonia-based cleaners on the door may cause it to crack. Allow the incubator to dry completely inside and outside, before reconnecting the power and using it again.

## Accessories

| Description                                | Quantity | Cat. No. |
|--------------------------------------------|----------|----------|
| Extra shelf, 286 x 285 mm, stainless steel | 1        | 390-0483 |
| Tube rack, 2 x 6 positions, Ø 32 mm        | 1        | 390-0386 |

## Technical service

### Web Resources

Visit the VWR website at [www.vwr.com](http://www.vwr.com) for:

- Complete technical service contact information.
- Access to VWR Online Catalogue, and information about accessories and related products.
- Additional product information and special offers.

### Contact us

For information or technical assistance contact your local VWR representative or visit [www.vwr.com](http://www.vwr.com).

## Warranty

**VWR International** warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

## Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

## Equipment disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle-end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

## Local VWR offices in Europe and Asia Pacific

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### India

VWR Lab Products Private Limited  
No.139. BDA Industrial Suburb,  
6th Main, Tumkur Road,  
Peenya Post,  
Bangalore, India – 560058  
Tel.: +91-80-28078400  
Email: vwr\_india@vwr.com

### Ireland / Northern Ireland

VWR International Ltd / VWR  
International (Northern Ireland) Ltd  
Orion Business Campus  
Northwest Business Park  
Ballycoolin  
Dublin 15  
Tel.: +353 01 88 22 222  
Email sales.ie@vwr.com

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VWR International AB
VWR International GmbH