VWR

Solar Battery Pack 

INSTRUCTION MANUAL

Model	European Catalogue Number
SBP10	630-1637

Version: 3
Issued: 19, 05, 2014

Legal Address of Manufacturer

Europe

VWR International bvba
Researchpark Haasrode 2020
Geldenaaksebaan 464
B-3001 Leuven
+ 32 16 385011
<http://be.vwr.com>

Country of origin: **ITALY**

Table of Contents

Safety information

Package contents

Symbols and conventions

Instructions for use

Product Specifications

Overview

Description of Buttons and Switches

Connection guide

Operation

Troubleshooting

Technical service

Warranty

Compliance with local laws and regulations

Disposal

Safety Information

 Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position.

Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

Package Contents

DESCRIPTION	QUANTITY
Solar battery pack unit	1
USB / Micro-USB cable	1
USB / Jack cable	1
Power supply OUTPUT 5Vdc	1

Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.


	CAUTION This symbol indicates a potential risk and alerts you to proceed with caution
---	---

Instructions for use

Product Specifications

Battery	polymer lithium-ion
Capacity	2600 mAh
Input (Micro-USB port)	5Vdc
Charging time (through USB)	about 5 hours
Output (USB port)	5.5Vdc / 1A
Solar panel power	0,7W max
Solar panel output current	80mA (at 1000W/m ² light)
Working temperature	-10°C ~ +55°C

Overview

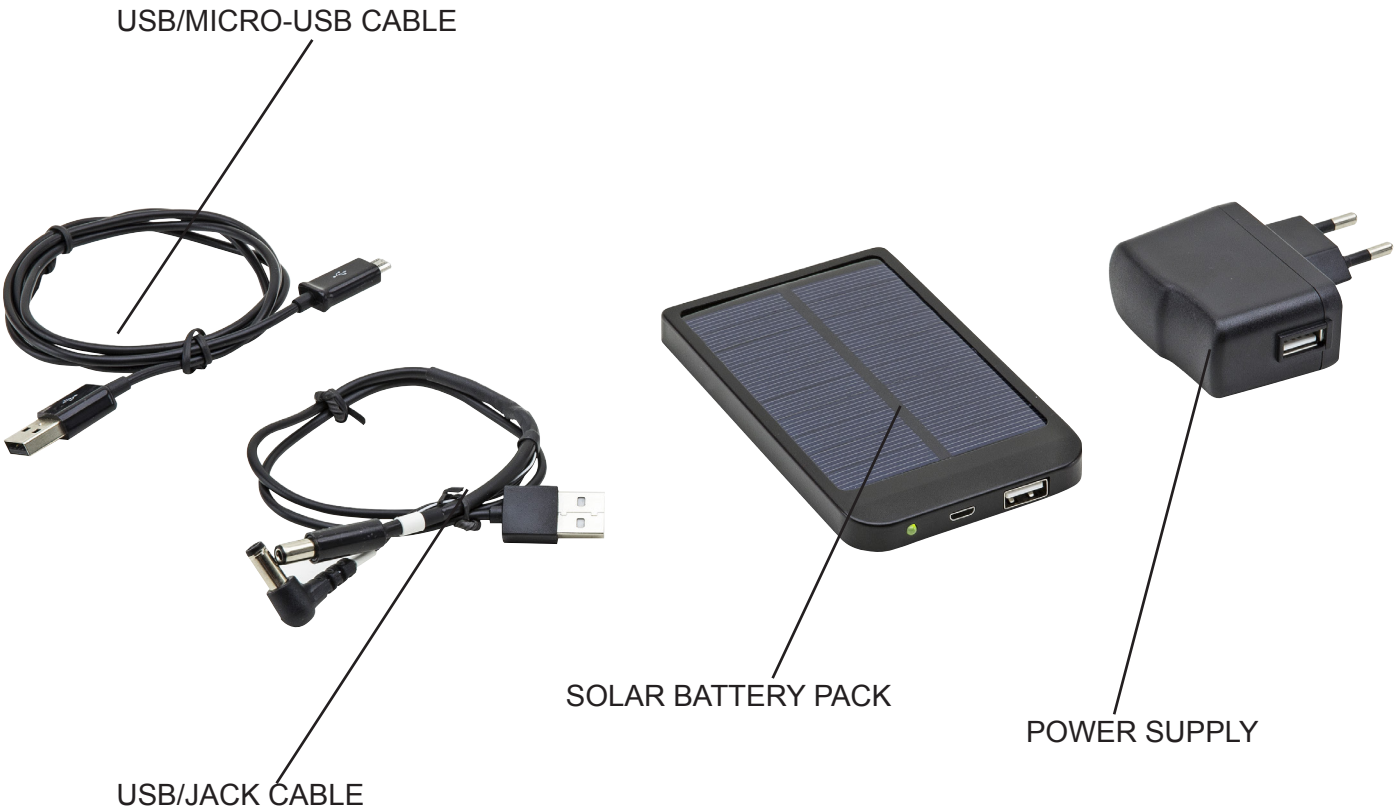Description of Buttons and Switches

Connection guide

Plug your microscope to the battery pack using the provided USB/Jack cable:

Jack 1 - Compatible models

Series 200 (ML214 - ML224 - BL214 -BL224)
Series 100
Series 100D
Series 100POL

Jack 2 - Compatible models

Series 300
Series 200 (STB250 - SZB250)
Series 800

Operation

Fully recharge the battery via USB cable and 5V power supply when using it for the first time.

Connect the USB/Jack cable to the USB output on the battery pack and to the input Jack of the microscope.

Recharging

a) Expose the front solar panel to sun, the LED indicator will blink.

b) Connect via USB/micro-USB cable to the provided power supply. c) Connect via USB/micro-USB cable to a PC USB port.

Troubleshooting

Review the information in the table below to troubleshoot operating problems.

PROBLEM	CAUSE	SOLUTION
AFTER CHARGING FOR LONG TIME, INDICATORS DON'T SHOW FULL CAPACITY.	Charging time is not enough	Charge it 8 hours continuously
	Power supply or battery pack is faulty	Refer to technical assistance
MICROSCOPE LIGHT IS NOT WORKING.	No power left	Recharge the battery
	Something wrong with cables and connections	Check that USB is well inserted into the output connector of the battery. Check that the jack connector is well inserted into the input of the microscope.
	Wrong settings of microscope switches	Check that the main power switch of the microscope is in ON position. Check that the brightness adjustment knob of the microscope is set to the desired value.

Technical service

Web Resources

Visit the VWR's website at **www.vwr.com** for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit. **www.vwr.com**.

Warranty

VWR International warrants that this product will be free from defects in material and workmanship for a period of 24 months from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

Disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it is your responsibility to correctly dispose of your equipment the end of its life cycle by handling it over to an authorized facility for separate collection and recycling. It is also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

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Ireland / Northern Ireland

VWR International Ltd /
VWR International (Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: 01 88 22 222
Fax: 01 88 22 333
E-mail: sales@ie.vwr.com

Italy

VWR International PBI S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: 02-3320311/02-487791
Fax: 800 152999/02-40090010
E-mail: info@it.vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: 020 4808 400
Fax: 020 4808 480
E-mail: info@nl.vwr.com

New Zealand

Global Science - A VWR Company
241 Bush Road
Albany 0632, Auckland
Tel.: 0800 734 100
Fax: 0800 999 002
E-mail: sales@globalscience.co.nz

Norway

VWR International AS
Haavard Martinsens vei 30
0978 Oslo
Tel.: 02290
Fax: 815 00 940
E-mail: info@no.vwr.com

Poland

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: 058 32 38 200 do 204
Fax: 058 32 38 205
E-mail: info@pl.vwr.com

Portugal

VWR International -
Material de Laboratório, Lda
Edifício Neopark
Av. Tomás Ribeiro, 43- 3 D
2790-221 Carnaxide
Tel.: 21 3600 770
Fax: 21 3600 798/9
E-mail: info@pt.vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel.: +65 6505 0760
Fax: +65 6264 3780
E-mail: sales@sg.vwr.com

Spain

VWR International Eurolab S.L.
C/ Tecnología 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: 902 222 897
Fax: 902 430 657
E-mail: info@es.vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: 08 621 34 00
Fax: 08 621 34 66
E-mail: kundservice@se.vwr.com

Switzerland

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: 044 745 13 13
Fax: 044 745 13 10
E-mail: info@ch.vwr.com

Turkey

Pro Lab Laboratuvar Teknolojileri Ltd.Şti.
a VWR International Company
Orta Mah. Cemal Gürsel Caddesi
Ördekcioglu İşmerkezi No.32/1
34896 Pendik - İstanbul
Tel.: +90216 598 2900
Fax: +90216 598 2907
Email: info@pro-lab.com.tr

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard - Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: 0800 22 33 44
Fax: 01455 55 85 86
E-mail: uksales@uk.vwr.com

**GO TO VWR.COM FOR THE
LATEST NEWS, SPECIAL OFFERS
AND DETAILS OF YOUR LOCAL
VWR DISTRIBUTOR**

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