

ProSieve™ Color Protein Markers

Volume: 500 µl Cat. No. 50550
Volume: 100 µl Cat. No. 50552

Store: -20°C

For best results, lyophilized material should be stored at -20°C and used within five years from date of manufacture (DOM).

Stable for three months once reconstituted. Do not freeze/thaw reconstituted vials of ProSieve™ Color Marker more than five times.

Recommended procedure

First Time Use

1. Add 500 µl water (Cat. No. 50550) or 100 µl water (Cat. No. 50552).
2. Warm briefly to 37°C and mix well.
3. Load 10 µl per well on standard polyacrylamide mini-gels.
4. Aliquot and store at -20°C.

Subsequent Use

Warming aliquots to 37°C is not required for subsequent use of reconstituted ProSieve™ Color Protein Marker.

For Research Use Only.

All trademarks herein are marks of the Lonza Group or its affiliates.

© 2007 Lonza Rockland, Inc.
All rights reserved.