ENGLISH

FRIDGE / FREEZER THERMOMETER

FEATURES

- Fridge / Freezer high and low temperature alarm setting
- Maximum / minimum temperature memory
- Fridge / Freezer or Room temperature display
- Waterproof fridge sensor
- °C / °F Exchange
- Magnetic attachment
- Velcro Magic Tape Fixing

SPECIFICATION

Measuring range	-10 ~ 50°C / 14 ~ 122°F (room temperature) -50 ~ 70°C / -58 ~ 158°F (fridge / freezer temperature)
Display accuracy	±1°C / ±1.8°F over full range
Display resolution	0.1°
Display reading update	10 seconds
Battery	1.5 volt, size A76 or equivalent x1 piece
Display size	26(W) x 13(H) mm
Battery Life	About 1 year
Fridge sensor cord length	1 meter
Product size	63(W) x 48(H) x 12 (D) mm
Accessories	1) Sensor holder x1 piece 2) Battery x1 piece 3) Velcro magic tape x1 piece

INSTALLATION

- Pull down battery cover.
- Install the battery with correct polarity positioning.
- The thermometer will have a full segment display for about 3 seconds and then display the fridge sensor temperature normally.
- Peel off the adhesive tape cover sheet on the sensor holder.
- Firmly attach the sensor holder onto the desired position inside the fridge / freezer.
- Insert the fridge sensor into the sensor holder.
- Attach thermometer onto the desired space outside the fridge / freezer by using the supplied Velcro magic tape or magnets on the back of thermometer.
- Peel off display protective sheet.

OPERATION

°C / °F Exchange

Press [°C/°F] button to select temperature unit.

Maximum/minimum reading memory

- Press [Max/Min] button to display maximum measured value, press the button once more to display minimum measured value, press the button again to return to normal display.
- Press and hold [Max/Min] button about 2 seconds to reset memory to current reading.

Room or fridge / freezer temperature display

Press [ROOM/FRIDGE] button to display room or fridge / freezer temperature.

Fridge/freezer High and low temperature alarm setting

- Press and hold [] button for about 2 seconds, the HIGH and FRIDGE icons will flash.
- Press [] button to set the alarm high limit.
- Press [] button again, the LOW and FRIDGE icons will flash.
- Press [] button to set the alarm low limit.
- Press [] button once more to finish setting.
 - During limits setting, holding [] button will advance the value automatically.
 - If the fridge sensor reading is out of the set limit then the alarm will sound.

Alarm on / off

Press [] button to switch alarm off or on.

NOTE

Do not put thermometer inside the fridge / freezer

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