## Sonda 50 10 T PT-BR

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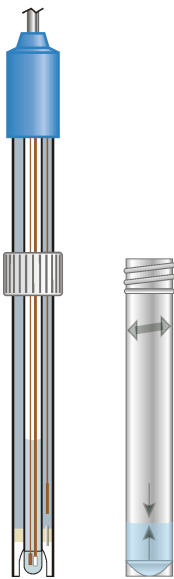
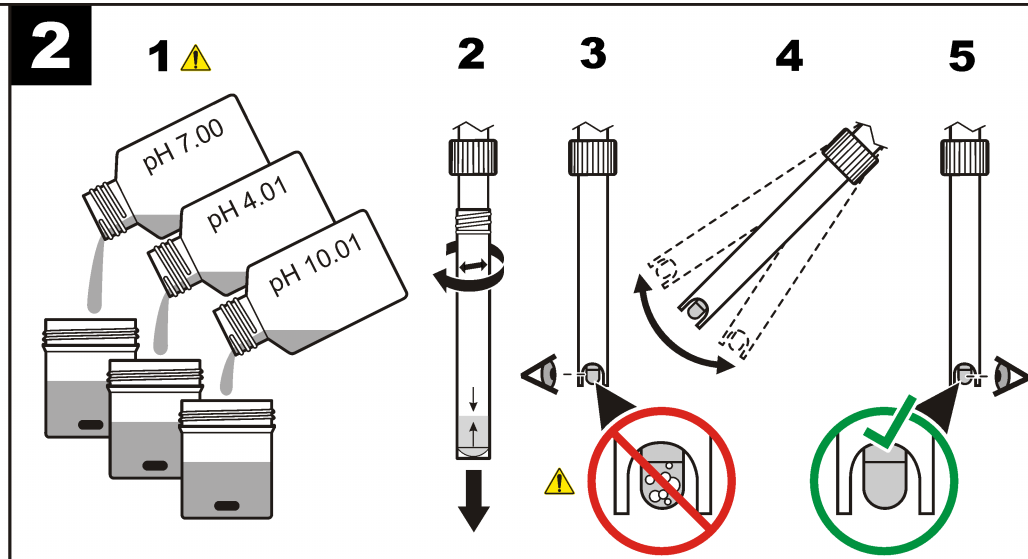
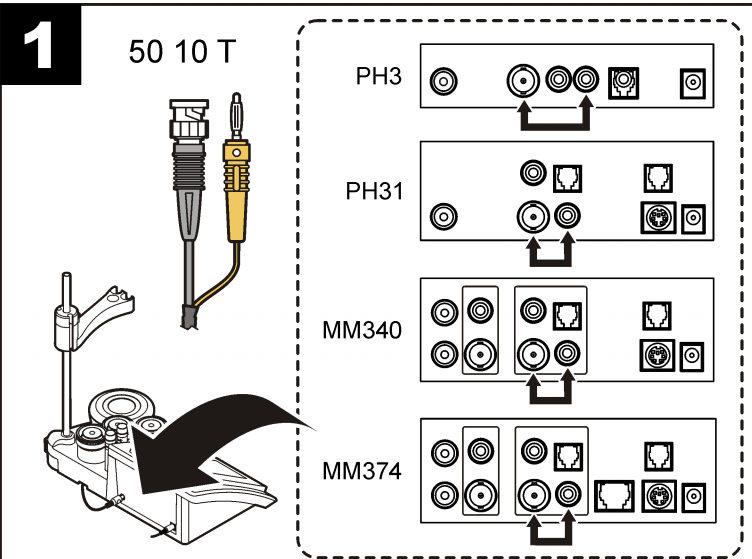
### 50 10 T probe

Refer to the meter manual for more information.

**Caution**  
 ⚠ *Personal injury hazard. Glass components can break. Handle with care to prevent cuts.*  
 ⚠ *Chemical exposure hazard. Refer to the current material safety data sheets (MSDS) for safety protocols.*

**Introduction**  
 This pH gel-filled electrode with temperature sensor is used in general water samples.

**Technical information**

Range	0 to 14 pH
Junction	Ceramic
Reference	Ag/AgCl
Temperature	Pt 1000
Minimum depth	12 mm (0.5 in.)
Operating temp	0 to 80 °C (32 to 176 °F)
Limits	Low conductivity, dirty or viscous solutions; colloids or solids in suspension; samples with sulfides

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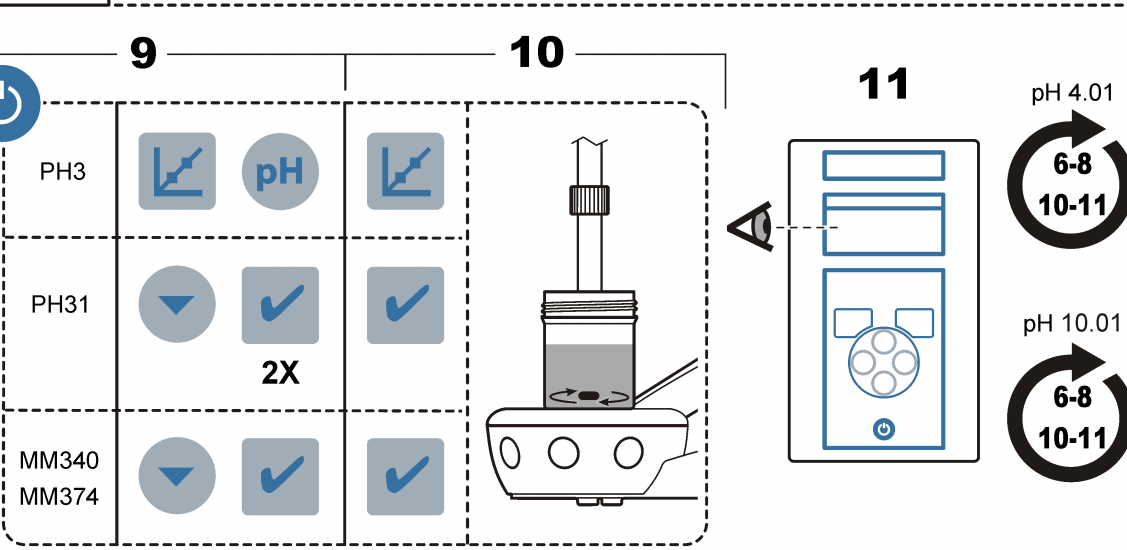
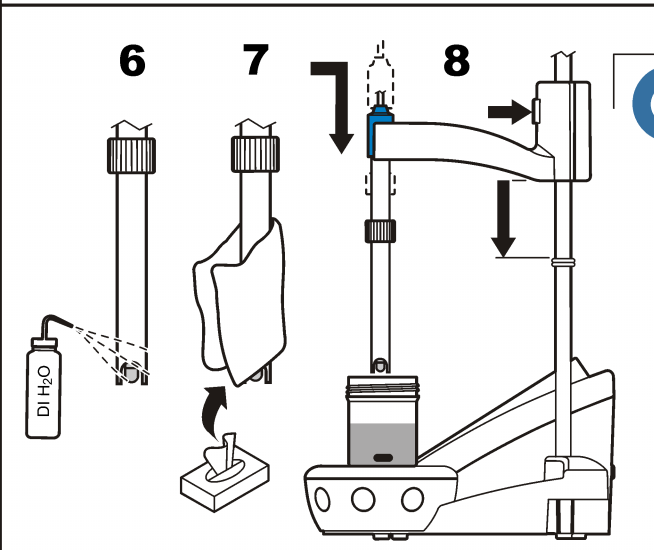
ES

**1 Connect the probe**  
**2 Calibration**  
 ⚠ *Air bubbles in or under the probe tip when submerged can cause slow stabilization or error in measurement.*  
**3 Measurement**

**Maintenance and storage**  
 Cleaning: soak the probe in a cleaning solution. Rinse with deionized water.

Contamination	Cleaning agent
Proteins	Pepsin cleaning solution
Grease, oils, fats	Electrode cleaning solution
Limescale	0.1 N HCl solution

Hydration: soak the dry probe in a dilute HCl solution for several hours.  
 Storage: put the probe in the protector with KCl 3 M solution.

**Troubleshooting**  
 Be sure that the samples are suitable for the application.

**Warranty**  
 6 months for manufacturing faults only. Damage from use is not covered.

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