



CE US Pro in action!

# Return to work

## Cleaning and disinfection with CE Duo™ Sponge and CE UltraSpeed Pro® for Controlled Non-Classified areas

### What to Clean/Disinfect?

#### High Touch Surfaces

##### Horizontal Surfaces:

- Tables, counter tops & work surfaces
- Chair seats, & floor mats
- Floors & ceilings

##### Vertical Surfaces:

- Walls, windows, softwalls

##### Uneven Surfaces:

- Misc. equipment
- Pipes, tanks, hoods
- Door handles, knobs, light switches, faucets and sinks



#### High Touch Devices

##### Work Devices:

- Computer components, telephones, and cell phones

### When to Clean?

Per SOP or



Start and end of the work day



After each shift


### What to Use?

#### Sponges

CE Duo™ CleanTech™ Sponge



Pre-saturate with Vileda's 6 Liter Bucket

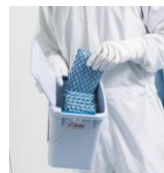

#### Mops

CE UltraSpeed Pro System with MicroIntensive™ L Mop



#### Disinfectant

Per SOP or visit EPA's site for recommended disinfectant to combat Covid-19


[Click here for the EPA List N](#)

### How to Clean/Disinfect?

#### Wiping with Sponges


Use one sponge per surface to prevent cross contamination.

Wiping with the sponges should be done in a pull & lift fashion with overlapping strokes. This can be done horizontally or vertical as shown below. Following this step is critical in removing contamination instead of spreading in circles.

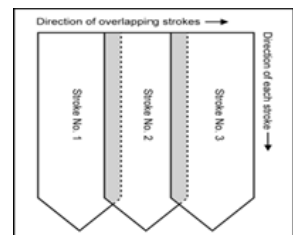

#### Mopping

With Mops use a modified figure 8 (S-Method). Right to left, then left to right, with overlapping strokes

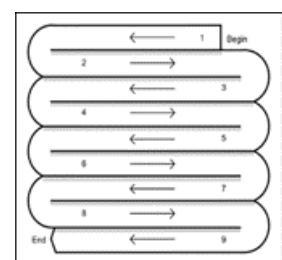

When the mop or sponge begins to dry out re-dip into the solution to re-saturate.



Accessories are optional\*

## CE Duo™ CleanTech™ Sponge

VWR Cat. No.	Description	Color	Size	Pieces/PK	PKS/CS
75871-654	CE Duo™ CleanTech™ Sponge	Blue/White	4 x 7 in	6	10
10153-180*	6l Bucket - accessory	Grey	11 x 8 x 7 in	NA	10
10153-182*	6l Bucket Lid - accessory	Grey	11 x 8 x 1 in	NA	12

## CE UltraSpeed Pro® System

VWR Cat. No.	Description	Color	Size	Pieces/PK	PKS/CS
76170-718	CE US Pro® Double Bucket Kit	Grey	11 & 26 qt	NA	1
76170-720	MicroIntensive® L Mop (single – use)	White	16 in	1	50
76170-720	CE US Pro® Frame	Grey	16 in	NA	10
89066-060	Vileda Aluminum Extendable Handle	Blue w/ coding	40 x 71 in	NA	10



800 932 5000 | VWR.COM

Prices, product, and/or services details are current when published and subject to change without notice. | Certain products or services may be limited by federal, state, provincial, or local regulations. | VWR, part of Avantor, makes no claims or warranties concerning sustainable/green products. Any claims concerning sustainable/green products are the sole claims of the manufacturer and not those of VWR International, LLC and/or Avantor, Inc. or affiliates. All prices are in US dollars unless otherwise noted. Offers valid in US, void where prohibited by law or company policy, while supplies last. | Trademarks are owned by Avantor, Inc. or its affiliates, unless otherwise noted. | Visit [vwr.com](http://vwr.com) to view our privacy policy, trademark owners, and additional disclaimers. © 2020 Avantor, Inc. All rights reserved.

0620 Lit. No. 06001W