
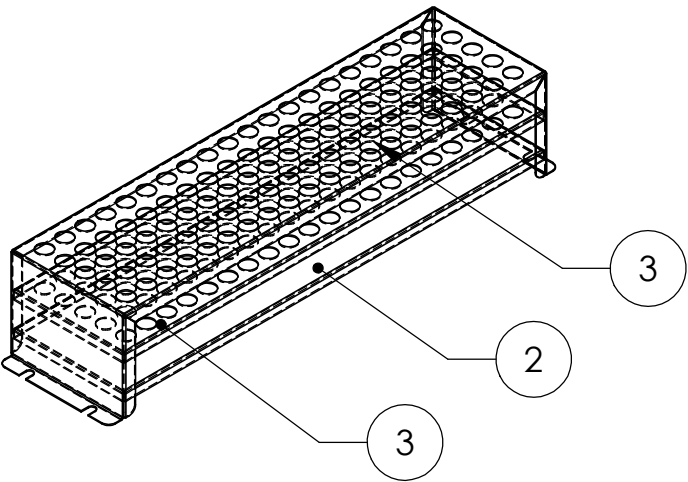


| | | | | |
|------------------------------|--|--|--|--|
| DWG BY: DKB DATE: 3/14/02 | TOLERANCE UNLESS OTHERWISE SPECIFIED .X = ± .063 .XX = ± .032 .XXX = ± .015 | SHELDON MANUFACTURING CORNELIUS, OREGON 10-13 MM SHAKER ASSEMBLY | SCALE: 1:4 PART#: 9751177 MATERIAL: ASSEMBLY | NOTE: BEND RADIUS .032" UNLESS OTHERWISE NOTED |
| | | | | |

| CHG. BY | CHG. # | REV. | DATE | DESCRIPTION |
|---------|--------|------|------|-------------|
| | | | | |

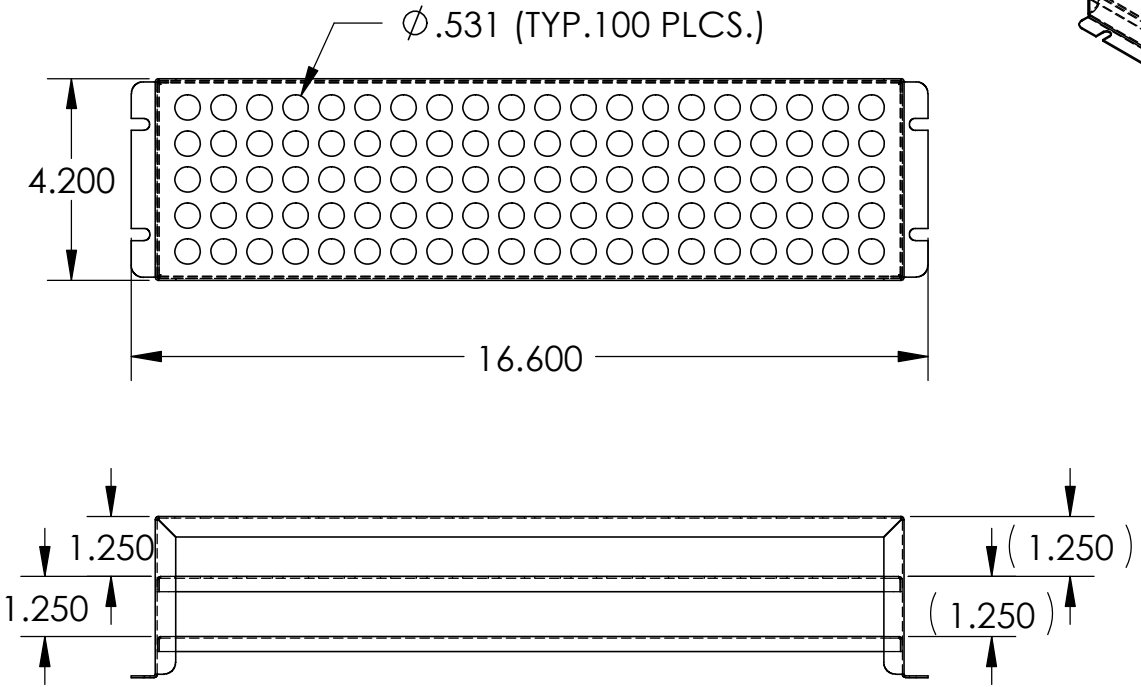

| ITEM NO. | QTY. | PART NO. | MATERIAL | DESCRIPTION |
|----------|------|----------|------------------|-----------------------|
| 1 | 1 | 5121163 | .036 SST. 304 2B | SHAKER TOP 10-13MM |
| 2 | 1 | 5121162 | 20 SST 304 2B | SHAKER SHELF |
| 3 | 1 | 5121164 | .036 SST. 304 2B | SHAKER SHELF 10-13 MM |