

Lentiviral Vector Production

Streamlined gene delivery systems.

PRODUCTION METHODS

Lentiviral Vector Construction:

1. Clone gene of interest into a modified Lentiviral vector.
2. Purify the constructed Lentiviral vector plasmid and packaging plasmids.

To Produce Lentiviral Particles by Transient Transfection In 10 mL Cell Culture Dish:

1. Day 0: One day prior to the transfection, plate 293T cells in 10 mL DMEM/10% FBS at a density of 2×10^6 per 100-mm tissue culture plate. Incubate.
2. Day 1: On the day of transfection, change culture medium with 10 mL fresh medium 1 hour prior to the transfection.
3. Mix DNAs used for Lentiviral particle production in a sterile 6 mL polypropylene tube.
 - I. Adjust the volume to 437 μ L with TE79/10.
 - II. Add 63 μ L of 2 M CaCl_2 and mix well.
 - III. Add 500 μ L of 2 \times HBS with constant agitation.
 - IV. Sit the mixture at room temperature for 30 minutes to allow calcium phosphate-DNA to precipitate.
4. Add the precipitate by drop into tissue culture plates in which 293T cells are at least 80–90% confluent.

Tips to titer: The cell density is critical for vector production. The best results are obtained when the plate is 90% confluent on the day of transfection. Vi-Cell for precise cell counting can help you reach the ideal cell confluence to increase titer.

5. After 6–8 hours, replace the culture medium with 6 mL fresh DMEM/10% FBS and continue the incubation.
6. Day 2–4: Collect the culture supernatant and replace by 6 mL fresh culture medium. Filter the collection through a sterile 0.4 μ m syringe filter, and store at ultra-low temperature (ULT).

SUPPORTING PRODUCTS

OPTIMA XPN-100
with NVT rotors (100, 90, 65 & 65.5)

MICROFUGE 16, GeXP

OPTIMA MAX-XP
with MLA rotors (150 & 130) and TLA rotors (120.1 & 120.2)

AVANTI J-26S XP
with JA rotors (10, 14) and JLA rotors (16.250, 10.500), JA rotors (17, 20 & 25.5)



ALLEGRA X-15R, ALLEGRA X-14, ALLEGRA X-30



VI-CELL

Polyethylene Glycol (PEG) Purifies and Concentrates Lentiviral Particles:

1. Mix thawed collection with 40% PEG solution to a final PEG concentration of 10%. Incubate the mixture in ice for 3–6 hours.
2. Spin at 2,000 × g for 30 minutes.
3. Discard the supernatant, disperse viral particle pellet by gentle pipetting in 1/20 of the original harvest volume of PBS (Phosphate Buffered Saline) or media of your choice.
4. Place the tubes into buckets. Weigh and balance them.
5. Spin at 100,000 × g (24,500 RPM) at 4°C in a SW 32 Ti rotor for 90 minutes, in a Beckman Optima X Series ultracentrifuge.
6. Remove the supernatant by inversion of the tubes or pipetting; be careful not to dislodge the viral pellet.
7. Re-suspend the pellet in PBS or the media of your choice.
8. Pipette up and down or shake for a few minutes, if necessary, to fully dissolve the pellet.
9. Aliquot and store at desired temperature; ultra-low temperature (ULT) storage is recommended for long term.

IMPROVED PROCESS:

Rotors	Tube	Part Number	Adapter	Process Advantages
SW 55 Ti	3.2 mL <i>g</i> -Max, konical and BioSafety with Quick-Seal	358647	355535 and 358153	Increased concentration, biosafety, reduced sample volume
SW 32.1 Ti	4.5 mL <i>g</i> -Max and BioSafety with Quick-Seal	356562	355579	Reduced sample volume, biosafety
	8.0 mL <i>g</i> -Max and BioSafety with Quick-Seal	344621	355579	Reduced sample volume, biosafety
SW 32 Ti	15 mL <i>g</i> -Max and BioSafety with Quick-Seal	343664	355536	Reduced sample volume, biosafety
	8.4 mL <i>g</i> -Max, konical and BioSafety with Quick-Seal	358652	355536 and 358156	Reduced sample volume, biosafety, increased concentration
TLS-55*	2.2 mL Ultra-Clear	347356	—	Miniaturization
MLS-50*	5.0 mL Ultra-Clear	344057	—	Miniaturization
SW 41 Ti	13.2 mL Ultra-Clear	344059	—	Reduced sample volume

*TLS and MLS rotors are used with the Optima MAX-XP tabletop ultracentrifuge.

References:

1. Hsin-Lung Lo and Jiing-Kuan Yee; Production of Pseudotype-Retroviral Vectors—Current Protocols in Human Genetics. John Wiley & Sons, Inc. pp. 12.71-11: 2007.
2. Jane C. Burns, Theodore Friedmann, Wolfgang Driever, Michelle Burrascano, and Jiing-Kuan Yee; Vesicular stomatitis virus G glycoprotein pseudotyped retroviral vectors: Concentration to very high titer and efficient gene transfer into mammalian and nonmammalian cells—Proc. Natl. Acad. Sci. Vol. 90: pp. 8033-8037.
3. Steven R. Bartz and Marie A. Vodicka; Production of High-Titer Human Immunodeficiency Virus Type 1 Pseudotyped with Vesicular Stomatitis Virus Glycoprotein—METHODS: A Companion to Methods in Enzymology. 12: pp. 337-342.
4. Jiing-Kuan Yee, Atsushi Miyanohara, Patricia Laporte, Kathy Bouic, Jane C. Burns, and Theodore Friedmann; A general method for the generation of high-titer, pantropic retroviral vectors: Highly efficient infection of primary hepatocytes—Proc. Natl. Acad. Sci. Vol. 91: pp. 9564-9568.
5. Richard A. Klinghoffer, Brian Roberts, James Annis, Jason Frazier, Patrick Lewis, Peter S. Linsley, and Michele A. Cleary; An Optimized Lentivirus-Mediated RNAi Screen Reveals Kinase Modulators of Kinesin-5 Inhibitor Sensitivity—ASSAY and Drug Development Technologies. Volume 6: pp. 105-119.
6. Jean-Michel Garcia, Anhui Gao, Pei-Lan He, Joyce Choi, Wei Tang, Roberto Bruzzone, Olivier Schwartz, Hugo Naya, Fa-Jun Nan, Jia Li, Ralf Altmeyer, and Jian-Ping Zuo; High-throughput screening using pseudotyped lentiviral particles: A strategy for the identification of HIV-1 inhibitors in a cell-based assay—Antiviral Research. 81: pp. 239-247.
7. H-L Lo, T Chang, P Yam, PM Marcovecchio, S Li, JA Zaia, and J-K Yee; Inhibition of HIV-1 replication with designed miRNAs expressed from RNA polymerase II promoters—Gene Therapy. 14: pp. 1503-1512.



Life Sciences

OptiSeal and Optima are trademarks of Beckman Coulter, Inc. Beckman Coulter, Quick-Seal, Allegra, Avanti, Vi-Cell, and the stylized logo are trademarks of Beckman Coulter, Inc. and are registered with the USPTO.

For Beckman Coulter's worldwide office locations and phone numbers, please visit "Contact Us" at www.beckmancoulter.com