07519-15 Pump Head with 07519-75 Large Cartridge
Tête de pompe 07519-15 à cartouche épaisse 07519-75
Pumpenkopf 07519-15 mit großer Schlauchkassette 07519-75
Cabeza de bomba 07519-15 con cartucho grande 07519-75
Testa pompante 07519-15 con cartuccia grande 07519-75

Model Nos. Modèles n°
Modellnummern Números de
modelo Modelli n°

MFLX07519-10
MFLX07519-15
MFLX07519-20
MFLX07519-25

OPERATING MANUAL

CARTRIDGE PUMP HEAD SYSTEM

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MFLX07519-75

**PUMP FOR LIQUIDS
ORIGINAL INSTRUCTIONS**

CONTENTS

Title	Page
SAFETY PRECAUTIONS	2
INTRODUCTION	3
(1) Application Data	3
(2) General Description	3
INSTALLATION	4
SETUP	5
(1) Select Cartridges	5
(2) Load Cartridges	5
(3) Cartridge Removal	7
(4) Bi-Directional Pumping	7
(5) Partial Bank Pumping	8
OPERATION	8
(1) Relative Pulsation Rates	8
(2) Select Tubing and Pump Heads	9
(3) Select Pump Speed	10
(4) Adjust Occlusion Settings on Cartridges	10
(5) Pump Overloading	11
MAINTENANCE	11
Service Parts	11
SPECIFICATIONS	12
WARRANTY	12
PRODUCT RETURN	12
TECHNICAL ASSISTANCE	12
APPENDIX A MASTERFLEX® L/S® Pump Drive Types	13
APPENDIX B Cartridge Loading Capacity	13

C-FLEX, PHARMED, NORPRENE, TYGON —Reg TM Saint-Gobain Performance Plastics Corp.
SANTOPRENE—Reg TM Exxon Mobil
Trademarks bearing the ® symbol in this publication are registered in the U.S. and in other countries.

US & Canada only
Toll Free 1-800-MASTERFLEX | 1-800-637-3739
Outside US & Canada
1-847-381-7050

***EN809 manufactured by:**
Masterflex LLC.
28092 W Commercial Avenue, Barrington, IL 60010
masterflex.tech@avantorsciences.com
www.avantorsciences.com/masterflex

SAFETY PRECAUTIONS

WARNINGS: *Tubing breakage may result in fluid being sprayed from pump. Use appropriate measures to protect operator and equipment.*

Turn off drive before removing or installing Cartridges. Safety guards are provided to minimize risk of fingers getting caught between the roller mechanism and the base of the module. However, be safe—Keep your fingers away from these areas.

CAUTIONS: *Excessive occlusion can cause high pump temperatures and early tubing and roller failure. Do not set occlusion tighter than what is recommended in Table 4.*

Use only tubing/cartridge/pump combinations defined by Tables 1, 2, and 3. Use of other combinations could cause malfunction or damage to the pump.

Operating the pumps at speeds in excess of 250 rpm may result in premature wear of the pump rollers and tubing.

Explanation of Symbols

CAUTION: *Risk of Danger. Consult Operator's manual for nature of hazard and corrective actions.*

CAUTION: *Risk of crushing. Keep fingers away from rotor while pump is in operation. Stop pump before loading or unloading tubing.*

WARNING: PRODUCT USE LIMITATION

These products are not designed for, nor intended for use in patient connected applications, including, but not limited to, medical and dental use and, accordingly, have not been submitted for FDA approval.

INTRODUCTION

The instructions in this manual are task-oriented for easy reference. You can go directly to a particular section and quickly find the answers. Appendix A and B list MASTERFLEX® L/S® pump drive types which can be used with this system.

Pump Heads

MODEL	TYPE
Model 07519-10	6-Roller, 2 Large or 4 Small Channel, No Speed Reduction
Model 07519-20	8-Roller, 2 Large or 4 Small Channel, No Speed Reduction
Model 07519-15	6-Roller, 6 Large or 12 Small Channel, 5:1 Speed Reduction
Model 07519-25	8-Roller, 6 Large or 12 Small Channel, 5:1 Speed Reduction

Cartridges

MODEL	TYPE
Model 07519-75	Large
Model 07519-85	Small

These Cartridge Pumps are designed to provide up to 12 simultaneously driven pump channels and the ability to provide nearly pulse-free flow. In addition, the pump systems incorporate the following features:

- 5:1 speed reduction on models 07519-15 and 07519-25, providing ultra-low flow rates.
- 12 cartridge capacity of models 07519-15 and 07519-25 permit dispensing simultaneously into 12 separate containers with one set-up and pump cycle.
- Six and eight roller designs provide reduced pulsation.
- Elimination of “suckback” and greatly reduced pulsation by pairing cartridges with offset occlusion.
- Low friction rollers for minimum abrasion of tubing.
- Auto-resetting limit clutch built in to protect the Pump and Drive in case of excessive overload on models 07519-15 and 07519-25.

(1) Application Data

The six- or eight-roller Cartridge Pump Systems are designed for multi-channel use where low pulsation and/or close occlusion control of a variety of different flows is required in low-flow applications. Tubing found in Table 3 can be used to deliver from .6 µL/min to 350 mL/min.

(2) General Description

Each Pump Head System (see Figure 1) is composed of a multi-roller Pump Head and two sizes of snap-in tubing cartridges, each with individually adjustable occlusion settings. The occlusion “wedge” design provides accurate and more balanced flow control.

The Pump Heads contain either six or eight rollers. All models can operate with a variety of tubing. The six roller models are compatible with the large cartridge and MASTERFLEX® L/S® thin wall tubing, and the eight roller models are compatible with the small cartridge and .9 mm wall Microbore Tubing Sets (see Tables 1, 2, and 3).


The Pump Head Systems have an off-center occlusion path that allows two oppositely mounted cartridges to be connected together for virtually pulseless flow with no suckback.

Both the six- and the eight-roller Pump Heads are available in two sizes. The short Pump Heads hold up to 2 large or 4 small cartridges and operate at the same speed as the drive. The long Pump Heads hold up to 6 large or 12 small cartridges and operate at one-fifth the speed of the drive.

Models 07519-15 and 07519-25 have an auto resetting limit clutch built in to protect the Pump and Drive in the event of an excessive overload. When an overload occurs, the clutch emits an audible clatter indicating the pump rotor is not rotating.

The Pump Head quickly mounts to most existing MASTERFLEX® L/S® Pump Drive types, (see Appendix A and B).

MODELS 07519-15 & 07519-25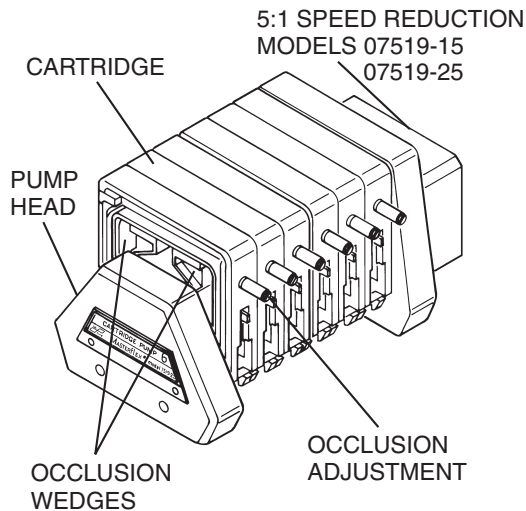

MODELS 07519-10 & 07519-20

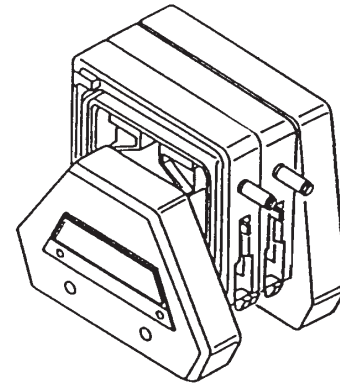
FIGURE 1 PUMP HEADS WITH CARTRIDGES

INSTALLATION

Tool Required: Long hex key (provided).

Mount the Pump Head to a MASTERFLEX® L/S® Pump Drive as follows:

1. Connect the Pump to Drive by aligning the tang on Pump Head (see Figure 2) with the slot in the motor Drive shaft.

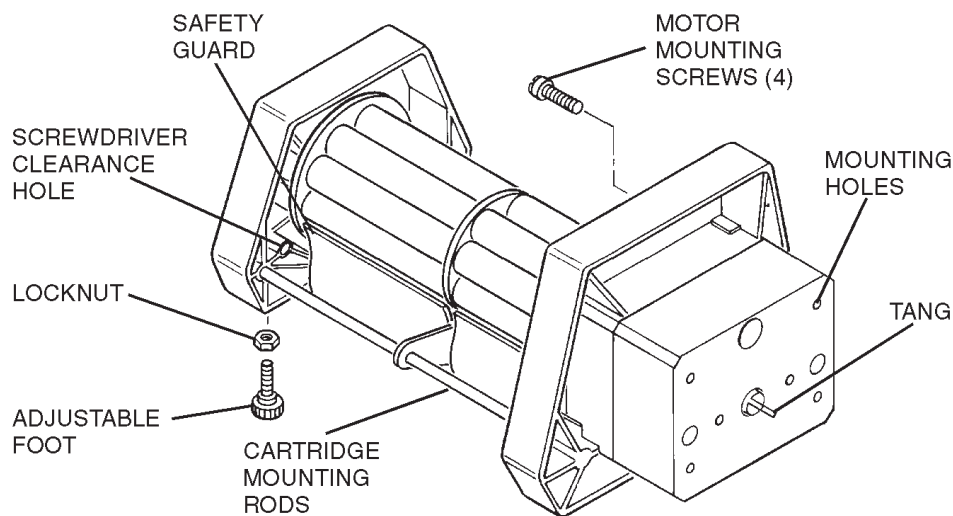
FIGURE 2 PUMP MOUNTING

2. Attach Pump Head to drive unit with four screws (provided) using the long hex key (also provided).
3. If the Pump Head requires support, install the appropriate length support screw with plastic foot and locknut (three provided). Level the Pump Head with the drive unit by adjusting the foot. Tighten the locknut against the bottom of the pump body.

SETUP

Two cartridge sizes (0.72 in and 0.36 in thick) (see Figure 3) accommodate a broad range of tubing sizes, as noted in Tables 1 and 2.

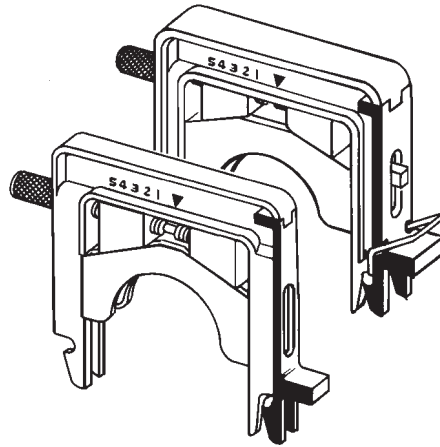
FIGURE 3 SMALL AND LARGE CARTRIDGES

(1) Select Cartridges

NOTE: These Pump Heads are designed to work with the Model 07519-75 and 07519-85 Cartridges. Do not use Model 07519-70 or 07519-80 Cartridges, which are designed for different Pump Heads.

Depending on which Pump Head is used, up to six large or twelve small tubing cartridges can be teamed up with the Pump Head. But there is a limit, depending on available drive power and pump load. For full tubing flow rates at 0 psi, Appendix B shows cartridge capacities according to tubing size and drive Type (see Appendix A for description of drive types). Cartridge sizes can be intermixed on the six-roller pump as long as the total load does not exceed the drive capability.

The small Cartridge is designed to operate with MASTERFLEX® Precision Tubing sizes L/S® 13 and L/S® 14 in either six-roller pump head and will also accommodate specially designed Microbore Tubing Sets (.9 mm wall) when used in one of the eight-roller pump heads.

The large Cartridge is designed to operate with MASTERFLEX® Precision Tubing sizes L/S® 14, L/S® 16, L/S® 25 and L/S® 17 in either six-roller pump head.

**Use only MASTERFLEX Precision Tubing and Microbore Tubing Sets with
MASTERFLEX pumps to ensure optimum performance.
Use of other tubing may void applicable warranties.**

(2) Load Cartridges

The Cartridges' occlusion surfaces are positioned asymmetrically around the pump roller assembly (see Figure 3), so orientation on the Pump Head may be important for certain applications.

To achieve parallel, synchronized pulsations and fluid "pillows" through all channels, orient Cartridges in the same direction.

To achieve nearly pulseless flow, install Cartridges in alternate directions and connect equal lengths of tubing from the discharge side of two alternate cartridges together through a Y connector as shown (see Figure 4). The asymmetric design of the cartridges provides overlapping pulses through the two tubings, which act to almost completely cancel pulsation in the combined flow channel.

NOTE: The joined flow rate will be greater than the single channel flow rate for any given tubing size and pump speed.

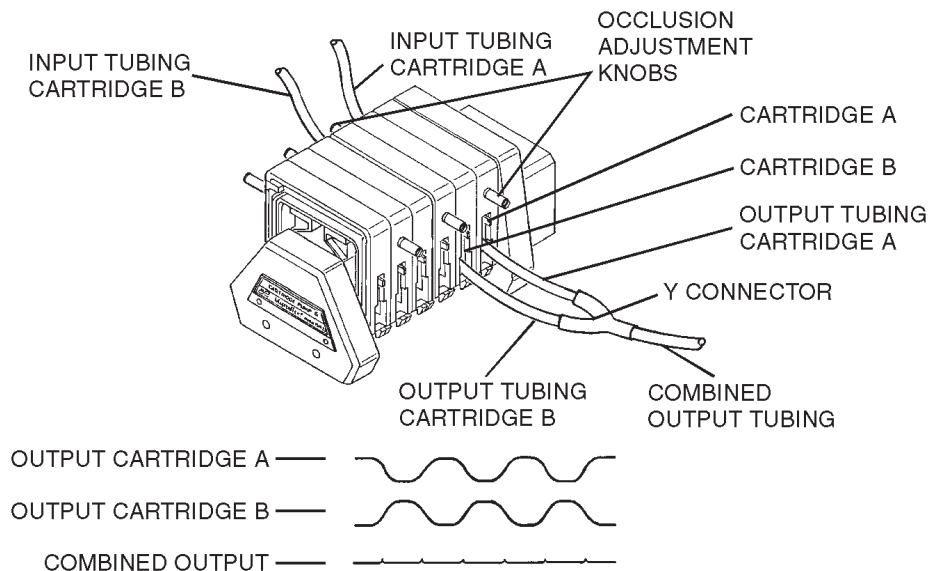
FIGURE 4 REDUCED PULSATION CONNECTION

WARNINGS: *Tubing breakage may result in fluid being sprayed from pump. Use appropriate measures to protect operator and equipment.*

Turn off drive before removing or installing Cartridges. Safety guards are provided to minimize risk of fingers getting caught between the roller mechanism and the base of the module. However, be safe—Keep your fingers away from these areas.

CAUTION: *Use only tubing/cartridge/pump combinations defined by Tables 1, 2, and 3. Use of other combinations could cause malfunction or damage to the pump.*

With the drive stopped, follow this procedure to install Cartridges.

1. Turn the Occlusion Adjustment Knob (see Figure 5) counterclockwise to maximum open. This step is not necessary when the occlusion is set at the factory setting or if the occlusion has already been set appropriately during previous pumping.
2. Set the Tubing Retainer according to tubing size (see Figure 7A) when using the large Cartridge. (See Table 4.) No retainer adjustment is required when using the small Cartridge. (**Note:** The retainer button located on the Occlusion Adjustment Knob-side of the large Cartridge is a loose component and may fall out. To reassemble, push it back into place while pressing in the button.)
3. Select tubing and place in Cartridge (see Figure 7B). If using a Microbore Tubing Set (see Figure 6), place the Collar up against the Tubing Retainer (see Figure 7C) on both sides of the Cartridge.

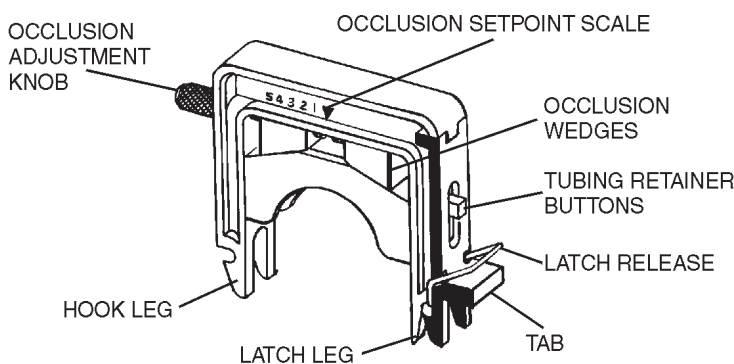
FIGURE 5 CARTRIDGE ASSEMBLY

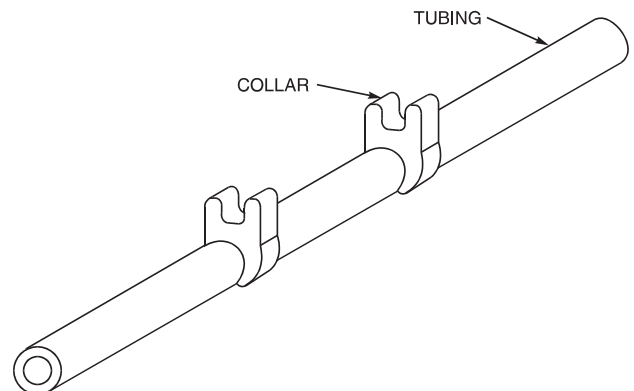
FIGURE 6 MICROBORE TUBING SET

4. Attach the Hook Leg of the Cartridge onto one of the Support Rods while holding the tubing between the Cartridge legs and against the Tubing Retainer (see Figure 7D).
5. Swing the Cartridge with the tubing over the rotor and push down on the Tab (see Figure 5) until the Latch snaps closed (see Figure 7E). **Note:** Depress the Cartridge as required in the location indicated in Figure 7E to facilitate latching and unlatching.
6. Lightly pull the tubing at the outlet and inlet to remove the slack when using continuous tubing. No slack removal is required when using a Microbore Tubing Set.
7. Adjust the occlusion setting (see OPERATION section). For a nominal setting with MASTERFLEX tubing, turn adjustment screw until inside edge of the wedge is aligned midway between #3 and #4 on the label.

(3) Cartridge Removal

Follow this procedure to unload the Cartridge.

1. Turn the Drive OFF.
2. Push down on latch side of Cartridge (see Figure 7E) and pull Latch Leg outward to unhook from rod. The Latch Release (see Figure 5) can be depressed to aid in this process as desired.
3. Pivot Cartridge upwards and unhook Hook Leg from rod and remove Cartridge.

(4) Bi-Directional Pumping

Fluids can be pumped into and back out of containers by reversing the drive. On drives that cannot be reversed, use longer tubing lengths and simply reverse the loaded Cartridge.

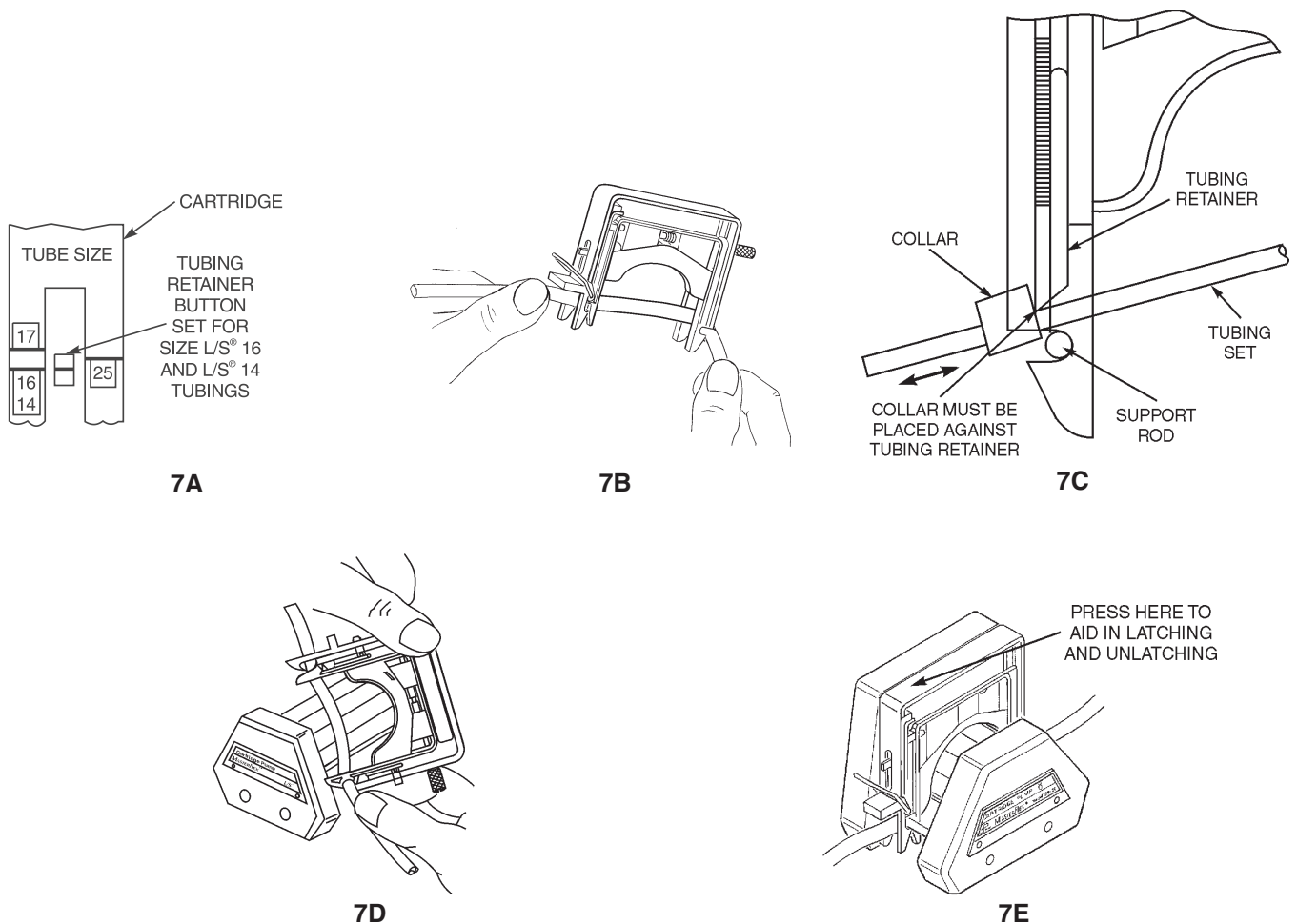
FIGURE 7 TUBING AND CARTRIDGE LOADING

(5) Partial Bank Pumping

The Pump can be operated with either a partial or a full bank of Cartridges. There is an elastic, FKM fluoroelastomer O-ring on each of the two metal rods of the pump (see Figure 8) that holds the Cartridges in place on the pump. These O-rings should be adjusted as follows. If the pump is partially loaded with Cartridges, slide the O-rings along the rods, up against the front of the outermost Cartridge, to keep it in place (note position A, for example, in Figure 8). If the pump is fully loaded with Cartridges, slide the O-rings to the front of the pump as far as possible (note position B in Figure 8). When using speed-reduced models (07519-15, -25) and the pump is less than half full of Cartridges, load the Cartridges into the front half of the pump so that the O-rings can be utilized.

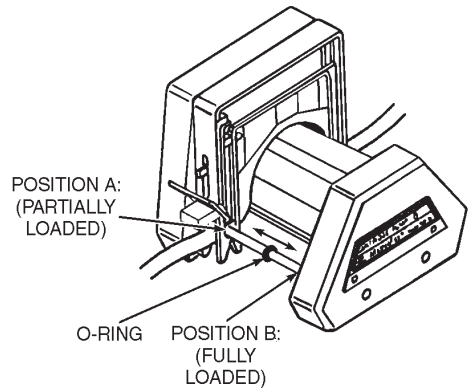
FIGURE 8 O-RING POSITION

OPERATION

This section describes the procedures for obtaining the desired performance.

(1) Relative Pulsation Rates

The relative pulsation rate in the pump output is dependent on several factors, including number of rollers on the pump rotor, cartridge size, and whether cartridges are paired.

Using the range of oscillation observed in a ball type flowmeter as a measure, approximate relative pulsation values for any given tubing size and flow rate are shown in the following chart. The chart is based on using a three-roller MASTERFLEX® L/S® pump as the benchmark with an arbitrary value of 10.

MASTERFLEX® L/S® Pump Type	Small Cartridge	Large Cartridge
6-Roller	4	4
8-Roller	3	—
6-Roller Paired Cartridges	1.5	>1
8-Roller Paired Cartridges	<1	—

(2) Select Tubing and Pump Heads

Use Tables 1, 2, and 3 to select the tubing, Cartridge, and Pump Head to provide the required single channel flow rate.

CAUTIONS: Use only tubing/cartridge/pump combinations defined by Tables 1, 2, and 3. Use of other combinations could cause malfunction or damage to the pump.

Operating the pumps at speeds in excess of 250 rpm may result in premature wear of the pump rollers and tubing.

TABLE 1 SMALL CARTRIDGE FLOW RATES—07519-85

Pump Head	Drive rpm	Pump rpm	Units	MICROBORE TUBING SET (mm ID)						MASTERFLEX® L/S® TUBING	
				.19	.25	.89	1.42	2.06	2.79	L/S® 13	L/S® 14
				FLOW RATES							
07519-25 8-Roller	1 600	.2 120	μL/min mL/min	.50 .30	1.0 .60	10 6.2	23 14	43 26	68 41	Not Compatible	
07519-15 6-Roller	1 600	.2 120	μL/min mL/min	Not Compatible						8.7 5.2	33 20
07519-20 8-Roller	1 250	1 250	μL/min mL/min	2.8 .70	5.2 1.3	52 13	120 29	220 54	340 85	Not Compatible	
07519-10 6-Roller	1 250	1 250	μL/min mL/min	Not Compatible						40 10	160 39

TABLE 2 LARGE CARTRIDGE FLOW RATES—07519-75

Pump Head	Drive rpm	Pump rpm	Units	MASTERFLEX® L/S® TUBING			
				L/S® 14	L/S® 16	L/S® 25	L/S® 17
				FLOW RATES			
07519-15 6-Roller	1 600	.2 120	μL/min mL/min	33 20	110 65	200 120	280 170
07519-10 6-Roller	1 250	1 250	μL/min mL/min	160 39	520 130	960 240	1400 350

- NOTES:** 1) Improper tubing stretch could result in reduced flow rate (see LOAD CARTRIDGE section).
 2) Models 07519-10 and 07519-20 operate at the same speed as the drive. Models 07519-15 and 07519-25 operate at 1/5 the speed of the drive.

TABLE 3 COMPATIBLE TUBING

Tubing Material	*MICROBORE TUBING SET (mm ID)			MASTERFLEX® L/S® TUBING	
	.19	.25	.89, 1.42, 2.06, 2.79	L/S® 13, 14, 16, 25	L/S® 17
Soft					
Silicone (platinum)			X	X	X
Silicone (peroxide)				X	X
TYGON® Silicone				X	X
C-FLEX® (50 A)				X	X
Stiff					
BPT (65)				X	
TYGON LFL				X	
NORPRENE® food (A 60 F)				X	
TYGON lab (R-3603)				X	
NORPRENE (A 60 G)				X	
TYGON fuel (F-4040-A)				X	
PVC	X	X	X		
FKM			X		
SANTOPRENE®		X	X		

*Use only specially designed MICROBORE TUBING SETS with Cartridges (Figure 6).

(3) Select Pump Speed

Select fixed speed drive or adjust speed of variable speed drive to provide desired nominal flow rate within the rotor rpm speed range shown for the pump model selected.

(4) Adjust Occlusion Settings on Cartridges

CAUTION: *Excessive occlusion can cause high pump temperatures and early tubing and roller failure. Do not set occlusion tighter than what is recommended in Table 4.*

NOTE: With large and small cartridges, the scale identified as “Mflex” provides nominal occlusion for MASTERFLEX® L/S® precision tubing at the #3 – #4 setting. With small cartridges only, the scale identified as “Micro” provides nominal occlusion for .9 mm (0.035 in) wall Microbore Tubing Sets at the #3 – #4 setting.

With a multi-channel cartridge system, flow rate, pressure sensitivity and tubing life can be fine-tuned—even while the pump is running. No lost time, and with operating speeds up to 600 rpm, you get higher maximum flows with the same size tubing.

Figure 9A (large and small cartridge) shows the wedges adjusted at #3 on the occlusion setpoint scale for the MASTERFLEX tubing.

Figure 9B (small cartridge only) shows the wedges adjusted at the #3 – #4 setting on the scale for the Microbore Tubing Sets.

The Occlusion Adjustment Knob, (see Figures 9A and 9B), controls a screw mechanism to move the occlusion wedges. Once a Cartridge is adjusted to a desired setting for a particular application, there is no need to reset. Or, you can record settings and quickly reproduce them at any later time.

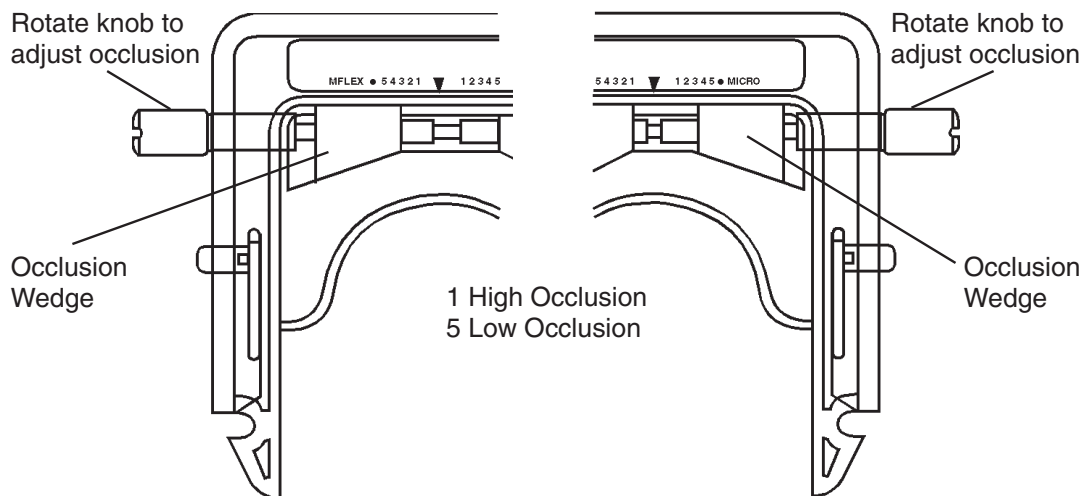
FIGURE 9A OCCLUSION INDICATIONS FOR THE MASTERFLEX TUBING

FIGURE 9B OCCLUSION INDICATIONS FOR THE MICROBORE TUBING SETS

(a) Normal Occlusion Settings

The following table shows the appropriate occlusion settings to provide satisfactory performance for various tubing sizes. (For optimum occlusion settings, see following section.)

TABLE 4 RECOMMENDED OCCLUSION SETTINGS

Tubing	Large Cartridge Mflex Scale	Small Cartridge	
		Mflex Scale	Micro Scale
Microbore Tubing Set	N/A	N/A	#3 Micro Scale
MASTERFLEX® L/S® Precision	#3 – #4 Mflex scale	#3 – #4 Mflex scale	N/A

(b) Occlusion Setting Procedure:

1. Select the recommended occlusion value from the table.
2. Turn the Occlusion Adjustment Knob to align the inside edge of the white Wedge with the scale number. (Clockwise rotation increases the occlusion.)

(c) Optimized Occlusion Settings

Some applications require additional fine-tuning of the occlusion setting to vary the flow rate for a particular tubing to reduce flow variations caused by changes in system pressure, or to increase tubing life.

1. Adjust the Occlusion Wedges as described in the preceding table.
2. Refine this setting, depending on your objectives:

NOTE: To Maximize Tubing Life

While running the pump, simply reduce the adjusted occlusion setting by turning the Occlusion Adjustment Knob counterclockwise to move the wedges to a higher setpoint scale number (toward 5). Continue turning the knob counterclockwise until the flow drop-off or the pressure sensitivity becomes unacceptable, then turn knob slightly clockwise.

NOTE: To Reduce or Eliminate Flow Drop-Off Rate with Pressure Fluctuations

While running the pump, alternate the backpressure between the expected high and low values and increase the adjusted occlusion setting. Turn the Occlusion Adjustment Knob clockwise to move the wedges to a lower setpoint scale number (toward 1) until the flow drop-off is minimized.

NOTE: To Fine-Adjust the Flow Rate

Increase or decrease the occlusion value to vary the flow.

(5) Pump Overloading

Models 07519-15 and 07519-25 have an auto-resetting limit clutch built in to protect the Pump and Drive in the event of excessive loading. When an overload occurs, the pump rotor will stop rotating and the clutch will emit an audible clatter.

The clutch capacity is sufficient to drive most combinations of tubing size and material at any pump speed with a full complement of channels operating. Some larger sizes of stiff tubings pumping through tight occlusion settings at higher speeds will cause an overload condition. This will limit the number of channels that can be run.

MAINTENANCE

No maintenance required. Wipe Pump with a clean cloth and a mild detergent. Never immerse or use excessive fluid.

Service Parts:

Hex Key (9/64 in Hex)

Models 07519-10 and 20:	A-3502
Models 07519-15 and 25:	A-4376

Mtg. Screws (8-32)

Models 07519-10 and 20:	B-1079-0407
Models 07519-15 and 25:	B-1079-0408

Leveling Feet

Models 07519-15 and 25 only:	A-4128-0002, (2 in)
	A-4128-0003, (1-1/4 in)
	A-4128-0004, (1-5/8 in)

SPECIFICATIONS

PUMP HEAD

Speed Range:	ROTOR	DRIVE
Models 07519-10 and 07519-20:	0–250 rpm	0–250 rpm
Models 07519-15 and 07519-25:	0–120 rpm	0–600 rpm
Operating Temperature Range:	0° to 40°C (32°F to 104°F)	
Storage Temperature Range:	–40°C to 60°C (40°F to 140°F)	
Noise level:	<70 dBA @ 1 meter	
Construction and Chemical Resistance:	Fiberglass-reinforced polysulfone housing; stainless steel rotor assembly with PTFE based plastic rollers, sealed stainless steel ball bearings, stainless steel structural parts and Buna N safety guard.	
Compliance:	EN809 (EU Machinery Directive)	

CARTRIDGES

Construction:	Durable polycarbonate frame with filled nylon and acetal components: filled polysulfone occlusion bed; anodized aluminum adjustment knob on stainless steel screw, FKM O-ring.
Dimensions:	
Large:	4 in H x 5 in W x 0.72 in Thick (102 mm x 127 mm x 18.3 mm)
Small:	4 in H x 5 in W x 0.36 in Thick (102 mm x 127 mm x 9.1 mm)

WARRANTY

Use only MASTERFLEX Precision Tubing and Microbore Tubing Sets with MASTERFLEX Pumps to ensure optimum performance. Use of other tubing may void applicable warranties.

The Manufacturer warrants this product to be free from significant deviations from published specifications. If repair or adjustment is necessary within the warranty period, the problem will be corrected at no charge if it is not due to misuse or abuse on your part as determined by the Manufacturer. Repair costs outside the warranty period, or those resulting from product misuse or abuse, may be invoiced to you.

The warranty period for this product is two (2) years from the date of purchase.

PRODUCT RETURN

To limit charges and delays, contact the seller or Manufacturer for authorization and shipping instructions before returning the product, either within or outside of the warranty period. When returning the product, please state the reason for the return. For your protection, pack the product carefully and insure it against possible damage or loss. Any damages resulting from improper packaging are your responsibility.

TECHNICAL ASSISTANCE

If you have any questions about the use of this product, contact the Manufacturer or authorized seller.

APPENDIX A
MASTERFLEX® L/S® PUMP DRIVE TYPES
FOR USE WITH CARTRIDGE PUMP SYSTEMS

Max rpm	Power (hp)	Torque oz-in	Drive Type
600	1/10	180	I
120	1/20	180	
600	1/20	90	II
100	1/10	360	III

NOTE: Check Max rpm and Power of the Drive or use the Torque rating to determine the Drive Type.

APPENDIX B
CARTRIDGE LOADING CAPACITY

Model 07519-20 (8-Roller, No Speed Reduction)

Microbore Tubing Set		Max. # of Cart. Drive Type		
		I	II	III
.19, .25 mm dia.	Soft	4	4	4
	Stiff	4	4	4
.89 mm dia.	Soft	4	4	4
	Stiff	4	3	4
1.42, 2.06 mm dia.	Soft	4	4	4
	Stiff	4	2	4
2.79 mm dia.	Soft	4	2	4
	Stiff	3	1	4

Model 07519-10 (6-Roller, No Speed Reduction)

MASTERFLEX® L/S® Tubing		Drive Type		
		I	II	III
13	Soft	4	4	4
	Stiff	4	3	4
14	Soft	4	4	4
	Stiff	4	2	4
16	Soft	2	2	2
	Stiff	2	2	2
25	Soft	2	1	2
	Stiff	2	1	2
17	Soft	2	1	2

Model 07519-25 (8-Roller, 5:1 Speed Reduction)

Microbore Tubing Set	Max. # of Cart. All Drives
.19, .25, .89, 1.42 mm dia.	Soft 12 Stiff 12
2.06 mm dia.	Soft 12 Stiff 10
2.79 mm dia.	Soft 12 Stiff 7

Model 07519-15 (6-Roller, 5:1 Speed Reduction)

MASTERFLEX® L/S® Tubing		All Drives
13	Soft	12
	Stiff	12
14	Soft	12
	Stiff	10
16	Soft	6
	Stiff	6
25	Soft	6
	Stiff	4
17	Soft	4

NOTE: See Tables 1, 2, and 3 for cartridge/tubing/pump compatibility.

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