## PHC745

### Basic User Manual

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



www.hach.com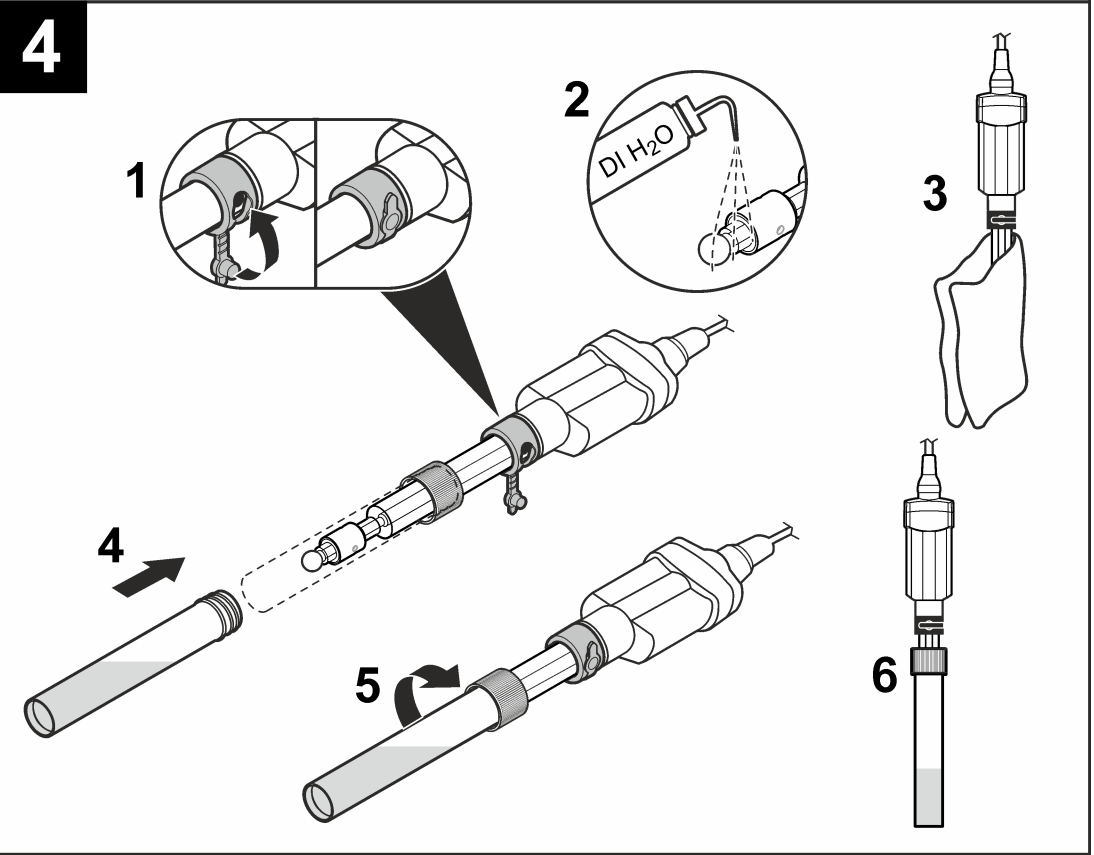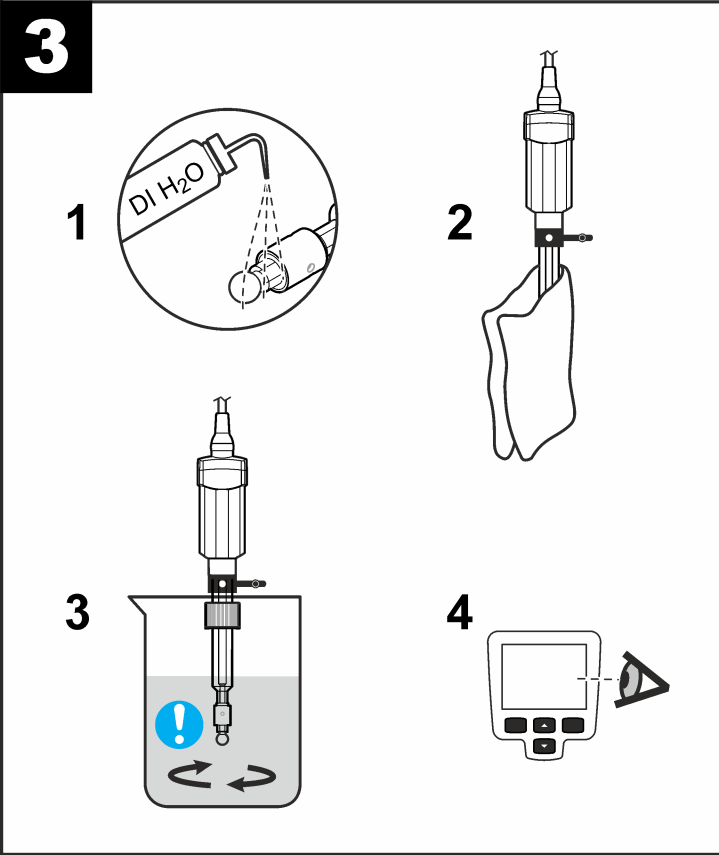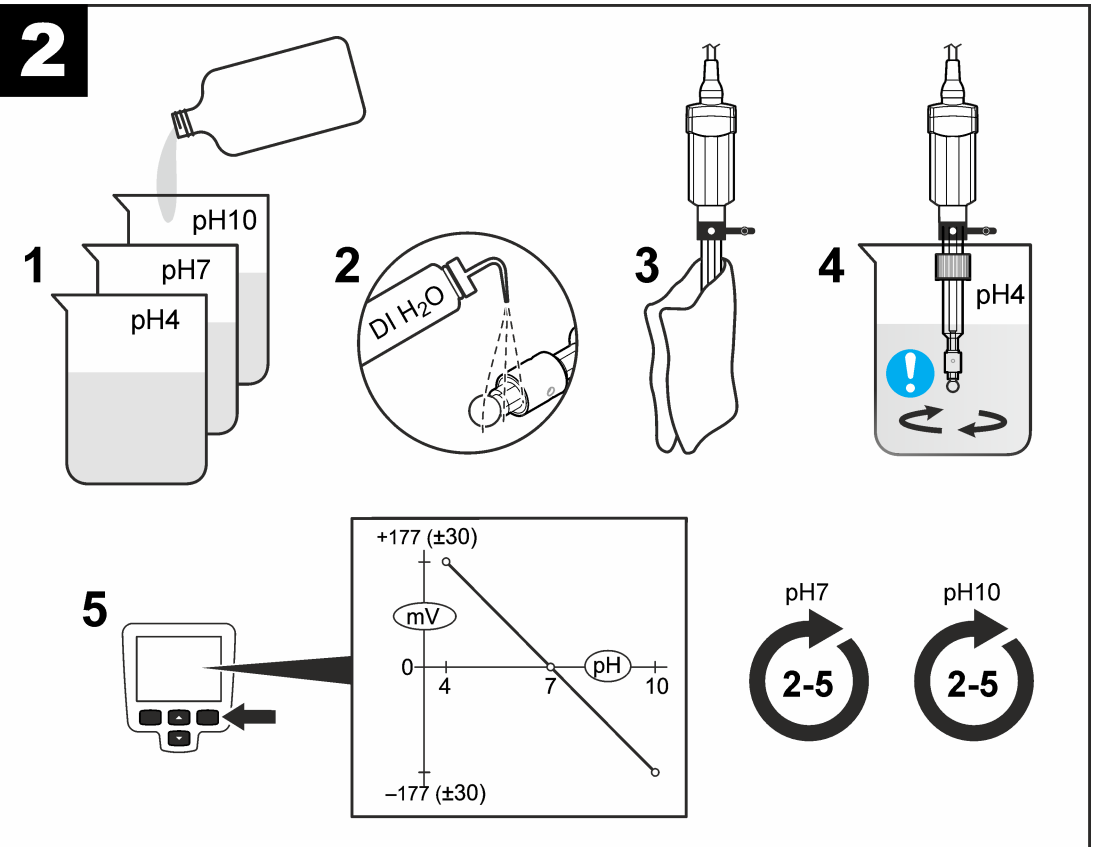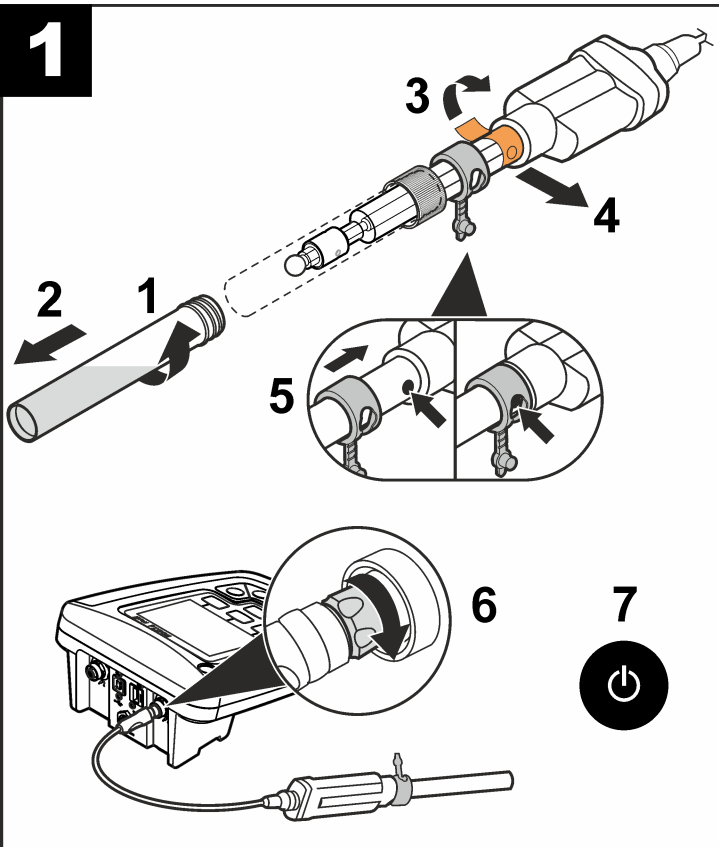

EN

**PHC805**  
Additional information is available on the manufacturer's website.

**Introduction**  
The probe is a pH electrode with refillable electrolyte, a glass body and an integrated temperature sensor. The probe has a glass sleeve junction that measures pH in wastewater, emulsions and viscous liquids.

**Warning**  
⚠ *Chemical exposure hazard. Obey laboratory safety procedures and wear all of the personal protective equipment appropriate to the chemicals that are handled. Refer to the current safety data sheets (MSDS/SDS) for safety protocols.*

**Items to collect:**  
HQd meter, pH standard solutions, beakers, DI water, no-lint cloth, filling solution

**1 Connect the probe**  
The sleeve junction must be open to let the electrolyte touch the sample.

**2 Calibration**  
⚠ *Air bubbles under the probe tip when submerged can cause slow stabilization or error in measurement. If bubbles are present, gently shake the probe until bubbles are removed.*

**3 Measurement**

**4 Storage**  
Always close the sleeve junction when the probe is not in use.

**Technical information**

pH range	pH 0 to 14
Slope	-59 mV/pH (85 to 115% at 25 °C)
Temperature	-10 to 100 °C (14 to 212 °F)
Minimum sample depth	18 mm (0.71 in.)
Filling solution	Saturated KCl
Storage solution	Saturated KCl

**Maintenance and storage**  
For maintenance and storage details, refer to the expanded version of this manual on the manufacturer's website.

**Warranty**  
1 year for manufacturing faults. This warranty does not cover improper use or wear.

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