

DWG. THA DATE 3/18/02
 BY: TOLERANCE UNLESS OTHERWISE SPECIFIED
 .X = ± .063 ANGULAR
 .XX = ± .032 HOLE DIA
 .XXX = ± .015 ± .005


SHEL-LAB
 SHELDON MANUFACTURING CORNELIUS, OREGON

18-20 MM SHAKER ASSY

SCALE: 1:6
 PART# 9751179
 MATERIAL ASSEMBLY
 REVISION A+

NOTE
 BEND RADIUS .032"
 UNLESS OTHERWISE NOTED

CHG. BY	CHG. #	REV.	DATE	DESCRIPTION

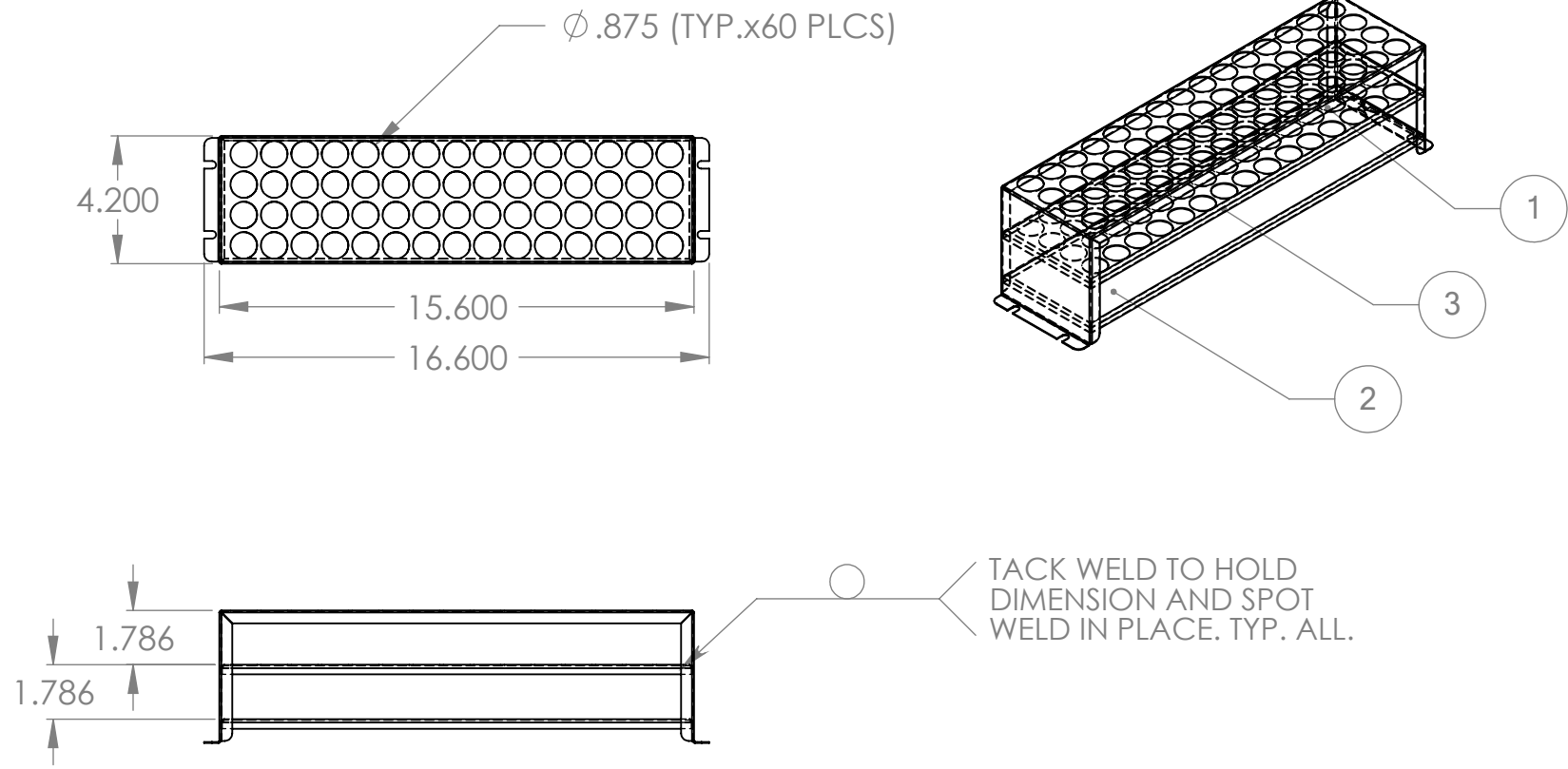

ITEM NO.	QTY.	PART NO.	MATERIAL	DESCRIPTION
1	1	5121167	.036 SST. 304-2B	SHAKER TOP 18-20 MM
2	1	5121162	20 SST 304 2B	SHAKER SHELF
3	1	5121168	20 SST 304 2B	SHAKER SHELF 18-20 MM