


STANDARD BAGS

OVER **15** OPTIONS AVAILABLE

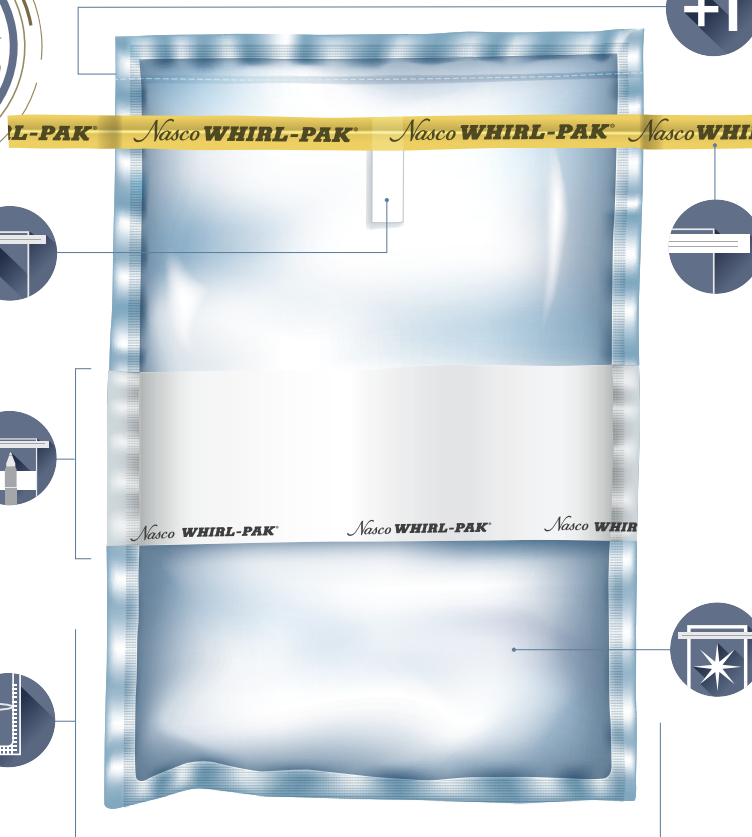

OPENING PULL TABS

- for easy opening
- non-slip textured Tyvek®

CONVENIENT WRITE-ON STRIP

- makes labeling quick and easy
- clear sample identification
- for organization & traceability

CONTINUOUS PERIMETER SEAL

- maximum protection against leaks

+1

POST-MANUFACTURING STERILIZATION

- increased sample integrity
- perforated closure to maintain sterility

TRUSTED, PUNCTURE-PROOF CLOSURE TABS

- tape extends beyond wire to eliminate sharp ends and provide a sturdy grip
- wire tabs hold the top open for easy access to add to or remove sample
- worry-free, no-leak closure for sample integrity

BLENDED, VIRGIN POLYETHYLENE FILM

- exceptional clarity
- superior strength
- sterilized after manufacturing
- FDA approved for food contact

premium quality with aseptic sampling assurance

smarter sustainability
LESS STORAGE SPACE • EASY TO SHIP • USER FRIENDLY • COST EFFECTIVE



ideal for use in food & bev, meat & poultry, dairy, water, cannabis, and many more industries.

Whirl-Pak® sterilized sampling bags are R-Nase, D-Nase, Pyrogen and BPH free and comply with regulatory and industrial requirements:





500 BAGS VS. 500 BOTTLES


GUARANTEED STERILE
Manufactured according to quality management systems ISO9001:2015

- o Uses less shelf space
- o Easier disposal
- o Durable
- o Lower shipping costs
- o Built-in closure
- o Lower total cost



SOLIDS OR LIQUIDS - LEAK-PROOF CLOSURE

STEP 1

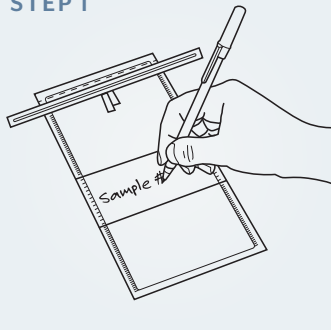

Label your sample if using a write-on bag.

STEP 2




To open, tear off the top along the perforation.

STEP 3



Use the Tyvek® pull tabs to open without contamination.

STEP 4

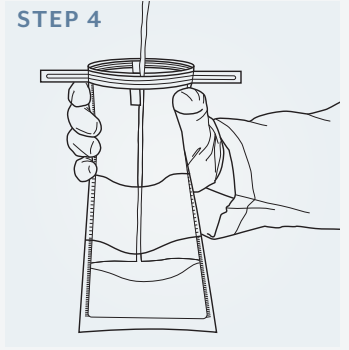


Place sample in bag. Perfect for liquids or solids.

STEP 5A




Hold bag by the tabs and whirl 3 to 4 times to make a seal.

STEP 5B

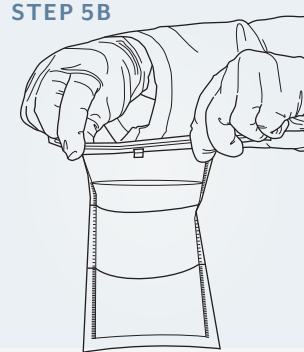


Or, fold the tab over tightly 3 to 4 times to close.

STEP 6A

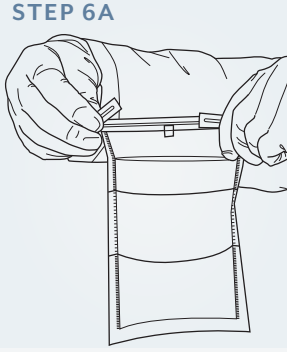

Secure closure, bend wire ends.

STEP 6B

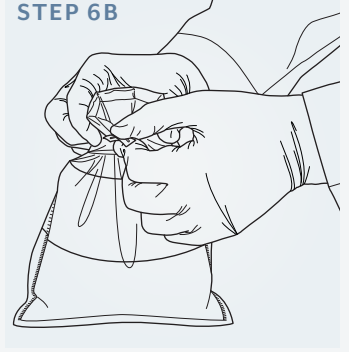

Or, bring the two wire ends together and twist tie them.



800 932 5000 | VWR.COM

Prices, product, and/or services details are current when published and subject to change without notice. | Certain products or services may be limited by federal, state, provincial, or local regulations. | VWR, part of Avantor, makes no claims or warranties concerning sustainable/green products. Any claims concerning sustainable/green products are the sole claims of the manufacturer and not those of VWR International, LLC and/or Avantor, Inc. or affiliates. All prices are in US dollars unless otherwise noted. Offers valid in US and Canada, void where prohibited by law or company policy, while supplies last. | Trademarks are owned by Avantor, Inc. or its affiliates, unless otherwise noted. | Visit vwr.com to view our privacy policy, trademark owners, and additional disclaimers. © 2020 Avantor, Inc. All rights reserved.