# Quick Start Guide

## VWR® PCR Thermal Cycler XT<sup>96</sup> / XTender<sup>96</sup>

	EU cat. no	US cat. no
XT96	732-1359	76452-151
XT96 Gradient	732-3428	76452-153
XTender96	732-3657	76452-155
XTender96 Gradient	732-3658	76452-157

## CONTENTS

1	Safety instructions	3
2	License Notice	3
3	System Overview	4
4	Installation	4
4.1	Content of delivery	4
4.2	Installing the thermal cycler	4
4.3	Operating the thermal cycler	4
5	Cycler Operation	5
5.1	Start screen = 'Run'	5
5.2	Running stored programs ('Run'):	5
5.3	Incubation mode	6
5.4	Create or edit PCR programs ('Programs')	6
5.5	Further dialogues: GLPs, Explorer, Diagnose and System	8
5.6	Further questions...?	10
6	Technical data	11
7	Maintenance and repair of the Thermal cycler	12
7.1	Cleaning, maintenance and repair	12
7.2	Temperature check up	12
7.3	Calibration	12

A more detailed instruction manual is available  
for download on our website [vwr.com](http://vwr.com).

## 1 SAFETY INSTRUCTIONS

Before the first use of the thermal cycler please read the entire instruction manual.

Special notice should be taken of the following:

Caution - dangerous voltage

Please ensure the voltage indicated on the device exactly matches your local electrical supply.

Caution - dangerous explosive material

Explosive or reactive material mustn't be heated in the thermal cycler.

Caution - liquids

Ensure that no liquids can enter the device.

Caution - hot surfaces

The thermal plate and the heated lid quickly reach temperatures above 50 °C. There is danger of burning! Keep the lid closed until the temperatures reach 30 °C or less.

Only use materials (tubes, tube caps, plates and sealing films) which are heat resistant to 120 °C.

Caution – environment

The devices ventilation system must not be covered.

Danger – line voltage

Incorrectly grounded products can be extremely hazardous in the event of a fault.

Use only a 3-conductor power cable (3x 0.75 mm<sup>2</sup>) with protective ground. The power connector may only be plugged into a socket with a protective ground. The protection must not be nullified by an extension cable without protective ground.

## 2 LICENSE NOTICE

This instrument is licensed for research and development and for uses other than human *in vitro* diagnostics under one or more of the following patents of Applera Corporation: U.S. Patent Nos. 5,656,493, 5,038,852, 5,333,675, 5,475,610 (claims 1-159 and 164-166), 6,703,236 (claims 7-10) and 7,238,517 or corresponding claims in their non U.S. counterparts. No right is conveyed expressly, by implication or by estoppels under any other patent of Applera, including but not limited to U.S. Patent No. 6,814,934 and its non U.S. counterparts, which describe and claim thermal cyclers capable of real time detection.

### 3 SYSTEM OVERVIEW

VWR thermal cyclers are the ideal PCR devices for all molecular biological and biochemical laboratories working in the field of basic research or routine diagnostic. The 250 °C HTR (High Temperature Range) Peltier technology allows quick changes in temperature up to 4 °C/sec (max). The individual monitoring and controlling of each of the eight Peltier elements per block equipped with 'Long Life Technology' ensure the tightest correlation of target and actual temperature across the entire thermal plate leading to highest reproducibility of the results. To avoid condensation each block is equipped with a heated lid.

The VWR thermal cyclers are equipped with a touch-sensitive, graphic coloured TFT display of the newest generation, allowing operation of the device by simply touching symbols and graphic elements on the screen. This leads to easy and intuitive handling and programming of the device. Additionally, a mouse can be connected via USB. Furthermore, an Ethernet port is available to connect the devices with Microsoft Windows® or Linux networks. Transfer of data can be done by using commercially available USB memory sticks enabling a nearly unlimited saving space for programs and GLP reports, which can be printed using a printer connected to the thermal cycler via the network.

**Note:** The USB ports support only standard cable mice and standard USB sticks.  
The maximum acceptable connecting cable length of the port is 3 m!

### 4 INSTALLATION

#### 4.1 Content of delivery

1x Thermal cycler

1x Power cable

1x Short manual

(A detailed instruction manual in English or German is available for download on our website [vwr.com](http://vwr.com))

1 PC software link: VWR PCR Cycler XT<sup>96</sup> software

Please check delivery for completeness and transport damage upon arrival. If any transport damage is noticed, contact your distributor or manufacturer immediately.

**NOTICE: IF THERE IS ANY DAMAGE TO THE DEVICE DO NOT USE!**

#### 4.2 Installing the thermal cycler

Check the packaging for any transport damage. Remove the packaging and place the thermal cycler onto a solid surface. The device should not be exposed to direct sunlight. There should be enough space to make sure that the ventilation grills on the front and back are not covered and sufficient ventilation is provided. The accessibility of the power switch must not be hindered. There should be a distance of at least 25 cm (10 inch) to the wall or the next device. Two thermal cyclers should not be placed back-to-back or back to front. Room temperature between 4 °C / 39 °F and 25 °C / 77°F is advised.

All original packaging should be stored in case a need arises to return the instrument to the supplier.

#### 4.3 Operating the thermal cycler

Before starting up the thermal cycler for the first time make sure that the voltage range of the device is suitable to the electricity supply. Plug in the mains plug of the thermal cycler to the electricity supply. The outlet is located at the back of the device. The serial number is written onto a typed label at the backside of the device.

The device should only be used by authorised staff. To ensure this, a user administration option with appointing rights for single users is available in the device.


## 5 CYCLER OPERATION

The operation of the thermal cycler can be run by the internal touch screen, either using the touch-sensitive surface of the display or by using a mouse connected to a free USB port.

If the user must enter numbers or text the corresponding field has to be touched on the touch screen or clicked with the mouse. An appropriate keyboard will appear on the screen according to the expected entry.


### 5.1 Start screen = 'Run'

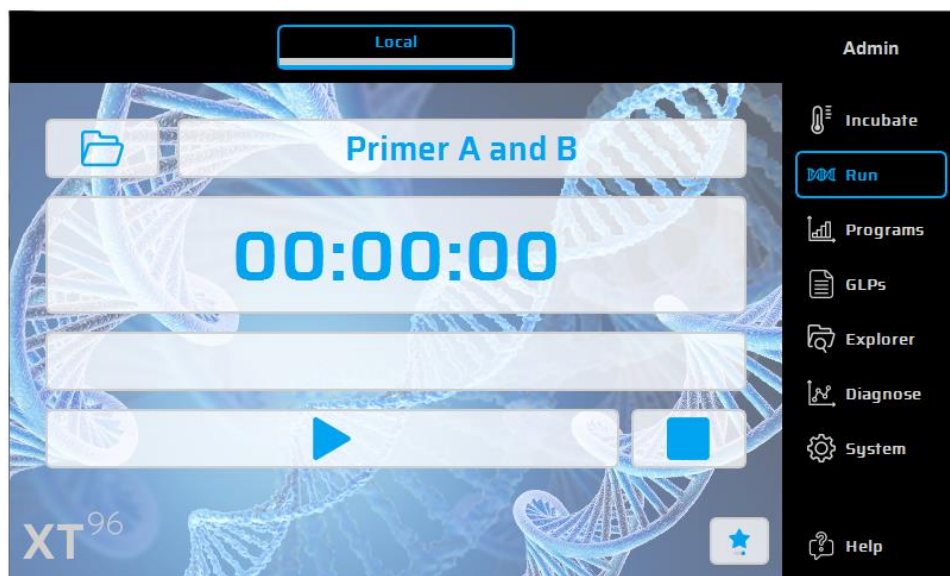The last PCR program run is displayed in the 'Run' dialogue and can directly be restarted with the 'Start PCR' button.

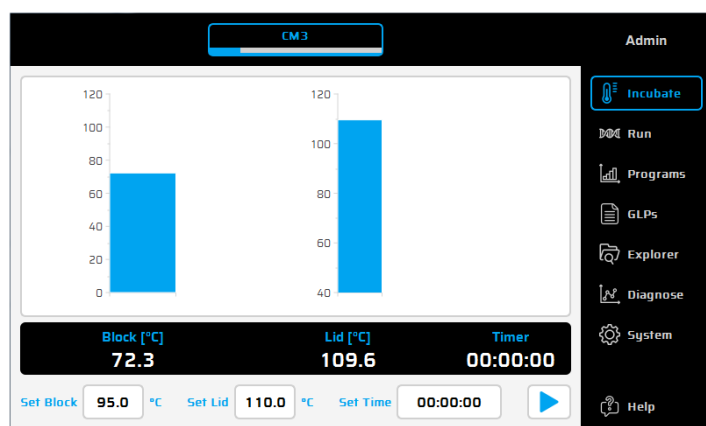
### 5.2 Running stored programs ('Run'):

1. Open the selection list in the 'Run' option using the  button.
2. Select the recently used program in the selection list and load the program with the 'Open Recent Program' button. Alternatively use the Program Wizard for new Programs or the File Explorer to load stored programs.
3. Start the loaded program by using the  button.

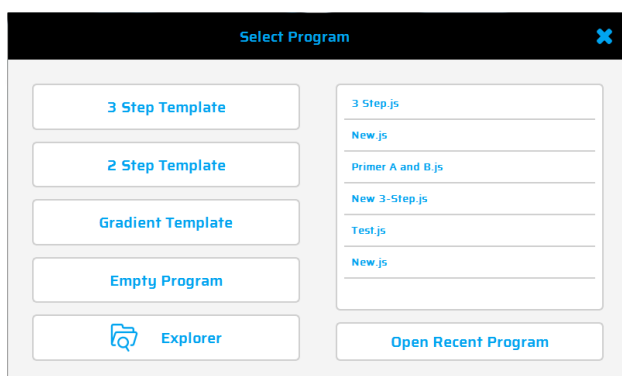
### 5.3 Incubation mode

Use the  **Incubate** button to set the device to incubation mode. The temperature for the thermal plate and lid can be set manually. The clock shows the time elapsed in the incubation mode or can be used as timer.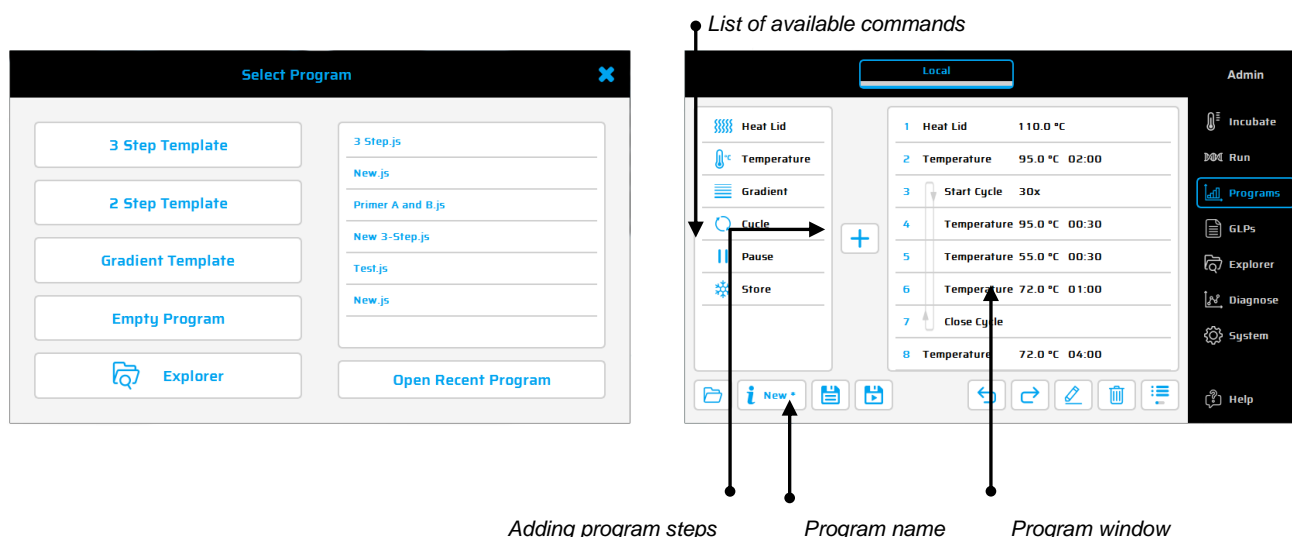

### 5.4 Create or edit PCR programs ('Programs')

1. To create or edit PCR programs select the 'Programs' option.
2. By touching the 'Programs' option for the first time during a session, the selection dialogue shows where a new program can be selected  or an existing program can be opened for editing.
3. On the left side of the program editor all available commands are displayed. On the right side of the editor a list of commands (Steps) of which the program is composed of is shown. After selecting a program command from the list in the left window it can be inserted into the desired place using the  button.
4. Using a double click or the 'Edit' button  parameters, single steps can be changed later.
5. Selected program steps can be deleted from the program by using the 'Delete' button .
6. Use the 'Save' button  to save a program.
7. Save and start programs in one step by simply using the 'Save&Run' button .
8. For even faster programming the VWR thermal cycler comes with a selection of template programs, which can easily be edited and saved under a new program name.

## Available commands:

 Heat lid	<b>Heat lid</b> Turns the heated lid on/off.	 Cycle	<b>Cycle</b> Combines Open Loop, Close Loop.
 Temperature	<b>Temperature</b> Temperature step for definable time, Temperature and Time Incr./Decr. function.	 Pause	<b>Pause</b> Interrupts the program temporarily at defined temperature settings.
 Gradient	<b>Gradient</b> Realises different temperatures at the same time for protocol optimisation.	 Store	<b>Store</b> Holds the defined temperature, turns off the heated lid automatically.

## Sample program

To create a new program, click on 'Empty Program', '3 Step Template', '2 Step Template' or 'Gradient Template'. Enter required values and confirm with **OK** to access the programming functions.

To enter a new command into a program the desired line has to be selected. After selecting the program command from the list in the left window it can be inserted above the desired place via the '+' button. Alternatively, the command can be inserted by double clicking on the program command. Once a command is added the corresponding parameters will need to be adapted. To enter a parameter the desired input field has to be selected

and click on  or by double click.

## Gradient Technology

The temperature of each of the eight rows of the block can be individually controlled!

The temperature range is defined by the 'Center' Temperature (55°C) and 'Gradient  $\pm$ ' Temperature (5°C). The gradient is always shown from the top to the bottom over the 8 rows of the block.

## Program options

When setting up a new program or when saving during programming (via the Info button)  program options can be set. The following options are available for the user:

### 1. Power Fail Denaturation

If a power failure happens while a program is running the thermal cycler will restart the program at the point it was stopped as soon as the power returns. If the 'Power Fail Denaturation' option is activated, the thermal block will heat to the stated temperature for the set time before continuing with the program.

### 2. Global Program Ramp

By using this function a global ramp rate in °C/s can be programmed. This global ramp rate will be used in all temperature and gradient steps of the PCR program but can be changed in a single temperature or gradient step manually by entering the desired ramp value at the step.


### 3. Tube Control

Using tube control gives you the ability to adjust temperature regulation for a temperature step according to sample volume. If necessary, the temperature regulation will perform short-term deliberate overshoots and undershoots depending on the entered sample volume during PCR run. Especially if combined with the emulation mode, this function can be helpful if the emulated cycler uses the tube control function before too.

## 5.5 Further dialogues: GLPs, Explorer, Diagnose and System

**GLPs:** Displaying and organising GLP (Good Laboratory Practice) reports created by the thermal cycler. This option is mainly for labs that need to save GLP reports for documentation purposes.


**Explorer:** The 'Explorer' dialogue is used for the data management with the thermal cycler.

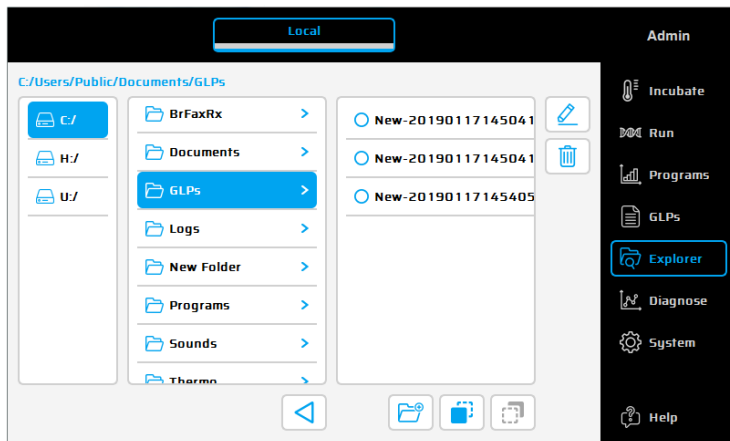**Diagnose:** Monitoring the actual temperatures of each of the eight Peltier elements during a run. In the event of an instrument malfunction have a look at the Peltiers' temperature profile.

**System:** Configuration of the system and the general changing device settings.

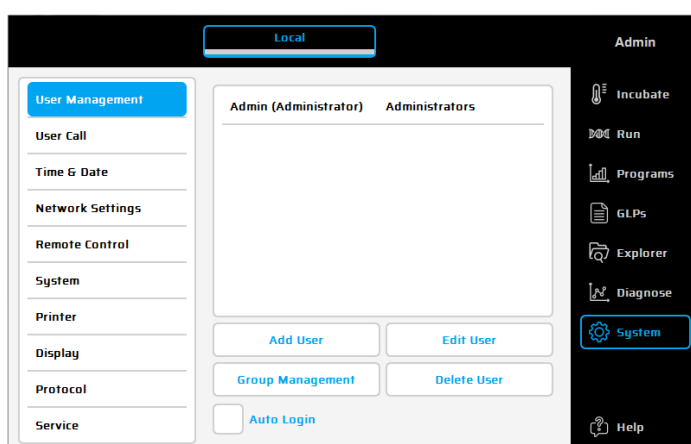## Function buttons

<b>User Management</b>	<b>User management</b> Defines user and user groups (plus their access rights). The default password for the user 'Administrator' is: 'admin'. Via 'Logout' button (upper right corner of the screen) the current user can be changed.	<b>Remote Control</b>	<b>Remote Control</b> For PC and Master/Slave control.
		<b>System</b>	<b>System</b> Shows the serial number of thermal cycler, the Software version and if the gradient function is loaded.
<b>User Call</b>	<b>User call</b> Setting of acoustic signals (Pause, Store, Program End).	<b>Printer</b>	<b>Printer</b> For Printer connection via network
<b>Time &amp; Date</b>	<b>Time / Date</b> Using this button, the time and date can be set.	<b>Display</b>	<b>Display</b> To adjust the backlight of the display
<b>Network Settings</b>	<b>Network Settings</b> Displays/ Changes network settings.	<b>Service</b>	<b>Service</b> Service function (only for authorised technical service)

## 5.6 Further questions...?

Direct help :

Use this button to activate (deactivate) the direct help mode. If the mouse pointer is visible a question mark appears alongside. By touching a function button or other dialogue element the corresponding help window will appear.

A more detailed instruction manual is available for download on our website [vwr.com](http://vwr.com).

## 6 TECHNICAL DATA

### General characteristics

- 8 Peltier elements with Long-Life-technology and 8 control circuits with Pt 1000 temperature sensors
- Maximal heating and cooling rate: 4 °C/s
- Block uniformity (at 72 °C):  $\pm 0.35$  °C
- Temperature range thermo block: 4 to 100 °C
- Control accuracy thermos block:  $\pm 0.2$  °C
- Variable ramping: 0.1 to 3.0 °C/s
- Increment/Decrement time: 0:00:01 to 0:59:59
- Increment/Decrement temperature: 0.1 to 3 °C
- Lid heating temperature range: 40 to 115 °C
- 2 x USB, 1 x Ethernet interfaces (MS Windows® or Linux)  
Note: USB interfaces only support standard USB sticks, cord mouse and keyboards!
- Dimensions (B x H x D): 22 x 19 x 33 cm
- Weight incl. block: 8.9 kg
- Pollution rate: 2
- Environmental temperature range: +10 °C to +30 °C
- Max. relative humidity: 70 %
- Maximum height above sea level: 2000 m
- Noise level: < 37 dB (A)

### Gradient feature (optional)

- Maximal gradient range over 8 columns, respectively: 30 °C ( $\pm 15$  °C)
- Temperature range Gradient: 35 to 100 °C
- Gradient accuracy:  $\pm 0.2$  °C

### User interface/functions

- Touch sensitive (capacitive) TFT-Display (800 x 480)
- Direct help function
- Internal flash memory for 500.000 typical PCR programs in free configurable folders/subfolders. Unlimited number of programs via network PC or USB stick
- Free PC software for the generation of PCR programs on the computer
- Unlimited number of steps per program
- Password protected user accounts with variable access rights
- GLP reports for complete recording of all runs
- Quick start function of the last used program
- Automatic restart after power failure, Power-Fail-Denaturation, Instant Incubation
- Remote control and monitoring of instruments via PC-software
- MP3 sounds free of choice
- Master-/Slave-control

### Block capacity

- 96 well Universal block with lid heating and automatic height adaption for 96 x 0.2 ml tubes and 96 well PCR plates

### Power supply

EU/UK Version: 220 - 240 V AC, 50/60 Hz, max. 860 VA

US Version: 120 V, 60 Hz, max. 900 VA

For indoor use, only!

## 7 MAINTENANCE AND REPAIR OF THE THERMAL CYCLER

### 7.1 Cleaning, maintenance and repair

For safety reasons the device must be switched off and the mains plug pulled out before cleaning is carried out! The surface of the thermal cyclers can be wiped off with a damp cloth or ethanol. Do not use aggressive or scouring cleaners or organic solvents for cleaning. The device should be protected from aggressive chemicals. Make sure that no liquid penetrates the interior of the device.

If sample liquids penetrate into the drill holes of the thermal cyclers clean them immediately with mild soap followed by distilled water. You can use mild disinfectants for decontamination.

Maintenance should be carried out regularly to detect possible technical failures.

The electric fuses are situated at the back of the device between the power switch and the mains plug. After pulling out the mains plug and pushing the safety catch upwards the fuses can be changed. Use only fuses with correct values (indications about the fuse type are situated at the back of the device).

The device should only be opened by a qualified specialist. Unauthorised work on the device voids the warranty. All repairs should only be carried out by authorised persons using original replacement parts. The replacement of single components of the device should not be done by the user but exclusively by authorised specialists. Please therefore send the device back to the manufacturer. The surface of the device should be decontaminated with a lint-free cloth soaked with 70% ethanol before sending. The drill holes of the thermo block should be cleaned with a 1,5% hypochlorite solution with the help of a cotton bud. Please fill in and sign the decontamination certificate and send it to the manufacturer together with the device.

### 7.2 Temperature check up

Maintenance of the temperature at regular intervals (e.g. annually) is recommended. The thermometer should be equipped with wired sensing devices (e.g. PT1000) and should have an accuracy of 1/10 class B ( $\Delta T = \pm 0,03\text{ }^{\circ}\text{C}$ ) measurements in the range of 0 to +100 °C. While processing the temperature control room temperature should not exceed +25 °C.

### 7.3 Calibration

If there is a reasonable suspicion that the device has deviated from its calibration range, because the variations of the temperature check up are considerably out of the acceptable tolerance, a new calibration must be carried out by the manufacturer. Please contact your local distributor.

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The screenshot shows the 'Gradient' control window. On the left, a vertical list of temperatures from 50.0 to 60.0 is visible. The main control area has three input fields: 'Temp (°C)' set to 55.0, 'Gradient (°C)' set to 5.0, and 'Ramp (°C/s)' set to 0.0. Below these are time units: 'Hours' (0), 'Minutes' (0), and 'Seconds' (30). A numeric keypad with digits 0-9, a decimal point, and a minus sign is on the right, along with a blue 'ok' button. A back arrow is also present next to the keypad.

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