

VWR® Automated cell counters

01. SIMPLICITY

02. ACCURACY

03. AFFORDABILITY

04. TIME SAVINGS



SAMPLE
PREPARATION
AND LOAD

INSERT
COUNTING
SLIDE

CELL
COUNTING

PROFILE
SETTINGS


SAVE
RESULTS

VWR® Automated cell counters

Save time and reduce variability with the VWR® automated cell counter and VWR® automated cell counter Fluo. Intelligent image recognition and advanced optical imaging technology ensure accurate cell analysis as well as cell concentration and size determination for a wide range of cell sample types. These machines pack a variety of features into a compact design made to fit virtually any workspace.



Simplicity:

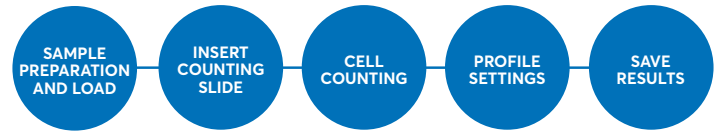


- **Accuracy:** Advanced algorithm enables fast recognition of clustered cells and flexible parameters setting include circularity, diameter and brightness will exclude debris and impurities
- **Affordability:** Brightfield Cell counter or Upgradable Fluo Cell counter compatible with affordable single use counting slides
- **Time savings:** Features like rapid capture and autosave quickly generate and save your results

SPECIFICATIONS

- Wide application: cell counting, cell viability, AO/PI staining, cell transfection efficiency, cell expression
- All-in-one machine with a compact design
- Dual-channel option for the VWR® automated cell counter Fluo
- Three interchangeable fluorescent cubes (green, blue, red)
- Cell counting slide consumables with two chambers
- Accurate cell identification in only 9 seconds
- User-friendly interface
- Adjustable parameters (focus, brightness, cell size range, roundness range)
- Data recording system can save more than 1000 counting reports

Specifications	VWR® Automated Cell Counter	VWR® Automated Cell Counter Fluo
Brightfield	Yes	Yes
Fluorescence option	No	Yes - upgradable (two interchangeable fluo cube channels)
Fluo cube colors	No	Green, Red, Blue
Counting time manual	5s'	5s'
Counting time auto focus	9s'	9s'
Imaging parameters	Auto focusing & Manual focusing / 5 megapixels, 2.5x Optical Magnification	Auto focusing & Manual focusing / 5 megapixels, 2.5x Optical Magnification
Compatible Dyes	Trypan blue	Trypan blue, AO, PI, GFP, DAPI
Application	Cell counting, Cell viability, Cell size - brightness and circularity	Cell counting, Cell Viability, Cell size - brightness and circularity, Transfection efficiency, AO/GFP, DAPI, PI cell RFU
Counting clumps, not debris	Yes	Yes
Counting range (cell density)	10 ⁴ -10 ⁷	10 ⁴ -10 ⁷
Cell size range (µm)	4-60	4-60
Sample volume (µl)	10	10
Pre-dilution calculator	Yes	Yes
Cell splitting calculator	Yes	Yes
Gating	Yes	Yes
Cell imaging and histogram	Yes	Yes
Disposable consumable	2 samples slide	2 samples slide
Counting PBMCs	Yes	Yes
CV (Hela Cell) - accuracy	< 5%	< 5%
Calibration	No calibration	No calibration
Display	7" Capacitive touch screen	7" Capacitive touch screen
IT compatibility	Double USB	Double USB
Output data	JPG, TIF, PDF, CSV, BMP, PNG	JPG, TIF, PDF, CSV, BMP, PNG
Built-in memory	1,000 counting reports	1,000 counting reports
Standalone	Yes	Yes
Ordering information	49862-2000	49893-2000


VWR® AUTOMATED CELL COUNTERS INTERFACE

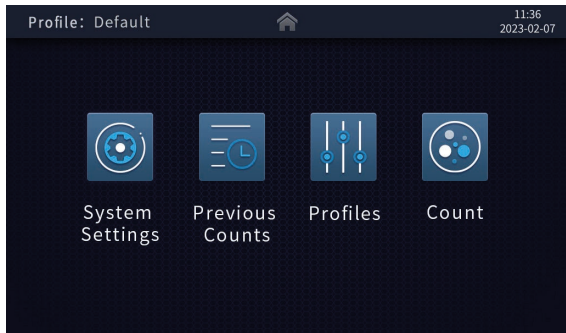

MAIN INTERFACE

VWR® automated cell counters feature a user-friendly interface including easily accessible icons for:

- System Settings
- Previous Counts
- Profiles
- Count

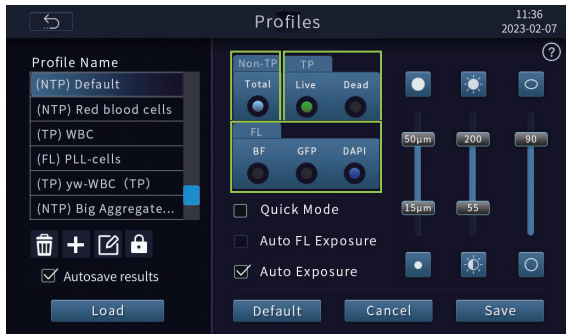

MODE SELECTION


VWR® automated cell counters have three different modes :

- **[Non-TP] mode:** Count total cell concentration, obtain the total cell size distribution histogram and average diameter
- **[TP] mode:** Count total cell concentration, dead/live cell concentration and proportion, obtain the size distribution histogram and average diameter of dead/live cells
- **[FL] mode:** Count total cell concentration, different fluorescent cell concentration and proportion, obtain the total cell size distribution histogram and RFU distribution




SETTINGS

Non-TP and TP mode

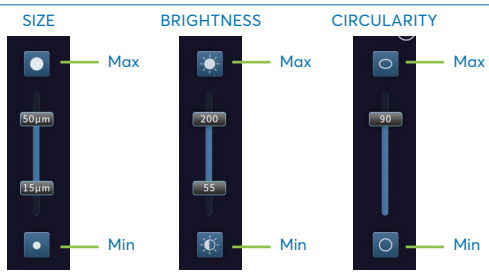

FL mode

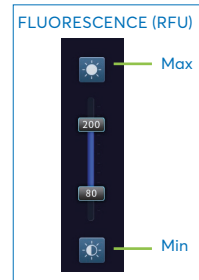

ORDERING INFORMATION

Description	Display	Pk	Cat. No.
Cell counters			
VWR® automated cell counter	7" capacitive touch screen, Brightfield	1	49862-2000
VWR® automated cell counter Fluo	7" capacitive touch screen, Brightfield, fluorescent	1	49893-2000
Consumables for VWR® automated cell counters			
VWR® cell counting slide (2 samples/slide)		50	10228-0050



INTERCHANGEABLE FLUO CUBES

VWR® automated cell counter Fluo dual channel:

detachable building block design enables easy installation and replacement of the fluorescent cubes. Fluorescent cubes sold separately.

Description	Compatibles Dyes	Excitation (nm)	Emission (nm)	Pk	Cat. No.
Fluo cubes					
VWR® Fluo cube blue	DAPI, Hoechst, BFP	375/28 nm	460/50 nm	1	49893-4950
VWR® Fluo cube green	Acridine orange(AO), Calcein-AM, Alexa Fluor 488, SYBR green, FITC, GFP	480/30 nm	535/40 nm	1	49893-4951
VWR® Fluo cube red	Propidium iodide(PI), DsRed, dTomato, RFP	540/25 nm	620/60 nm	1	49893-4952



Setting science in motion to create a better world



AUSTRIA

VWR International GmbH
Graumannsgasse 7
1150 Wien
Tel.: +43 1 97 002 0
info.at@vwr.com

BELGIUM

VWR International bv
Researchpark Haasrode 2020
Geldenaaksebaan 464
3001 Leuven
Tel.: +32 (0) 16 385 011
vwr.be@vwr.com

CZECH REPUBLIC

VWR International s. r. o.
Veetee Business Park
Pražská 442
CZ - 281 67 Stříbrná Skalice
Tel.: +420 321 570 321
info.cz@vwr.com

DENMARK

VWR International A/S
Tobaksvejen 21
2860 Søborg
Tel.: +45 43 86 87 88
info.dk@vwr.com

FINLAND

VWR International Oy
Valimatie 9
00380 Helsinki
Tel.: +358 (0) 9 80 45 51
info.fi@vwr.com

FRANCE

VWR International S.A.S.
Le Périgares – Bâtiment B
201, rue Carnot
94126 Fontenay-sous-Bois cedex
Tel.: 0 825 02 30 30* (national)
Tel.: +33 (0) 1 45 14 85 00 (international)
info.fr@vwr.com

* 0,18 € TTC/min + prix appel

GERMANY

VWR International GmbH
Hilpertstraße 20a
D - 64295 Darmstadt
Tel.: 0800 702 00 07* (national)
Tel.: +49 (0) 6151 3972 0 (international)
info.de@vwr.com
*Freecall

HUNGARY

VWR International Kft.
Simon László u. 4.
4034 Debrecen
Tel.: +36 52 521130
info.hu@vwr.com

IRELAND

VWR International Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 (0) 1 88 22 222
sales.ie@vwr.com

ITALY

VWR International S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: +39 02 3320311
info.it@vwr.com

THE NETHERLANDS

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: +31 (0) 20 4808 400
info.nl@vwr.com

NORWAY

VWR International AS
Brynsalleen 4,
0667 Oslo
Tel.: +47 22 90 00 00
info.no@vwr.com

POLAND

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: +48 58 32 38 200
info.pl@vwr.com

PORTUGAL

VWR International - Material de
Laboratório, Lda
Centro Empresarial de Alfragide
Rua da Indústria, nº 6
2610-088 Amadora
Tel.: +351 21 3600 770
info.pt@vwr.com

SPAIN

VWR International Eurolab S.L.U.
C/ Tecnología 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: +34 902 222 897
info.es@vwr.com

SWEDEN

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: +46 (0) 8 621 34 00
kundservice.se@vwr.com

SWITZERLAND

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: +41 (0) 44 745 13 13
info.ch@vwr.com

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard - Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: +44 (0) 800 22 33 44
uksales@vwr.com

CHINA

VWR (Shanghai) Co., Ltd
Bld.No.1, No.3728 Jinke Rd,
Pudong New District
Shanghai, 201203- China
Tel.: 400 821 8006
info_china@vwr.com

ÍNDIA

Avantor Performance Materials India
Limited
17th Floor, Building No. 5, Tower C
DLF Cyber City Phase - III
Gurgaon - 122002, Haryana
Tel.: +91-1244-65-6700
help@avantoric.com

KOREA

VWR International -
17, Daehak 4-ro, Yeongtong-gu,
Suwon-si, Gyeonggi-do
Tel.: +82 31 645 7256
saleskorea@avantorsciences.com

MIDDLE EAST & AFRICA

VWR International FZ-LLC
Office 203, DSP Lab Complex,
Dubai Science Park,
Dubai, United Arab Emirates
Tel.: +971 4 5573271
Info.mea@vwr.com

SINGAPORE

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel.: +65 6505 0760
sales.sg@vwr.com

GO TO [VWR.COM](https://www.vwr.com) FOR THE LATEST NEWS, SPECIAL OFFERS AND DETAILS FROM YOUR LOCAL VWR SUPPORT TEAM