



## Protocol for Plasmid DNA Purification from Cultured *Escherichia coli* Cells Using Pall Nucleic Acid Binding Nanosep<sup>®</sup> Centrifugal Device

### 1. Consumables and Reagents

**Table 1***Consumables for pDNA Purification*

Supplier	Product Description	Part Number
Pall Laboratory	Nucleic Acid Binding (NAB) Nanosep Centrifugal Device	76360-454, 76360-456
VWR	Ethanol (not denatured)	71001-866
VWR	Spectrophotometer Cuvettes ~100 $\mu$ L (260/280 nm)	47743-840
VWR	Tubes 15 mL (RNase-DNase free)	21008-103
VWR	Microcentrifuge Tubes 1.5 mL (RNase-DNase free)	76005-210

**Table 2***Reagents for pDNA Purification*

Supplier	Product Description	Part Number
Various	Resuspension Buffer	N/A
Various	Lysis Buffer	N/A
Various	Neutralization Buffer	N/A
Various	Binding Buffer	N/A
Various	Wash Buffer	N/A
Various	RNase A (100 mg/mL)	N/A
Various	Nuclease-free Water	N/A
VWR	Tris Buffer pH 7.0 (1M)	97062-936

### 2. Instruments

- Microcentrifuge
- Spectrophotometer
- Vortex

### 3. Important Points Before Starting

- Add the provided RNase A solution to resuspension buffer prior to use. Mix and store at 2 – 8 °C
- Add Ethanol (not denatured) to the wash buffer prior to use. See bottle label for instructions.
- Check commercially supplied buffers before use. If precipitate is observed, heat to 37 °C and gently agitate – Do not vigorously shake the lysis buffer bottle.
- Ensure that the lid of the lysis buffer is closed immediately after use to avoid acidification of buffer from CO<sub>2</sub> in the air.
- All centrifugation steps are performed at room temperature at 10,000 – 14,000 x g.
- For each NAB Nanosep device insert there are three receiver tubes. This is enough to complete the below process. Use only the receiver tubes provided with the NAB Nanosep device.
- All buffers should be allowed to equilibrate to room temperature before use.
- Briefly centrifuge tubes after vortexing to remove drops from inside the lid.
- Centrifugation at full speed will not affect yield or purity of the DNA. If, after centrifugation the sample has not completely passed through the membrane, centrifuge again until all the solution has passed through.

### 4. Protocol

*This protocol is designed for purification of plasmid DNA from 1 – 5 mL (overnight) cultures of E. coli in rich-medium*

1. Harvest the bacterial cells by centrifugation at 3,000 x g in a table-top microcentrifuge for 10 minutes at room temperature.
2. Discard the supernatant and resuspend pelleted bacterial cells in 250 µL of resuspension buffer (with added RNase A) without creating cell clumps. Transfer to a microcentrifuge tube.
3. Add 250 µL lysis buffer and invert the tube 4 – 6 times to mix. Avoid vortexing as this can shear the DNA. If necessary, continue inverting until the solution becomes viscous and slightly clear. Do not allow the lysis to continue for more than 5 minutes.
4. Add 350 µL of neutralization buffer and mix immediately by inverting 4-6 times. Avoid localized precipitation by mixing thoroughly, immediately after the addition of Buffer N3.
5. Centrifuge for 5 – 10 minutes at 10,000 – 14,000 x g.
6. Transfer 500 µL of the supernatant to the NAB Nanosep device at a time and centrifuge for 45-60 seconds at 10,000 – 14,000 x g discarding the flow through. Retain the receiver tube and repeat this step with the remaining supernatant.
7. Add 500 µL of binding buffer to wash the NAB Nanosep device, centrifuging for 60 seconds at 10,000 – 14,000 x g. Discard the flow-through and retain the receiver tube for the next step.
8. Perform a second washing step by adding 500 µL of wash buffer and centrifuging for 60 seconds at 10,000 – 14,000 x g.
9. Discard the flow-through and centrifuge again at full speed for an additional 1 minute to remove any residual buffer.

10. Transfer the NAB Nanosep device to a clean receiver tube (provided) and add 50  $\mu\text{L}$  of nuclease-free water (or elution buffer) to the center of the membrane of each NAB Nanosep device. Leave to stand for 1 minute.
11. Centrifuge for 60 seconds at 10,000 – 14,000 x g and recover the purified plasmid DNA in the receiver tube.

### Storage of DNA

Purified plasmid DNA can be stored in nuclease-free water at  $-20\text{ }^{\circ}\text{C}$  for several months.

### Quantification of DNA

DNA concentration can be determined by measuring the absorbance at 260 nm ( $A_{260}$ ) in a spectrophotometer (see details below). For lower concentrations of DNA however, it can be difficult to determine these amounts photometrically. Smaller quantities of DNA can be accurately quantified using fluorometric quantification.


### Spectrophotometric quantification of DNA

$A_{260}$  readings should be greater than 0.10 and lower than 1.0 to ensure significance. An absorbance reading of 1.0 at 260 nm corresponds to 50  $\mu\text{g}$  of DNA per mL. This is only valid for measurements at neutral pH however. As a result, if it is necessary to dilute the DNA sample, ensure that the dilution buffer is of neutral pH.




Laboratory

Ordering:  
vwr.com

© 2020 Pall Corporation. Pall , and Nanosep are trademarks of Pall Corporation. ® indicates a trademark registered in the USA.  
**Filtration. Separation. Solution.** is a service mark of Pall Corporation.

4/20, PDF, GN19.1003

**Filtration. Separation. Solution.<sup>SM</sup>**

 **VWR**<sup>™</sup>  
part of avantor

800 932 5000 | [VWR.COM](http://VWR.COM)

Prices and product details are current when published and subject to change without notice. | Certain products may be limited by federal, state, provincial, or local regulations. | VWR, part of Avantor, makes no claims or warranties concerning sustainable/green products. Any claims concerning sustainable/green products are the sole claims of the manufacturer and not those of VWR International, LLC and/or Avantor, Inc. or affiliates. All prices are in US dollars unless otherwise noted. Offers valid in US and Canada, void where prohibited by law or company policy, while supplies last. | Trademarks are owned by Avantor, Inc. or its affiliates, unless otherwise noted. | Visit [vwr.com](http://vwr.com) to view our privacy policy, trademark owners, and additional disclaimers. © 2019 Avantor, Inc. All rights reserved.

0420 Lit. No. 160158W