
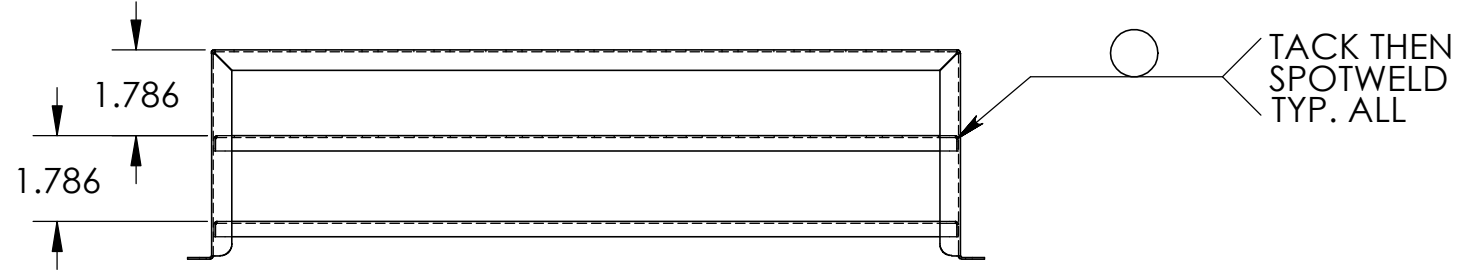
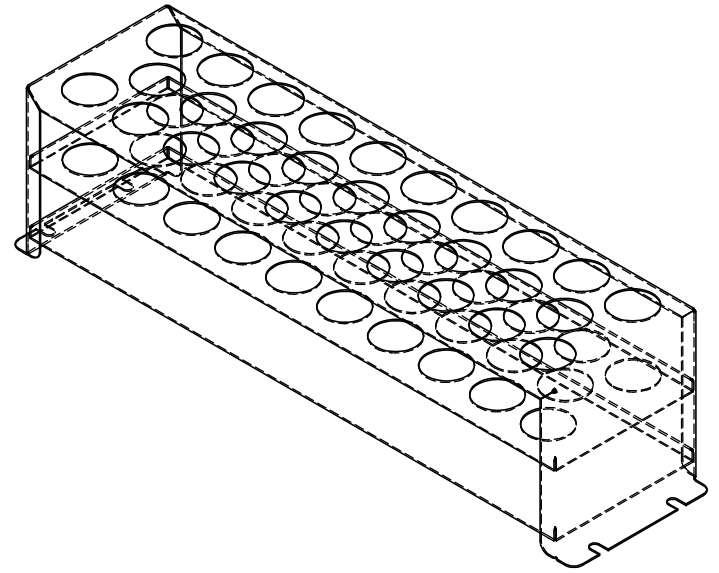


DWG BY: THA DATE: 7-18-02 TOLERANCE UNLESS OTHERWISE SPECIFIED .X = ± .063 ANGULAR .XX = ± .032 HOLE DIA .XXX = ± .015 ± .005	SHEL-LAB SHELDON MANUFACTURING CORNELIUS, OREGON 50ML SHAKER ASSY.	SCALE: 1:4 PART#: 9751181 MATERIAL: ASSY	NOTE BEND RADIUS .032" UNLESS OTHERWISE NOTED
--	--	--	---

CHG. BY	CHG. #	REV.	DATE	DESCRIPTION

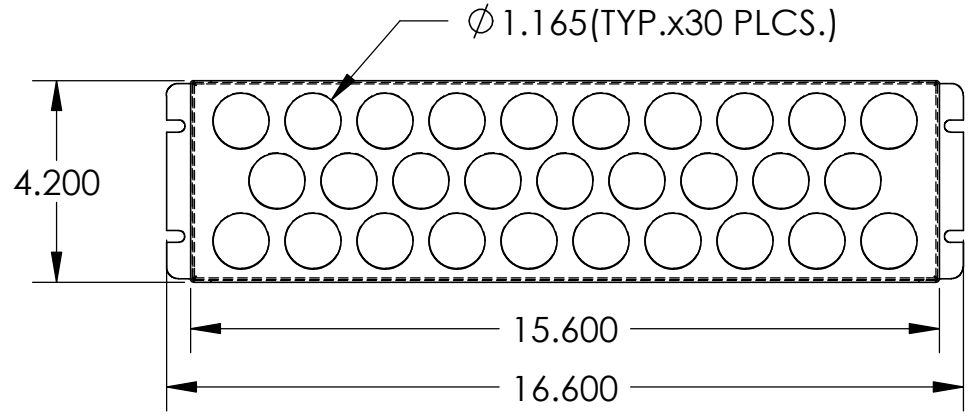

ITEM NO.	QTY.	PART NO.	MATERIAL	DESCRIPTION
1	1	5121184	20 GA 304-2B	SHAKER TOP 50ML
2	1	5121183	.036 SST. 304 2B	SHAKER SHELF 50ML
3	1	5121162	20 SST 304 2B	SHAKER SHELF