


GUIDE TO DONNING AND DOFFING OF NON-STERILE CLEANROOM GLOVES


DONNING GLOVES

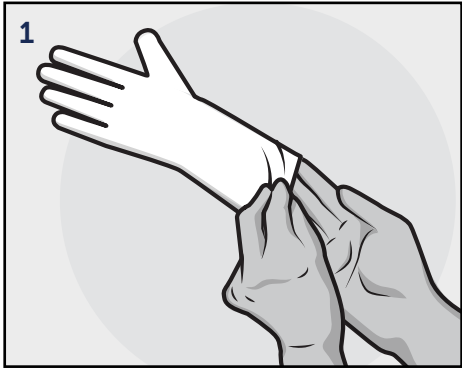


1 Grasp the beaded cuff of one glove and carefully pull over the opposite hand.



2 Gently ease the glove up over the hand and wrist.

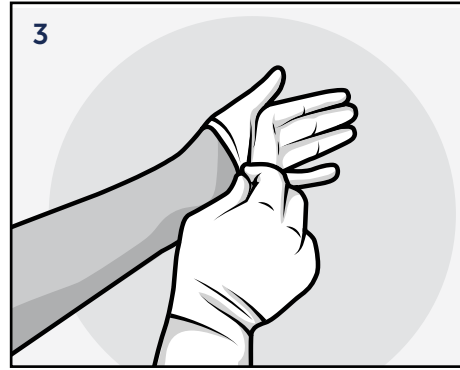

3 Grasp the second glove at the edge of the beaded cuff.

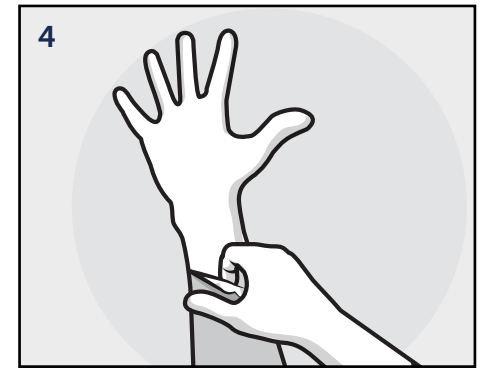

4 Gently ease the glove up over the hand and wrist.




5 You are now ready to perform your task.


DOFFING GLOVES



1 Grasp the exterior of one glove with your other gloved hand.

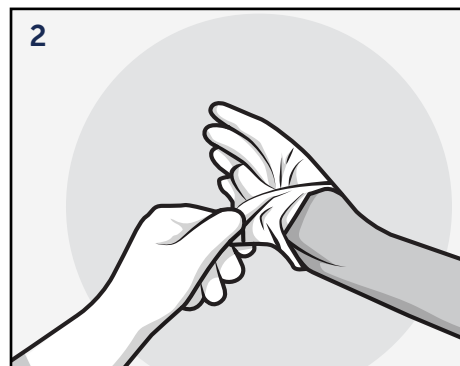


2 Carefully pull the glove off your hand, turning it inside-out.

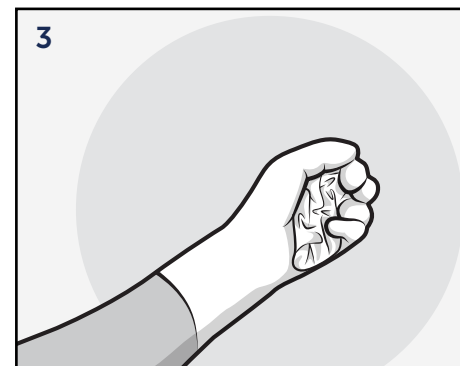


3 Ball the glove up and hold in your other gloved hand.



4 Slide your ungloved fingers into the opening of the other glove.

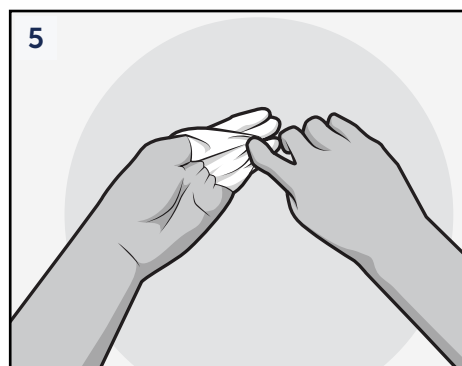

5 Carefully pull the glove off your hand, turning it inside-out and wrapping it around the doffed glove in your palm.

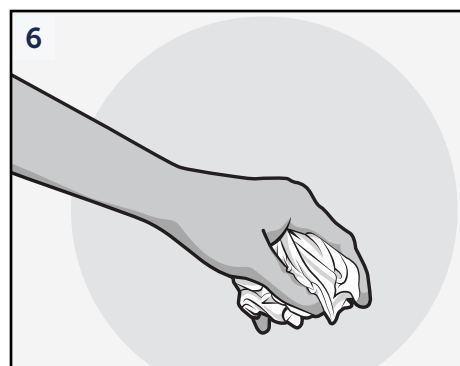

6 Discard appropriately.

avantor™ delivered by VWR™ VWR.COM

Prices, product, and/or services details are current when published and subject to change without notice. | Certain products or services may be limited by federal, state, provincial, or local regulations. | VWR, part of Avantor, makes no claims or warranties concerning sustainable/green products. Any claims concerning sustainable/green products are the sole claims of the manufacturer and not those of VWR International, LLC and/or Avantor, Inc. or affiliates. All prices are in Canadian dollars unless otherwise noted. Offers valid in Canada, void where prohibited by law or company policy, while supplies last. | Trademarks are owned by Avantor, Inc. or its affiliates, unless otherwise noted. | Visit vwr.com to view our privacy policy, trademark owners, and additional disclaimers. © 2022 Avantor, Inc. All rights reserved.

0722 Lit. No. 080098W

For more information, contact your HALYARD* representative or distributor, or visit www.purezerogloves.com