

# IMPORTANT! Please Read

## Conventional serological pipetting

When you first receive your ali-Q™ 2, press the purple button to wake it out of shipping mode.



Note:

1. The motor will activate at seemingly random times when pressing the aspirate or dispense buttons. Don't worry. It's normal!
2. ali-Q 2 will go into sleep mode after 3 minutes of non-use but will wake with any movement of the unit.

For more information on ali-Q's intelligent measuring system, see Operating Manual at [www.vistalab.com/support](http://www.vistalab.com/support)

## Aliquoting/repeat pipetting

speed dial



**ali-Q 2 VS model only:**

Set desired aliquoting speed using speed dial

aliquot dials




ali-Q 2 & ali-Q 2 VS



ali-Q 2 LS

**All ali-Q 2 models:**

1. Set aliquot volume on aliquot dial.
2. Aspirate adequate volume for total number of desired aliquots.
3. Press and HOLD the purple  aliquot button until the aliquot is dispensed. You will hear and feel the motor come on. Repeat press and hold for each aliquot. For maximum precision, discard first aliquot.

aliquot button



Note: if you don't hold the button long enough for the full aliquot volume, you will hear an "alert" sound and red light will flash.