
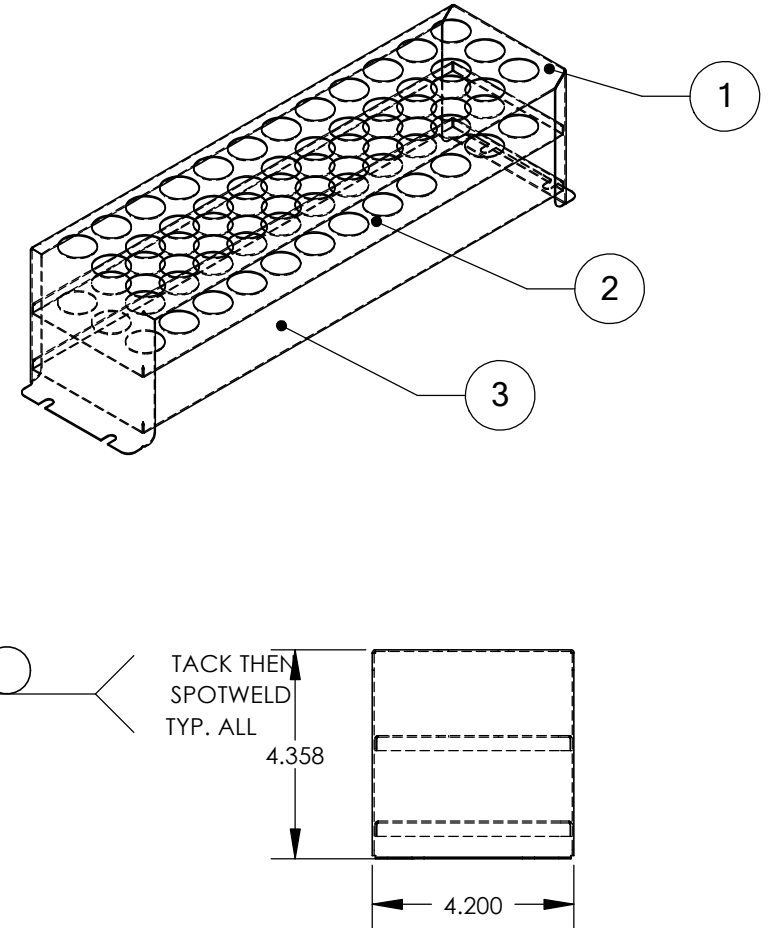


DWG BY: DKB	DATE: 3/18/02	TOLERANCE UNLESS OTHERWISE SPECIFIED: X = ± .063 ANGULAR .XX = ± .032 HOLE DIA .XXX = ± .015 ± .005	SHELDON MANUFACTURING CORNELIUS, OREGON	SCALE: 1:4	REVISION	NOTE BEND RADIUS .032" UNLESS OTHERWISE NOTED
			22-25 MM SHAKER ASSY	PART#: 9751180	MATERIAL: ASSEMBLY	

CHG. BY	CHG. #	REV.	DATE	DESCRIPTION
THA		A1	7-18-02	RECREATED CORRUPT DRAWING

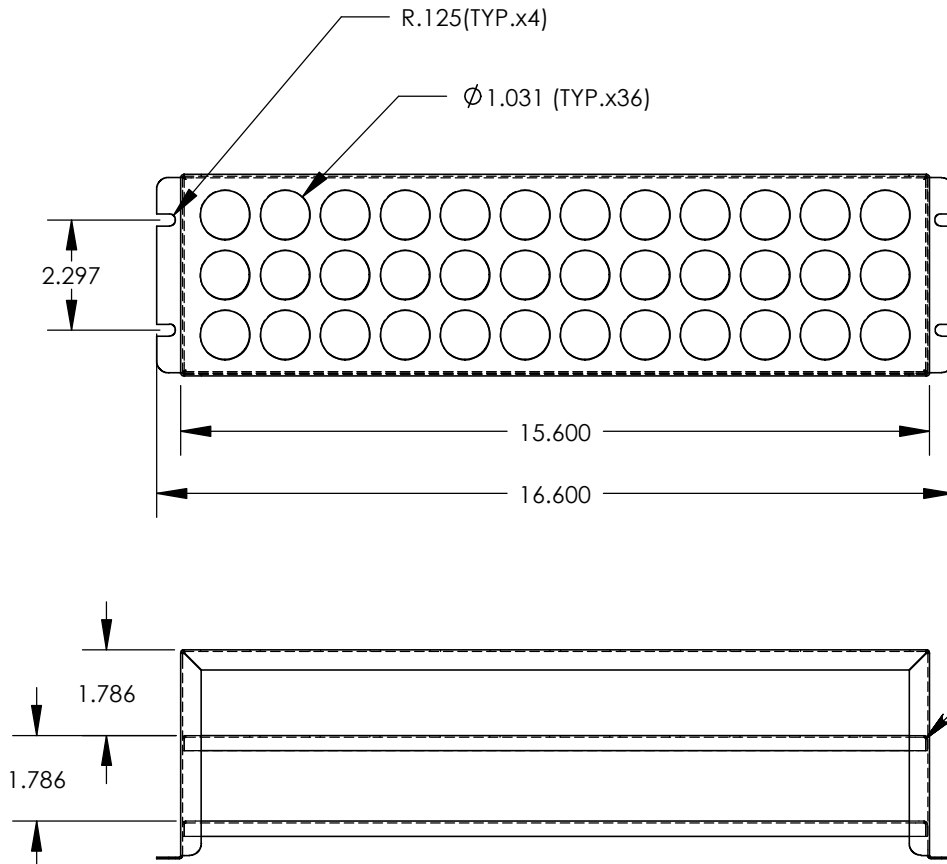

ITEM 1 NOT SHOWN,
RESIDES IN ASSEMBLY
AS PLACE HOLDER.

ITEM NO.	QTY.	PART NO.	MATERIAL	DESCRIPTION
1	1	5121169	.036 SST. 304 2B	SHAKER TOP 22-25 MM
2	1	5121170	20 GA 304-2B	SHAKER SHELF 22-25 MM
3	1	5121162	20 SST 304 2B	SHAKER SHELF