


USER MANUAL

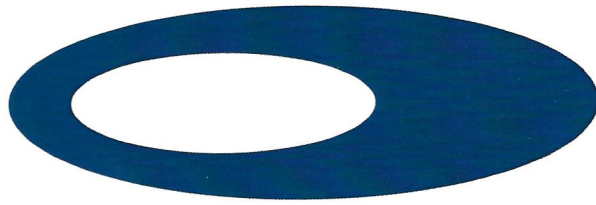

BootieButler[®]

Automatic Shoe Cover Systems



KBLG
Faster, Cleaner,
Safer!



How to properly load dispenser



Remove plastic bag.


Self Adjustable Track (with Allen wrench provided) to accommodate the size of the shoe cover cartridge.



Align the blue "clips" with the slots of the adjustable tracks (Note arrow pointing downward). Make sure ALL CLIPS enter the slots keeping the round tip on the outside of the track (as seen on picture).



Enlarged to show detail.

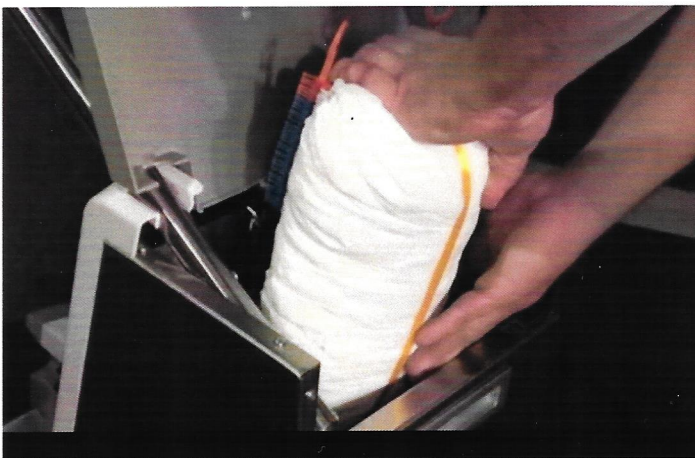

Push downward and forward (keeping the cartridge at a 90 degree angle), allowing the clips to enter easily in the slots. Note the use of both hands to ensure all clips enter the tracks.




Note the cartridge's position should be 3-4 inches BELOW top of the tracks. Remove the yellow ribbon and cardboard.




Place one hand on top of the cartridge. While holding down the shoe covers, remove the orange tabs, one at a time.



Push cartridge down until it stops at the underside of the resistance "tongue".

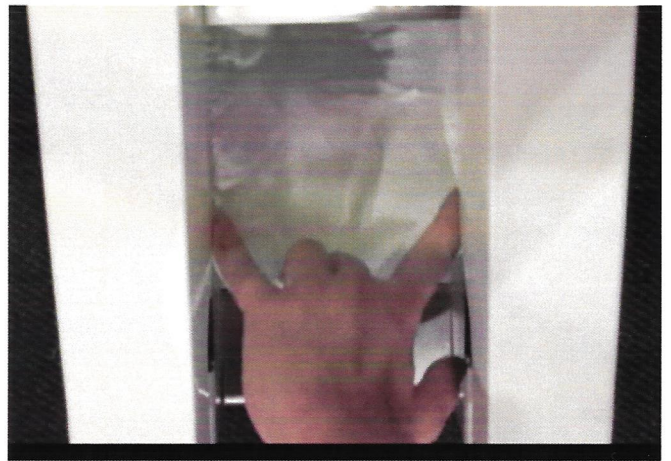

Prime the dispenser by facing the front and pulling out the front part of the exposed shoe cover. Then with your index and pinky fingers, pull gently until the first shoe cover "locks" into place on the holding bars on each side of the dispenser.

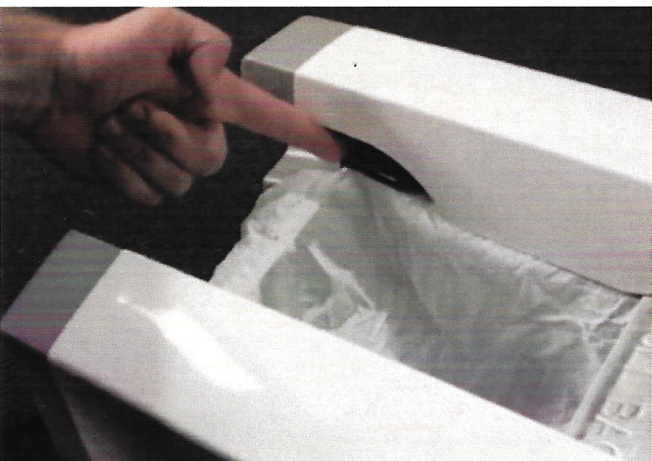

Properly loaded and ready for use!!

Now, putting shoe covers on... **EASY AS 1, 2, 3!**



Heel first.
Keep elastic level with top of shoe.

Don't press down to the floor with your heel, OR AT ALL.
Keep shoe 3-4" off of, and parallel to the ground.



SLOWLY pull back until cover stretches beyond the tip of your shoe.
Then... tilt your toe down below the elastic of the shoe cover.



Pull back until shoe cover is released.
Repeat on other foot.

Packaging Specification:

Machine Measurements: 29 9/64 x 12 19/32 x 31.5 Inches
Case Measurement (for shipping purposes): 32 x 16 x 32 Inches
Case Weight (for shipping purposes): 52 Lbs.

Packaging List:

Shoe Cover Dispenser: 1 piece
Handle: 1 piece
Screws: 4 pieces
Allen Wrench: 1 piece
User Manual: 1 piece

Customer Service:

BootieButler
3925 Papermill Dr.
Knoxville, TN 37909
info@bootiebutler.com
orders@bootiebutler.com
800.710.9863



For more information, visit our Video Section on our web page
www.BootieButler.com or call us at (800) 710-9863