



GUIDE TO DONNING AND DOFFING OF STERILE CLEANROOM GLOVES

DONNING GLOVES

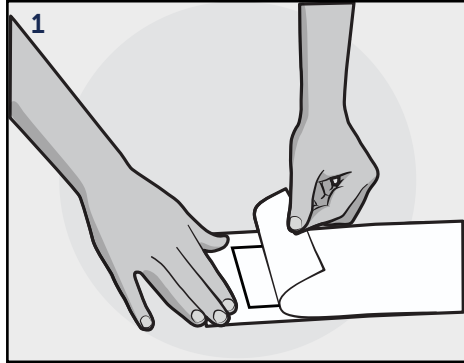


1 Peel open the pouch which keeps the product sterile. Open clear wallet.

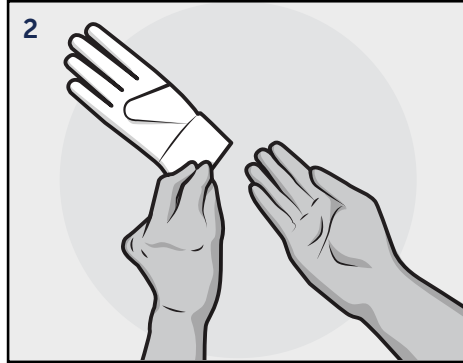


2 Grasp the interior of one glove and carefully pull over the opposite hand.




3 Gently ease the glove up over the hand and wrist, leaving the cuff folded.

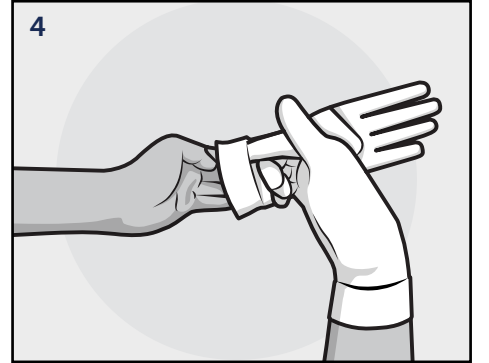


4 Lift the second glove under the folded cuff using gloved fingers.

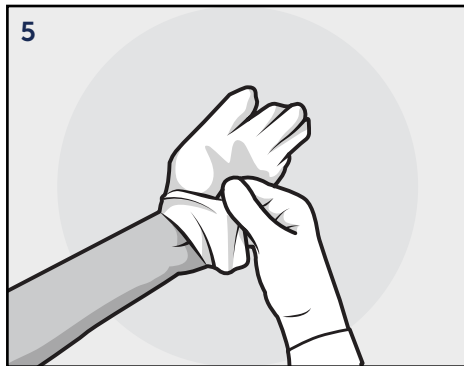

5 Slide your second hand into the glove by holding it under the folded cuff, then unfold the cuff over the wrist and arm.

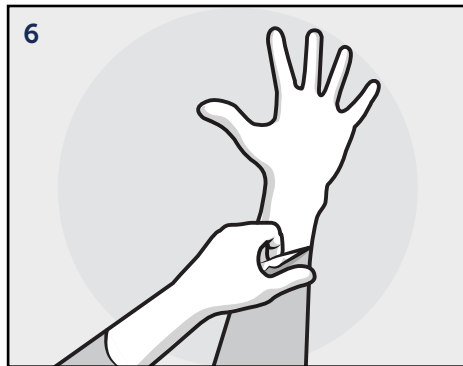


6 Using gloved fingers, unfold the second cuff over your wrist and arm.

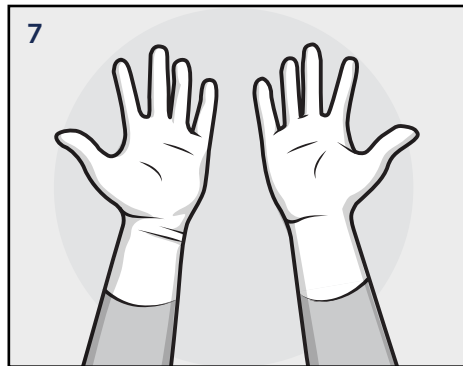


7 To remain sterile, do not touch anything with your gloved hands.


DOFFING GLOVES

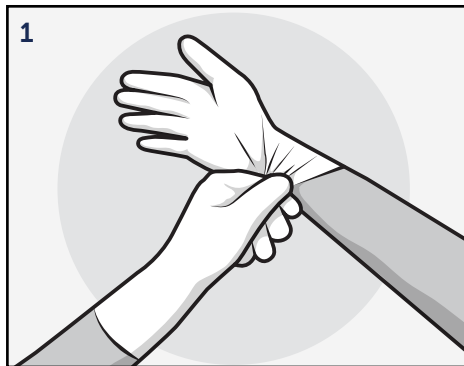


1 Grasp the exterior of one glove with your other gloved hand.

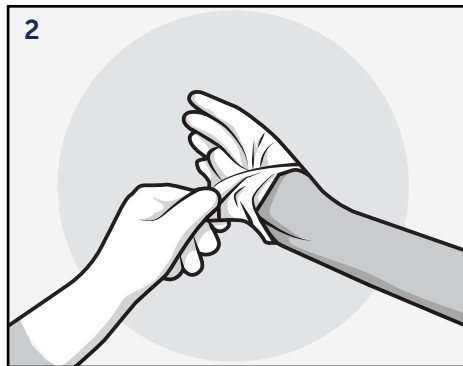


2 Carefully pull the glove off your hand, turning it inside-out.

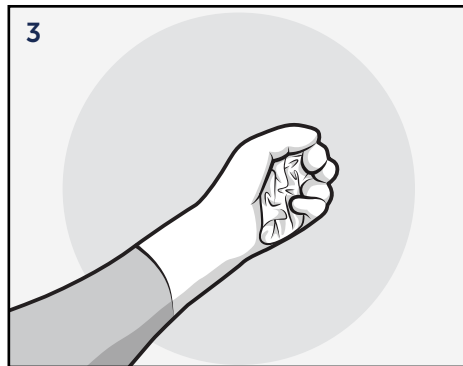


3 Ball the glove up and hold in your other gloved hand.

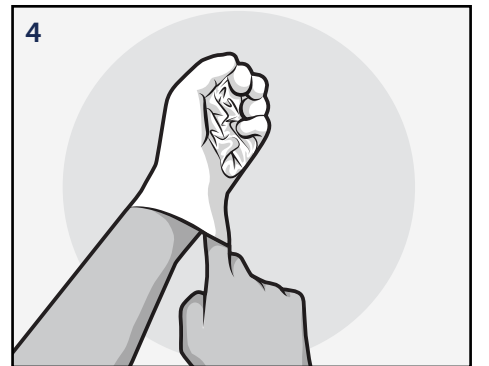

4 Slide your unglved fingers into the opening of the other glove.

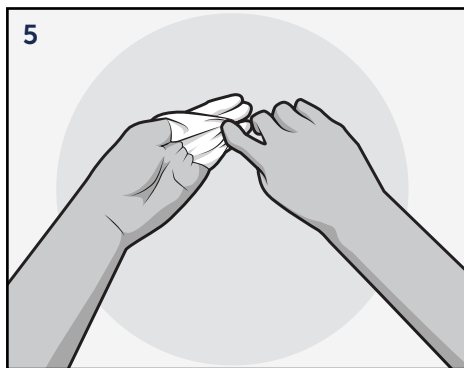

5 Carefully pull the glove off your hand, turning it inside-out and wrapping it around the doffed glove in your palm.

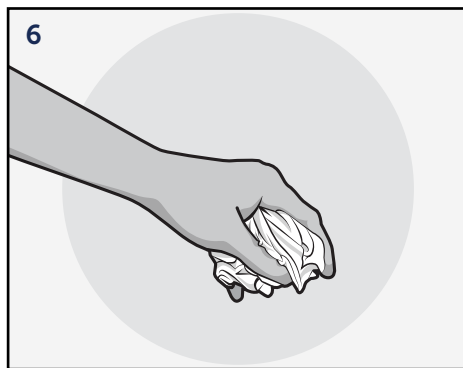

6 Discard appropriately.

HALYARD* recommends double donning/gloving. Don the gloves so that the sleeve cuffs are between the two pairs of gloves.

For more information, contact your HALYARD* representative or distributor, or visit www.purezerogloves.com

avantor™ delivered by VWR™ VWR.COM

Prices, product, and/or services details are current when published and subject to change without notice. | Certain products or services may be limited by federal, state, provincial, or local regulations. | VWR, part of Avantor, makes no claims or warranties concerning sustainable/green products. Any claims concerning sustainable/green products are the sole claims of the manufacturer and not those of VWR International, LLC and/or Avantor, Inc. or affiliates. All prices are in Canadian dollars unless otherwise noted. Offers valid in Canada, void where prohibited by law or company policy, while supplies last. | Trademarks are owned by Avantor, Inc. or its affiliates, unless otherwise noted. | Visit vwr.com to view our privacy policy, trademark owners, and additional disclaimers. © 2022 Avantor, Inc. All rights reserved.

0722 Lit. No. 080098W