



## Pall Nucleic Acid Binding Nanosep<sup>®</sup> Centrifugal Device

### Description

The Pall Nanosep device for nucleic acid binding (NAB) incorporates a dual layer silica-based quartz glass fiber media to allow for efficient binding of DNA and RNA, while providing smooth flow and rapid processing of samples. This media offers researchers the flexibility to purify plasmid DNA, genomic DNA or total RNA from a variety of starting materials: a single column for multiple applications. The new NAB Nanosep device offers flexibility without sacrificing on quantity or quality of the nucleic acid. The Pall NAB Nanosep device is a multipurpose centrifugal device providing flexibility in applications, supported by comprehensive protocols for sample processing.

### Maximum yields and quality of nucleic acid purification

- Rapid processing of samples
- Silica-based quartz glass fiber media that allows efficient binding of DNA and RNA
- Constructed of low-binding polypropylene
- Ultrasonically welded seals prevent bypass or seal failure
- Fits standard centrifuge rotors that accept 1.5 mL tubes

### Downstream Applications

#### Purified Plasmid DNA

- Restriction digest
- Cloning
- Sanger sequencing

#### Purified Genomic DNA

- qPCR
- Next Generation Sequencing

#### Purified Total RNA

- RT-PCR
- cDNA library construction



### Specifications

#### Materials of Construction

Dual Filter Media: Binder-free, silica-based glass fiber  
Housing: Polypropylene

#### Effective Filtration Area

0.28 cm<sup>2</sup>

#### Sample Volume

Up to 500 µL

#### Dimensions

Overall length (fully assembled): 45 mm (1.8 inches)  
Column insert height: 23 mm (0.9 inches)


#### Operating specs

Operating Temperature Range: 15 – 30 °C (59 – 86 °F)  
Max. Centrifugal Force: 14,000 x g  
Minimum Elution Vol.: 40 µL  
Base Pair Binding from 50 bp  
Binding Capacity, up to: 160 µg pure RNA,  
69 µg pure genomic DNA, 11.5 µg pure plasmid DNA

## Performance

**Figure 1**


Genomic DNA isolated from freshly harvested *E. coli* cells, CHO cells and plant leaves compared to the competitor commercial product with commercial buffers for genomic DNA isolation.

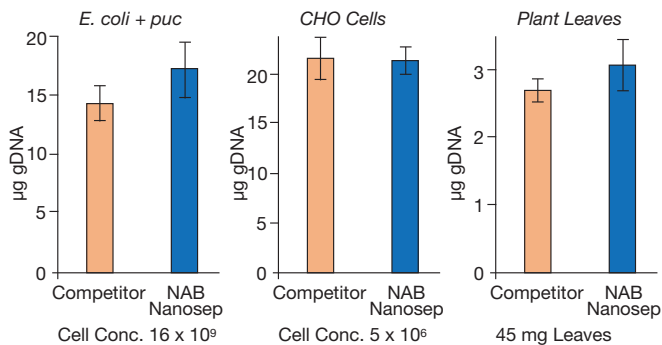

Purity ( $A_{260}/A_{280}$ ) *E. coli*  $1.77 \pm .05$ , CHO  $1.79 \pm .06$ , Basil  $1.84 \pm .04$

**Figure 2**


Total RNA isolated from freshly harvested *E. coli* cells, CHO cells and plant leaves compared to the competitor commercial product with commercial buffers for RNA isolation.

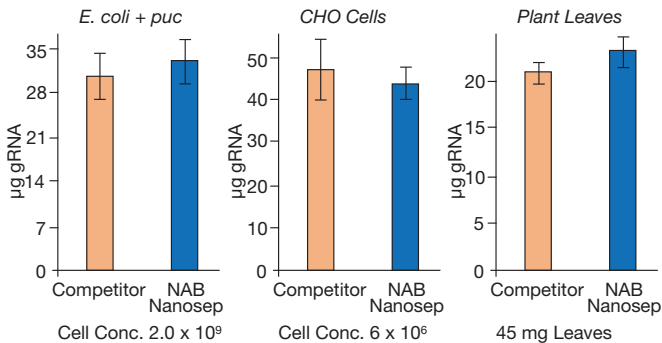

Purity ( $A_{260}/A_{280}$ ) *E. coli*  $2.05 \pm .04$ , CHO  $2.04 \pm .01$ , Basil  $2.07 \pm .01$

**Figure 3**

Plasmid DNA isolated from freshly harvested *E. coli* cells compared to the competitor commercial product with commercial buffers.

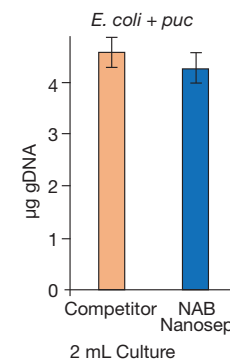

Purity ( $A_{260}/A_{280}$ )  $1.79 \pm .03$

Yield, purity and reproducibility of nucleic acid isolated from a variety of starting materials are similar for the NAB Nanosep (blue bars) device when compared to the three different devices provided with three different kits from a single competitor (Orange bars). This demonstrates the cross functional capabilities of the NAB Nanosep device; a single device, capable of replacing several competitor devices for all nucleic acid applications without compromising on yield or purity. This is a single competitive product that has been compared to a competitor's products following the manufacturers protocols.

## Ordering Information

### NAB Nanosep Device

VWR Cat. No.	Pall Part No.	Description	Pkg
76360-454	ODNABC33	NAB Nanosep Device	24/pkg*
76360-456	ODNABC34	NAB Nanosep Device	100/pkg*

\*Both pack sizes come with 2 additional filtrate tubes for each device

### Accessories and Replacement Parts

VWR Cat. No.	Pall Part No.	Description	Pkg
29300-650	FDX001X34	Nanosep Filtrate Tube	100/pkg

### Related Products

VWR Cat. No.	Pall Part No.	Description	Pkg
10158-728	8133	AcroPrep™ Advance 96 Well Long Tip Filter Plate for Nucleic Acid Binding	5/pkg



Ordering:  
vwr.com

© 2020 Pall Corporation. Pall, PALL, AcroPrep, and Nanosep are trademarks of Pall Corporation. ® indicates a trademark registered in the USA. Filtration. Separation. Solution. is a service mark of Pall Corporation.

1/20, PDF, GN19.1024

Filtration. Separation. Solution.™

800 932 5000 | VWR.COM

Prices and product details are current when published and subject to change without notice. | Certain products may be limited by federal, state, provincial, or local regulations. | VWR, part of Avantor, makes no claims or warranties concerning sustainable/green products. Any claims concerning sustainable/green products are the sole claims of the manufacturer and not those of VWR International, LLC and/or Avantor, Inc. or affiliates. All prices are in US dollars unless otherwise noted. Offers valid in US and Canada, void where prohibited by law or company policy, while supplies last. | Trademarks are owned by Avantor, Inc. or its affiliates, unless otherwise noted. | Visit vwr.com to view our privacy policy, trademark owners, and additional disclaimers. © 2019 Avantor, Inc. All rights reserved.