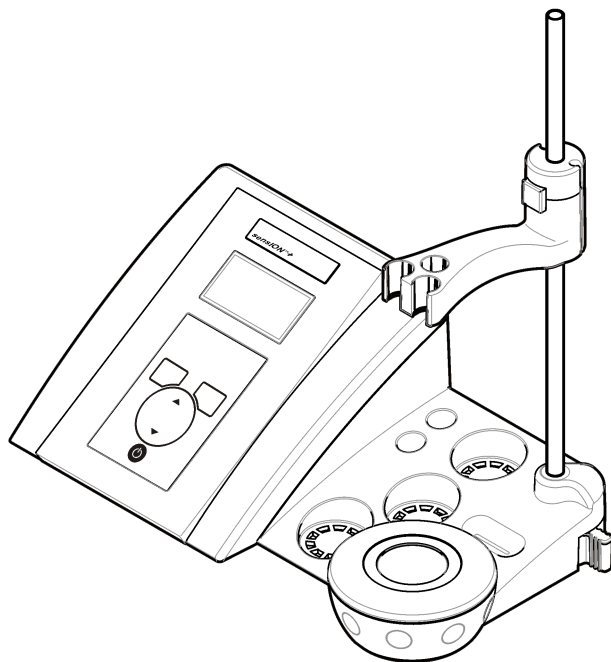
DOC022.97.90250

senSIOn™ + MM340

04/2011, Edition 1

User Manual
Manuel d'utilisation
Manual del usuario
Manual do usuário
用户手册
取扱説明書
사용 설명서
คู่มือผู้ใช้

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Specifications

Specifications are subject to change without notice.

Specification	Details
Dimensions	35 x 20 x 11 cm (13.78 x 7.87 x 4.33 in.)
Weight	1100 g (2.43 lb)
Meter enclosure	IP42
Power requirements (external)	100–240 V, 0.4 A, 47–63 Hz
Meter protection class	Class II
Storage temperature	–15 to +65 °C (5 to +149 °F)
Operating temperature	0 to 40 °C (41 to 104 °F)
Operating humidity	< 80% (non-condensing)
Resolution	pH: 0.1/0.01/0.001, ORP: 0.1/1 mV, ISE: programmable, temperature: 0.1 °C (0.18 °F)
Measuring error (± 1 digit)	pH: ≤ 0.002, ORP: ≤ 0.2 mV, temperature: ≤ 0.2 °C (≤ 0.36 °F)
Reproducibility (± 1 digit)	pH: ± 0.001, ORP: ± 0.1 mV, temperature: ± 0.1 °C (± 0.18 °F)
Data storage	330 results and last 9 calibrations
Connections	2 Combined or indicator probe: BNC connector (Imp. >10 ¹² Ω); 2 Reference electrodes: banana connector; A.T.C. type Pt 1000: banana or telephonic connector magnetic stirrer: RCA connector; 2 magnetic stirrers: RCA connector
Temperature correction	Manual, Pt 1000 temperature probe (A.T.C.), isopotential pH programmable, standard value 7.00
Measurement display lock	Continuous measurement, by stability
Display	Liquid crystal, backlit, 128 x 64 dots

Specification	Details
Keyboard	PET with protective treatment
Certification	CE

General information

Revised editions are found on the manufacturer's website.

Safety information

Please read this entire manual before unpacking, setting up or operating this equipment. Pay attention to all danger and caution statements. Failure to do so could result in serious injury to the operator or damage to the equipment.

Make sure that the protection provided by this equipment is not impaired, do not use or install this equipment in any manner other than that specified in this manual.

Use of hazard information

▲ DANGER

Indicates a potentially or imminently hazardous situation which, if not avoided, will result in death or serious injury.

▲ WARNING

Indicates a potentially or imminently hazardous situation which, if not avoided, could result in death or serious injury.

▲ CAUTION

Indicates a potentially hazardous situation that may result in minor or moderate injury.

NOTICE

Indicates a situation which, if not avoided, may cause damage to the instrument. Information that requires special emphasis.

Precautionary labels

Read all labels and tags attached to the instrument. Personal injury or damage to the instrument could occur if not observed. A symbol, if noted

on the instrument, will be included with a danger or caution statement in the manual.

	<p>This symbol, if noted on the instrument, references the instruction manual for operation and/or safety information.</p>
	<p>Electrical equipment marked with this symbol may not be disposed of in European public disposal systems after 12 August of 2005. In conformity with European local and national regulations (EU Directive 2002/98/EC), European electrical equipment users must now return old or end-of-life equipment to the Producer for disposal at no charge to the user.</p> <p><i>Note: For return for recycling, please contact the equipment producer or supplier for instructions on how to return end-of-life equipment, producer-supplied electrical accessories, and all auxiliary items for proper disposal.</i></p>

Product overview


The sensION™+ meters are used with probes to measure various parameters in water.

The sensION™+ MM340 meter has two measurement channels to measure pH, ORP (mV) or ISE (concentration) with a selective probe. Measurement data can be stored and transferred to a printer or PC.

Product components

Refer to [Figure 1](#) to make sure that all components have been received. If any items are missing or damaged, contact the manufacturer or a sales representative immediately.

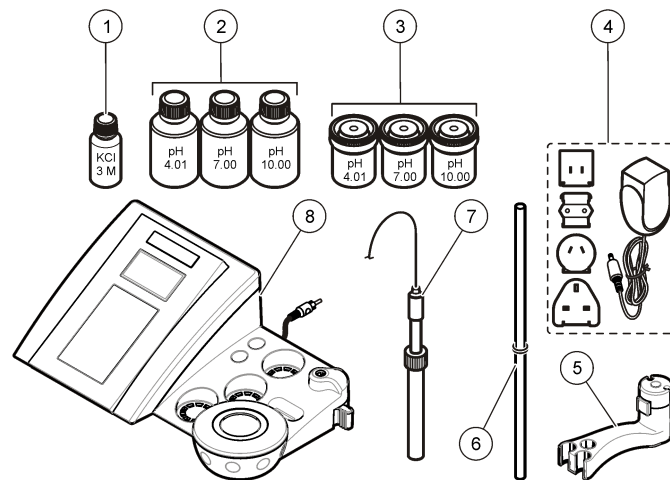Figure 1 Meter components





<p>1 Electrolyte for the probe</p>	<p>5 Probe holder</p>
<p>2 Buffer solutions (pH 4.01, pH 7.00 and pH 10.00)</p>	<p>6 Rod with o-ring</p>
<p>3 Calibration beakers (with magnetic bar inside)</p>	<p>7 Probe (included with kits only)</p>
<p>4 Power supply</p>	<p>8 Meter</p>

Installation

Assemble the probe holder

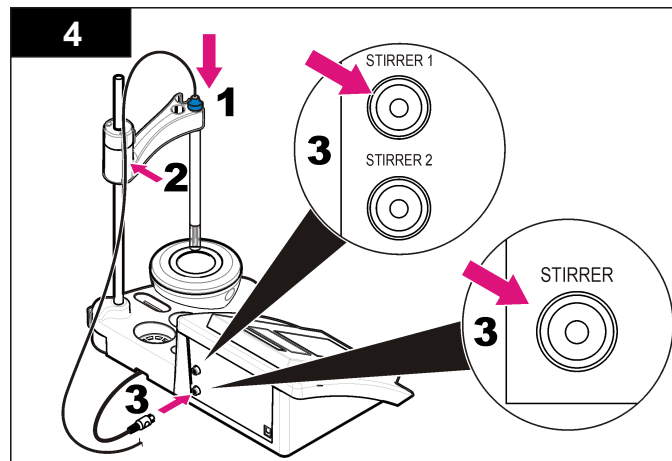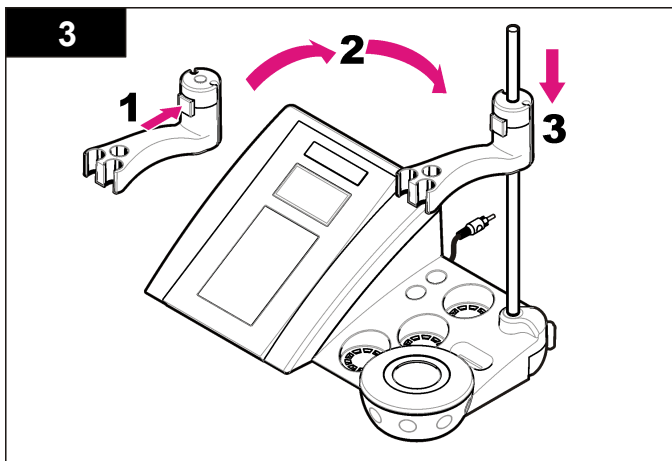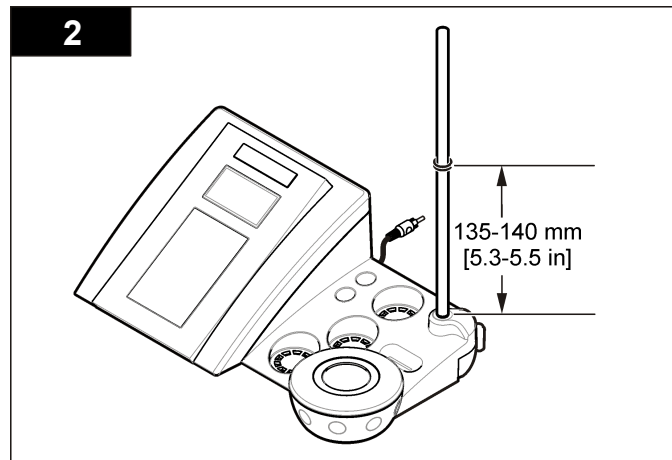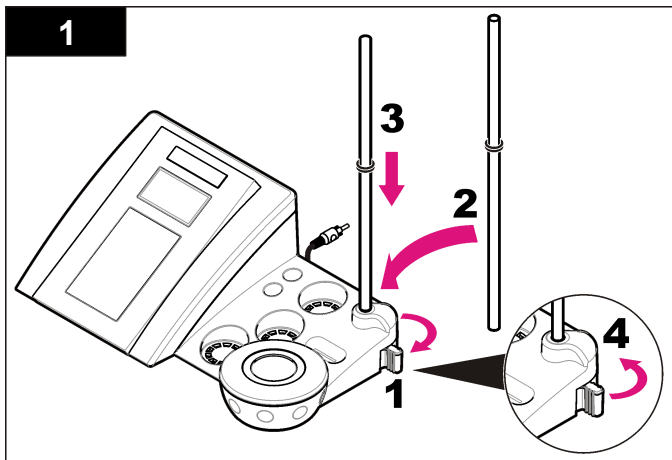Follow the numbered steps to assemble the probe holder and to connect the magnetic stirrer.

Connect to AC power

⚠ DANGER


Electrocution hazard. If this equipment is used outdoors or in potentially wet locations, a Ground Fault Circuit Interrupt (GFCI/GFI) device must be used to connect the equipment to its main power source.

The meter can be powered by AC power with the universal power adapter.


1. Select the correct adapter plug for the power outlet from the adapter kit.
2. Connect the universal power adapter to the meter ([Figure 2](#)).
3. Connect the universal power adapter to an AC receptacle ([Figure 3](#)).
4. Turn the meter on.

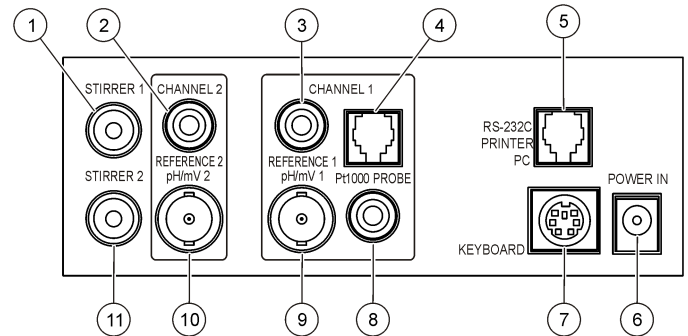Figure 2 Connector panel

1 Magnetic stirrer 1 connector, Channel 1	7 PC keyboard, mini DIN connector
2 Reference electrode (separated electrodes) connector, Channel 1	8 Temperature probe connector, Channel 2
3 Reference electrode (separated electrodes) connector, Channel 2	9 Combined pH electrode (or indicator) connector, Channel 2
4 Separated temperature probe connector, Channel 2	10 Combined pH electrode (or indicator) connector, Channel 1
5 RS-232C for printer or PC connector	11 Magnetic stirrer 2 connector, Channel 2
6 Power supply	

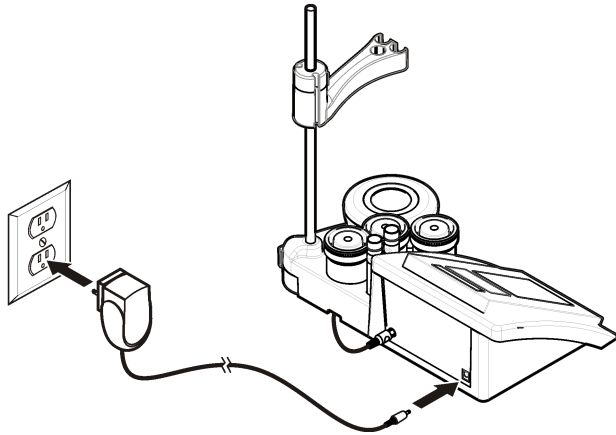
Figure 3 AC power connection

User interface and navigation

User interface

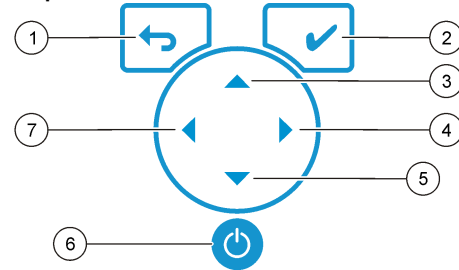Keypad description


1 RETURN key: cancel or exit the current menu screen to the previous menu screen	5 DOWN key: scroll to other options, change a value
2 MEASUREMENT key: confirm the selected option	6 ON/OFF: turn on or turn off the meter
3 UP key: scroll to other options, change a value	7 LEFT key: change between Channel 1 and Channel 2, enter numbers and letters
4 RIGHT key: change between Channel 1 and Channel 2, enter numbers and letters	

Display description

The meter display shows the concentration, units, temperature, calibration status, operator ID, sample ID, date and time.

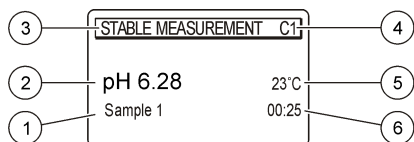Figure 4 Single screen display

1 Sample ID	4 Sample temperature (°C or °F)
2 Measurement unit and value (pH, ORP (mV) or ISE)	5 Visual measurement timer
3 Measurement mode or time and date	

Navigation

Use the to return to the previous menu. Use the measure key to take a sample measurement or to confirm options. Use the arrow keys to scroll to other options or to change a value. To change the parameters use the arrow keys and . Refer to each task for specific instructions.

Start-up

Turn the meter on and off

NOTICE

Make sure that the probe is connected to the meter before the meter is turned on.

Push to turn on or turn off the meter. If the meter does not turn on, make sure that the AC power supply is properly connected to an electrical outlet.

Change the language

The display language is selected when the meter is powered on for the first time.

1. Use the or to select a language from a list.
2. Confirm with . The measurement screen shows DATA OUTPUT.
3. Select Deactivated if no printer or PC is connected and confirm. Refer to [Select the data output](#) on page 10 for more information about Data Output.

Standard operations

Calibration

⚠ WARNING

Chemical exposure hazard. Obey laboratory safety procedures and wear all of the the personal protective equipment appropriate to the chemicals that are handled. Refer to the current material safety data sheets (MSDS) for safety protocols.

Calibration settings

The calibration settings contain Calibration type, Calibration frequency and Display options.

1. From the main menu use the or to select CALIBRATION. Confirm.
2. Use the to enter the calibration menu.
3. Use the or to select the following options:

Option	Description
Stability C.:	Criterion by stability—select Fast, Standard or Strict.
Calibration type	Calibration type—select Technical buffers, DIN19266 Buffers, User Buffers, To a X value, data introduction or Theoretical calibration. Refer to Calibration types for more information.
Cal. frequency	Calibration reminder—can be set between 0–7 days (default daily). The display shows the remaining time to the new calibration. Refer to Set the calibration reminder on page 9 for more information.
Display mV	Display mV—select YES or NO to show mV.
Standard 220 mV	Standard 220 mV—calibrate with a 220 mV standard solution
To a X value	To a X value—use a specific standard solution. Adjust the value during the calibration.
Data introduction	Data introduction Compensation—enter the compensation value
Factory adjust	Factory adjust—enter the compensation value

Calibration types

Different calibration types can be selected.

1. From the main menu use the ▲ or ▼ to select CALIBRATION. Confirm.
2. Use the ▲ to enter the calibration menu.
3. Use the ▲ or ▼ to select Calibration type.

Option	Description
Technical buffers	pH 2.00, 4.01, 7.00, 9.21 and 10.90 at 25 °C (77 °F)
DIN19266 Buffers	pH 1.679, 4.006, 6.865, 7.000, 9.180, 10.012 and 12.454
User Buffers	Selected when the technical or DIN19266 buffers are not used. Refer to Technical buffer solutions (DIN 19267) on page 17 for pH values of specific buffer sets at varying temperatures.
Calibration to a X value	To adjust manually any scale value of the measured pH.
Data introduction	Manual probe constant introduction.
Theoretical calibration	The probe calibration data is replaced at 25 °C (77 °F).

Calibration procedure

This procedure is for general use with liquid calibration solutions. Refer to the documents that are included with each probe for additional information.

Note: Solutions should be stirred during calibration. For more information about the stirring settings, refer to [Change the stirring settings](#) on page 11.

1. Pour the buffers or calibration solutions into the labeled calibration beakers.
2. From the main menu use the ▲ or ▼ and ◀ and ▶ to select the CALIBRATION parameter. Confirm.
3. If required select the Operator ID (1 to 10) and confirm.
4. Rinse the probe with deionized water and put the probe into the first calibration beaker. Be sure that there are no air bubbles in the membrane.

5. Push ✓ to Start calibration.
6. Push ✓ to measure the first calibration solution. The next calibration solution is shown.
7. Rinse the probe with deionized water and put the probe into the first calibration beaker. Be sure that there are no air bubbles in the membrane.
8. Push ✓ to measure the second calibration solution. The next calibration solution is shown.
9. Rinse the probe with deionized water and put the probe into the first calibration beaker. Be sure that there are no air bubbles in the membrane.
10. Push ✓ to measure the third calibration solution. When the calibration is good, the display briefly shows Calibration OK and then returns to the main menu.
Note: When a printer is connected the print menu opens and the result can be printed.

View the calibration data

Data from the most recent calibration can be shown.

1. From the main menu use the ▲ or ▼ to select DATA LOGGER. Confirm.
2. Select Display data.
3. Select Calibration data and confirm with ✓. The last calibration data is shown.
 - pH-the slope and offset values are shown alternating with the deviation (in %) and calibration temperature.
 - ORP-the measured mV value and calibration temperature are shown.
 - Conductivity-the cell constant and calibration temperature for each standard are shown.

Set the calibration reminder

The calibration reminder can be set between 0 to 23 hours or 1-7 days (default 0 days). The display shows the remaining time to the new calibration.

Note: When 0 days is selected, the calibration reminder is turned off.

1. From the main menu use the ▲ or ▼ to select CALIBRATION. Confirm.
2. Use the ▲ to enter the calibration menu.
3. Use the ▲ or ▼ to select Cal. frequency and confirm.
4. Use the ◀ and ▶ to advance to the next step and use the ▲ or ▼ to change a value. Confirm.
Push ✓ to start the calibration.

Sample measurements

Each probe has specific preparation steps and procedures for taking sample measurements.

1. From the main menu use the ▲ or ▼ and ◀ and ▶ to select MEASURE. Confirm.
2. Use the ▲ to change the following settings. Confirm every entry.

Option	Description
Resolution	Select the resolution: 1, 0.1, 0.01 (default) or 0.001
Measure	Stability—select By stability Criterion: Fast (variation < 0.02 pH in 6 s), Standard (variation < 0.01 pH in 6 s) or Strict (variation < 0.002 pH in 6 s). In continuous—enter the time interval for the In continuous Acquis. interval (data storage or printing data). By time—enter the time interval for data storage or printing data.
Display mV	Display mV—select YES or NO to show mV.
Limits	Limits—select YES or NO. YES: Enter the upper and lower limits. An acoustic warning appears when the measurement is out of limit. The report output shows an A next to the measured value when the measurement was out of limit.
Isopotential	Isopotential—change the Isopotential pH value in Data introduction. Select Calculate to calibrate the probe again.

3. Push ✓ to start the measurement.

Note: If the measurement is not stabilizing after 120 seconds, the meter turns automatically into the continuous measurement mode.

Advanced operations

Change the measurement units

The measurement units can be changed individually for each channel.

1. From the main menu use the ▲ or ▼ to select SYSTEM. Confirm.
2. Use the ▲ or ▼ to select Measurement units and confirm.
3. Select Channel 1 or Channel 2 and confirm.
4. Select ORP (mV), pH or ISE and confirm.

Use a sample ID

The sample ID tag is used to associate readings with a particular sample location. If assigned, stored data will include this ID.

1. From the main menu use the ▲ or ▼ to select SYSTEM. Confirm.
2. Use the ▲ or ▼ to select Sample ID and confirm.
3. Use the ▲ or ▼ to select

Option	Description
Automatic	A consecutive number will be automatically assigned to every sample.
Manual	A keyboard or a barcode scanner is required to enter the sample ID name before taking a measurement (maximum 15 characters).

Select the data output

Data can be stored or transferred to a printer or to a PC.

1. From the main menu use the ▲ or ▼ to select SYSTEM. Confirm.
2. Use the ▲ or ▼ to select Data Output and confirm.
3. Use the ▲ or ▼ to select

Option	Description
Deactivated	Select Deactivated if no printer or PC is connected.

Option	Description
For Printer	Select Dot matrix printer or Thermal printer.
For Computer	Select Terminal, ComLabo or ComLabo Easy. The ComLabo Software controls several modules, pH and conductivity meters, automatic burettes, Samplers and so on from a computer. The ComLabo Easy software gets pH and conductivity data from a PC.

Change the date and time

The date and time can be changed from the Date / Time menu.

1. From the main menu use the ▲ or ▼ to select SYSTEM. Confirm.
2. Use the ▲ or ▼ to select Date / Time and confirm.
3. Use the ◀ and ▶ to advance to the next step and use the ▲ or ▼ to change a value. Confirm.
The current date and time will be shown on the display.

Adjust the display contrast

1. From the main menu use the ▲ or ▼ to select SYSTEM. Confirm.
2. Use the ▲ or ▼ to select Display contrast and confirm.
3. Use the ◀ and ▶ to adjust the contrast of the display and confirm.

Adjust the temperature

The temperature measurement can be adjusted at 25 °C (77 °F) and/or 85 °C (185 °F) to increase accuracy.

1. Put the probe and a reference thermometer in a container of water at approximately 25 °C and allow the temperature to stabilize.
2. Compare the temperature read by the meter with that of the reference thermometer. The difference is the adjustment value for the meter.
Example: reference thermometer: 24.5 °C; meter: 24.3 °C. Adjustment value: 0.2 °C.
3. Enter the adjustment value for the 25 °C reading:
 - a. From the main menu use the ▲ or ▼ to select SYSTEM. Confirm.

- b. Use the ▲ or ▼ to select Readjust temp. and confirm.
 - c. Select Channel 1 or Channel 2 and confirm. The temperature for channel 1 is measured by the conductivity cell and channel 2 is measured by the pH probe. If no Conductivity cell is connected, the temperature value must be entered manually or the temperature measured in channel 2 can be applied to the meter.
 - d. Use the ▲ or ▼ to select 25 °C and confirm.
 - e. Use the arrow keys to enter the adjustment value for 25 °C. Confirm.
4. Put the probe and a reference thermometer in a container of water at approximately 85 °C and allow the temperature to stabilize.
 5. Compare the temperature from the meter with that of the reference thermometer. The difference is the adjustment value for the meter.
 - a. Use the ▲ or ▼ to select 85 °C and confirm.
 - b. Use the arrow keys to enter the adjustment value for 85 °C. Confirm.
 - c. Select Save changes and confirm.

Change the stirring settings

The magnetic stirrer can be turned on and the stirring speed can be changed in the Stirring menu.

1. From the main menu use the ▲ or ▼ to select SYSTEM. Confirm.
2. Use the ▲ or ▼ to select Stirring and confirm.
3. To turn on/off the stirrer push ✓.
4. When the stirrer is turned on, use the ▲ or ▼ to change the stirring speed in %.

Note: Use the ▲ or ▼ to change the stirring speed during calibration and during a measurement.

Turn the stirrer on or off

The stirrer 1 works with Channel 1 and 2 (Stirrer 1). A second stirrer can be connected to channel 2 (Stirrer 2). For the stirrer 2 activation refer to the following steps.

1. From the main menu use the ▲ or ▼ to select SYSTEM. Confirm with ✓.
2. Use the ▲ or ▼ to select Stirrer N.2 and confirm with ✓.
3. Use the ▲ or ▼ to select YES to turn the stirrer 2 on.
Note: Select NO to turn the stirrer 2 off.

Change the temperature units

The temperature units can be changed to Celsius or Fahrenheit.

1. From the main menu use the ▲ or ▼ to select SYSTEM. Confirm.
2. Use the ▲ or ▼ to select Temperature units and confirm.
3. Use the ▲ or ▼ to select between Celsius or Fahrenheit and confirm.

Data logger

Display data

The Display data log contains Measurement data, Electrode report and Calibration data. The stored data can be send to a printer or to a PC. When the data log becomes full (400 data points), the oldest data point is deleted when a new data point is added.

1. Use the ◀ and ▶ to select Channel 1 or Channel 2 and confirm.
2. From the main menu use the ▲ or ▼ to select SYSTEM. Confirm.
3. Use the ▲ or ▼ to select DATA LOGGER and confirm.
4. Use the ▲ or ▼ to select Display data and confirm.
5. Use the ▲ or ▼ to select

Option	Description
Measurement data	Measurement data—stores automatically each time a sample is measured
Electrode report	Electrode report—stores automatically the electrode history and measurement conditions
Calibration data	Calibration data—stores automatically the current calibration

Delete data

The entire measurement data or electrode report log can be deleted to remove data that has already been sent to a printer or PC.

1. From the main menu use the ▲ or ▼ to select SYSTEM. Confirm.
2. Use the ▲ or ▼ to select DATA LOGGER and confirm.
3. Use the ▲ or ▼ to select Erase and confirm.
4. Use the ▲ or ▼ to select Measurement data or Electrode report and confirm. Confirm again to delete the data.
The entire log is deleted at once.

Send data to a printer or to a computer

NOTICE

The data output (printer or PC) needs to be selected first, so that the Print menu is available (refer to [Select the data output](#) on page 10).

Note: Refer to [Report output](#) on page 12 to select the report output type.

1. From the main menu use the ▲ or ▼ to select SYSTEM. Confirm.
2. Use the ▲ or ▼ to select DATA LOGGER and confirm.
3. Use the ▲ or ▼ to select Print and confirm. Select one of the following options and confirm with ✓ to print the data: Measurement data, Electrode data, Calibration data, Calibration report or Instrument condit.

Report output

NOTICE

The data output (printer or PC) needs to be selected first, so that the Type of report menu is available (refer to [Select the data output](#) on page 10).

Different report output types can be selected when a printer or a PC is connected.

1. From the main menu use the ▲ or ▼ to select SYSTEM. Confirm.
2. Use the ▲ or ▼ to select Type of report and confirm.

3. When a printer or a computer and Terminal is connected, use the ▲ or ▼ to select

Option	Description
Reduced	Select Several or One sample as an output format
Standard	Select Several or One sample as an output format. Select several: Users: The user name appears on the printed report (17 characters). Header: The company name can be added as a header (40 characters) and appears on the print report. Identify sensor: The sensor model and the sensor serial number can be added and appears on the print report.
GLP	Select Several or One sample as an output format. Select several: Users: The user name appears on the printed report (17 characters). Header: The company name can be added as a header (40 characters) and appears on the print report. Identify sensor: The sensor model and the sensor serial number can be added and appears on the print report.

4. When a computer is connected and ComLabo Easy (refer to for more information) is selected, use the ▲ or ▼ to select

Option	Description
Users	The user name appears on the printed report (17 characters).
Identify sensor	The sensor model and the sensor serial number can be added and appears on the print report.

Maintenance

▲ WARNING

Multiple hazards. Do not disassemble the instrument for maintenance. If the internal components must be cleaned or repaired, contact the manufacturer.

▲ CAUTION

Personal injury hazard. Only qualified personnel should conduct the tasks described in this section of the manual.

Clean the instrument

NOTICE

Never use cleaning agents such as turpentine, acetone or similar products to clean the instrument including the display and accessories.

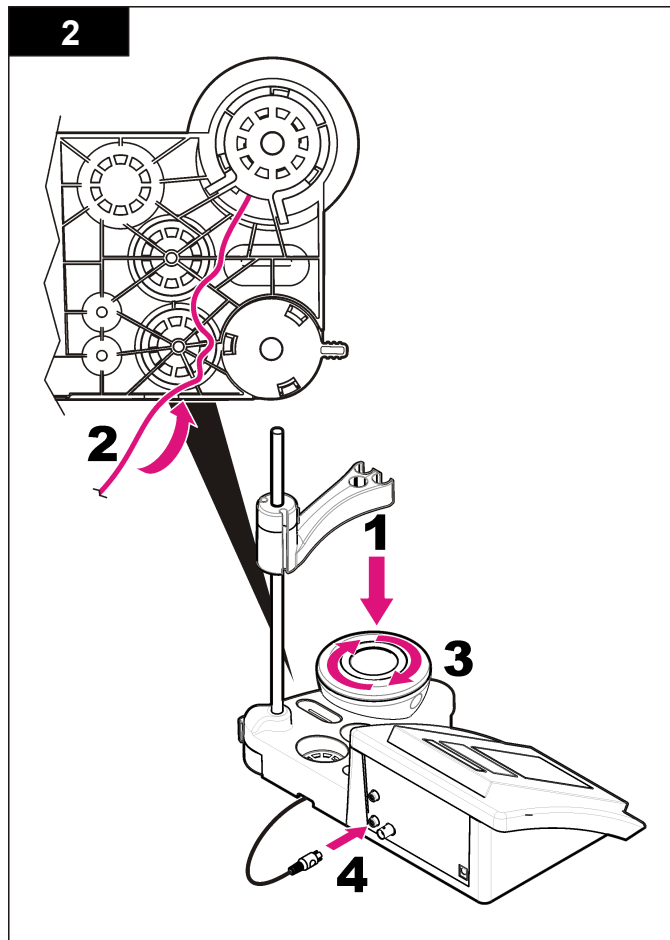
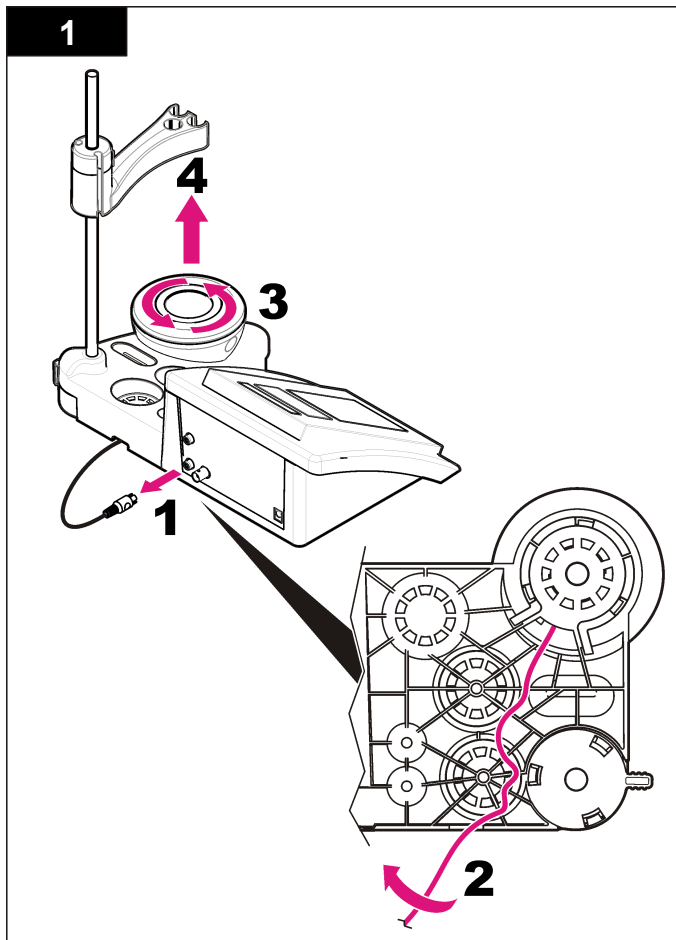
Clean the exterior of the instrument with a moist cloth and a mild soap solution.

Clean the probe

Clean the probe as needed. Refer to [Troubleshooting](#) on page 15 for more information about cleaning. Refer to the probe documentation for information about the probe maintenance.

Replace the magnetic stirrer

If the magnetic stirrer does not start, follow the numbered steps to replace the magnetic stirrer.

Troubleshooting

Refer to the following table for common problem messages or symptoms, possible causes and corrective actions.

Table 1 Calibration warnings and errors

Error/Warning	Solution
Asymmetry > 58 mV	Slope out of range (accepted values ± 58 mV). Calibrate again. Connect a new probe.
Buffer or electrode in poor conditions.	Examine the probe: Clean the probe (refer to Clean the probe on page 13 for more information); make sure that there are no air bubbles in the membrane. Shake the probe like a thermometer; connect a different probe to verify if problem is with probe or meter. Examine the buffer solution: Make sure that the buffer used matches the buffer specified in configuration; make sure of the temperature specification in configuration; use new buffer solution.
BUFFER NOT RECOGNIZED	Calibrate again. Examine the probe: Clean the probe (refer to Clean the probe on page 13 for more information); make sure that there are no air bubbles in the membrane. Shake the probe like a thermometer; connect a different probe to verify if problem is with probe or meter.
Sens. (a) < 70%	Examine the buffer solution: Make sure that the buffer used matches the buffer specified in configuration; make sure of the temperature specification in configuration; use new buffer solution.
UNSTABLE READING Time t > 100 s	Calibrate again. Examine the probe: Clean the probe (refer to Clean the probe on page 13 for more information); make sure that there are no air bubbles in the membrane. Shake the probe like a thermometer; connect a different probe to verify if problem is with probe or meter. Make sure that the membrane and the diaphragm are properly immersed in the sample.

Table 1 Calibration warnings and errors (continued)

Error/Warning	Solution
Check temperature. Check electrode	Make sure of the temperature specification in configuration. Examine the probe: Clean the probe (refer to Clean the probe on page 13 for more information); make sure that there are no air bubbles in the membrane. Shake the probe like a thermometer; connect a different probe to verify if problem is with probe or meter.
Electrode in poor conditions.	Examine the probe: Clean the probe (refer to Clean the probe on page 13 for more information); make sure that there are no air bubbles in the membrane. Shake the probe like a thermometer; connect a different probe to verify if problem is with probe or meter.
Check the electrode.	Examine the probe: Clean the probe (refer to Clean the probe on page 13 for more information); make sure that there are no air bubbles in the membrane. Shake the probe like a thermometer; connect a different probe to verify if problem is with probe or meter.
SAME BUFFERS	Calibrate again. Examine the probe: Clean the probe (refer to Clean the probe on page 13 for more information); make sure that there are no air bubbles in the membrane. Shake the probe like a thermometer; connect a different probe to verify if problem is with probe or meter. Examine the buffer solution: Use new buffer solution.

Table 2 Measurement warnings and errors

Error/Warning	Solution
pH 12.78 19°C	Calibrate again
pH out of range	Examine the probe: Clean the probe (refer to Clean the probe on page 13 for more information); make sure that there are no air bubbles in the membrane. Shake the probe like a thermometer; connect a different probe to verify if problem is with probe or meter.
Temp out of range °C	Examine the temperature sensor. Connect a different probe to verify if problem is with probe or meter.

Table 2 Measurement warnings and errors (continued)

Error/Warning	Solution
Time > 60 s	Make sure that the membrane and the diaphragm are properly immersed in the sample.
Time > 150 s	
Time > 300 s	Examine the temperature.
	Examine the probe: Clean the probe (refer to Clean the probe on page 13 for more information); make sure that there are no air bubbles in the membrane. Shake the probe like a thermometer; connect a different probe to verify if problem is with probe or meter.
10.389 mg/L >>>>>	ISE measurement: The measured value is superior to the highest standard used in calibration. Measure again.
00012 00:13	
0.886 mg/L <<<<<<	ISE measurement: The measured value is inferior to the lowest standard used in calibration. Measure again.
00018 00:11	

Replacement parts and accessories

Note: Product and Article numbers may vary for some selling regions. Contact the appropriate distributor or refer to the company website for contact information.

Replacement parts

Description	Item no.
sensION+ PH3 Lab pH-meter with accessories, without probe	LPV2000.97.0002
sensION+ PH31 Lab pH-meter, GLP, with accessories, without probe	LPV2100.97.0002
sensION+ MM340 Lab pH & Ion-meter, GLP, 2 channels, with accessories, without probe	LPV2200.97.0002
sensION+ EC7 Lab conductivity meter, with accessories, without probe	LPV3010.97.0002

Replacement parts and accessories (continued)

Description	Item no.
sensION+ EC71 Lab conductivity meter, GLP, with accessories, without probe	LPV3110.97.0002
sensION+ MM374, 2 channel Lab meter, GLP, accessories, without probes	LPV4110.97.0002

Consumables

Description	Item no.
pH buffer solution 4.01, with certificate, 125 mL	LZW9460.99
pH buffer solution 7.00, with certificate, 125 mL	LZW9461.97
pH buffer solution 10.00, with certificate, 125 mL	LZW9470.99
pH buffer solution 4.01, with certificate, 250 mL	LZW9463.99
pH buffer solution 7.00, with certificate, 250 mL	LZW9464.97
pH buffer solution 10.00, with certificate, 250 mL	LZW9471.99
pH buffer solution 4.01, with certificate, 1000 mL	LZW9466.99
pH buffer solution 7.00, with certificate, 1000 mL	LZW9467.97
pH buffer solution 10.00, with certificate, 1000 mL	LZW9472.99
Electrolytic solution (KCl 3M), 125 mL	LZW9510.99
Electrolytic solution (KCl 3M), 250 mL	LZW9500.99
Electrolytic solution (KCl 3M), 50 mL	LZW9509.99
Electrolytic solution 0.1 M, 125 mL	LZW9901.99
Enzyme solution	2964349
Pepsin Cleaning Solution	2964349
Electrode cleaning solution	2965249

Replacement parts and accessories (continued)

Description	Item no.
0.1 N HCl solution	1481253
Ethanol, 95% (Grease, oils, fats)	2378900

Accessories

Description	Item no.
Thermal printer, RS232, for sensION+ benchtop	LZW8201.99
Thermal paper for printer LZW8201, bag with 5 rolls	LZW9117.99
Dot-impact printer, RS232, for sensION+ instruments	LZW8200.99
Standard paper for printer LZW8200, bag with 10 rolls	LZW9000.99
Ribbon for printer LZW8200, 3 units	LZW9001.99
LabCom Easy PC SW, for sensION+ GLP, CD, cable, USB adapter	LZW8997.99
LabCom PC SW, for sensION+ GLP, CD, cable, USB adapter	LZW8999.99
RS232 cable, for sensION+ benchtop instruments	LZW9135.99
RS232 cable for sensION+ benchtop, with USB adapter	LZW9135USB.99
Magnetic stirrer with sensor holder, for sensION+ MM benchtop	LZW9319.99
3x50 mL printed beakers for benchtop pH calibration	LZW9110.97
Three-sensor holder, for sensION+ benchtop instruments	LZW9321.99
Holder and clamp for three sensors	LZW9155.99
Pyrex glass chamber, continuous flow measurements	LZW9118.99
PP protector, electrode storage	LZW9161.99

Standard solutions

Technical buffer solutions (DIN 19267)

Refer to [Table 3](#) pH and ORP (mV) values of specific buffer sets at varying temperatures.

Table 3 pH, ORP (mV) and temperature values

Temperature		pH					mV
°C	°F						
0	32	2.01	4.01	7.12	9.52	10.30	—
10	50	2.01	4.00	7.06	9.38	10.17	245
20	68	2.00	4.00	7.02	9.26	10.06	228
25	77	2.00	4.01	7.00	9.21	10.01	220
30	86	2.00	4.01	6.99	9.16	9.96	212
40	104	2.00	4.03	6.97	9.06	9.88	195
50	122	2.00	4.06	6.97	8.99	9.82	178
60	140	2.00	4.10	6.98	8.93	9.76	160
70	158	2.01	4.16	7.00	8.88	—	—
80	176	2.01	4.22	7.04	8.83	—	—
90	194	2.01	4.30	7.09	8.79	—	—

Buffer solutions (DIN 19266)

Refer to [Table 4](#) for pH and ORP (mV) values of specific buffer sets at varying temperatures.

Table 4 pH, ORP (mV) and temperature values

Temperature		pH					mV	
°C	°F							
5	32	1.668	4.004	6.951	7.087	9.395	10.245	13.207
10	50	1.670	4.000	6.923	7.059	9.332	10.179	13.003

Table 4 pH, ORP (mV) and temperature values (continued)

Temperature		pH					mV	
°C	°F							
20	68	1.675	4.001	6.881	7.016	9.225	10.062	12.627
25	77	1.679	4.006	6.865	7.000	9.180	10.012	12.454
30	86	1.683	4.012	6.853	6.987	9.139	9.966	12.289
40	104	1.694	4.031	6.838	6.970	9.068	9.889	11.984
50	122	1.707	4.057	6.833	6.964	9.011	9.828	11.705
60	140	1.723	4.085	6.836	6.968	8.962	—	11.449
70	158	1.743	4.126	6.845	6.982	8.921	—	—
80	176	1.766	4.164	6.859	7.004	8.885	—	—
90	194	1.792	4.205	6.877	7.034	8.850	—	—

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P.O. Box 389, Loveland, CO 80539-0389 U.S.A.
(800) 227-4224 (U.S.A. only)