

Think. Inside. The. Box.

CENTRA®-R120*



Pre-treatment

LC175/LC177

Lifetime expectancy: 6 months

What is purified from the water?

Chlorine and particulates

Impact of not changing:

Chlorine, chloramine and organics remain increasing the burden on downstream consumables which will need more frequent and expensive changes



Reverse Osmosis

LC180

Lifetime expectancy: 24-36 months

What is purified from the water?

Majority of organics and inorganics

Impact of not changing:

High levels of organics and inorganics remain in the water reducing the functional life of the DI pack



DI Pack

LC187

Lifetime expectancy: 6 months

What is purified from the water?

Ions

Impact of not changing:

Weakly bound ions are released back into the water reducing resistivity



UV light

LC118

Lifetime expectancy: 12 months

What is purified from the water?

TOC, inactivates DNA & RNA

Impact of not changing:

Uncontrolled bacterial contamination
TOC levels not reduced



Composite Vent Filter

LC136M2

Lifetime expectancy: 6 months

What protection does it provide?

Reduces CO₂ and airborne organics from entering the sealed system

Impact of not changing:

Increase in airborne contaminants and potential recontamination of purified water



Ultra-Micron Filter

LC109

Lifetime expectancy: 6 months

What is purified from the water?

Particles & Microorganisms

Impact of not changing:

Reduced flow-rates and bacterial specification will not be maintained

*A combination of treatments and technologies produce the final water specification. Optional degassing membrane available