## ESPAÑOL

### Instrucciones

#### INSTALACIÓN

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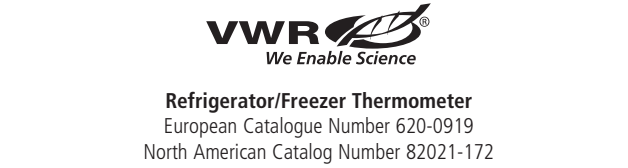
### Instructions

#### INSTALLATION

Pull down the battery cover, install the battery with correct polarity, and close the battery compartment. The thermometer will have a full segment display for about 3 seconds and then display the fridge sensor temperature normally. Place probe in test condition, record your reading when the reading becomes stable. Attach the digital display portion of thermometer onto the desired space outside the refrigerator/freezer by using the supplied tape or magnets on the back of thermometer. Peel off the protective film on the display.

#### OPERATION

**°C/°F Exchange:**

Press the [°C/°F] to select temperature unit.

#### Maximum/Minimum Memory:

- Press and hold [MAX/MIN] for about 2 seconds, the HI and FRIDGE icons and numeric display will flash.
- Press [MAX/MIN] once more to show the minimum reading.
- Press [MAX/MIN] again to return to normal display.
- Press and hold [MAX/MIN] about 2 seconds to reset memory to current reading.

#### Room or Fridge/Freezer Temperature Display:

Press [ROOM/FRIDGE] button to display room or refrigerator/freezer temperature.

#### Fridge/Freezer High and Low Temperature Alarm Setting:

- Press and hold [alarm] for about 2 seconds, the HI and FRIDGE icons and numeric display will flash.
- Press [up arrow] to set the alarm high limit. Arrow is on the back of the unit above the battery cover.
- Press [alarm] again, the LO and FRIDGE icons and numeric display will flash.
- Press [up arrow] to set the alarm low limit.
- Press [alarm] once more to finish setting.

\* During alarm limit setting, holding down [up arrow] will advance the value automatically. If the fridge sensor reading is out of the set limit, the alarm will sound.

#### Alarm On/Off:

Press [alarm] to switch alarm off or on.

#### NOTE:

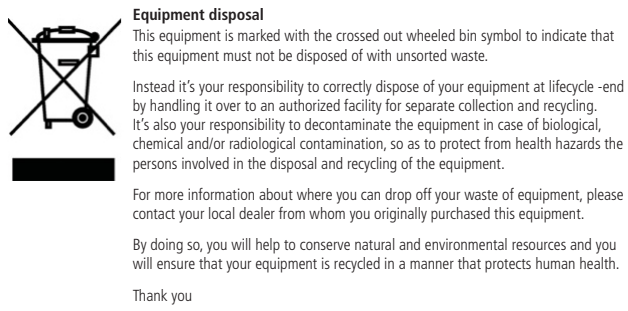
- If the display becomes dim, replace the battery with same type of battery.
- After each measurement, clean the probe immediately to avoid cross contamination.
- Do not put the digital display thermometer unit inside the freezer, refrigerator, incubator or oven.

#### ERROR SYMBOLS

Symbol	Description	Action required
X	Low battery voltage	Replace the batteries
LLL	1) Sensor open circuit 2) The reading is out of low range (-50°C)	1) Return the thermometer for repair 2) Keep the measurement above -50°C
HHH	1) Sensor short circuit 2) The reading is out of high range (70°C)	1) Return the thermometer for repair 2) Keep the measurement below 70°C

#### Battery Handling/Packaging

- Remove batteries from equipment immediately after they fail to operate the equipment.
  - If equipment uses two or more batteries, always replace batteries in complete sets.
  - Do not attempt to recharge alkaline batteries, they are not rechargeable.
  - Do not handle hot or warm batteries.
  - Do not heat, incinerate, crush, puncture or mutilate batteries.
  - Wear personal protective equipment if batteries show signs of leakage, bulging, swelling or deformity.
- Warning:** If potassium hydroxide electrolyte from an alkaline battery comes in contact with the skin, do not try to neutralize the electrolyte with vinegar or any other acidic solutions. Neutralization may trap electrolyte on the skin. Flush the affected skin area with copious amounts of water. If the battery electrolyte gets into your eyes, it can cause severe damage and or blindness.

#### Technical service

#### Web Resources

Visit the VWR website at **www.com** for:

- Complete technical service contact information
- Access to the VWR Online Catalogue, and information about accessories and related products
- Additional product information and special offers

**Contact us** For information or technical assistance contact your local VWR representative or visit **www.com**.

#### Warranty

VWR warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

#### Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

<b>Manufactured for</b>		
<b>Europe</b> VWR International bvba Researchpark Haasrode 2020 Geldenaaksebaan 464 B-3001 Leuven + 32 16 385011 http://www.vwr.com	<b>United States</b> VWR International 1310 Goshen Parkway West Chester, PA 19380 800-932-5000 http://www.vwr.com	