



MCO-80ICL-PA | 10046-928

CytoGrow ReachIn

CO₂ Incubator

30.1 cu.ft. | 851 L



The ReachIn CO₂ incubator is designed to save lab space while offering the flexibility to grow your culturing needs. The control system provides a precise and repeatable chamber environment.

Horizontal Laminar Airflow

The cross shelf, horizontal laminar airflow system promotes optimum temperature uniformity throughout the chamber which contributes to quick temperature recovery after door openings. Side plenum walls are perforated for positive, even airflow. The horizontal airflow allows for full use of solid shelves providing maximum culturing capacity.

Contamination Control

Incubator interior, shelves and plenum system incorporate inCu-saFe[®] copper enriched, stainless steel alloy which offers germicidal properties of elemental copper. InCu-saFe prevents corrosion and discoloration of the incubator interior.


Control and Monitoring

The incubator functions are managed by a fully integrated microprocessor controller with a range of setpoints and alarms for temperature and CO₂. Two levels of elevated humidity can be selected. All functions are programmable through the control panel. Independent LED displays for CO₂ and temperature are standard.



CO₂ Control

IR sensor coupled with the PID microprocessor controller is used to provide superior CO₂ control and exceptionally fast recovery after door openings. Gas system design minimizes CO₂ consumption.

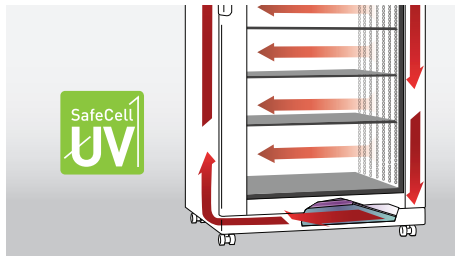

Optional UV Protection

Located safely below the interior base, the optional patented SafeCell™ UV lamp destroys airborne contaminants as they pass over the humidity reservoir surface. Pathogens introduced during door openings are ultimately removed.



User Convenience

Softkey eight button menu for intuitive programming. LED indicators monitor incubator functions. Remote alarm contacts provided standard.


Large Capacity CO₂ Incubator

The MCO-80ICL-PA | 10046-928 is ideal for culturing large volumes of biological samples, performing microbiological studies and working with large volume cell apparatus.

Humidity Control

Humidity reservoir heaters are located outside the walls of the reservoir. They are not as susceptible to corrosion and scaling from water as competitive systems are. The unit can be set to normal (80% RH or above) or high (90% RH or above) humidity mode.


Rapid CO₂ Density Recovery

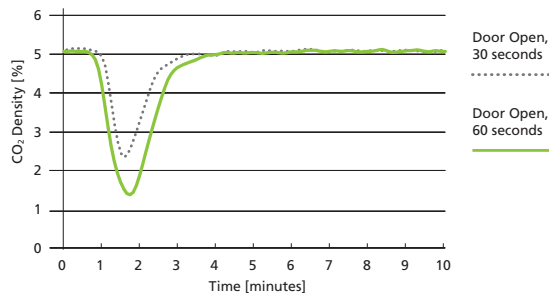

IR sensor achieves ultrafast CO₂ recovery without overshoot, even following multiple door openings.

Dimensions

Unit : inch | mm

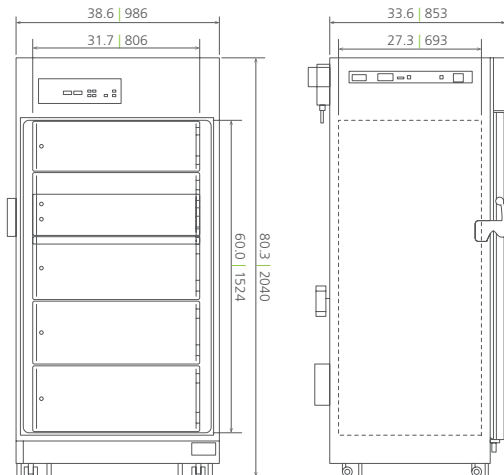

1.800.932.5000 | VWR.COM

Prices and product details are current when published and subject to change without notice. | Certain products may be limited by federal, state, provincial, or local regulations. | VWR, part of Avantor, makes no claims or warranties concerning sustainable/green products. Any claims concerning sustainable/green products are the sole claims of the manufacturer and not those of VWR International, LLC and/or Avantor, Inc. or affiliates. All prices are in US dollars unless otherwise noted. Offers valid in US and Canada, void where prohibited by law or company policy, while supplies last. | Trademarks are owned by Avantor, Inc. or its affiliates, unless otherwise noted. | Visit vwr.com to view our privacy policy, trademark owners, and additional disclaimers. © 2019 Avantor, Inc. All rights reserved.

Model Number	MCO-80ICL-PA 10046-928	With Optional UV Decontamination
External Dimensions (W x D x H) ¹⁾	inches mm	38.6 x 37.2 x 80.3 986 x 945 x 2040
Internal Dimensions (W x D x H)	inches mm	31.7 x 27.3 x 60.0 806 x 693 x 1524
Volume	cu.ft. liters	30.1 851
Net Weight	lbs kg	606 275
Performance		
Warranty ²⁾		3 years parts and labor
Temperature Control Range and Fluctuation ³⁾	°C	+5 above ambient to +50, ± 0.1
Temperature Uniformity ³⁾	°C	± 0.5 (9 point measurement)
CO ₂ Control Range and Fluctuation ³⁾	%	0 to 20, ± 0.15
Humidity Level & Fluctuation	% RH	Normal mode: over 80 (high mode: over 90)
Control		
Controller		Microprocessor
Temperature Sensor		Thermistor
Display	qty	2; LED (1 for temperature and 1 for CO ₂) readable to 0.1 increments
Sensor	CO ₂	IR
Construction		
Exterior Material		Painted steel
Interior Material		Stainless steel copper enriched alloy
Outer Door	qty	1; Dual pane heated glass with latch (provision for padlock)
Shelves	qty	5 (stainless steel copper enriched alloy)
Shelf Dimension (W x D x H)	inches mm	30.6 x 25.9 x 0.4 776 x 659 x 10
Max. Load per Shelf - Distributed Evenly	lbs kg	66.1 30
Max. Total Load	lbs kg	330.0 150
Max. Shelf Capacity	qty	18
Access Port / Position	qty	2; right side and left side
Access Port Diameter	inches mm	1.6 40 (with silicone (non-VOC) stopper)
Casters	qty	4; Dual wheel (front 2 swivel, back 2 fixed)
Leveling Feet	qty	4
Humidity Reservoir Drain	qty	Drain valve – lower side front (drainage tray provided)
Humidity Reservoir Material		Stainless steel
Decontamination Control		
InCu-safe Chamber, Air Plenum and Shelves	passive	Included (stainless steel copper enriched alloy)
SafeCell UV Light System	passive/active	Optional Included
Alarms (V=Visual Alarm, Buzzer Alarm, R=Remote Alarm)		
Power Failure		R
Temperature Deviation	high	V-B-R
Gas Deviation	CO ₂	V-B-R
Supply Empty	CO ₂	V-B-R
Door Open		V-B
Low Water in Humidity Reservoir		V
Electrical and Noise Level		
Power Supply		115V, 1Ø, 60Hz, NEMA 5-20P requires NEMA 5-20R receptacle
Noise Level ⁴⁾	dB(A)	33
Outlets		
Chamber Duplex – Vapor Proof Cover	qty	1; 115v 3 amps max rating
Cabinet Outlet	qty	1; 115v 1 amps max rating
Options		
SafeCell UV Light System		MCO-80UVS-PA Included
Humidity Reservoir–Auto Fill System		MCO-80AS-PW
CO ₂ Gas Pressure Regulator	psi	0 – 15; MCO-100L 10119-838
Automatic CO ₂ Cylinder Changeover System		MCO-80GC-PW 10119-834
Inner Door Kit		MCO-80ID-PW
InCu-safe Solid Shelf with Brackets		MCO-80ST-PW
InCu-safe Solid Shelf Reinforced with Brackets		MCO-80AICRSLF (Max 3 Shelves)
Cell Roller Mounting Ramp Kit		MCO-80RBS-PW
4-20mA Analog Output		MCO-420MA-PW 10119-842
Optional Communication System		
Wireless, Cloud-based, Automatic Data Management		LabAlert® Monitoring System

¹⁾ Exterior dimensions of cabinet excluding handle, rear stand-off brackets and other external projections. Consult sales rep for doorway entry instructions, less than 37.2".

²⁾ Current warranty offered at time of printing and may be subject to change; US and Canada only

³⁾ Ambient temperature (in a +20°C to +35°C), setting 37°C, CO₂ 5%, no load, air temperature measured at incubator center

⁴⁾ Nominal value – Background noise 20dB(A)

Note: Additional options available.

Specifications are subject to change without notice.

Performance data herein is based on independent testing at time of publication.