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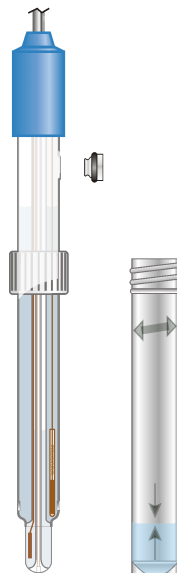
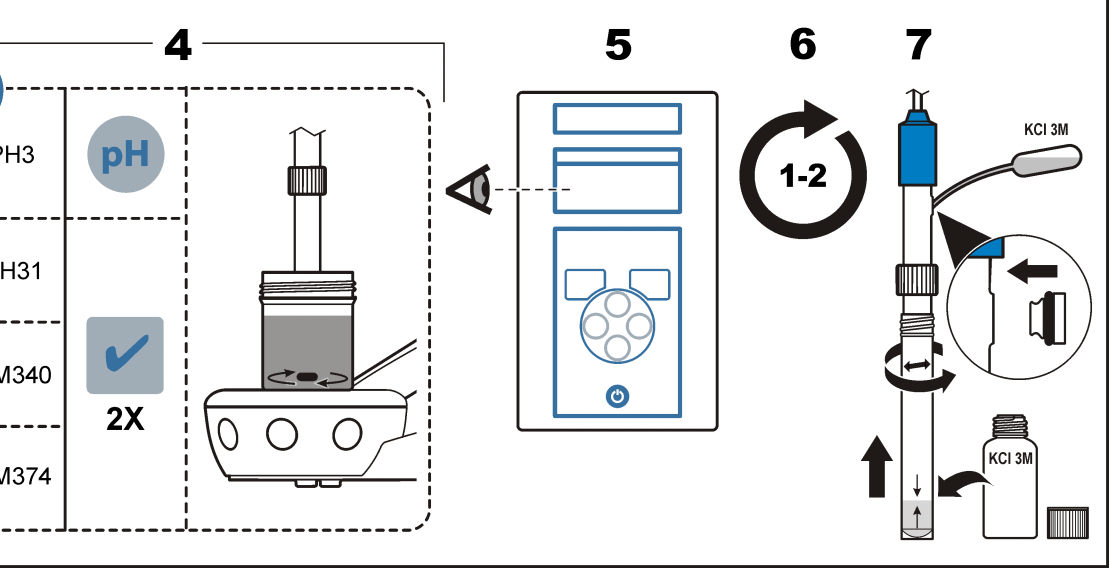
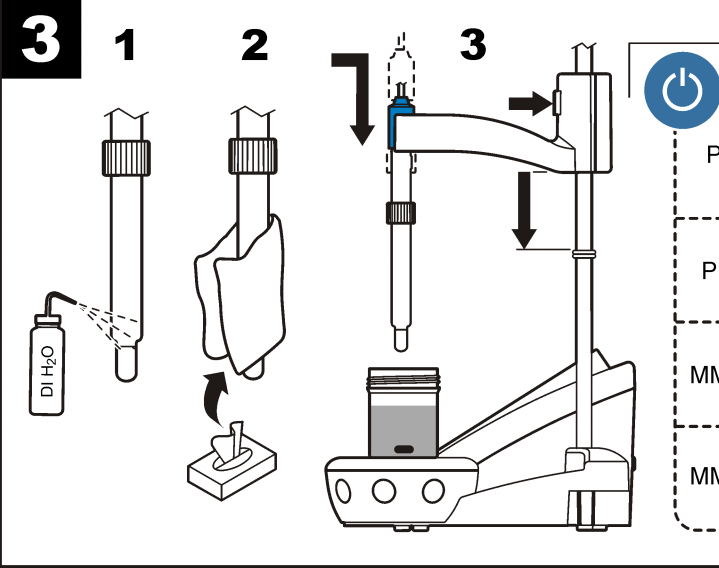
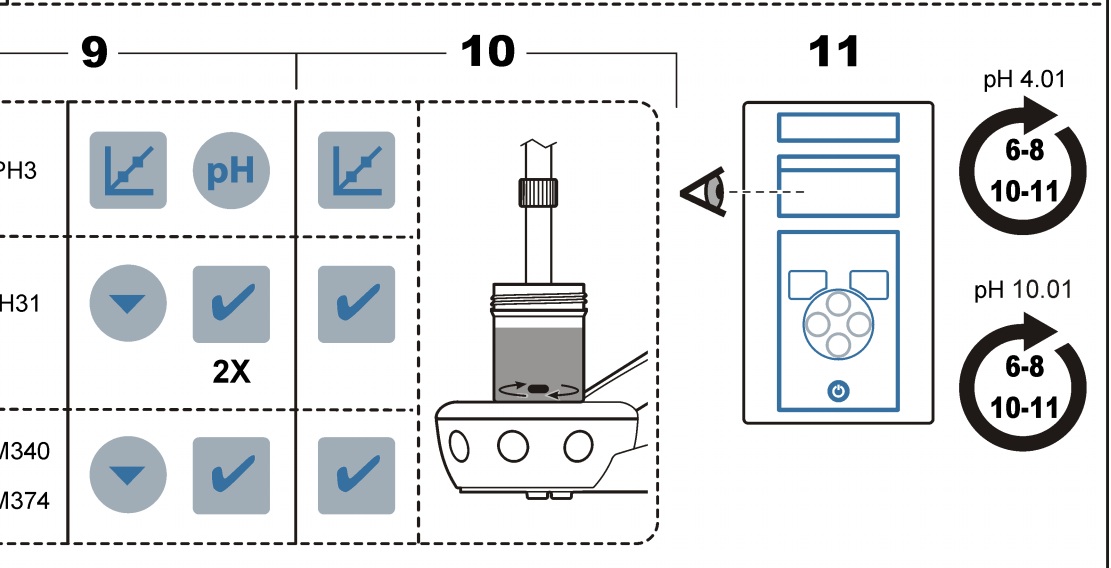
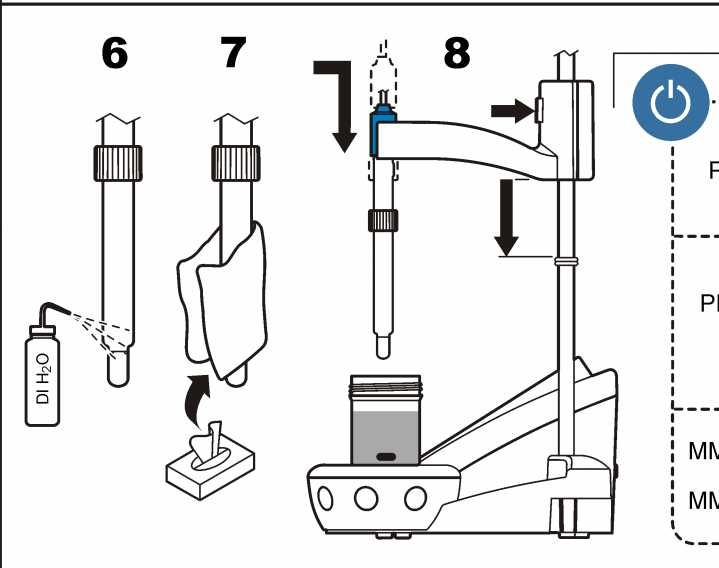
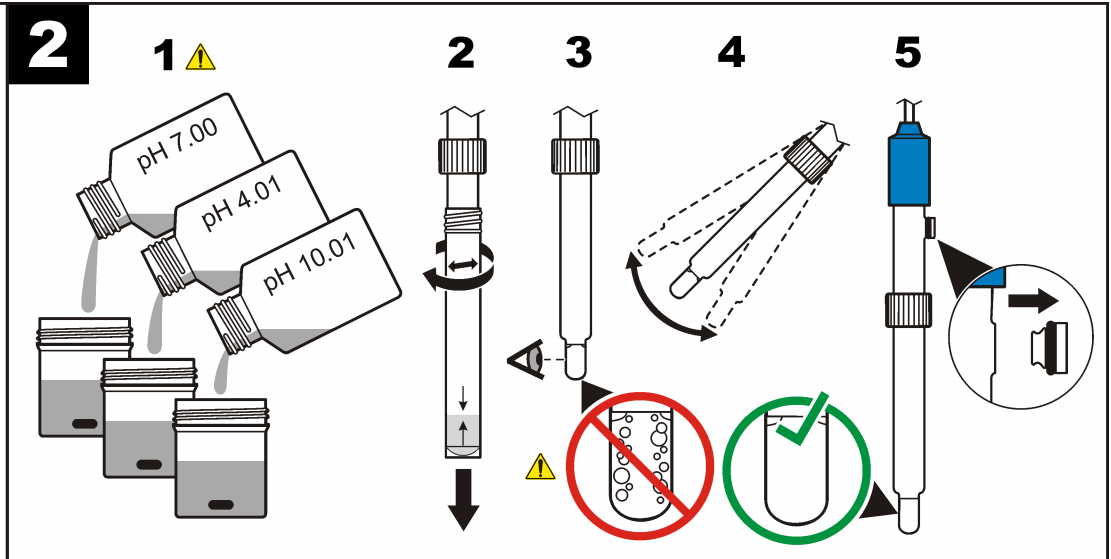
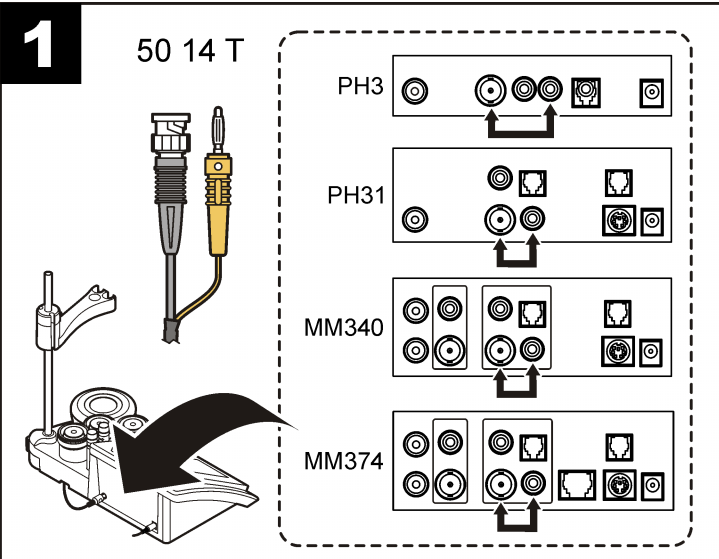
**50 14** EN

Refer to the meter manual for more information.

**Caution**  
 ⚠ Personal injury hazard. Glass components can break. Handle with care to prevent cuts.  
 ⚠ Chemical exposure hazard. Refer to the current material safety data sheets (MSDS) for safety protocols.

**Introduction**  
 This pH refillable probe with temperature sensor is used in general water samples, samples with sulfides and samples with reduction substances.

**Technical information**

Range	2 to 14 pH
Junction	Ceramic (2)
Electrolyte	KCl, 3 M
Reference	Ag ion barrier cartridge
Temperature	Pt 1000
Minimum depth	20 mm (0.8 in.)
Operating temp	-10 to 100 °C (14 to 212 °F)
Limits	Viscous solutions, distilled water, samples with colloids

Range	2 to 14 pH
Junction	Ceramic (2)
Electrolyte	KCl, 3 M
Reference	Ag ion barrier cartridge
Temperature	Pt 1000
Minimum depth	20 mm (0.8 in.)
Operating temp	-10 to 100 °C (14 to 212 °F)
Limits	Viscous solutions, distilled water, samples with colloids

**1 Connect the probe**  
**2 Calibration**  
 ⚠ Air bubbles in or under the probe tip when submerged can cause slow stabilization or error in measurement.

**3 Measurement**  
**Maintenance and storage**  
 Cleaning: soak the probe in a cleaning solution. Rinse with deionized water.

Contamination	Cleaning agent
Proteins	Pepsin cleaning solution
Grease, oils, fats	Electrode cleaning solution
Limescale	0.1 N HCl solution

Hydration: soak the dry probe in a dilute HCl solution for several hours.

Storage: put the probe in the protector with KCl 3 M solution.

**Troubleshooting**  
 Be sure that the samples are suitable for the application.

**Warranty**  
 1 year for manufacturing faults only. Damage from use is not covered.

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