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INSTRUCTIONS FOR
MASTERFLEX® E/S® AUTOMOTIVE POWER ADAPTER
MODEL NO. 07571-50

 **WARNING: PRODUCT USE LIMITATION** 

This product is not designed for, nor intended for use in, patient-connected applications, including, but not limited to, medical and dental use and, accordingly, has not been submitted for FDA approval.

INTRODUCTION AND GENERAL DESCRIPTION

The MASTERFLEX Automotive Power Adapter is intended to provide 12V DC power to the Portable Sampling Drives 07571-00 and 07571-05 from an automotive electrical system through the cigarette lighter present in most foreign and domestic vehicles.

The Automotive Power Adapter is current limited by the customer replaceable fuse. The fuse is a fast acting type F (3AG) rated @ 5.0A 250V.

FUSE REPLACEMENT

1. Disconnect the adapter cable from the cigarette lighter in the vehicle. Disconnect the other end from the Portable Sampling Drive.
2. To access and replace the fuse, unscrew the **ADAPTER TIP** from the **ADAPTER BODY**, the **FUSE** will slide out and may be replaced. The MASTERFLEX fuse replacement part number is 77500-03.

 **WARNING:** Replace the fuse only with same type and rating as the original. Use of a fuse type of rating other than specified can result in damage to equipment or injury to the operator.

3. To install, reverse the replacement process.

ORIGINAL INSTURCTIONS

A-1299-0829B
Edition 08

Trademarks bearing ® symbol in this publication are registered in the U.S. and in other countries.

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Masterflex LLC