
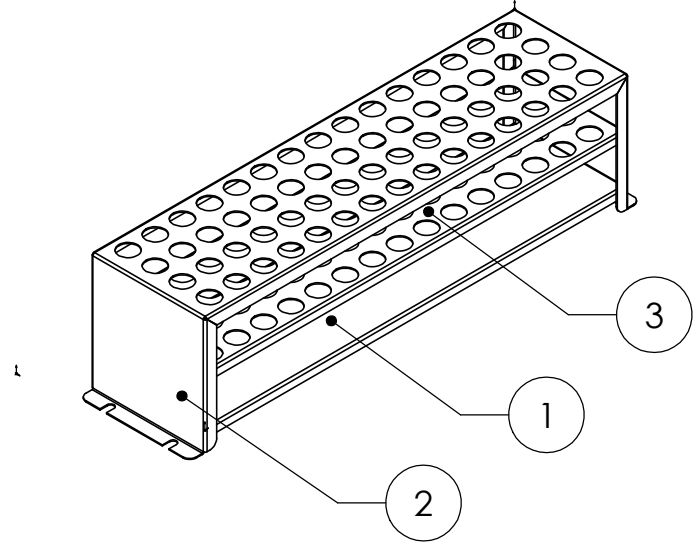


DWG BY: THA DATE: 3/18/02 TOLERANCE UNLESS OTHERWISE SPECIFIED .X = ± .063 ANGULAR .XX = ± .032 HOLE DIA .XXX = ± .015 ± .005	SHEL-LAB SHELDON MANUFACTURING CORNELIUS, OREGON 14-16 MM SHAKER ASSY	SCALE: 1:4 PART#: 9751178 MATERIAL: ASSEMBLY	NOTE BEND RADIUS .032" UNLESS OTHERWISE NOTED
--	---	--	---

CHG. BY	CHG. #	REV.	DATE	DESCRIPTION

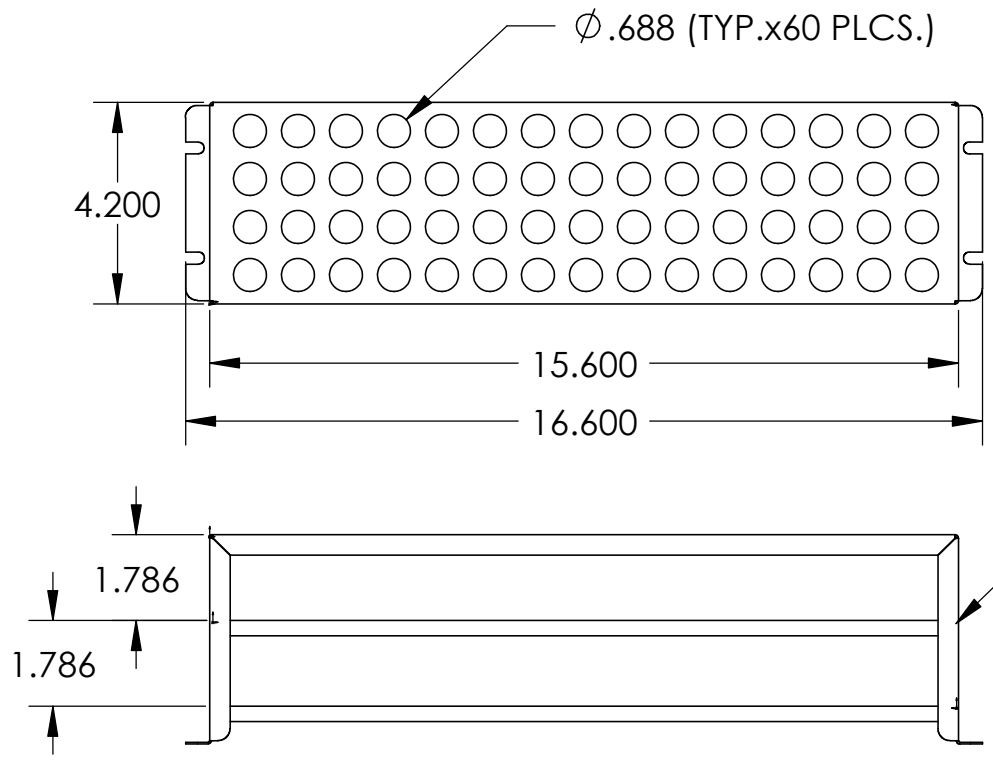

ITEM NO.	QTY.	PART NO.	MATERIAL	DESCRIPTION
1	1	5121162	20 SST 304 2B	SHAKER SHELF
2	1	5121165	.036 SST. 304 2B	SHAKER TOP 14-16MM
3	1	5121166	.036 SST. 304 2B	SHAKER SHELF 14-16 MM