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


www.hach.com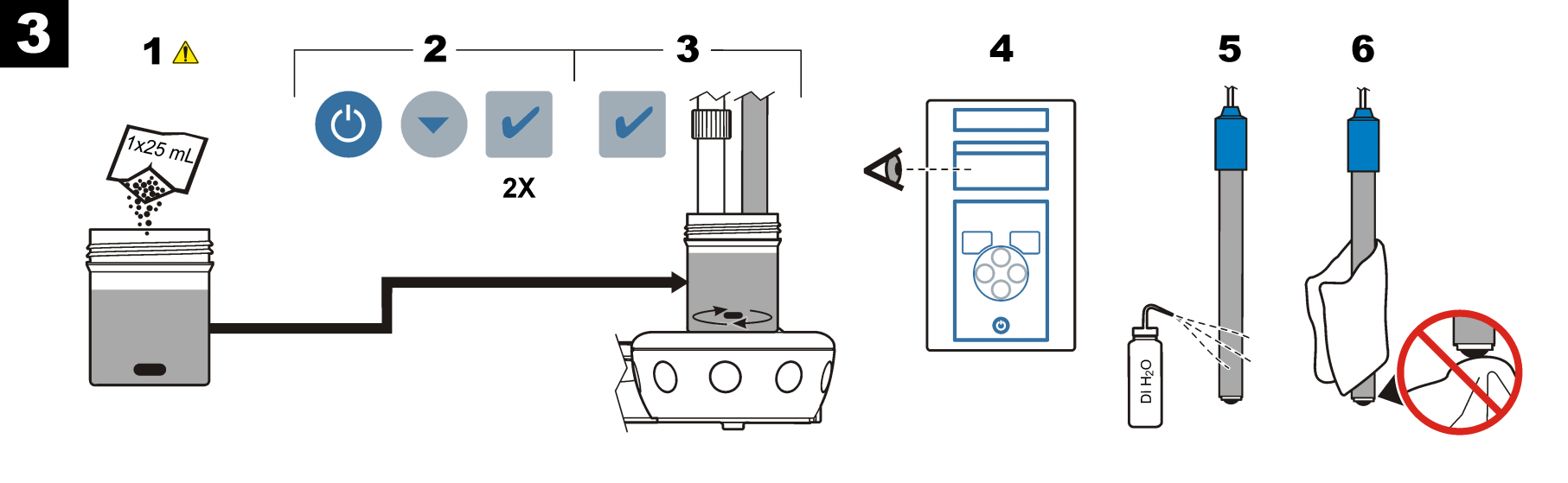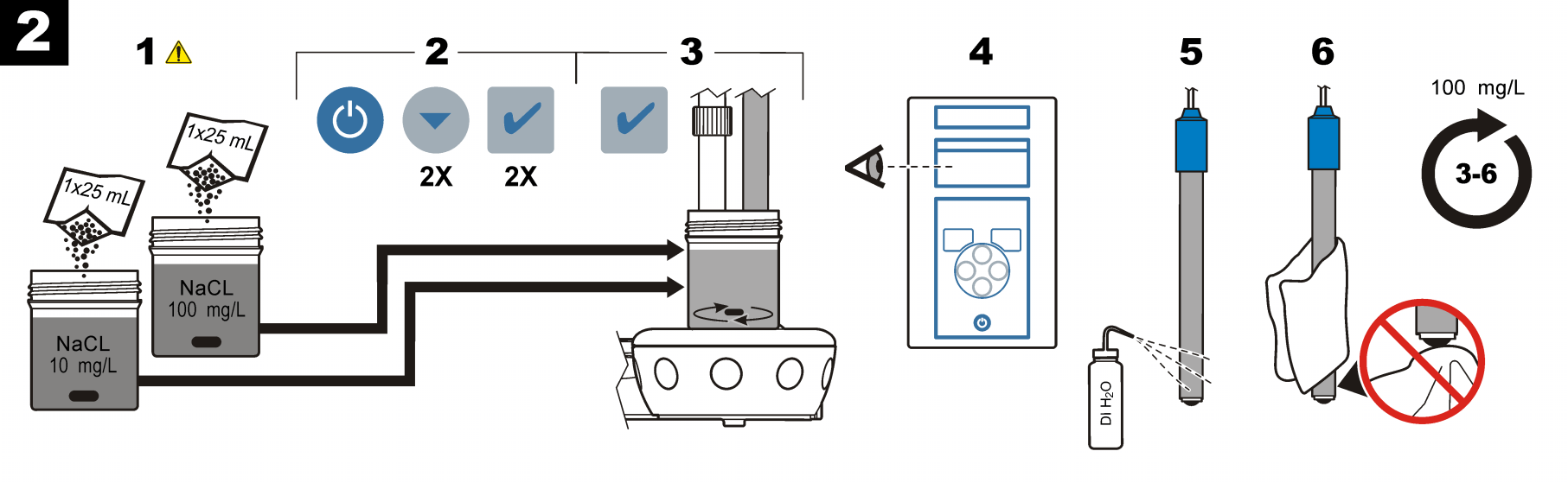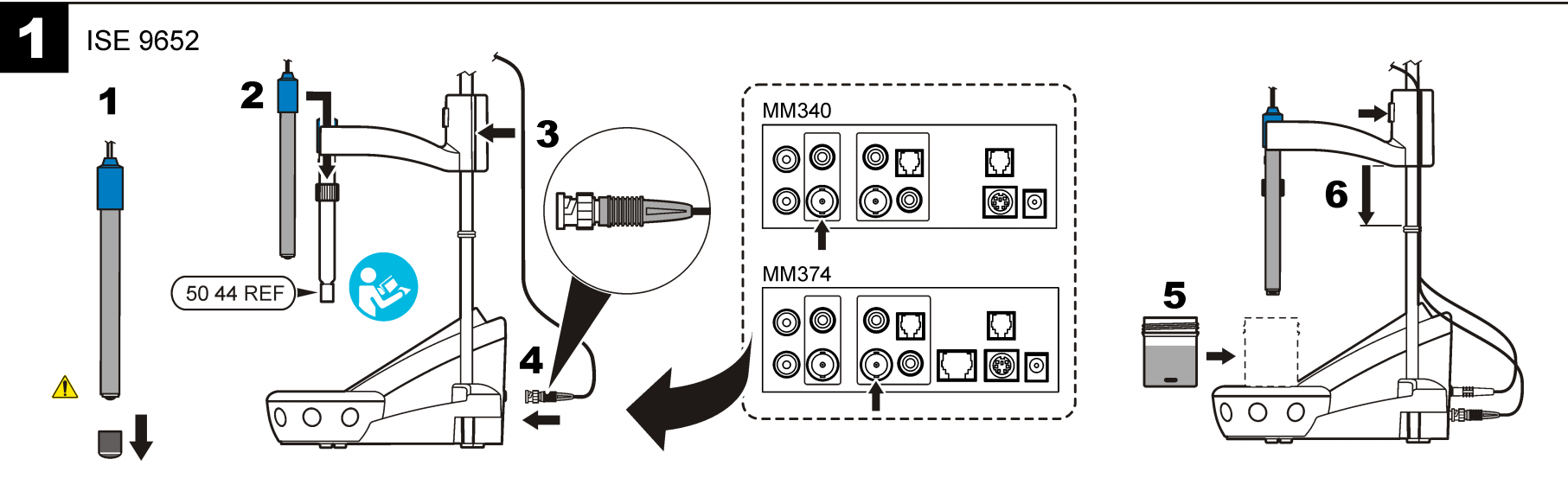

## ISE Cl<sup>-</sup> 9652 EN

Refer to the meter manual for more information.

**Caution**  
 ⚠ Personal injury hazard. Glass components can break. Handle with care to prevent cuts.  
 ⚠ Chemical exposure hazard. Refer to the current material safety data sheets (MSDS) for safety protocols.

**Introduction**  
 This probe measures chloride concentration in drinking and subterranean water, wastewater and food. Use with a reference electrode.

**Technical information**

Range	1 mg/L to 3.5 g/L Cl <sup>-</sup>
pH range	3 to 9
Linear range	3.5 mg/L to 3.5 g/L
Slope	58 mV/decade, 90 to 110% at 25 °C (77 °F)
Ionic strength adjuster (ISA)	Hach ISA 2318069
Operating temp	5 to 50 °C (41 to 122 °F)
Interferences	I <sup>-</sup> , Br <sup>-</sup> , CN <sup>-</sup> , Ag <sup>+</sup> , S <sup>2-</sup> ; remove with 1:1 NaBrO <sub>3</sub>

**Preparation**  
 Soak the electrode in a 0.01 M NaCl solution for 20 to 30 minutes.

**1 Connect the probe**

**2 Calibration**  
 Add 1 ISA pillow per 25 mL of standard.

**3 Measurement**  
 Add 1 ISA pillow per 25 mL of sample.

**Maintenance and storage**  
 Cleaning: rinse with distilled water between each standard and each sample.  
 Storage: clean and dry the probe for storage.

Damage or deposits on the membrane change the response of the electrode.

**Troubleshooting**  
 The ionic strength and temperature of samples and standards must be the same. Make sure that the stir speed is the same for all standards and samples.

**Warranty**  
 6 months for manufacturing faults only. Damage from use is not covered.

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