

## Instruction Manual

Product	Size	Catalog Number
Human Cardiac Myocytes (HCM)	500,000 cryopreserved cells	C-12810
	500,000 proliferating cells	C-12811

### Product Description

Cardiac muscle, a unique type of striated muscle in the human heart, is composed of single Cardiac Myocytes (HCM). Because they are responsible for the myogenic contraction of the whole cardiac muscle, Cardiac Myocytes are the most physically energetic cells in the body. PromoCell Human Cardiac Myocytes are produced at PromoCell's cell culture facility from normal human ventricle tissue of the adult heart. Unlike freshly isolated rod-shaped myocytes, PromoCell HCM are suitable for long-term experiments, as they are prepared according to a special protocol.

Shortly after isolation, all PromoCell Human Cardiac Myocytes are cryopreserved at passage 2 (P2) using PromoCell's proprietary serum-free freezing medium, Cryo-SFM. Each cryo vial contains more than 500,000 viable cells after thawing.

Proliferating cell cultures are made from

500,000 cryopreserved cells that have been thawed and cultured for three days at PromoCell.

### Quality Control

Rigid quality control tests are performed for each lot of PromoCell Human Cardiac Myocytes.

They are tested for cell morphology, adherence rate, and cell viability. Furthermore, flow cytometric analyses for cell-type specific markers, e.g. sarcomeric  $\alpha$ -actinin and slow muscle myosin, are carried out for each lot (see page 5). Growth performance is tested through multiple passages up to 15 population doublings (PD) under culture conditions without antibiotics and antimycotics.

In addition, all cells have been tested for the absence of HIV-1, HIV-2, HBV, HCV, HTLV-1, HTLV-2 and microbial contaminants (fungi, bacteria, and mycoplasma).

A detailed certificate of analysis (CoA) for each lot can be downloaded at: [www.promocell.com/coa](http://www.promocell.com/coa)

### Intended Use

PromoCell Human Cardiac Myocytes are for *in vitro* research use only and not for diagnostic or therapeutic procedures.

### Warning

Although tested negative for HIV-1, HIV-2, HBV, HCV, HTLV-1 and HTLV-2, the cells – like all products of human origin – should be handled as potentially infectious. No test procedure can completely guarantee the absence of infectious agents.

### Follow appropriate safety precautions!

*After delivery, start immediately with the protocol for cryopreserved cells (see page 2) or the protocol for proliferating cells (see page 3).*

*Start immediately after delivery.  
Use aseptic techniques and a laminar flow bench.*

### Protocol for Cryopreserved Cells

**Straight after arrival, store the cryopreserved cells in liquid nitrogen or seed them immediately.**

**Note:** Storage at  $-80^{\circ}\text{C}$  is not sufficient for cell preservation and causes irreversible cell damage.

#### 1. Prepare the medium

Calculate the needed culture surface area according to the plating density (see page 5) and the lot-specific cell numbers stated on the certificate of analysis. Fill the appropriate volume of PromoCell Growth Medium (at least 9 ml per vial of cells) in cell culture vessels. Place the vessels in an incubator ( $37^{\circ}\text{C}$ , 5%  $\text{CO}_2$ ) for 30 minutes.




#### 2. Thaw the cells

Remove the cryovial from the liquid nitrogen container and immediately place it on dry ice – even for short transportation. Under a laminar flow bench, briefly twist the cap a quarter turn to relieve pressure, then re-tighten. Immerse the vial into a water bath ( $37^{\circ}\text{C}$ ) just up to the screw cap for 2 minutes. Ensure that no water enters the thread of the screw cap.



#### 3. Disinfect the vial and seed the cells

Thoroughly rinse the cryovial with 70% ethanol under a laminar flow bench. Then, aspirate the excess ethanol from the thread area of the screw cap. Open the vial and transfer the cells to a cell culture vessel containing the prewarmed medium from step 1.

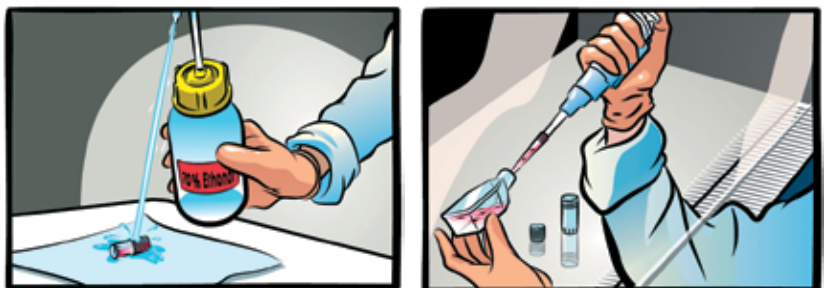

#### 4. Incubate the cells

Place the vessel in an incubator ( $37^{\circ}\text{C}$ , 5%  $\text{CO}_2$ ) for cell attachment. Replace the medium after 16-24 hours. The cells should be subcultured according to the subcultivation protocol (see page 4) once they have reached 70–90% confluency.



*Start immediately after delivery.  
Use aseptic techniques and a laminar flow bench.*

### Protocol for Proliferating Cells

#### 1. Incubate the cells

Unpack the culture vessel, do not open the lid, and immediately place it in an incubator (37°C, 5% CO<sub>2</sub>) for 3 hours to allow the cells to recover from the transportation.




#### 2. Replace the transport medium

Carefully open the vessel, rinse the inner side of the lid with 70% ethanol, and let air dry. Aspirate the transport medium from the vessel. Add 10 ml of the appropriate PromoCell Cell Growth Medium.



#### 3. Check and incubate the cells

Check the cell density. Open the lid half a turn and place the vessel in an incubator (37°C, 5% CO<sub>2</sub>). The cells should be subcultured according to the subcultivation protocol (see page 4) once they have reached 70–90% confluency.

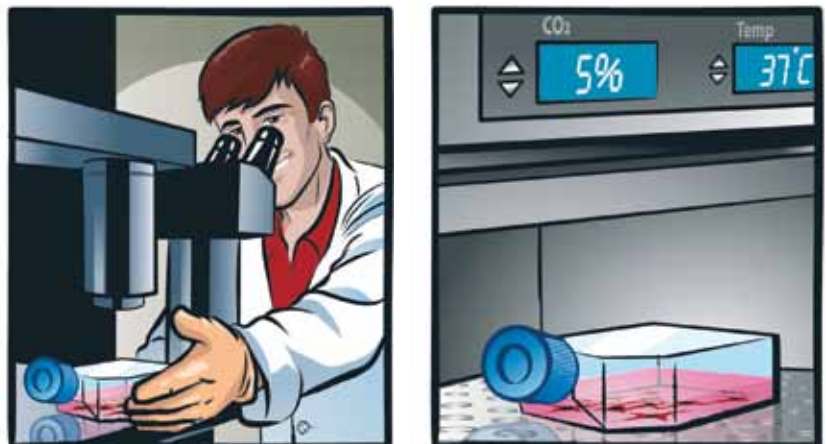

*Use aseptic techniques and a laminar flow bench.*

## Subcultivation Protocol

### 1. Prepare the reagents and wash the cells

Place the PromoCell DetachKit at room temperature for at least 30 minutes to adjust the temperature of the reagents. Carefully aspirate the medium from the culture vessel. Add 100  $\mu$ l Hepes BSS Solution per  $\text{cm}^2$  of vessel surface to wash the cells and agitate the vessel carefully for 15 seconds.



### 2. Detach the cells

Carefully aspirate the Hepes BSS from the culture vessel. Add 100  $\mu$ l Trypsin/EDTA Solution per  $\text{cm}^2$  of vessel surface. Note: We recommend detaching the cells at room temperature. Close the vessel and examine the cells under a microscope. When the cells start to detach, gently tap the side of the vessel to loosen the remaining cells.




### 3. Neutralize the trypsin and harvest the cells

Add 100  $\mu$ l Trypsin Neutralization Solution per  $\text{cm}^2$  of vessel surface and gently agitate. Carefully aspirate the cell suspension and transfer it to a centrifugation tube. Spin down the cells for 3 minutes at 220 x g.



### 4. Incubate the cells

Discard the supernatant (step 1), add 1 ml of the appropriate PromoCell Cell Growth Medium (step 2), and resuspend the cells by carefully pipetting up and down. Plate the cells according to the recommended seeding density in new cell culture vessels containing PromoCell Cell Growth Medium pre-warmed to 37°C. Place the vessels in an incubator (37°C, 5%  $\text{CO}_2$ ).

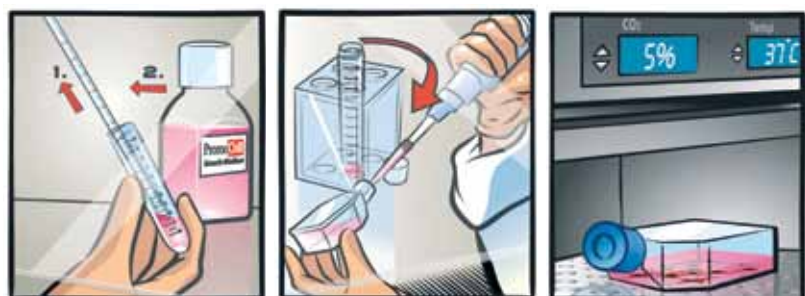

## Specifications

Product	Recommended Culture Media*	Plating Density	Passage after Thawing	Marker	Population Doublings
Human Cardiac Myocytes (HCM)	C-22070	10,000 - 15,000 cells per cm <sup>2</sup>	P2	Sarcomeric $\alpha$ -Actinin <sup>+</sup> Slow muscle myosin <sup>+</sup>	> 15

## Related Products

Product	Size	Catalog Number
Myocyte Growth Medium (Ready-to-use)	500 ml	C-22070
Myocyte Growth Medium Kit	500 ml	C-22170
Myocyte Basal Medium	500 ml	C-22270
Myocyte Basal Medium, phenol red-free	500 ml	C-22275
Myocyte Growth Medium SupplementMix	for 500 ml	C-39275
Myocyte Growth Medium SupplementPack	for 500 ml	C-39270
DetachKit	30 ml	C-41200
	125 ml	C-41210
	250 ml	C-41220
Cryo-SFM	30 ml	C-29910
	125 ml	C-29912
HCM Pellet	1 million cells per pellet	C-14080

\*The catalog numbers in this table are for media in ready-to-use packaging.

**PromoCell GmbH**

Sickingenstr. 63/65  
69126 Heidelberg  
Germany

Email: [info@promocell.com](mailto:info@promocell.com)  
[www.promocell.com](http://www.promocell.com)

**USA/Canada**

Phone: 1 – 866 – 251 – 2860 (toll free)  
Fax: 1 – 866 – 827 – 9219 (toll free)

**Deutschland**

Telefon: 0800 – 776 66 23 (gebührenfrei)  
Fax: 0800 – 100 83 06 (gebührenfrei)

**France**

Téléphone: 0800 90 93 32 (ligne verte)  
Téléfax: 0800 90 27 36 (ligne verte)

**United Kingdom**

Phone: 0800 – 96 03 33 (toll free)  
Fax: 0800 – 169 85 54 (toll free)

**Other Countries**

Phone: +49 6221 – 649 34 0  
Fax: +49 6221 – 649 34 40