VWR[®] Thermometer

INSTRUCTION MANUAL

**North American Catalog Numbers: 61161-364, 61161-336,
89094-778**

European Catalogue Numbers: 620-1661, 620-2023

Version: 1
Issued: 5th January 2015

Legal Address of Manufacturer

United States

VWR International
Radnor Corporate Center
Building One, Suite 200
100 Matsonford Road, P.O. Box 6660
Radnor, PA 19087
800-932-5000
<http://www.vwr.com>
Country of origin - USA

Europe

VWR International bvba
Researchpark Haasrode 2020
Geldenaaksebaan 464
B-3001 Leuven
+ 32 16 385011
<http://be.vwr.com>

Intended use

This item is intended for general laboratory use to monitor temperature in refrigerators and freezers.

Product Specifications

Range: -58 to 158°F and -50 to 70°C

Resolution: 1°

Accuracy: ±1°C

Dimensions: 70 x 108 x 19 mm.

Weight: 114 g.

Description of Buttons and Switches / Operational keys

Key	Functions	Location
Memory Clear	Clears the stored memory	Front
Mode	To view either the alarm set, or normal display	Front
Alarm on/off	Activates the alarm	Front

Operation

With probe sensor attached the unit simultaneously displays current probe temperature and probe minimum/maximum temperatures.

With the probe sensor removed the unit simultaneously displays current ambient temperature and ambient minimum/maximum temperatures.

Resetting the Unit

Use a pointed object to push the RESET button on the back of the unit. It is necessary to reset the unit whenever a change is made to C/F temperature, FAST/NORMAL sampling rate, or the probe sensor is plugged in or removed. HI/LO temperature alarm settings and MIN/MAX memories will be cleared after resetting the unit.

NOTE: Reset the unit whenever the probe sensor is removed or inserted.

Display Modes

The unit has two display modes: NORMAL DISPLAY MODE and ALARM DISPLAY MODE. To change from one mode to another press the MODE button once.

NORMAL DISPLAY MODE - is indicated by the display of current temperature and the MIN (minimum) and MAX (maximum) temperatures. These temperatures relate to external (probe) temperature if the probe is plugged in and to internal (ambient) temperature if the sensor probe is removed.

ALARM DISPLAY MODE - is indicated by the display of current temperature and the HI and LO alarm settings. These temperatures relate to external (probe) temperature if the probe is plugged in and to internal (ambient) temperature if the sensor probe is removed.

Resetting Min/Max Memories

There are two ways to reset the memories:

1. To reset both minimum and maximum memories - press MEMORY CLEAR button.
2. To reset individual minimum and maximum memories - press either the MIN or MAX button (on back of unit) once to reset the respective memory. The other memory will remain unaffected.

NOTE: You must be in Normal Display Mode to reset the memories.

To display the temperature readings in Fahrenheit (F) or Celsius (C) slide the switch on the back of the unit to F or C.

NOTE: The unit must be reset after changing the temperature unit.

Selecting the Sampling Rate

To select normal sampling rate (60 seconds) or fast sampling rate (10 seconds) slide the switch on back of unit to NORMAL or FAST.

NOTE: The unit must be reset after changing the sampling rate. Using the NORMAL sampling rate will extend battery life.

Setting the High/Lo Alarm

1. Press the MODE button until unit is in the ALARM DISPLAY MODE indicated by the display of current temperature and HI and LO alarm settings.
2. To set the high temperature alarm limit press the HI button, to advance the display until the desired high temperature is displayed. Press and hold the HI button to rapidly advance the display.
3. To set the low temperature alarm limit press the LO button, to advance the display until the desired low temperature is displayed. Press and hold the LO button to rapidly advance the display.

Alarm

The unit will sound the alarm in both NORMAL DISPLAY MODE and ALARM DISPLAY MODE.

Slide the ALARM ON/OFF switch to ON. When the current temperature display rises above the HI set point, or falls below the LO set point, the alarm will sound for 1 minute and issue a 3 second repeater beep every minute thereafter, for up to 12 hours. The alarm will continue to sound even if the current temperature display returns to an "in range" condition.

To temporarily silence the alarm press either the HI or LO button the back of the unit. The unit is still active

and the alarm will sound again if the temperature goes out of range. To permanently silence the alarm slide the ALARM ON/OFF switch to OFF position.

Troubleshooting

If this product does not function properly for any reason, please replace the battery with a new high quality battery (see "Battery Replacement" section). Low battery power can occasionally cause any number of "apparent" operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

Technical service

Web Resources

Visit the VWR's website at www.vwr.com for:

1. Complete technical service contact information
2. Access to VWR's Online Catalogue, and information about accessories and related products
3. Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit.

www.vwr.com.

Warranty VWR International warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of purchase. If a defect is present, VWR will, at its option, repair, replace, or refund the purchase price of this product at no charge to you, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear.

For your protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Equipment disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste. Instead it's your responsibility to correctly dispose of your equipment at lifecycle - end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment. For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment. By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

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Local VWR offices in Europe and Asia Pacific

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India

VWR Lab Products Private Limited
135/12, Brigade Towers, 2nd Floor
Front wing, Brigade Road,
Bengaluru, India – 560 025
Tel: +91-2522-647911/922 (Mumbai)
Tel: +91-80-41117125/26 (Bangalore)
Fax +91-80-41117120
E-mail: vwr_india@vwr.com

Ireland / Northern Ireland

VWR International Ltd / VWR International
(Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: 01 88 22 222
Fax: 01 88 22 333
E-mail sales@ie.vwr.com

Italy

VWR International PBI S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: 02-3320311/02-487791
Fax: 02-332031307/02-40090010
E-mail: info@it.vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: 020 4808 400
Fax: 020 4808 480
E-mail: info@nl.vwr.com

New Zealand

VWR International LP
241 Bush Road
Albany
0632 New Zealand

Norway

VWR International AS
Haavard Martinsens vei 30
0978 Oslo
Tel.: 02290
Fax: 815 00 940
E-mail: info@no.vwr.com

Poland

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: 058 32 38 210
Fax. 058 32 38 205
E-mail: labart@pl.vwr.com

Portugal

VWR International - Material de
Laboratório, Lda
Edifício Neopark
Av. Tomás Ribeiro, 43- 3 D
2790-221 Carnaxide
Tel.: 21 3600 770
Fax: 21 3600 798/9
E-mail: info@pt.vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel: +65 6505 0760
Fax: +65 6264 3780
E-mail: sales@sg.vwr.com

Spain

VWR International Eurolab S.L.
C/ Tecnología 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: 902 222 897
Fax: 902 430 657
E-mail: info@es.vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: 08 621 34 00
Fax: 08 621 34 66
E-mail: info@se.vwr.com

Switzerland

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: 044 745 13 13
Fax: 044 745 13 10
E-mail: info@ch.vwr.com

Turkey

Pro Lab Laboratuvar Teknolojileri Ltd.Şti.
a VWR International Company
Orta Mah. Cemal Gürsel Caddesi
Ördekcioglu İşmerkezi No.32/1
34896 Pendik - Istanbul
Tel.: +90216 598 2900
Fax: +90216 598 2907
Email: info@pro-lab.com.tr

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard
Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: 0800 22 33 44
Fax: 01455 55 85 86
E-mail: uksales@uk.vwr.com