

LavaLAMP DNA and RNA Isothermal Amplification Kits

Rapid and sensitive loop-mediated amplification assays available in convenient master mix and flexible component kit formats

LavaLAMP™ DNA Master Mix

- The LavaLAMP DNA Master Mix greatly simplifies reaction optimisation by limiting optimisation to target specific components and conditions, such as LAMP primer design, target concentration and reaction temperature. This master mix is heat stable at 90 °C for ≤5 minutes, which enables the addition of a reaction preheating step which may, depending on the target, increase assay sensitivity and decrease time to results (TTR).

LavaLAMP DNA Component Kit

- The LavaLAMP DNA Component Kit, designed for LAMP assay development, contains a the same unique, highly processive DNA polymerase enzyme as the LavaLAMP DNA Master Mix and individual component reagents for complete control and optimisation for your LAMP assays.
- This kit is an exact deconstruction of the LavaLAMP DNA Master Mix. When set up with the initial conditions described in its User Manual, the reactions will match the Master Mix-based reaction composition.

• Loop-mediated isothermal amplification (LAMP):

Allows researchers to run assays outside of high-tech laboratories using low complexity instrumentation.

- **Single enzyme:** All LavaLAMP assays are built around a proprietary DNA polymerase with strand displacement activity that works on both RNA and DNA templates. Master mix formulations have been specifically optimised for each template type.

- **Lyophilisation-ready reagents:** Specially formulated for room temperature stability in the absence of glycerol and betaine. Liquid formulations can be seamlessly freeze-dried for long term storage at ambient temperatures. Ideal for point of care (POC) applications.

Faster detection with LavaLAMP DNA Master Mix

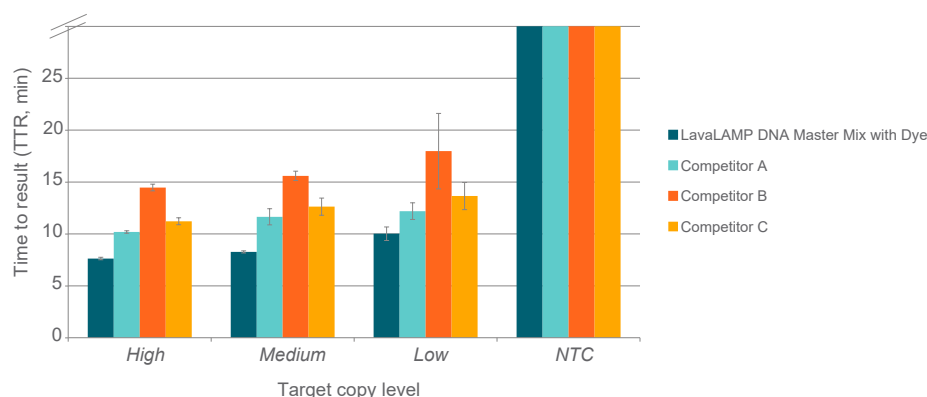


Figure 1. Loop-mediated isothermal amplification (LAMP) with real-time fluorescent detection of amplified products. LAMP reactions were set up using the indicated kits according to manufacturer's recommendations. Target DNA (*C. difficile*) at varying input amounts, *tcdA* target LAMP primers, and Green Fluorescent Dye (LavaLAMP Kit) were included in all reactions. Reactions were run on a CFX96 Thermal Cycler (Bio-Rad) at the following temperatures: LavaLAMP; 68 °C; other kits at the recommended 65 °C and fluorescence was measured in real-time to determine the TTR. NTC denotes No Target Control.

LavaLAMP RNA Master Mix

- The LavaLAMP RNA Master Mix is built around a proprietary DNA polymerase with strand displacement activity that works on RNA and DNA templates, making LAMP possible for RNA targets/templates. Just one enzyme is needed to convert the RNA target into cDNA to then act as template for additional LAMP. By combining this enzyme with an optimised master mix formulation, we've made an easy-to-use system that enables fast, sensitive detection of specific RNA targets. A "pre-optimised reaction formulation" lets you shift focus to optimising the most important parameters: RNA LAMP primer design and reaction temperature.


LavaLAMP RNA Component Kit



- The LavaLAMP RNA Component Kit employs the same DNA polymerase used in the LavaLAMP RNA Master Mix, compatible with both RNA and DNA templates, making RNA LAMP possible using a single DNA polymerase to convert RNA into cDNA before amplification into a detectable, target-specific dsDNA. Because this kit is a deconstruction of the LavaLAMP RNA Master Mix, reactions set up using initial conditions outlined in its User Manual will match the reaction compositions of the Master Mix reaction. The individual reagent format enables complete control to optimise reaction conditions for your target and primer set.

Performance comparison: LavaLAMP RNA Component Kit vs. competitor kits

Figure 2. (A) RNA LAMP reactions were set up according to kit manufacturer's recommendations. Target MS2 RNA at varying input amounts, MS2 target RNA LAMP primers, and Green Fluorescent Dye (LavaLAMP Kit) were included in all reactions. Reactions were run on a CFX96 Thermal Cycler (Bio-Rad) at 68 °C for LavaLAMP kits and at 65 °C for other polymerases and fluorescence was measured over 60 minutes to determine the TTR.

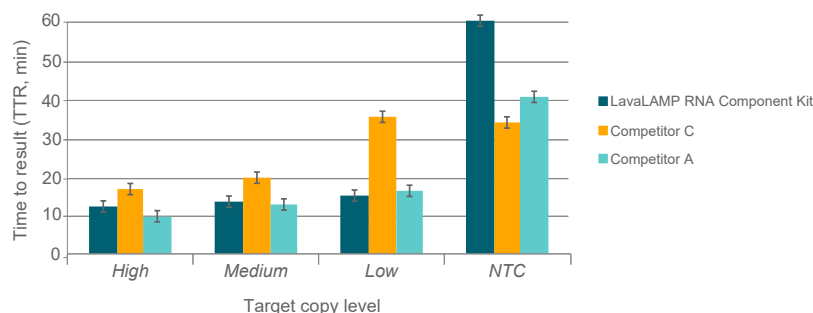

Ordering information

Cat no.	Size	Description
75996-746*	200 rxn	LavaLAMP DNA Master Mix
75996-748*	200 rxn	LavaLAMP DNA Master Mix with Dye
76203-026	200 rxn	LavaLAMP RNA Master Mix
76203-028	200 rxn	LavaLAMP RNA Master Mix with Dye
76180-468	500 rxn	LavaLAMP DNA Component Kit
76180-470	500 rxn	LavaLAMP DNA Component Kit with Dye
76210-506	500 rxn	LavaLAMP RNA Component Kit
76210-508	500 rxn	LavaLAMP RNA Component Kit with Dye
76203-024	200 rxn	Green Fluorescent Dye
76180-472	500 rxn	Green Fluorescent Dye

* These products are coming soon to Canada. Please contact your VWR sales rep to learn about availability.

Master Mix kits contain: LavaLAMP DNA or RNA Master Mix, DNA or RNA Positive Control LAMP Primer Mix, and DNA or RNA Positive Control as appropriate for each DNA or RNA kit. The LavaLAMP DNA or RNA Master Mix with Dye also contains Green Fluorescent Dye for fluorescent detection of amplified DNA.

Component kits contain: 10X LavaLAMP DNA or RNA Buffer, LavaLAMP DNA or RNA Enzyme, Magnesium Sulfate, 100 mM, DNA or RNA Positive Control LAMP Primer Mix and DNA or RNA Positive Control as appropriate for each DNA or RNA kit. The LavaLAMP DNA or RNA Component Kit with Dye also contains Green Fluorescent Dye for fluorescent detection of amplified DNA.

Note: Standard LavaLAMP reactions are 25 µL.

BIOSEARCH™
TECHNOLOGIES
GENOMIC ANALYSIS BY LGC

**Integrated tools.
Accelerated science.**



For Research Use Only. Not for use in diagnostic procedures.

All trademarks and registered trademarks mentioned herein are the property of their respective owners. All other trademarks and registered trademarks are the property of LGC and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details. No part of this publication may be reproduced or transmitted in any form or by any means, electronic or mechanical, including photocopying, recording or any retrieval system, without the written permission of the copyright holder. © LGC Limited, 2020. All rights reserved. GEN/626/SW/1020



800 932 5000 | VWR.COM

Prices, product, and/or services details are current when published and subject to change without notice. | Certain products or services may be limited by federal, state, provincial, or local regulations. | VWR, part of Avantor, makes no claims or warranties concerning sustainable/green products. Any claims concerning sustainable/green products are the sole claims of the manufacturer and not those of VWR International, LLC and/or Avantor, Inc. or affiliates. All prices are in US dollars unless otherwise noted. Offers valid in US and Canada, void where prohibited by law or company policy, while supplies last. | Trademarks are owned by Avantor, Inc. or its affiliates, unless otherwise noted. | Visit vwr.com to view our privacy policy, trademark owners, and additional disclaimers. © 2020 Avantor, Inc. All rights reserved.