

Preserve Sample Volume And Avoid **FACS***/ **FLOW** Clogs With...

FLOWMI™

Cell Strainers
40µm and 70µm Porosity



actual
size

Ideal for Small Volume Straining

vwr.com/flowmi

These Unique Strainers are Perfect For Fast and Efficient Filtering of Small Volume Samples (up to 1000 μ l) Prior to FLOW or FACS Analysis.



Preserve your sample volume; FLOWMI strainers maintain sample volume and desired cells when straining small sample sizes



Avoid wasted time and costly delays; fast and efficient filtering using FLOWMI strainers decreases the potential for clogging of FLOW or FACS instruments



Improves your work flow; FLOWMI strainers are oriented in a compact tray for direct press and fit attachment to a P1000 pipette tip



Quality assurance; Sterile, FLOWMI strainers are housed in a compact, covered tray with a re-sealable bag to maintain sterility. Tray features a sliding cover for easy, one-handed use



Use your favorite tips; FLOWMI strainers are compatible with a wide variety of P1000 tips including VWR brand, Axygen, Nichiryo and Eppendorf

Because a Sample is a Terrible Thing to Waste Smart. Effective. Innovative.



Aspirate
sample
into
1000µl
tip



Press
FLOWMI
onto
pipette tip




Strain
sample
into
tube



Eject to
dispose
of tip and
FLOWMI


FLOWMI™ Delivers Results You Can Count On

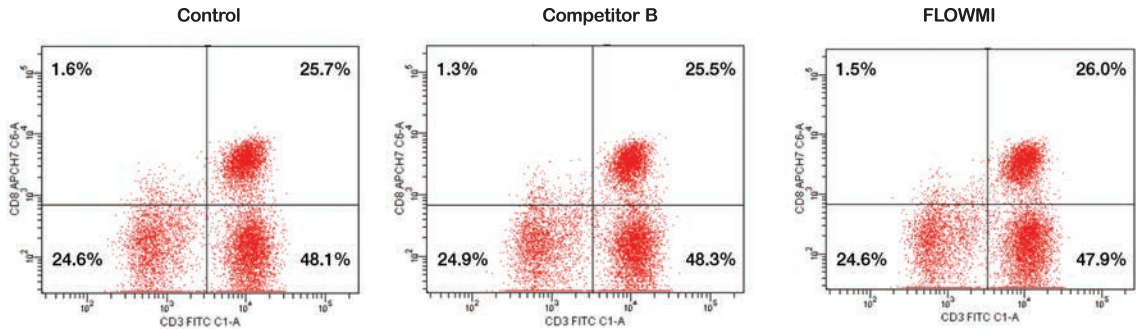

Cytograms of CD-Chex™ Plus cells from left to right – Control (non-filtered); Sample filtered using Competitor; and Sample filtered using FLOWMI. Cytograms and “Percent of Parent” data demonstrate no loss of specific cell types using FLOWMI vs. competing filter.

Flowmi is recommended for use with samples having a maximum concentration of 2 million cells/ml (40µm) or 4 million cells/ml (70µm).

FLOWMI™ Ranks Best in Sample Protection

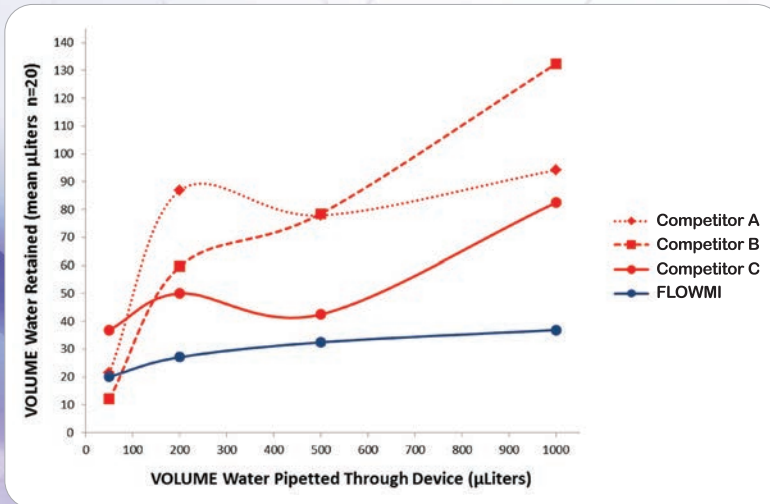

FLOWMI™ and three competing strainers were examined at four pipetted volumes (50, 200, 500, and 1000 µl). The Volume of Water Retained (µl) by each strainer, at each pipetted volume, is represented by the mean difference in weight before and after pipetting water through the strainer. (n=20 per pipetted volume; each trial using a new strainer).

Results demonstrate minimal loss of sample volume due to fluid retention inside FLOWMI™ Cell Strainer.

To see FLOWMI™ in action
and for additional technical data
please visit vwr.com/flowmi



Now Available!
70µm Porosity



VWR Cat. No.	Porosity
10032-802	40 micron
10204-924	70 micron

Patent Pending **50 Strainers/pack**

Flowmi is recommended for use with samples having a maximum concentration of 2 million cells/ml (40µm) or 4 million cells/ml (70µm).

avantor™
delivered by **VWR™**

VWR.COM

Prices, product, and/or services details are current when published and subject to change without notice. | Certain products or services may be limited by federal, state, provincial, or local regulations. | VWR, part of Avantor, makes no claims or warranties concerning sustainable/green products. Any claims concerning sustainable/green products are the sole claims of the manufacturer and not those of VWR International, LLC and/or Avantor, Inc. or affiliates. All prices are in US dollars unless otherwise noted. Offers valid in US and Canada unless otherwise noted, void where prohibited by law or company policy, while supplies last. | Trademarks are owned by Avantor, Inc. or its affiliates, unless otherwise noted. | Visit vwr.com to view our privacy policy, trademark owners, and additional disclaimers. © 2022 Avantor, Inc. All rights reserved.

0522 Lit. No. 190173WREV