

Safety: When melting wax and sectioning tissue, students should be supervised by a knowledgeable, trained instructor. Students should wear protective gloves of heavyweight cotton Kelnit, safety goggles, and aprons or smocks when handling melted wax.

- 1. Melt the paraffin granules in a waterbath of approximately 52°C (126°F). Caution: Do not overheat! Overheated wax will produce smoke and will catch fire.
- 2. Prepare a section of material to be cut, approximately 1" long and $\frac{1}{4}$ " diameter.
- 3. Place the section in the microtome so that a small portion protrudes from the top.
- 4. Carefully pour the melted wax onto the specimen, filling the space around the section with melted wax (see Figure 1). Caution: Avoid contact with skin; melted wax can cause second- and third-degree burns.
- 5. Let wax cool; do not move specimen while wax solidifies.
- 6. Using the enclosed safety razor, shave the protruding section until it is flush with the top of the microtome (see Figure 2). Safety: Always move the razor blade away from you.
- 7. Twist the threaded bolt slightly so the section rises to just above the top of the microtome.
- 8. Shave as thin a specimen as possible from the section.

 Note: To obtain a suitable specimen, you may need to practice this technique a number of times.
- 9. Using a pipet, apply a drop of water to the center of a clean microscope slide. Using fine-tip forceps, place the thin section on top of the water; the water will help flatten the thin section.




WARD'S Natural Science 5100 West Henrietta Road P.O. Box 92912 Rochester, New York 14692-9012

Fax: **1-800-635-8439**Phone: **1-800-962-2660**Website: **www.wardsci.com**

E-mail: customer_service@wardsci.com

©1997 WARD'S Natural Science Establishment, Inc. All Rights Reserved.

250-1407

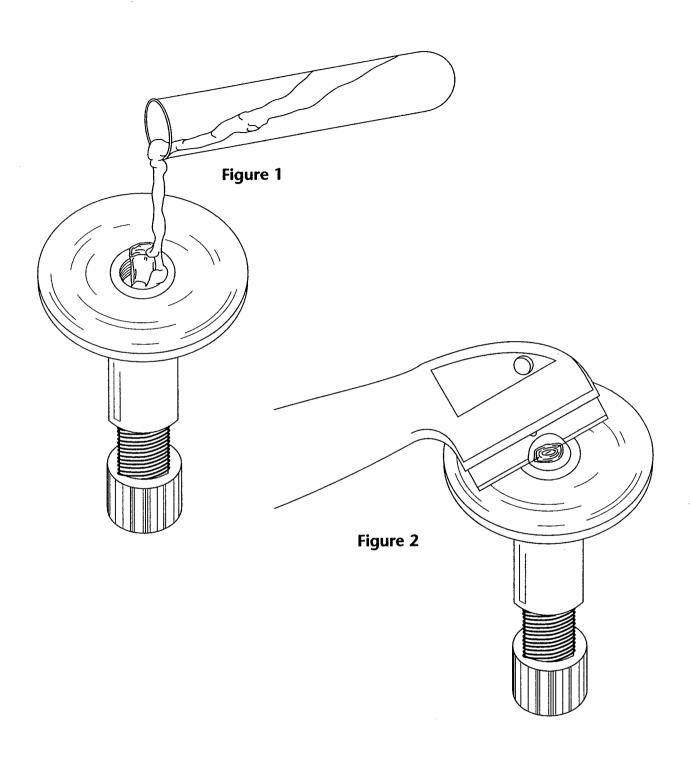

WARD'S

WARD'S Natural Science 812 Fiero Lane P.O. Box 5010 San Luis Obispo, California 93403-5010

Fax: 805-781-2704 Phone: 805-781-2700 Toll-Free: 1-800-872-7289 WARD'S Natural Science Ltd. 397 Vansickle Road St. Catharines, Ontario L2S 3T5

> Fax: 905-984-5952 Phone: 905-984-8900 Toll-Free: 1-800-387-7822