

HiTrap rProtein A FF, HiTrap Protein A HP and HiTrap Protein G HP are part of the range of prepacked, ready to use columns for preparative affinity chromatography. Fast, simple and easy separations are provided by the combination of easy to use HiTrap column and the various affinity media.

HiTrap rProtein A FF, HiTrap Protein A HP and HiTrap Protein G HP 1 ml and 5 ml columns allow convenient purification of polyclonal and monoclonal antibodies from cell culture supernatants, serum and ascites.

- Rapid and convenient preparative purification of polyclonal and monoclonal antibodies
- Very high purity in one step
- High binding capacities
- Simple and proven method giving reproducible results
- Simple operation with a syringe, a pump, an ÄKTA™ system, or other chromatography systems.

The basis for antibody purification is the high affinity and specificity of protein A and protein G for the Fc-region of IgG from a variety of species. Protein A and protein G have been immobilized to several matrices resulting in excellent purification of IgG and IgG subclasses from ascites fluid, cell culture supernatants, and serum.

The degree to which protein A and protein G bind to IgG varies with respect to both origin and antibody subclass and may even vary within a single subclass. The binding capacity of protein A and protein G for IgG depends on the source species of the particular immunoglobulin. The capacity



Fig 1. HiTrap rProtein A FF, HiTrap Protein A HP and HiTrap Protein G HP are designed for puification of monoclonal and polyclonal IgG.

depends also upon several other factors such as flow rate during sample application, and sample concentration.

The specificity of the recombinant protein A for the Fc-region of IgG is similar to native protein A and provides excellent purification in one step, see Table 2 for more details.

Column characteristics

HiTrap column is made of polypropylene, a material which is biocompatible and does not interact with biomolecules. The column is delivered with a stopper on the inlet and a snapoff end on the outlet. Both ends have 1/16" fittings for easy connection to ÄKTA systems.

Media characteristics

rProtein A Sepharose Fast Flow

rProtein A and protein A share similar specificity for the Fc-region of IaG, but recombinant protein A (rProtein A) offers several potential advantages. Since rProtein A has been engineered to include a C-terminal cysteine, controlled epoxy chemistry is used to favor single point oriented immobilization via thioether coupling and results in enhanced binding capacity for IgG. Furthermore, rProtein A is produced in E. coli and no human IgG affinity step is used during validated fermentation and purification processes, minimizing risk om human IgG contamination.

Recombinant (E. coli) protein A ligand is immobilized to Sepharose[™] Fast Flow, a robust highly cross-linked agarose with spherical 90 um beads.

Protein A Sepharose High Performance and Protein G Sepharose High Performance

Sepharose High Performance is the base matrix for HiTrap Protein A HP and HiTrap Protein G HP. The carbohydrate nature of the agarose base provides a hydrophilic and chemically favourable environment for coupling, while the highly cross-linked structure of the 34 µm spherical beads ensures excellent chromatographic properties. The protein A and protein G ligands are coupled to Sepharose High Performance by the N-hydroxysuccinimide activation method.

Protein A is a 42 000 molecular weight protein derived from a strain of Staphylococcus aureus. It consists of six regions, five of which bind IgG. As an affinity ligand, protein A is immobilized to the matrix so that these regions are free to bind. One molecule of immobilized protein A can bind at least two molecules of IgG.

Protein G, a cell surface protein from Group G Streptococci, is a type III Fc receptor and binds IgG with a non-immune mechanism similar to that of protein A. Here a recombinant form of the protein produced in E. coli, from which the albumin-binding region of the native protein has been genetically deleted, is used. Recombinant protein G contains two Fc-binding regions.

Fast kinetics with high dynamic capacities are properties of all HiTrap affinity columns. The binding capacity of rProtein A, protein A and protein G for IgG depends on the source species of the particular immunoglobulin. The total capacity depends also upon several other factors, such as flow rate during sample application and sample concentration. As a reference, the binding capacity for human IgG is approximately 20 mg IgG/ml medium for HiTrap Protein A HP, approximately 25 mg IgG/ml medium for HiTrap Protein G HP and approximately 50 mg/ml medium for HiTrap rProtein A FF.

Table 1 lists the main characteristics of HiTrap rProtein A FF, HiTrap Protein A HP and HiTrap Protein G HP.

Operation

All HiTrap columns are quick and easy to use. Instructions and connectors are included with each pack of columns. In general, the separation can be easily achieved with a syringe. using the luer adapter provided. Figure 2 illustrates this technique. Alternatively, the column can be operated using a laboratory pump or a chromatography system, for example when linear gradients are required or when large sample volumes are loaded. Two or more columns can be connected in series by screwing the end of one into the top of the next (back pressure will increase). The columns can not be opened or repacked.

Table 1. Main characteristics of HiTrap rProtein A FF, HiTrap Protein A HP and HiTrap Protein G HP

| Column volume | 1 ml and 5 ml |
|--|---|
| Column dimensions | 0.7 × 2.5 cm (1 ml) 1.6 × 2.5 cm (5 ml) |
| Ligand | Recombinant protein A (E. coli), protein A or protein G |
| Ligand concentration (Approx.) | 6 mg rProtein A/ml medium (HiTrap rProtein A FF) 3 mg protein A/ml medium (HiTrap Protein A HP) 2 mg protein G/ml medium (HiTrap Protein G HP) |
| Binding capacity (Approx.) | 50 mg human IgG/ml medium (HiTrap rProtein A FF) 20 mg human IgG (HiTrap Protein A HP) 25 mg human IgG (HiTrap Protein G HP) |
| Dynamic binding capacities* (HiTrap rProtein A FF) | 23 mg mouse monoclonal IgG _{2a} /ml medium 12 mg mouse monoclonal IgG ₁ /ml medium 11 mg monoclonal humanized IgG ₄ /ml medium |
| Average particle size | 90 µm (HiTrap rProtein A FF) 34 µm (HiTrap Protein A HP and HiTrap Protein G HP) |
| Bead structure | Highly cross-linked spherical agarose |
| Recommended flow rate | 1 and 5 ml/min for 1 and 5 ml columns respectively |
| Maximum flow rate [†] | 4 and 20 ml/min for 1 and 5 ml columns respectively |
| Column hardware pressure limit | 5 bar (0.5 MPa, 70 psi) |
| pH stability [‡] Working | 3 to 10 (HiTrap rProtein A FF) 3 to 9 (HiTrap Protein A HP and HiTrap Protein G HP) |
| Cleaning | 2 to 11 (HiTrap rProtein A FF) 2 to 10 (HiTrap Protein A HP) 2 to 9 (HiTrap Protein G HP) |
| Storage | +4°C to +8°C in 20% ethanol |

Running conditions for determining the dynamic binding capacity of HiTrap rProtein A FF:

Binding buffer: 20 mM sodium phosphate (+3 M NaCl for IgG1), pH 7.0

Elution buffer: 0.1 M sodium citrate, pH 3.0 Column: HiTrap rProtein A FF 1 ml

Flow rate: 1 ml/min

Monoclonal cell culture supernatants

Room temperature, aqueous buffers

The ranges given are estimates based on our knowledge and experience. Please note the following: pH stability, working refers to the pH interval where the medium is stable over a long period of time without adverse effects on its subsequent chromatographic performance. pH stability, cleaning refers to the pH interval for regeneration.







Fig 2. Using HiTrap rProtein A FF, HiTrap Protein A HP or HiTrap Protein G HP with a syringe.

(A) Prepare buffers and sample. Re-move the column's top cap and snap off the end. Wash and equilibrate. (B) Load the sample and begin collecting fractions.

(C) Elute and continue collecting fractions.

Applications

Protein A and protein G have different IgG binding specificities, depending on the origin of the IgG. Compared with protein A, protein G binds more strongly to polyclonal IgG, for example, from cow, sheep and horse. Furthermore, unlike protein A, protein G binds rat IgG, human IgG $_3$ and mouse IgG $_1$. Table 2 lists the relative binding strengths of polyclonal IgG from various species to protein G and protein A. Binding was measured in a competitive ELISA test. The amount of IgG required to give a 50% inhibition of binding of rabbit IgG conjugated with alkaline phosphatase was determined.

For more information, please refer to Antibody Purification Handbook, see ordering information.

Scale-up

The easiest way to scale-up is to go from a 1 ml HiTrap column to a 5 ml column. Alternatively, scale-up of small scale purifications can be done by coupling the columns in series (back pressure will increase).

Further scale-up can be done with bulk packages using nProtein A Sepharose Fast Flow, rProtein A Sepharose Fast Flow or Protein G Sepharose Fast Flow.

Storage

Recommended storage conditions for HiTrap rProtein A FF, HiTrap Protein A HP and HiTrap Protein G HP is in 20% ethanol at $+4^{\circ}$ C to $+8^{\circ}$ C.

Table 2. Relative binding strengths of protein A and protein G

| Species | Subclass | Protein A binding | Protein G binding |
|-----------------|---|-----------------------------------|-------------------------------------|
| Human | IgA IgD IgE IgG1 IgG2 IgG3 IgG4 IgM* | varible ++++ ++++ - ++++ variable | - - - ++++ ++++ ++++ |
| Avian egg yolk | IgY [†] | _ | - |
| Cow | | ++ | ++++ |
| Dog | | ++ | + |
| Goat | | _ | ++ |
| Guinea pig | IgG_1 | ++++ | ++ |
| | IgG ₂ | ++++ | ++ |
| Hamster | | + | ++ |
| Horse | | ++ | ++++ |
| Koala | | _ | + |
| Llama | | - | + |
| Monkey (rhesus) | | ++++ | ++++ |
| Mouse | IgG_1 | + | ++++ |
| | IgG_{2a} | ++++ | ++++ |
| | IgG_{2b} | +++ | +++ |
| | IgG_3 | ++ | +++ |
| | IgM* | variable | - |
| Pig | | +++ | +++ |
| Rabbit | no distinction | ++++ | +++ |
| Rat | IgG ₁ | - | + |
| | IgG_{2a} | - | ++++ |
| | IgG_{2b} | _ | ++ |
| | IgG ₃ | + | ++ |
| Sheep | | +/- | ++ |

^{*} Purified using HiTrap IgM Purification HP columns.

[†] Purified using HiTrap IgY Purification HP columns.

^{++++ =} strong binding

^{++ =} medium binding

^{+/-} = weak or no binding

HiTrap rProtein A FF

Purification of monoclonal mouse IgG_{2b} from ascites


Mouse IgG_{2b} was purified on HiTrap rProtein A FF 1 ml column operated with a syringe. The eluted pool contained 1 mg IgG_{2b} . The silver-stained SDS-PAGE confirmed that the eluted antibody was over 95 % pure, (Fig 3).


Sample: 1 ml mouse ascites containing IgG_{2b} , filtered through a 0.45 μm filter.

The sample was a kind gift from Dr. N. Linde, EC Diagnostics, Sweden.

Column: HiTrap rProtein A FF 1 ml
Binding buffer: 0.02 M sodium phosphate, pH 7.0
Elution buffer: 0.1 M sodium citrate, pH 3.0

Flow rate: approx. 1 ml/min
Instrumentation: Syringe

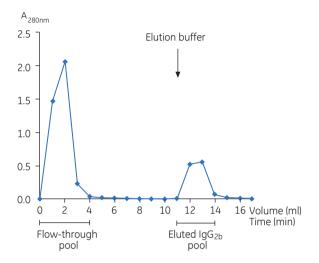

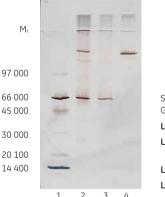

SDS-PAGE on PhastSystem™ using PhastGel™ Gradient 10–15, silver-staining

Lane 1: LMW markers

Lane 2: Starting material, mouse ascites,

diluted 1:10

Lane 3: Flow-through pool
Lane 4: Eluted IgG_{2b}

Figure 3. Purification of mouse IgG_{2b} from ascites on HiTrap rProtein A FF 1 ml column using a syringe.

Purification of monoclonal mouse IgG₁ from cell culture supernatant

Mouse IgG_1 was purified from 150 ml cell culture supernatant on HiTrap rProtein A FF 5 ml column.

The eluted pool contained 28 mg lgG₁.

The eluted IgG_1 was over 95 % pure according to SDS-PAGE with silver-staining, (Fig 4).


Sample: 150 ml of cell culture supernatant containing IgG, filtered through a


0.45 µm filter

Column: HiTrap rProtein A FF 5 ml

Binding buffer: 0.02 M sodium phosphate, 3 M NaCl, pH 7.0

Elution buffer: 0.1 M sodium citrate, pH 3.0 Flow rate: 5 ml/min (150 cm/h)
Instrumentation: FPLC System

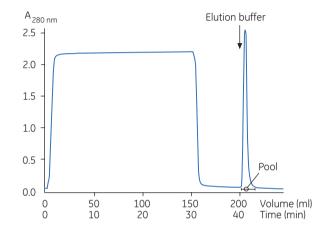

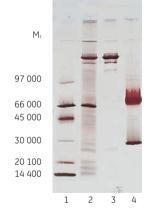

SDS-PAGE on PhastSystem using PhastGel Gradient 10–15, silver-staining

Lane 1: LMW markers

Lane 2: Starting material, cell culture

supernatant

Lane 3: Eluted IgG₁ pool, diluted 1:5,

non-reduced condition

Lane 4: Eluted IgG₁ pool, diluted 1:5, reduced condition

Fig 4. Purification of mouse IgG_1 from cell culture supernatant on HiTrap rProtein A FF 5 ml column.

HiTrap Protein A HP and HiTrap Protein G HP

Purification of monoclonal mouse IgG_{2b}

Mouse IgG_{2b} from hybridoma cell culture fluid was purified on HiTrap Protein A HP.

The purity was checked with SDS-PAGE, (Fig 5).

Sample: 10 ml mouse IgG_{2b} hybridoma cell culture fluid

Column: HiTrap Protein A HP 1 ml

Binding buffer: 0.02 M sodium phosphate, pH 7.0

Elution buffer: 0.1 M citric acid-NaOH, pH 3.0

Chromatographic

procedure: 2 ml binding buffer, 10 ml sample, 10 ml binding buffer

5 ml elution buffer, 5 ml binding buffer. The eluted fractions were


neutralized with 1 M Tris-HCl, pH 9.0


Electrophoresis: SDS-PAGE, PhastSystem, PhastGel Gradient 10-15, 1 µl sample,


silver stained

 $Immunodiffusion: 1\%\ Agarose\ A\ in\ 0.75\ M\ Tris,\ 0.25\ M\ 5,5-diethylbarbituric\ acid,$

5 mM Ca-lactate, 0.02% sodium azide, pH 8.6

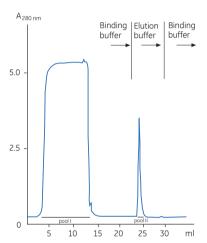

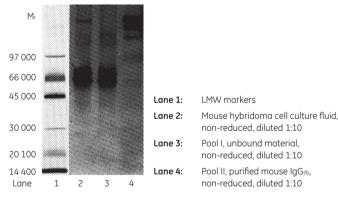

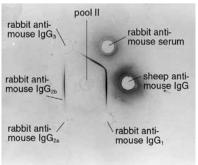

Immunodiffusion

Purification of mouse monoclonal IgG₁ from cell culture supernatant

Mouse monoclonal cell supernatant IgG_1 , anti-transferrin, was purified on HiTrap Protein G HP using syringe operation and pump operation.

The purity was checked with SDS-PAGE, (Fig 6).


Sample: 10 ml mouse monoclonal cell supernatant, IgG1, anti-transferrin

Column: HiTrap Protein G HP 1 ml
Binding buffer: 0.02 M sodium phosphate, pH 7.0
Elution buffer: 0.1 M glycine-HCl, pH 2.7


Electrophoresis: SDS-PAGE, PhastSystem, PhastGel Gradient 10–15, 1 µl sample,


silver stained

A) Syringe operation, approximately 60 drops/min

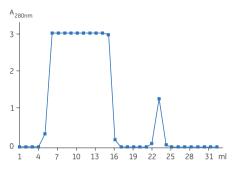

B) Pump operation, flow rate 2 ml/min

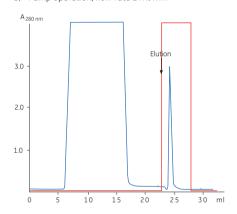

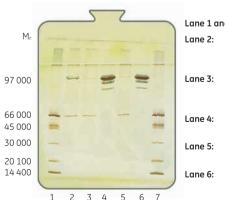

Lane 1 and 7: LMW markers

Crude cell culture supernatant, mouse IgG₁, diluted 1:10 Flow through, using peristaltic pump, diluted 1:10

Eluted mouse IgG1, using a peristaltic pump Flow through, using a

syringe, diluted 1:10 Eluted mouse IgG₁, using

a syringe

Fig 6. Purification of mouse monoclonal lgG_1 from cell culture supernatant (A) with syringe operation (B) with pump operation SDS-PAGE on PhastSystem using PhastGel 10-15, non-reduced condition, and silver staining.

Fig 5. Purification of monoclonal mouse IgG_{2b} on HiTrap Protein A HP.

Purification of monoclonal mouse IgG₁ from hybridoma cell culture

Mouse IgG₁ hybridoma cell culture fluid was purified on HiTrap Protein G HP. The purity was checked with SDS-PAGE, (Fig 7).

Sample: 12 ml mouse IgG₁ hybridoma cell culture fluid

HiTrap Protein G HP 1 ml Column:

1.0 ml/min Flow rate:

Binding buffer: 0.02 M sodium phosphate, pH 7.0 Elution buffer: 0.1 M glycine-HCl, pH 2.7

Chromatographic

5 ml binding buffer, 12 ml sample, 10 ml binding buffer Procedure:


6 ml elution buffer, 7 ml binding buffer. The eluted fractions were


neutralized with 1 M Tris-HCl, pH 9.0

Electrophoresis: SDS-PAGE, PhastSystem, PhastGel Gradient 10-15, 1 µl sample,

Immunodiffusion: 1% Agarose A in 0.75 M Tris, 0.25 M 5,5-diethylbarbituric acid,

5 mM Ca-lactate, 0.02% sodium azide, pH 8.6

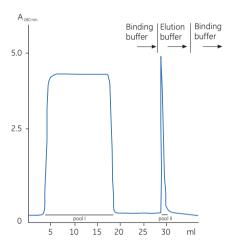

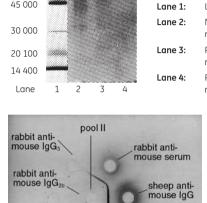

rabbit anti-

mouse IgG,

M,

97 000

66 000

45 000

rabbit anti-

mouse IgG_{2a}

LMW markers

Mouse hybridoma cell culture fluid,

non-reduced, diluted 1:10

Pool I, unbound material, non-reduced, diluted 1:10

Pool II, purified mouse IgG₁, non-reduced, diluted 1:10

Immunodiffusion

Fig 7. Purification of monoclonal mouse IgG1 on HiTrap Protein G HP, 1 ml.

Tables 3 and 4 list physio-chemical data for human and mouse immunoglobulins.

Table 3. Physio-chemical properties of human immunoglobulins

| Immunoglobulin | Heavy chain | Light chain | Sedimentation coefficient | M_r | M _r heavy chain | Carbohydrate content (%) | A _{280nm} | pl |
|------------------|-------------------------|----------------|---------------------------|---------|-------------------------------|--------------------------|--------------------|---------|
| IgG ₁ | λ_1 | κ, λ | 7S | 146 000 | 50 000 | 2-3 | 13.8 | 5.0-9.5 |
| IgG_2 | λ_1 | κ, λ | 7S | 146 000 | 50 000 | 2–3 | | 5.0-8.5 |
| IgG₃ | λ_1 | κ, λ | 7S | 170 000 | 60 000 | 2–3 | | 8.2-9.0 |
| IgG ₄ | λ_1 | κ, λ | 7S | 146 000 | 50 000 | 2-3 | | 5.0-6.0 |
| IgM | μ | κ, λ | 195 | 900 000 | 68 000 | 12 | 12.5 | 5.1-7.8 |
| IgA_1 | α_1 | κ, λ | 7S | 160 000 | 56 000 | 7-11 | 13.4 | 5.2-6.6 |
| IgA_2 | α_2 | κ, λ | 7S | 160 000 | 52 000 | 7–11 | | 5.2-6.6 |
| IgA _s | α_1 , α_2 | κ, λ | 11S | 370 000 | 52-56 000 | 11 | - | 4.7-6.2 |
| IgD | δ | κ, λ | 7S | 184 000 | 68 000 | 12 | 17.0 | - |
| IgE | 3 | κ, λ | 8S | 190 000 | 72 000 | 12 | 15.3 | _ |

Table 4. Physio-chemical properties of mouse immunoglobulins

| Immunoglobulin | Heavy chain | Light chain | Sedimentation coefficient | M_r | M _r heavy chain | Carbohydrate content (%) | pl |
|-------------------|---------------------|----------------|---------------------------|---------|-------------------------------|--------------------------|---------|
| IgG ₁ | λ_1 | κ, λ | 7S | 150 000 | 50 000 | 2–3 | 7.0-8.5 |
| IgG_{2a} | $\lambda_{2\alpha}$ | κ, λ | 7S | 150 000 | 50 000 | 2–3 | 6.5-7.5 |
| IgG _{2b} | λ_{2b} | κ, λ | 7S | 150 000 | 50 000 | 2–3 | 5.5-7.0 |
| IgG₃ | λ_3 | κ, λ | 7S | 150 000 | 50 000 | 2–3 | _ |
| IgM | μ | κ, λ | 195 | 900 000 | 80 000 | 12 | 4.5-7.0 |
| IgA | α | κ, λ | 7S | 170 000 | 70 000 | 7–11 | 4.0-7.0 |
| IgD | δ | κ, λ | 7S | 180 000 | 68 000 | 12-14 | _ |
| IgE | 3 | κ, λ | 8S | 190 000 | 80 000 | 12 | _ |

Ordering information

| Product | Quantity | Code number |
|----------------------|-------------------------|-------------|
| HiTrap Protein A HP | 5 × 1 ml | 17-0402-01 |
| HiTrap Protein A HP | $2 \times 1 \text{ ml}$ | 17-0402-03 |
| HiTrap Protein A HP | $1 \times 1 \text{ ml}$ | 29-0485-76 |
| HiTrap Protein A HP | $1 \times 5 \text{ ml}$ | 17-0403-01 |
| HiTrap Protein A HP | $5 \times 5 \text{ ml}$ | 17-0403-03 |
| HiTrap Protein G HP | $5 \times 1 \text{ ml}$ | 17-0404-01 |
| HiTrap Protein G HP | $2 \times 1 \text{ ml}$ | 17-0404-03 |
| HiTrap Protein G HP | $1 \times 1 \text{ ml}$ | 29-0485-81 |
| HiTrap Protein G HP | $1 \times 5 \text{ m}$ | 17-0405-01 |
| HiTrap Protein G HP | $5 \times 5 \text{ ml}$ | 17-0405-03 |
| HiTrap rProtein A FF | $5 \times 1 \text{ ml}$ | 17-5079-01 |
| HiTrap rProtein A FF | $2 \times 1 \text{ ml}$ | 17-5079-02 |
| HiTrap rProtein A FF | $1 \times 5 \text{ ml}$ | 17-5080-01 |
| HiTrap rProtein A FF | 5 × 5 ml | 17-5080-02 |

| Related products | Quantity | Code number |
|----------------------------------|-------------|-------------|
| HiTrap Desalting | 5 × 5 ml | 17-1408-01 |
| HiTrap Desalting | 1× 5 ml | 29-0486-84 |
| HiTrap Desalting | 100 × 5 ml* | 11-0003-29 |
| HiPrep™ 26/10 Desalting | 1 × 53 ml | 17-5087-01 |
| HiPrep 26/10 Desalting | 4 × 53 ml | 17-5087-02 |
| MAbTrap™ Kit | 1 kit | 17-1128-01 |
| nProtein A Sepharose 4 Fast Flow | 5 ml | 17-5280-01 |
| nProtein A Sepharose 4 Fast Flow | 25 ml | 17-5280-04 |
| rProtein A Sepharose 4 Fast Flow | 5 ml | 17-1279-01 |
| rProtein A Sepharose 4 Fast Flow | 25 ml | 17-1279-02 |
| Protein G Sepharose 4 Fast Flow | 5 ml | 17-0618-01 |
| Protein G Sepharose 4 Fast Flow | 25 ml | 17-0618-02 |

^{*} Pack size available by special order.

| Related products | Quantity | Code number |
|------------------------|-------------------------|-------------|
| HiTrap MabSelect SuRe™ | 5 × 1 ml | 11-0034-93 |
| HiTrap MabSelect SuRe | $1 \times 1 \text{ ml}$ | 29-0491-04 |
| HiTrap MabSelect SuRe | $1 \times 5 \text{ ml}$ | 11-0034-94 |
| HiTrap MabSelect SuRe | $5 \times 5 \text{ ml}$ | 11-0034-95 |
| MabSelect SuRe | 25 ml | 17-5438-01 |
| MabSelect Xtra™ | 25 ml | 17-5269-07 |
| MabSelect™ | 25 ml | 17-5199-01 |

| Accessories | Quantity | Code number |
|---|----------|-------------|
| 1/16" male/luer female* | 2 | 18-1112-51 |
| Tubing connector flangeless/M6 female | 2 | 18-1003-68 |
| Tubing connector flangeless/M6 male | 2 | 18-1017-98 |
| Union 1/16" female/M6 male | 6 | 18-1112-57 |
| Union M6 female /1/16" male | 5 | 18-3858-01 |
| Union luerlock female/M6 female | 2 | 18-1027-12 |
| HiTrap/HiPrep, 1/16" male connector for ÄKTA design | 8 | 28-4010-81 |
| Stop plug female, 1/16"† | 5 | 11-0004-64 |
| Fingertight stop plug, 1/16"‡ | 5 | 11-0003-55 |
| | | |

| Related literature | Code number |
|--|-------------|
| Antibody Purification Handbook | 18-1037-46 |
| Affinity Chromatography Handbook, Principle and Methods | 18-1022-29 |
| Affinity Chromatography Columns and Media Product Profile | 18-1121-86 |
| Convenient Protein Purification, HiTrap Column Guide | 18-1129-81 |

^{*} One connector included in each HiTrap package.

† Two, five, or seven stop plugs female included in HiTrap packages depending on products.

† One fingertight stop plug is connected to the top of each HiTrap column at delivery.

For local office contact information, visit www.gelifesciences.com/contact

www.gelifesciences.com/hitrap www.gelifesciences.com/protein-purification

GE Healthcare Bio-Sciences AB Björkgatan 30 751 84 Uppsala Sweden



 $\ensuremath{\mathsf{GE}}$ and $\ensuremath{\mathsf{GE}}$ monogram are trademarks of General Electric Company.

HiTrap, Sepharose, MabSelect SuRe, MabSelect, MabSelect Xtra, ÄKTA, PhastSystem, PhastGel, MAbTrap and HiPrep are trademarks of General Electric Company or one of its subsidiaries.

 $\ \odot$ 2005–2014 General Electric Company—All rights reserved. First published 2005

All goods and services are sold subject to the terms and conditions of sole of the company within GE Healthcare which supplies them. A copy of these terms and conditions is available on request. Contact your local GE Healthcare representative for the most current information.

GE Healthcare UK Limited Amersham Place, Little Chalfont Buckinghamshire, HP7 9NA

GE Healthcare Europe, GmbH Munzinger Strasse 5 D-79111 Freiburg Germany

GE Healthcare Bio-Sciences Corp. 800 Centennial Avenue, P.O. Box 1327 Piscataway, NJ 08855-1327 USA

GE Healthcare Japan Corporation Sanken Bldg., 3-25-1, Hyakunincho Shinjuku-ku, Tokyo 169-0073 Japan