

GE Healthcare

Amersham
Interleukin-10 [(h)IL-10]
human, Biotrak ELISA
system

96 wells

Product Booklet

Code: RPN2755



Page finder

1. Legal	3
2. Handling	4
2.1. Safety warnings and precautions	4
2.2. Storage	4
2.3. Expiry	4
3. Contents of the assay system	5
4. Description	6
5. Introduction	7
6. Summary of the assay	9
7. Assay methodology	10
7.1. Materials and equipment required	10
7.2. Sample preparation	10
7.3. Critical parameters	11
7.4. Assay procedure	12
7.5. Calculation of results	18
8. Additional information	20
8.1. Specificity	20
8.2. Calibration	20
8.3. Reproducibility	20
8.4. Sensitivity	20
8.5. Recovery	21
8.6. Expected values	21
9. References	22
10. Related products	23

1. Legal

GE, imagination at work and GE monogram are trademarks of General Electric Company.

Amersham and Biotrak are trademarks of GE Healthcare companies.

© 1995–2009 General Electric Company – All rights reserved.

First published 1995.

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them.

A copy of these terms and conditions is available on request.

Contact your local GE Healthcare representative for the most current information.

<http://www.gelifesciences.com>

GE Healthcare UK Limited.

Amersham Place,

Little Chalfont, Buckinghamshire,

HP7 9NA UK

2. Handling

2.1. Safety warnings and precautions

Warning: For research use

only. Not recommended or intended for diagnosis of disease in humans or animals. Do not use internally or externally in humans or animals.

All chemicals should be considered as potentially hazardous. We therefore recommend that this product is handled only by those persons who have been trained in laboratory techniques and that it is used in accordance with the principles of good laboratory practice. Wear suitable protective clothing such as laboratory overalls, safety glasses and gloves. Care should be taken to avoid contact with skin or eyes. In the case of contact with skin or eyes wash immediately with water. See material safety data sheet(s) and/or safety statement(s) for specific advice.

2.2. Storage

Store at 2–8°C

2.3. Expiry

The expiry date is stated on the package and will normally be at least 4 weeks from the date of despatch

3. Contents of the assay system

This pack contains the following assay components, sufficient material for 96 wells.

All reagents are stored refrigerated at 2–8°C. Refer to the expiry date on the kit box.

(h)IL-10 microplate - 96 well polystyrene microplate (12 strips of 8 wells) coated with an antibody against (h)IL-10.

Biotinylated antibody reagent - antibody against (h)IL-10 conjugated to biotin, with preservative, 8 ml.

(h)IL-10 standard - 2 vials of prediluted recombinant human IL-10 lyophilized.

Streptavidin-HRP concentrate - streptavidin conjugated to HRP, with preservative, 50 µl.

Streptavidin-HRP dilution buffer - with preservative, 14 ml.

Standard diluent - with preservative, 12 ml.

Wash buffer concentrate - 30-fold concentrated solution, with preservative, 50 ml.

Pre-mixed TMB substrate reagent - 12–13 ml.

Stop solution - <1% sulfuric acid, 13–15 ml.

Plate covers - 4 adhesive strips.

4. Description

The Biotrak™ human interleukin-10 ELISA system from GE Healthcare provides a simple, specific, reliable and precise quantitative determination of (h)IL-10 in cell culture supernatants, plasma, serum and urine.

The assay system is based on a solid phase ELISA, which utilizes an antibody for (h)IL-10 bound to the wells of a microplate together with a biotinylated antibody to (h)IL-10 and streptavidin conjugated to horseradish peroxidase. Although the Biotrak (h)IL-10 immunoassay contains recombinant (h)IL-10 and antibodies raised against recombinant (h)IL-10 it has been shown to quantitate accurately both natural (h)IL-10 and recombinant (h)IL-10.

(h)IL-10 can be measured in the approximate range 15–600 pg/ml (0.75–30 pg/well) in less than 4 hours using the protocol provided with the kit. Each pack contains sufficient material for 96 wells. If one standard curve is constructed, 42 unknowns can be measured in duplicate.

- High sensitivity - <3 pg/ml (0.15 pg/well)
- Same day protocol
- Pre-coated plate
- Specific for (h)IL-10

5. Introduction

IL-10 was recently identified as the cytokine produced by the T_H2 subpopulation of T cells that inhibit the synthesis of immunostimulatory cytokines by the T_H1 cells. Human IL-10 can be produced by B and T cells, activated mast cells, macrophages monocytes and keratinocytes and is expressed as a noncovalently linked homodimer. Human IL-10 is a 178 amino acid protein with a molecular weight of 18 kDa with two N-linked glycosylation sites.

IL-10 has been shown to exhibit profound inhibitory effects on monocytes including down-regulation of MHC class II antigen expression and suppression of IL-1 α and β , IL-6, IL-8, GM-CSF, G-CSF and TNF α production. IL-10 has been shown to synergize with IL-2 and IL-4 to promote the proliferation of thymocytes. IL-10 has also been shown to act in conjunction with IL-3 and IL-4 to enhance the survival of mast cells. As yet the IL-10 receptor has not been described.

Although IL-10 has been demonstrated to play an important role in several *in vitro* phenomena, the detection of IL-10 *in vivo* using bioassays or immunoassays has been difficult. Treatment of mice with anti-IL-10 antibodies resulted in the reversible depletion of Ly-1⁺ B cells, reduced serum IgM and IgA levels and attenuated antibody responses to hapten antigens. These mice also exhibited elevated levels of circulating IgG_{2a} and IgG_{2b} antibodies and TNF α . Mice lacking the IL-10 gene locus did not display any of the above effects, suggesting that immune complexes may have mediated the changes observed in anti-IL-10 treated mice. IL-10 production can be induced by mitogenic lectins and LPS. Agents that inhibit IL-10 production include IL-4 and IFN γ .

The suppressive effects of IL-10 on monocytes and T_H1 cytokine synthesis suggests that IL-10 may have utility as a general suppressor of immune function. The targets for such

immunosuppressive drugs include infectious disease, transplantation, induction of tolerance and possibly cancer. IL-10 is currently in preclinical studies to evaluate its potential in various disease states. In addition, IL-10 may be useful in promoting T_H2 -mediated (humoral) immunity where desired. It has been shown that as disease progresses, HIV infected patients switch from a T_H1 to a T_H2 state. Since IL-10 is the T_H2 cytokine that down regulates T_H1 responses, it could be responsible for the shift.

6. Summary of the assay

This assay employs the quantitative 'sandwich' enzyme immunoassay technique. An antibody specific for (h)IL-10 has been coated on the microplate provided in the kit. Samples are pipetted into the wells along with biotinylated antibody reagent, and the (h)IL-10, if present, is bound by both the immobilized antibody and biotinylated antibody. After washing away any unbound sample proteins and biotinylated antibody, a streptavidin-HRP conjugate is added to the wells. Any (h) IL-10 which was bound by both the immobilized and the biotinylated antibody during the first incubation will be bound by the streptavidin conjugate. Following a wash to remove any unbound conjugate, a substrate solution is added to the wells and color develops in proportion to the amount of (h)IL-10 bound in the initial step.

In addition to the samples to be tested, a series of wells is prepared using known concentrations of the Baculovirus-derived human IL-10 standard. A curve, plotting the optical density versus the concentration of the standard well, is prepared. By comparing the optical density of the samples to this standard curve, the concentration of the (h)IL-10 in the unknown samples is then determined.

7. Assay methodology

Users are recommended to read this entire section before starting work.

7.1. Materials and equipment required

The following materials and equipment are required:

Pipettes or pipetting equipment with disposable tips (50 µl, 100 µl and 1.00 ml)

Disposable polypropylene test tubes - do not use polystyrene, polycarbonate or glass

Measuring cylinder 2 L

Distilled or deionized water

A centrifuge for preparing streptavidin-HRP solution

Plate reader capable of reading at 450 nm

Optional equipment

Assays may be performed with commercially available microplate washers to aid convenience and assay throughput.

7.2. Sample preparation

Serum, plasma, urine and cell culture supernatants

Serum, plasma, urine and culture supernatant samples that are to be assayed within 24 hours should be stored at 2–8°C. Specimens to be stored for longer periods of time should be frozen at -70°C to avoid loss of biologically active cytokine. Avoid freezing and thawing samples more than once. Test samples should be assayed in duplicate each time the ELISA is performed, 50 µl of sample per well is required in this way.

Dilution of test samples

If you suspect that the (h)IL-10 concentration of a sample exceeds the highest point of the standard curve, prepare one or more five fold

dilutions of the test sample. Mix thoroughly between dilutions and before assaying

7.3. Critical parameters

1. First select the number of strips to be run and allow samples and all reagents to reach room temperature prior to performing the assays. Do not use water baths to thaw samples or reagents.
2. Mix samples and all reagents thoroughly before use.
3. Avoid excessive foaming of reagents. Also avoid exposure of reagents to excessive heat or light during storage and incubation.
4. Avoid handling the tops of the wells both before and after filling.
5. Standards and samples should be assayed in duplicate.
6. Run a separate standard curve for each assay.
7. The total dispensing time for each plate should not exceed 20 minutes.
8. Use only coated wells from the same reagent batch for each assay. Also do not mix reagents from different kit lots.
9. It is important that the wells are washed thoroughly and uniformly. If using an automatic washer check operation of heads before starting. If washing by hand ensure that all wells are completely filled at each wash.
10. A small amount of precipitate may be present in some vials. It will not affect assay performance and should be ignored.

7.4. Assay procedure

Reagent preparation

Wash buffer concentrate

Any precipitate formed during storage will redissolve upon dilution. Dilute 30-fold to prepare 1500 ml of wash buffer. Store at 2–8°C until the expiry date of the kit. Do not use wash buffer if it becomes visibly contaminated on storage.

Streptavidin-HRP solution

Prepare the exact amount of streptavidin-HRP solution no more than 15 minutes prior to use.

The streptavidin-HRP concentrate may require spinning down to force the contents to the bottom of the vial. Add 30 µl of streptavidin-HRP concentrate per 12 ml of streptavidin-HRP dilution buffer in a plastic 15 ml tube and mix gently.

If running partial plates use 2.5 µl of streptavidin-HRP concentrate and 1 ml of streptavidin-HRP dilution buffer per strip being run.

Preparation of working standards

- a. Two vials of lyophilized standards are provided with this kit.
Reconstitute and use one vial per partial plate.
- b. Prepare standards shortly before use. Use within one hour of reconstitution.
Do not store reconstituted standards.
- c. When running culture supernatant samples, reconstitute standard in distilled or deionized water. Reconstitution volume is stated on the standard vial label. The standard will take approximately 1 minute to dissolve. Mix by gently inverting the vial. Use your culture medium to prepare the dilutions of the standard curve, go to step e) below for further instructions. If running a partial plate, refer to step a) above.

- d. If running serum, plasma or urine samples, reconstitute standard with distilled or deionized water. Reconstitution volume is stated on the standard vial label. The standard will take approximately 1 minute to dissolve. Mix by gently inverting the vial. Use the standard diluent provided to prepare the dilutions of the standard curve. If running a partial plate, refer to step a) above.
- e. Label 6 tubes, one for each standard curve point: 600 pg/ml, 240 pg/ml, 96 pg/ml, 38.4 pg/ml, 15.36 pg/ml and 0 pg/ml. Then prepare 1:2.5 serial dilutions for the standard curve as follows:
- f. Pipette 240 µl of appropriate diluent (see steps c) and d) above) into each tube.
- g. Pipette 160 µl of the reconstituted standard into the first tube, 600 pg/ml and mix.
- h. Pipette 160 µl of this dilution into the second tube labelled 240 pg/ml and mix.
- i. Repeat serial dilutions three more times. These concentrations, 600 pg/ml, 240 pg/ml, 96 pg/ml, 38.4 pg/ml, 15.36 pg/ml and 0 pg/ml are your standard curve.

Running partial plates

This ELISA provides the flexibility to run two partial plates on separate occasions. Decide the number of strips you wish to run, leaving the strips to be used in the frame. Remove the unnecessary strips and store them in the foil pouch with the desiccant provided at 2–8°C, making sure the foil pouch is sealed tightly.

When adding the TMB substrate reagent, pour out from the bottle **only** the amount needed to run the first half plate. Do not combine left over substrate with that reserved for the second half of the plate. Care must be taken to ensure that the remaining TMB substrate reagent is not contaminated. **If the substrate reagent is bright blue prior to use, it has been contaminated. DO NOT USE.**

Assay protocol

1. Prepare assay reagents and working standards as described in the previous sections.
2. Set up the microplate with sufficient wells to enable the running of all standards and samples as required (see figure 1).
3. Remove excess microplate strips from the frame and store in the resealable foil bag.
4. Add 50 μ l of standard or sample per well.
5. Add 50 μ l of biotinylated antibody reagent (has a cloudy white appearance) per well. Cover with adhesive strip and incubate for 2 hours at room temperature (20–25°C).
6. Aspirate or decant each well and wash, repeating the process twice for a total of three washes. Wash vigorously by filling each well with wash buffer using a washbottle, pipette or manifold dispenser. Complete removal of liquid at each step is essential for good performance. After the last wash, remove any remaining wash buffer by inverting the plate and blotting it against clean paper towelling.
7. Add 100 μ l of pre-diluted streptavidin-HRP conjugate. Cover with a new adhesive strip and incubate for 30 minutes at room temperature (20–25°C).
8. Repeat the aspiration/wash step as in step 6.
9. Add 100 μ l of TMB substrate solution into each well, incubate for 30 minutes at room temperature (20–25°C). If the substrate reagent is bright blue prior to use, do not use. **THE PLATE SHOULD BE DEVELOPED IN THE DARK.** Do not cover the plate with aluminium foil or an adhesive strip.
10. Add 100 μ l of stop solution to each well.
11. Determine the optical density of each well within 30 minutes, using a spectrophotometer set to 450 nm.


	1	2	3	4	5	6	7	8	9	10	11	12
A	0	0	S	S	S	S	S	S	S	S	S	S
B	15.36	15.36	S	S	S	S	S	S	S	S	S	S
C	38.4	38.4	S	S	S	S	S	S	S	S	S	S
D	96	96	S	S	S	S	S	S	S	S	S	S
E	240	240	S	S	S	S	S	S	S	S	S	S
F	600	600	S	S	S	S	S	S	S	S	S	S
G	S	S	S	S	S	S	S	S	S	S	S	S
H	S	S	S	S	S	S	S	S	S	S	S	S

Figure 1. Recommended positioning of standard (0–600 pg/ml) and sample wells (S).

Table 1. Assay protocol (all volumes are in microlitres)

	Zero Standard (B₀)	Standards	Samples
Standard	-	50	-
Standard diluent or cell culture media	50	-	-
Sample	-	-	50
Biotinylated antibody reagent	50	50	50
Cover plate, incubate at room temperature for 2 hours.			
Aspirate/decant and wash vigorously all wells three times with wash buffer.			
Streptavidin-HRP conjugate	100	100	100
Cover plate, incubate at room temperature for 30 minutes.			
Aspirate/decant and wash vigorously all wells three times with wash buffer.			
Substrate	100	100	100
Incubate at room temperature for 30 minutes in the dark.			
Stop solution	100	100	100
Determine optical density at 450 nm within 30 minutes.			

Summary of assay protocol

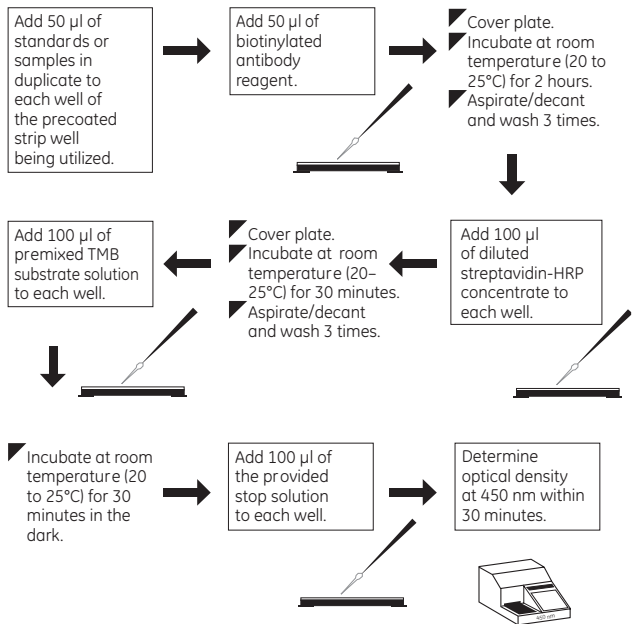

7.5. Calculation of results

Average the duplicate readings for each standard, control and sample and subtract the zero standard optical density.

Plot the optical density for the standards versus the concentration of the standards and draw the best curve. The data can be linearized by using a log/log plot and regression analysis can be applied to the log transformation.

Figure 2 shows such a plot of the data from table 2. The standard curve is provided for illustration only. A standard curve should be generated for each set of samples to be assayed. This allows for the measurement of 42 unknowns in duplicate.

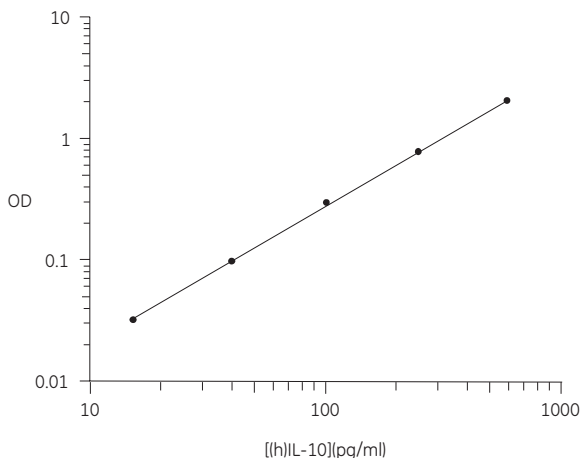


Figure 2. Standard curve

Typical assay data

The following data (table 2) were obtained for a standard curve using the protocol provided.

Table 2. Typical assay data

Tube	Optical density	Zero standard subtracted
Zero standard	0.020	-
15.36 pg/ml standard	0.053	0.033
38.4 pg/ml standard	0.124	0.104
96 pg/ml standard	0.305	0.285
240 pg/ml standard	0.776	0.756
600 pg/ml standard	1.989	1.969

8. Additional information

8.1. Specificity

This assay recognizes both natural and recombinant (h)IL-10. It does not cross-react with human IL-1 α , IL-2, IL-3, IL-4, IL-6, IL-7, IL-8, TNF α , TNF β or IFN γ . There is a 2% reactivity with recombinant mouse IL-10.

8.2. Calibration

The standard in this ELISA is calibrated to the NIBSC reference lot 93/722.

One (1)pg of Biotrak standard = 1.4 NIBSC pg.

8.3. Reproducibility

Within-assay precision

The within-assay coefficient of variation of the ELISA has been determined to be <10%.

Between-assay precision

The between-assay coefficient of variation of the ELISA has been determined to be <10%.

8.4. Sensitivity

The minimum detectable dose of (h)IL-10 was determined to be <3 pg/ml (0.15 pg/well), by adding two standard deviations to the optical density value of zero and calculating the corresponding concentration from the standard curve.

8.5. Recovery

Serum recovery in the ELISA has been determined by spiking recombinant cytokine into human serum, plasma, and urine samples and comparing it with spiked standard diluent control. The values below are typical recoveries:

Standard diluent spike	Recovery from serum
32 pg/ml	89%
110 pg/ml	93%
276 pg/ml	80%

Standard diluent spike	Recovery from plasma
20 pg/ml	94%
86 pg/ml	88%
264 pg/ml	84%

Standard diluent spike	Recovery from urine
38 pg/ml	77%
123 pg/ml	78%
358 pg/ml	75%

8.6. Expected values

Normal average levels of (h)IL-10 in serum samples are
 3.5 ± 3.7 pg/ml.

Range 0–14.1 pg/ml.

Normal average levels of (h)IL-10 in plasma samples are
 4.5 ± 4.57 pg/ml.

Range 1.25–15.56 pg/ml.

Normal average levels of (h)IL-10 in urine samples are
 0.7 ± 0.9 pg/ml.

Range 0–2.71 pg/ml.

9. References

1. BENJAMIN, D. *et al.*, *Blood*, **81**(4), p.1106, 1993.
2. CASSATELLA, M.A. *et al.*, *J. Exp. Med.*, **178** p.2207, 1993.
3. FLUCKIGER, A.C. *et al.*, *J. Exp. Med.*, **178**, p. 1473, 1993.
4. JINQUAN, T. *et al.*, *J. Immunology*, **151**, p.4545, 1993.
5. MOORE, K.W. *et al.*, *Annu. Rev. Immunol.*, **11**, p.165, 1993.
6. GOTLIEB, W.H. *et al.*, *Cytokine*, **4**(5), p.385, 1992.

10. Related products

Biotrak range of Easy ELISA systems

Interferon-alpha [(h)IFN α]	RPN5960
Interferon-gamma [(h)IFN γ]	RPN5961
Interleukin 1 β [(h)IL-1 β]	RPN5971
Interleukin-2 [(h)IL-2]	RPN5965
Interleukin-2 [(m)IL-2]	RPN5966
Interleukin-6 [(h)IL-6]	RPN5968
Interleukin-8/Neutrophil-activating peptide-1-1 [(h)IL-8/NAP-1]	RPN5969
Interleukin-10 [(m)IL-10]	RPN5963
Interleukin-13 [(h)IL-13]	RPN5972
Monocyte chemoattractant protein-1 [(h)MCP-1]	RPN5964
Transforming growth factor beta-1 [(h)TGF β_1]	RPN5970
Tumor necrosis factor-alpha [(h)TNF α]	RPN5967

Biotrak range of human cytokine and growth factor ELISA systems

Interleukin-1 α [(h)IL-1 α]	RPN2750
Interleukin-1 β [(h)IL-1 β]	RPN2751
Interleukin-6 [(h)IL-6]	RPN2754
Interleukin-8 [(h)IL-8]	RPN2764
Interleukin-12 p40/70 [(h)IL-12]	RPN2765
Interferon-gamma [(h)IFN γ]	RPN2757
Tumor necrosis factor-alpha [(h)TNF α]	RPN2758
Vascular necrosis factor-alpha [(h)TNF α]	RPN2779

Biotrak range of high sensitivity human cytokine and growth factor ELISA systems

Interferon-gamma [(h)IFN γ]	RPN2787
Interleukin-6 [(h)IL-6]	RPN2784
Interleukin-10 [(h)IL-10]	RPN2785

Tumor necrosis factor-alpha [(h)TNFα]	RPN2788
---------------------------------------	---------

Biotrak range of mouse cytokine and growth factor ELISA systems

Interferon-gamma [(m)IFNγ]	RPN2707
----------------------------	---------

Interleukin 1β [(m)IL-1β]	RPN2720
---------------------------	---------

Interleukin-4 [(m)IL-4]	RPN2712
-------------------------	---------

Interleukin-6 [(m)IL-6]	RPN2708
-------------------------	---------

Interleukin-12 [(m)IL-12]	RPN2702
---------------------------	---------

Tumor necrosis factor-alpha [(m)TNFα]	RPN2718
---------------------------------------	---------

Biotrak range of rat cytokine and growth factor ELISA systems

Cytokine-induced neutrophil chemoattractant

[(r)GRO/C1NC-1]	RPN2730
-----------------	---------

Interleukin-1β [(r)IL-1β]	RPN2743
---------------------------	---------

Interleukin-6 [(r)IL-6]	RPN2742
-------------------------	---------

Monocyte chemoattractant protein-1 [(r)MCP-1]	RPN2740
---	---------

Tumor necrosis factor-alpha [(r)TNFα]	RPN2744
---------------------------------------	---------

Cell Proliferation assay system and reagents

Cell Proliferation Biotrak ELISA	RPN250
----------------------------------	--------

Cell proliferation kit for immunocytochemistry	RPN20
--	-------

Monoclonal anti-bromodeoxyuridine	RPN202
-----------------------------------	--------

Cell Proliferation Labeling Reagent	RPN201
-------------------------------------	--------

GE Healthcare offices:

GE Healthcare Bio-Sciences AB
Björkgatan 30 751 84 Uppsala,
Sweden

GE Healthcare Europe GmbH
Munzinger Strasse 5 D-79111 Freiburg,
Germany

GE Healthcare Bio-Sciences Corp.
800 Centennial Avenue, P.O. Box 1327,
Piscataway, NJ 08855-1327,
USA

GE Healthcare Bio-Sciences KK
Sanken Bldg. 3-25-1, Hyakunincho,
Shinjuku-ku, Tokyo 169-0073,
Japan

For contact information for your local office,
please visit: www.gelifesciences.com/contact

GE Healthcare UK Limited
Amersham Place
Little Chalfont, Buckinghamshire,
HP7 9NA, UK

<http://www.gelifesciences.com>



imagination at work