

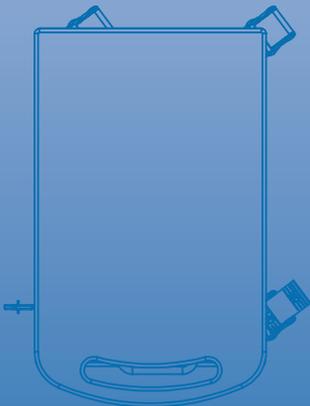
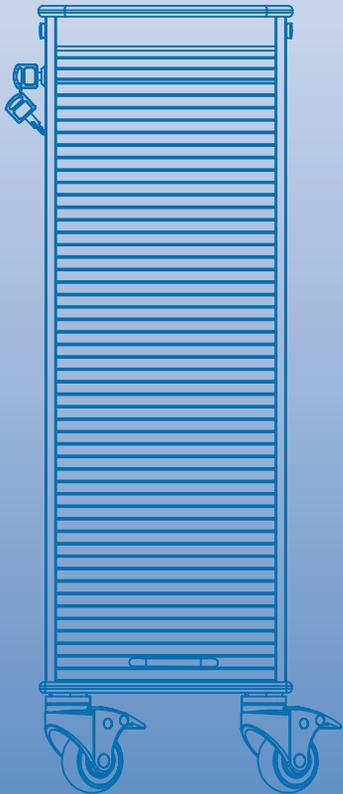


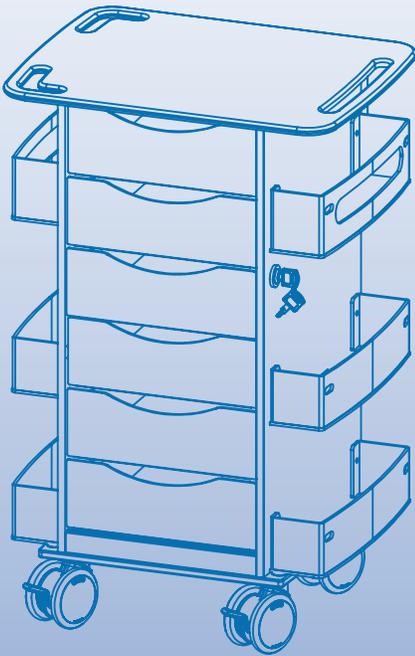

Cart Use and Care Manual

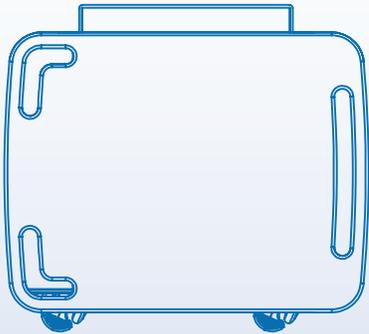

TrippNT's Lab/Healthcare Carts are designed to improve today for a better tomorrow. Built with the highest quality and backed by a total lifetime warranty, TrippNT carts are made for you.*

Thank You for Your Purchase



The cart is being used for its intended purpose by the original user,
For additional warranty information please call us toll-free at (800) 874-7768

Warnings



Important Warnings

The above symbol represents safety warnings that require significant attention when seen on the cart(s) or in the user manual. Failure to do so could result in minor injury, major injury, or even death.



Do not transport the cart up or down stairs or curbs



Service and Replacement

Do not attempt to service or replace any part of your TrippNT Cart unless directed to do so through TrippNT approved documentation (i.e., this User Manual or other instructions). Only TrippNT or a TrippNT-certified entity may service or replace the cart components. If any component on the cart is missing or damaged, the cart must not be used. Contact TrippNT immediately to request service.



Safe Working Loads

Safe working loads labels must be abided by for all elements of TrippNT



Lock casters when cart is not in transit



Carts must be lifted by 2+ persons



Remove all unsecured items before moving cart, all doors and drawers must be locked during transit



Do not sit or stand on carts. Do not stand or lean on the base of the cart



Carts must be pushed by its handle(s)

Statement of Use:

TrippNT Carts are designed to make tomorrow a better day for you. Our goal is to ensure that everyday is better than the next through exacting standards, quality customer service, and durable years of use. Our carts are constructed from corrosion-resistant materials and backed with a total replacement lifetime warranty.

Recommended Shipping and Operating Conditions:

- Temperature Range: - 4 F/ -20 C to 160 F/ 71 C

Terms of Use:

Removing or damaging the serial number and barcode label will void all cart warranties. Returns are not accepted as all TrippNT carts are made to order.

*For additional information or customization options please call
TrippNT toll-free at (800) 874-7768.*

Contact Us

Customer Service

TrippNT

10991 N Airworld Dr.

Kansas City, MO 64153

Customerservice@TrippNT.com

(800) 874-7768

Recommended Cleaning Methods

When cleaning your TrippNT product we recommend the following cleaning solutions:

- Isopropyl Alcohol
- Mild soap and water solution
- Any sanitizing cloth found in healthcare facilities such as Cavi Wipes, Sani-Cloth, or Opti-Cide

Materials:

- HDPE (High Density Polyethylene) is high-impact resistant and has a high melting point of 120 C (225 F).
- Acrylic is a shatter resistant transparent thermoplastic with a melting point of 160 C (320 F)
- ABS (Acrylonitrile Butadiene Styrene) is a thermoplastic polymer. ABS has no true melting point but should be stored in temperatures below 105 C (221 F)
- ACP (Aluminum Composite Panel) is a lightweight and sturdy sandwich panel with a melting point of 101 C (214 F)



Visually inspect parts for obvious wear or damage and remove any debris build up on castor or bearings monthly.