

GE Healthcare



illustra
MicroSpin
G-25 Columns

For the the purification of oligonucleotides and small DNA fragments

Product booklet

Code: 27-5325-01 (50 purifications)



Page finder

1. Legal	3
2. Handling	4
2.1. Safety warnings and precautions	4
2.2. Storage	4
2.3. Expiry	4
3. Components	5
3.1. Kit contents	5
3.2. Materials to be supplied by user	5
3.3. Equipment needed	5
4. Description	6
4.1. Introduction	6
4.2. The basic principle	7
4.3. Product specifications	8
4.4. When to use an illustra MicroSpin G-25 Column	9
5. Protocol	11
5.1. Protocol for purification of a range of sample types	12
6. Appendices	14
6.1. RPM calculation from RCF	14
6.2. Guidelines for use of an illustra MicroSpin G-25 column	14
6.3. Trouble shooting guide	16
6.4. Related products	17
 Quick Reference Protocol Card	Back Cover
Tear off sheet containing protocols for the experienced user purifying oligonucleotides or small DNA fragments	

1. Legal

Product use restriction

The **illustra™ MicroSpin™ G-25 Columns** and components have been designed, developed and sold **for research purposes only**. They are suitable **for *in vitro* use only**. No claim or representation is intended for their use to identify any specific organism or for clinical use (diagnostic, prognostic, therapeutic, or blood banking).

It is the responsibility of the user to verify the use of the **illustra MicroSpin G-25 Columns** for a specific application, as the performance characteristics of this product have not been verified for any specific organism.

GE, imagination at work and GE monogram are trademarks of General Electric Company.

illustra, GFX, PuReTaq, Hybond, Sephadex, Redivue, Rediprime, NICK, Megaprime, Rapid-Hyb, Hyperfilm, FideliTaq, MicroSpin, Hypercassette, ProbeQuant, AutoSeq, NAP and Ready-To-Go are trademarks of GE Healthcare companies.

All third party trademarks are the property of their respective owners.

© 2006–2010 General Electric Company-All rights reserved.
Previously published 2006.

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them. A copy of these terms and conditions is available upon request. Contact your GE Healthcare representative for the most current information (see back cover for contact details).

<http://www.gelifesciences.com>

GE Healthcare UK Limited.
Amersham Place, Little Chalfont,
Buckinghamshire. HP7 9NA UK.

2. Handling

2.1. Safety warnings and precautions

Warning: For research use only.

Not recommended or intended for diagnosis of disease in humans or animals. Do not use internally or externally in humans or animals.

All chemicals should be considered as potentially hazardous. Only persons trained in laboratory techniques and familiar with the principles of good laboratory practice should handle these products. Suitable protective clothing such as laboratory overalls, safety glasses and gloves should be worn. Care should be taken to avoid contact with skin or eyes; if contact should occur, wash immediately with water (See Material Safety Data Sheet(s) and/or Safety Statement(s) for specific recommendations).

2.2. Storage

All kit components should be stored at room temperature (20–25°C). **Do not freeze.**

2.3. Expiry

For expiry date please refer to outer packaging label.

3. Components

3.1. Kit contents

Identification	Pack size Cat. No.	50 purifications 27-5325-01
	illustra™ MicroSpin™ G-25 micro columns	50
	Collection tubes	50

Refer to the Certificate of Analysis for a complete list of kit components.

3.2. Materials to be supplied by user

Disposables:

1.5 ml DNase-free microcentrifuge tubes

3.3. Equipment needed

Microcentrifuge that accommodates 1.5 ml microcentrifuge tubes

Vortex mixer (optional)

4. Description


4.1. Introduction

illustra MicroSpin G-25 columns contain Sephadex™ G-25 DNA grade F resin. They allow DNA purification by the process of gel filtration. Molecules larger than the largest pores in the Sephadex are excluded from the gel and elute first. Intermediate size molecules penetrate the matrix to varying extents, depending on their size and the resin used. Penetration of the matrix retards progress through the column; very small molecules elute last. The volume required to elute these small molecules is dependent on the volume available both inside and outside the pores i.e. the bed volume.

Gel filtration resins do not exhibit a fixed exclusion limit when used in a spin-column format. Exclusion limits of gel filtration resins are only meaningful in continuous flow processes where the molecules being purified have sufficient time to reach equilibrium between the time spent in the gel filtration medium and the time spent in the eluent stream. In spin-column chromatography, the observed exclusion properties that allow the product to pass through the gel while the smaller impurities are retained depends on experimental factors, such as: the resin used, sample volume, product size, and the g forces used in the purification process.

4.2. The basic principle

Use of **illustra MicroSpin G-25 Columns** involves the following steps:

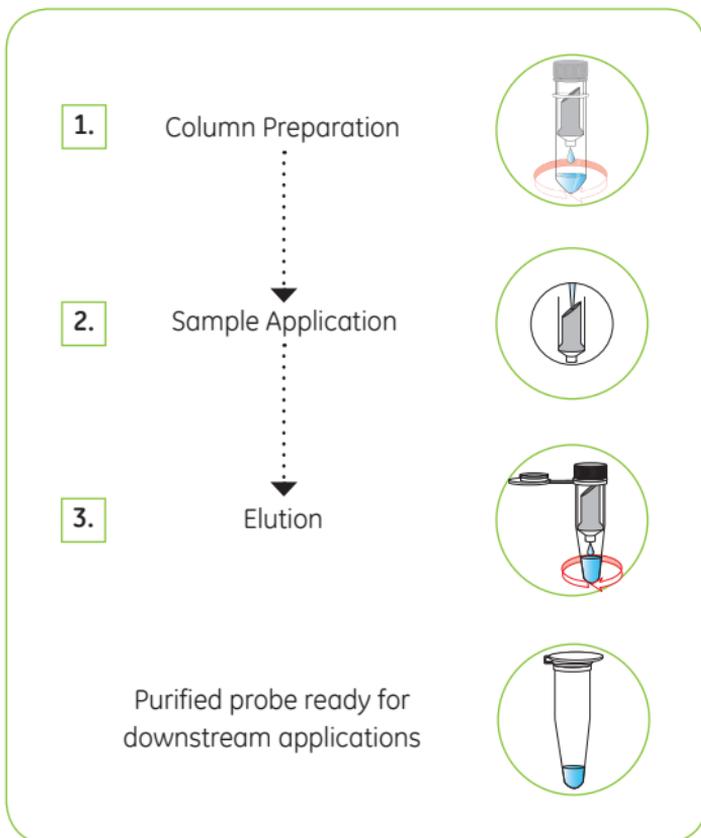

Step	Comments	Component
1. Column Preparation	The resin is re-suspended and excess storage buffer removed by centrifugation.	illustra MicroSpin G-25 column 
2. Sample Application	The sample is applied to the column.	
3. Elution	Purified sample is eluted by centrifugation.	

4.3. Product specifications

Table 1. illustra MicroSpin G-25 column specifications

Sample Type:	Oligonucleotide or small DNA fragment
Principle	Gel filtration.
Column matrix	Sephadex G-25 DNA grade F
Column storage buffer	Double distilled water containing 0.05% Kathon™ CG/ICP Biocide as preservative.
Input sample volume	100–150 µl when desalting oligonucleotides following de-protection in Ammonia. 25–50 µl when removing unincorporated labeled nucleotides from a oligonucleotide labeling reactions. 50 µl for buffer exchange or de-salting of DNA.

Sample Type:	Oligonucleotide or small DNA fragment
Percent sample recovery	Variable—depends on input sample.
Length of labeled DNA recovered	> 10 bases (N.B. there is no maximum length of probe that can be purified).
Nuclease Testing	Column components are tested in nickase, single and double-stranded exonuclease and RNase assays.
Major subsequent applications	Dependent on input sample, but includes blotting, PCR and sequencing applications.

4.4. When to use an illustra MicroSpin G-25 Column

The illustra MicroSpin G-25 column is designed for the rapid purification of DNA for use in a wide range of applications, including desalting, buffer exchange, and removal of labeled nucleotides from DNA labeling reactions. It is suitable for any DNA greater than 10 bases in length and is therefore ideal for the purification of oligonucleotides or very small DNA fragments following synthesis or a labeling reaction. It will not remove or denature enzyme. For guidelines to consider for use of illustra MicroSpin G-25 columns, please see section 6.2.

GE Healthcare provides a wide range of nucleic acid purification products, some of which may be better suited to your application. These products and the application for which they have been optimized are summarized in table 1 below. illustra AutoSeq™ G-50 columns and illustra ProbeQuant™ G-50 micro columns are provided pre-equilibrated in the optimal buffer for the application for which they are designed.

Application	Product	Product code	Pack size
PCR reaction and enzymatic DNA reaction purification 50 bp-10 kbp size range. Extraction of DNA from agarose gels.	illustra GFX™ PCR DNA and Gel Band Purification Kit	28-9034-70	100 purifications
		28-9034-71	250 purifications
Dye terminator removal from automated sequencing reactions.	illustra AutoSeq G-50	27-5340-01	50 purifications
		27-5340-02	250 purifications
		27-5340-03	1 000 purifications
Unincorporated labeled nucleotide removal from a DNA labeling reaction (> 20 mers).	illustra ProbeQuant G-50 Micro-Columns	28-9034-08	50 purifications
Purification of oligonucleotides following synthesis, buffer exchange and de-salting. Gravity format, 500 µl loading volume.	illustra NAP™-5 Columns	17-0853-01	20 purifications

5. Protocol

Note: Columns are NOT transferable between GE Healthcare kits, e.g., the composition of the MicroSpin G-25 columns is not the same as the composition of the MicroSpin G-50 columns.

Use of icons

The Key below describes the purpose of the icons used throughout the protocol booklet.



This icon is used to highlight particularly critical steps within the protocol that must be adhered to. If this advice is not followed it will have a detrimental impact on results.



This icon is used to highlight technical tips that will enhance the description of the step. These tips may indicate areas of flexibility in the protocol or give a recommendation to obtain optimum performance of the kit.

See section 3.2 and 3.3 for Materials & Equipment to be supplied by user.

See section 6.2 for guidelines to consider when purifying different sample types.

5.1. Protocol for purification of range samples

1. Column Preparation

- Re-suspend the resin in the column by vortexing.
- Loosen the cap one-quarter turn and twist off the bottom closure.
- Place the column in the supplied Collection tube for support.
- Spin for 1 minute at $735 \times g$.



Note: See section 6.1 for RPM calculation from RCF.



Note: Use columns immediately after preparation to avoid drying out of the resin. If the column resin appears dry, displaced or cracked after the first spin, this is usually indicative of over-centrifugation (too fast or too long). Re-hydrate the column with 250 μ l of double distilled water, vortex and re-centrifuge, checking the settings. Spin speed can be reduced by 20% if necessary.

Do not use the pulse button on the microcentrifuge as this may over-ride the speed setting.

- Proceed immediately to step 2 below

2. Sample Application

- Place the column into a fresh DNase-free 1.5 ml microcentrifuge tube (user supplied).



Loosen cap
twist off
bottom



1 minute
 $735 \times g$



- b. Slowly apply appropriate volume of sample to the top-center of the resin, being careful not to disturb the resin bed (see Table 2 below).



Table 2: Loading volumes for different sample types


Sample type	Loading volume (μ l)
Oligonucleotides following de-protection in Ammonia	100-150
Oligonucleotide labeling reaction	25-50
De-salting or buffer exchange	50

Where sample volume exceeds that shown in table, use multiple columns, or consider use of NAP-5 Columns. Where sample volume is less than the minimum shown in the table, dilute the sample to improve product recovery (note concentration will be reduced).

 **Note:** The resin will have come away from the column slightly to form a pillar. It is essential that the sample being purified is applied slowly and is not allowed to run down the sides of the resin bed. Avoid touching the resin bed with the pipette tip.

4. Elution

- a. Spin for 2 minutes at $735 \times g$.
The purified sample is collected in the bottom of the 1.5 ml microcentrifuge tube.
- b. Store the purified probe at -20°C .

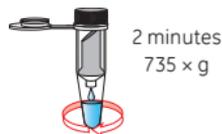

6. Appendices

6.1. RPM calculation from RCF

The appropriate centrifugation speed for a specific rotor can be calculated from the following formula:

$$\text{RPM} = 1\,000 \times \sqrt{(\text{RCF}/1.12r)}$$

Where RCF = relative centrifugal force, r = radius in mm measured from the center of the spindle to the bottom of the rotor bucket, and RPM = revolutions per minute.

For example, if an RCF of $735 \times g$ is required using a rotor with a radius of 73 mm, the corresponding RPM would be 3 000.

Table 1 below shows appropriate RPM for various microcentrifuges.

Table 1: Appropriate RPM for an RCF of $735 \times g$

Microcentrifuge	Appropriate RPM for an RCF of $735 \times g$
Heraeus Biofuge 15	2 800
Beckman GS15R	2 100
Hettich Mikro 24-48	2 630
Hettich Mikro EBA12	2 700
Eppendorf Centrifuge 5415C	3 000
Eppendorf Centrifuge 5417C	2 700

6.2. Guidelines for use of an illustra MicroSpin G-25 column

illustra MicroSpin G-25 Columns can be used for a wide variety of DNA purification applications. The DNA to be purified must be at least 10 bases in length. When using these columns, consider the following guidelines:

20x rule The best results will be obtained when the product being purified is at least 20 times larger than the largest impurity. If the

difference in size is less than 20-fold, either purity or yield may be compromised.

Purity versus yield In general, purity is inversely proportional to yield. Larger sample volumes will provide higher yield but lower purity, and vice-versa.

Non-specific binding The non-specific binding exhibited by the illustra MicroSpin G-25 column is relatively insignificant, allowing purification of samples in the nanogram range. There will be a uniform proportional loss of sample which is due to the nature of spin column chromatography.

Retention For a given sample volume, product retention is relative to molecular size. As the size of the product increases, its relative retention decreases.

Enzyme removal

For purification of DNA fragments 100 bp–10 kbp in length, following an enzymatic reaction, we recommend using the illustra GFX PCR DNA and Gel Band Purification Kit, as the enzyme will be removed during the spin column process. If using an illustra MicroSpin G-25 column, you must Phenol Chloroform extract prior to loading onto the column to ensure enzyme removal.

6.3. Trouble shooting guide

This guide may be helpful in the first instance, however if problems persist or for further information please contact GE Healthcare technical services. Telephone numbers are on the back page. Alternatively log onto <http://www.gelifesciences.com/illustra>

Problem: Poor sample purity.

Possible causes	Suggestions
	<ul style="list-style-type: none">• Ensure the sample volume was within acceptable range prior to loading (see section 6.2).• Ensure sample is CAREFULLY pipetted into centre of resin. Do not disturb the column. Do not allow the sample to run into the sides of the resin bed.• Use the column immediately after completing Column Preparation step. Do not allow the resin to become dried out or cracked.



6.4. Related products

A full range of Molecular Biology reagents can be found in the GE Healthcare catalog and on the web site

<http://www.gelifesciences.com/illustra>

A full range of Detection Products and available pack sizes can be found in the GE Healthcare catalog and on the web site

<http://www.gelifesciences.com/newhyperfilm>

If you need further information, GE technical services are happy to assist (world-wide phone numbers can be found on the back cover).

Application	Product	Product code	Pack size
Removal of protein or enzyme	Hybond™-N+ (82 mm)	RPN82B	50 discs
	Hybond-N+ (15 × 20 cm)	RPN1520B	10 sheets
	Hybond-NX (82 mm)	RPN82T	50 discs
	Hybond-NX (15 × 20 cm)	RPN1520T	10 sheets
	Hybond-N (82 mm)	RPN82N	50 discs
	Hybond-N (15 × 20 cm)	RPN1520N	10 sheets
	Hybond-XL (82 mm)	RPN82S	50 discs
	Hybond-XL (15 × 20 cm)	RPN1520S	10 sheets

Application	Product	Product code	Pack size
Removal of protein or enzyme <i>Continued</i>	Hybond blotting paper (20 × 20 cm)	RPN6101M	100 sheets
Radioactive labeling	Rediprime™ II DNA Labeling System	RPN1633	30 reactions
	Ready-To-Go™ DNA Labeling Beads (-dCTP)	27-9240-01	1 kit
	Megaprime™ DNA Labeling System, dNTP	RPN1604	30 reactions
	Megaprime DNA Labeling System, dCTP	RPN1606	30 reactions
	Nick Translation Kit, dNTP	N5500	20 reactions
	Nick Translation Kit, dCTP	N5000	20 reactions
	5'-End Labeling Kit	RPN1509	20 reactions
	Rapid-Hyb™ Buffer	RPN1635	125 ml
Detection	Hyperfilm™ MP (18 × 24 cm)	28-9068-43	50 sheets

Application	Product	Product code	Pack size
Detection <i>Continued</i>	Hyperfilm MP Enveloped (18 × 24 cm)	28-9068-50	50 sheets
	Hypercassette™	RPN11642	1
Purification of DNA probes and oligonucleotides	illustra ProbeQuant G-50 Micro Columns	28-9034-08	50 purifications
	illustra NICK™ Columns	17-0855-02	50 purifications
	illustra NAP-5™ Columns	17-0853-02	50 purifications
Purification of DNA from PCR, agarose gel bands and enzymes	illustra GFX PCR DNA & Gel Band Purification Kit	28-9034-70	100 purifications
	illustra GFX 96 PCR Purification Kit	28-9034-45	10 × 96 well plates
Preparation of PCR products for automated sequencing	ExoSAP-IT™	US78200	100 purifications
Dye terminator removal from automated sequencing reactions	illustra AutoSeq G-50 Columns	27-5340-01	50 purifications

Application	Product	Product code	Pack size
Kits containing ready-to-use mix for PCR amplification	illustra Hot Start Master Mix	25-1500-01	100 purifications
	illustra PuReTaq™ Ready-To-Go PCR Beads	27-9557-01	96 reactions in 0.2 ml tubes/plate
	illustra PuReTaq Ready-To-Go PCR Beads	27-9557-02	5 × 96 reactions in 0.2 ml tubes/plate
	FideliTaq™ PCR Master Mix Plus (2 ×)	E71182	100 reactions
	FideliTaq Master Mix Plus	E71183	100 reactions
Premixed nucleotides for PCR amplification	illustra DNA Polymerization Mix dNTP Set (A,C,G,T) 20 mM each	28-4065-57	10 µmol
	illustra DNA Polymerization Mix dNTP Set (A,C,G,T) 20 mM each	28-4065-58	40 µmol (4 × 10 µmol)
	illustra PCR Nucleotide Mix dNTP Set (A,C,G,T) 25 mM each	28-4065-60	500 µl

Application	Product	Product code	Pack size
Premixed nucleotides for PCR amplification	illustra PCR Nucleotide Mix dNTP Set (A,C,G,T) 2 mM each	28-4065-62	1 ml

Continued

GE Healthcare offices:

GE Healthcare Bio-Sciences AB
Björkgatan 30 751 84
Uppsala
Sweden

GE Healthcare Europe GmbH
Munzinger Strasse 5 D-79111
Freiburg
Germany

GE Healthcare UK Limited
Amersham Place
Little Chalfont
Buckinghamshire
HP7 9NA
UK

GE Healthcare Bio-Sciences Corp
800 Centennial Avenue
P.O. Box 1327
Piscataway
NJ 08855-1327
USA

GE Healthcare
Japan Corporation
Sanken Bldg. 3-25-1
Hyakunincho Shinjuku-ku
Tokyo 169-0073
Japan

For your local office contact information, visit
<http://www.gelifesciences.com/contact>

GE Healthcare UK Limited
Amersham Place,
Little Chalfont, Buckinghamshire,
HP7 9NA, UK

<http://www.gelifesciences.com>




imagination at work

27-5325-01PL AC 03/2010

28977285

The next two pages are a
protocol card.
If required please add to the
back page as a tear off addition

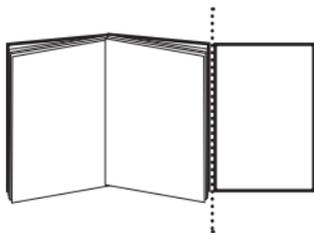

If not then delete these three
pages.

Quick Reference Protocol Card

illustra™ MicroSpin™ G-25 Columns

28-9034-08 (50 purifications)

Protocol for purification of a range of sample types

 : Add  : Spin


1. Column preparation

- Re-suspend the resin in the column by vortexing
 - Loosen the cap one-quarter turn and twist off the bottom closure
 - Place the column in the supplied Collection tube
-  1 minute 735 x g



2. Sample application

- Place the column into a fresh DNase-free 1.5 ml microcentrifuge tube (user supplied)
- Sample to the top-center of the resin with care (see Table below)

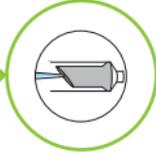

3. Elution

-  2 minutes 735 x g
- Retain eluate
- Store the purified probe at -20°C



Table: Loading volumes for different sample types

Sample type	Loading volume (µl)
Oligonucleotides following de-protection in Ammonia	100–150
Oligonucleotide labeling reaction	25–50
De-salt or buffer exchange	50



GE, imagination at work and GE monogram are trademarks of General Electric Company.
illustra and MicroSpin are trademarks of GE Healthcare Companies.
© 2006–2010 General Electric Company-All rights reserved.
Previously published 2006 July.

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them. A copy of these terms and conditions is available upon request.
Contact your GE Healthcare representative for the most current information.

<http://www.gelifesciences.com>

GE Healthcare UK Limited,
Amersham Place, Little Chalfont,
Buckinghamshire, HP7 9NA UK