



CE Duo in action!

# Return to work

Cleaning and disinfection with CE Duo™ CleanTech™ for Controlled Environments

## What to Clean/Disinfect?

### High Touch Surfaces

#### Horizontal Surfaces:

- All work surfaces
- Floors & ceilings

#### Vertical Surfaces:

- Walls, windows, softwalls

#### Uneven Surfaces:

- Misc. equipment
- Pipes, tanks, hoods
- Door handles, knobs, light switches



### High Touch Devices

#### Work Devices:

- Equipment & computer controllers

## When to Clean?

Per SOP or

 Start and end of the work day

 After each shift

## What to Use ?

### Wipes & Sponges

EvoControl® 400 Cleanroom Wipes



CE Duo™ CleanTech™ Sponge



Pre-saturate with Vileda's 6 Liter Bucket

### Mops

CE Duo™ CleanTech™ Mops with Vileda's Pre-Prepared System



### Disinfectant


Per SOP

## How to Clean/Disinfect?

### Wiping


Use one wipe or sponge per surface to prevent cross contamination.

Wiping should be done in a pull & lift fashion with overlapping strokes. This can be done horizontally or vertical as shown below. Following this step is critical in removing contamination instead of spreading in circles.

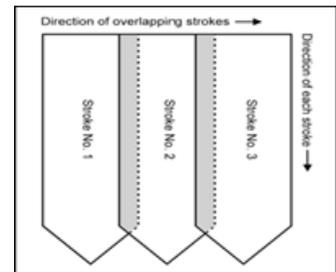

### Mopping

With Mops a modified figure 8 (S-Method), Right to left, then left to right, with overlapping strokes

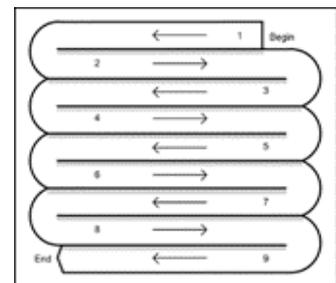

When the software begins to dry out it's time to replace with a new one.



Accessories are optional\*

## EvoControl<sup>®</sup> Wipes & CE Duo<sup>™</sup> CleanTech<sup>™</sup> Sponge

VWR Cat. No.	Description	Color	Size	Pieces/PK	PKS/CS
10030-966	EvoControl <sup>®</sup> Wipe	White	9 x 9 in	150	10
10030-968	EvoControl <sup>®</sup> Wipe	White	12 x 12 in	150	10
75871-654	CE Duo <sup>™</sup> CleanTech <sup>™</sup> Sponge	Blue/White	4 x 7 in	6	10
10153-180*	6l Bucket - accessory	Grey	11 x 8 x 7 in	NA	10
10153-182*	6l Bucket Lid - accessory	Grey	11 x 8 x 1 in	NA	12

## CE Duo<sup>™</sup> CleanTech<sup>™</sup> System

VWR. Cat. No.	Description	Color	Size	Pieces/PK	PKS/CS
89413-086	CE Duo <sup>™</sup> CleanTech <sup>™</sup> Microfiber Mop	Blue/White	19.7 in	1	50
76204-546	Swep Duo <sup>™</sup> Plus Frame w/Limiter	Grey	19.7 in	1	10
89066-060	Vileda Aluminum Extendable Handle	Blue w/coding	40 to 71 in	NA	10
10153-168*	CE Pre-Prepared Trolley - accessory	Stainless Steel	30 x 28 x 37 in	NA	1
10153-180*/ 10153-182*	Pre-prepared bucket & lid - accessory	Grey	15 x 11 x 8 in 15 x 11 x 1 in	NA	5